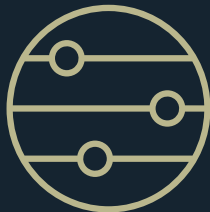


Mutation Surveyor



User Manual



For more information,
visit [SoftGenetics.com](https://www.softgenetics.com)



Release Information

<i>Document Version Number</i>	<i>MutationSurveyor-5.0-UG001</i>
<i>Software Version</i>	<i>5.2</i>
<i>Document Status</i>	<i>Final</i>
<i>Document Release Date</i>	<i>November, 2021</i>

Copyright

© 2021. SoftGenetics, LLC, All rights reserved.

The information contained herein is proprietary and confidential and is the exclusive property of SoftGenetics, LLC. It may not be copied, disclosed, used, distributed, modified, or reproduced, in whole or in part, without the express written permission of SoftGenetics, LLC.

Limit of Liability

SoftGenetics, LLC has used their best effort in preparing this guide.

SoftGenetics, LLC makes no representations or warranties with respect to the accuracy or completeness of the contents of this guide and specifically disclaims any implied warranties of merchantability or fitness for a particular purpose. Information in this document is subject to change without notice and does not represent a commitment on the part of SoftGenetics, LLC or any of its affiliates. The accuracy and completeness of the information contained herein and the opinions stated herein are not guaranteed or warranted to produce any particular results, and the advice and strategies contained herein may not be suitable for every user.

The software described herein is furnished under a license agreement or a non-disclosure agreement. The software may be copied or used only in accordance with the terms of the agreement. It is against the law to copy the software on any medium except as specifically allowed in the license or the non-disclosure agreement.

Trademarks

The name “*SoftGenetics*,” the SoftGenetics logo, and *Mutation Surveyor* are trademarks or registered trademarks of SoftGenetics, LLC. All other products and company names mentioned herein might be trademarks or registered trademarks of their respective owners.

Customer Support

Customer support is available to organizations that purchase Mutation Surveyor and that have an annual support agreement. Contact SoftGenetics, LLC at:

SoftGenetics, LLC
100 Oakwood Ave, Suite 350
State College, PA 16803
(814) 237-9340
(888) 791-1270 (US Only)
tech_support@softgenetics.com
www.softgenetics.com

Table of Contents

Chapter 1: Getting Started with Mutation Surveyor	19
Mutation Surveyor System Requirements	21
Installing Mutation Surveyor.....	22
To install Mutation Surveyor	22
Starting Mutation Surveyor	23
Mutation Surveyor Conventions.....	24
Genomic terms	24
Mutation terms	24
Reports and graphical output.....	25
The Mutation Surveyor Main Window	26
Title bar	26
Main menu	26
Main window toolbar	27
Browser pane.....	27
To navigate the Browser pane.....	28
Viewing and Reporting pane.....	30
Configuring User Management.....	31
To configure user management.....	31
To turn on user management	34
To turn off user management	36
Managing Groups in Mutation Surveyor	38
To manage groups in Mutation Surveyor.....	38
To add a new group.....	40
To edit a group	40
To delete a group	41
Managing Users in Mutation Surveyor.....	43
To manage users in Mutation Surveyor.....	43

To add a user	45
To edit a user	46
To delete a user	47
Chapter 2: Project Setup.....	49
Project File Types.....	50
Sample file	50
Reference file	50
GenBank file.....	50
Automatic GenBank download.....	51
Modifying the Default Mutation Surveyor Project Settings	52
To modify the default Mutation Surveyor project settings	53
Input tab	54
Contig tab.....	55
Mutation tab	59
Output tab	62
Display tab	64
2 Directions tab	65
Others tab	67
Database tab.....	69
Setting up a Single Mutation Surveyor Project.....	70
To set up a Mutation Surveyor project	70
Contig directory for a Mutation Surveyor project.....	72
Chapter 3: Graphical Analysis Display.....	75
Graphical Analysis Display Layout.....	77
Nucleotide Sequence frame.....	78
Amino Acid frame	79
Tracks frame	79

Sample Traces frame	79
Mutation trace legend	80
Mutation trace peak identification	82
Mutation table	83
GAD Toolbar	84
Search GenBank sequence	86
Variation Tracks Settings dialog box	87
Graphical Analysis Display Navigation and Analysis Options.....	90
GAD viewing and navigation options	90
GAD analysis options	91
Miscellaneous GAD options.....	93
Exporting the GAD	94
Chapter 4: Mutation Surveyor Graphical Reports	97
Project Reviewer Report	99
Project Reviewer report layout.....	99
Project Reviewer report toolbar	102
Project Reviewer Report Settings dialog box	104
Mutation tab	104
Color tab	105
Nomenclature tab	106
Display tab	107
Graphic Display Report.....	108
Graphic Display report layout	109
Graphic Display report toolbar	110
Project Display subreport	111
Project Display subreport toolbar	113
GAD Report	114
GAD Report toolbar	115

Clinical Report	116
Nucleotide Text Display Report.....	120
Nucleotide Text Display report layout	121
Nucleotide Text Display report toolbar	122
Chapter 5: Mutation Surveyor Standard Reports	125
Mutation Report.....	127
Mutation Surveyor toolbar	127
Mutation report fields.....	128
Mutation report codes	129
Substitution mutation code.....	129
Color codes	130
Insertion, duplication, and deletion mutation codes	131
Mutation report viewing and analysis options	132
Two Direction Report.....	134
Two Direction report layout	134
Two Direction report toolbar	135
Two Direction report mutation codes	136
Two Direction report viewing and analysis options	136
Two Direction report settings	136
Two Direction report context menu	137
Advanced Two Direction Report.....	140
Advanced Two Direction report toolbar	142
Advanced Two Direction report viewing and analysis options	144
Project Report	146
Project report toolbar	148
Project Sample Assembly subreport.....	150
Consensus Sequence Text function	151
Consensus Sequence Text window toolbar	151

Sequence Text Comparison function	152
Sequence Text Comparison window toolbar	153
Amino Acid Text Comparison function	154
Amino Acid Text Comparison window toolbar	155
Project report viewing and analysis options.....	155
HGVS Report.....	157
HGVS report toolbar	160
HGVS report viewing and analysis options.....	161
Chapter 6: Mutation Surveyor Custom Report.....	163
Custom Report.....	164
To create a Custom Report template.....	164
Format tab	166
Filters tab	170
Display tab.....	172
Mutation tab.....	173
Color tab	176
Nomenclature tab	177
Other tab.....	179
Custom Report toolbar.....	180
Custom Report viewing and analysis options	181
Chapter 7: Mutation Surveyor Customer Reports	183
JHU report.....	185
JHU Report toolbar	186
JHU report viewing and analysis options	187
JHU report settings.....	187
JHU report context menu.....	188
Tassel report.....	190

Tassel report toolbar	191
Tassel report viewing and analysis options.....	192
Emory report.....	193
Emory report toolbar	193
Emory report viewing and analysis options.....	194
Bentley report and MSM_Song report.....	196
Bentley/MSM_Song report toolbar	197
Bentley/MSM_Song reports viewing and analysis options.....	197
Bentley/MSM_Song report settings	198
Bentley/MSM_Song reports context menu	198
PrettyBase report	200
PrettyBase report toolbar	200
PrettyBase report viewing and analysis options.....	201
PrettyBase report settings.....	201
PrettyBase report context menu	202
Methylation report.....	203
Methylation report toolbar	204
Graphic Display of Methylation report.....	205
Methylation report context menu	206
Vertex report.....	207
Vertex2 report.....	209
Vertex2 report toolbar	211
Vertex2 report viewing and analysis options.....	211
Vertex3 report.....	212
Vertex3 report layout.....	213
Vertex3 report toolbar	214
Vertex3 report viewing and analysis options.....	214
Full Gene report	215

Full Gene Report Settings dialog box, Format tab.....	216
Full Gene Report Settings, Header tab.....	217
Full Gene Report Settings, Output tab.....	217
Full Gene report toolbar.....	219
Full Gene viewing and analysis options.....	220
FullBase report.....	222
FullBase report layout.....	223
FullBase report toolbar	223
FullBase viewing and analysis options	224
Chapter 8: GenBank File Editors.....	225
Overview of GenBank Files	227
To download a GenBank file.....	227
Advanced GBK File Editor Tool	231
To load a GenBank file into the Advanced GBK File Editor tool.....	231
Advanced GBK File Editor tool toolbar	232
Advanced GBK File Editor tool GenBank Tree File pane	233
Advanced GBK File Editor tool Tabs pane	235
Sequence tab	235
Basic Information tab.....	237
Allele Frequency tab.....	240
Modifying and adding variations.....	241
Advanced GBK File Editor tool functions.....	243
To search the loaded GenBank file for a specific sequence	243
To download a GenBank file with the same sequence.....	244
To adjust the number of the first coding base or the sequence start.....	244
To add ROIs that are based on CDSs.....	245
To reverse complement the currently loaded GenBank file.....	245
SEQ File Editor Tool	246

To load a SEQ file into the SEQ File Editor tool.....	246
SEQ File Editor tool layout	247
SEQ File Editor tool functions	250
To search the loaded SEQ file for a specific sequence	250
To adjust the number of the first coding base or the sequence start	251

Chapter 9: Mutation Surveyor Tools.....253

Filename Editor Tool	257
To programmatically change filenames	257
To manually change filenames	259
2D Filename Match Editor tool	261
To use the 2D Filename Match Editor tool	261
ND Filename Match Editor Tool	264
To use the ND Filename Match Editor	264
Print Header Editor Tool.....	267
To use the Print Header Editor tool	267
Create SCF Tool	271
Synthetic traces versus pseudo traces	271
To use the Create SCF tool	272
Convert to SCF tool.....	274
To use the Convert to SCF tool.....	274
Extract SGP Tool.....	275
To use the Extract SGP tool.....	275
SGP Comparison Tool	279
To use the SGP Comparison tool	279
Mutation Assembly Tool.....	284
To use the Mutation Assembly tool	284
Sequence Assembly Tool.....	287

To use the Sequence Assembly tool	287
Mask Vector Sequence Tool.....	291
To use the Mask Vector Sequence tool.....	291
Output Trace Data Tool	293
To use the Output Trace Data tool	293
Signal Factor Intensity and Signal Factor Dev	294
BasePatch Tool.....	295
To use the BatchPatch tool.....	295
The HLA Tool.....	297
To use the HLA tool	297
Chapter 10: Mutation Surveyor Specialized Applications	299
Methylation Detection	301
Mutation Quantifier Function.....	303
To run the Mutation Quantifier function	303
Mutation Quantifier report, Standardized Allele Ratio selected	305
Mutation Quantifier report, Simplified Allele Ratio selected.....	307
Mutation Quantifier toolbar	308
Quantification Group Editor dialog box	309
Mutation Quantifier report settings dialog box.....	310
Chapter 11: Multiple Projects	313
Modifying Mutation Surveyor Projects	315
The Log File Wizard in the Mutation AutoRun Tool	317
To open the Log File Wizard.....	317
To set up a new log file.....	318
To modify a log file.....	322
To modify an existing job in the log file.....	322
To add a new job to a log file.....	323

To remove a job from a log file	323
The Mutation AutoRun Tool	324
To use the Mutation AutoRun tool.....	324
Chapter 12: Reviewing and Editing Mutation Projects	327
Adding Mutation Calls	329
To add a substitution.....	329
To add a homozygous indel call.....	330
To add or edit a heterozygous indel (The Het-Indel tool).....	331
Heterozygous Indel Detection window toolbar.....	334
Editing Mutation Calls.....	335
To edit substitutions	335
To edit homozygous indels	336
To edit heterozygous indels	337
To edit Del/Ins	339
To edit a homozygous Del/Ins	339
To edit a heterozygous Del/Ins	339
Annotating Reported Variations	341
To annotate a variation	341
Creating a User Knowledge Database for Known Deletions.....	342
To create a custom variant database for known deletions.....	342
Viewing the Edit History (Audit Trail) for a Mutation.....	344
Index	345

Preface

Welcome to the *Mutation Surveyor User's Manual*. The purpose of the *Mutation Surveyor User's Manual* is to answer your questions and guide you through the procedures necessary to use the Mutation Surveyor application efficiently and effectively.

Using the manual

You will find the *Mutation Surveyor User's Manual* easy to use. You can simply look up the topic that you need in the table of contents or the index. Later, in this Preface, you will find a brief discussion of each chapter to further assist you in locating the information that you need.

Special information about the manual

The *Mutation Surveyor User's Manual* has a dual purpose design. It can be distributed electronically and then printed on an as-needed basis, or it can be viewed online in its fully interactive capacity. If you print the document, for best results, it is recommended that you print it on a duplex printer; however, single-sided printing will also work. If you view the document online, a standard set of bookmarks appears in a frame on the left side of the document window for navigation through the document. For better viewing, decrease the size of the bookmark frame and use the magnification box to increase the magnification of the document to your viewing preference.



If you do print the document using a single-sided printer, you might see a single blank page at the end of some chapters. This blank page has been added solely to ensure that the next chapter begins on an odd-numbered page. This blank page in no way indicates that your book is missing information.

Conventions used in the manual

The *Mutation Surveyor User's Manual* uses the following conventions:

- Information that can vary in a command—variable information—is indicated by alphanumeric characters enclosed in angle brackets; for example, <Mutation ID>. Do not type the angle brackets when you specify the variable information.
- A new term, or term that must be emphasized for clarity of procedures, is *italicized*.
- Mutation Surveyor provides multiple ways of accessing the same Mutation Surveyor functions (for example, a main menu option and a toolbar button). For brevity and ease of

use, this manual references only one method—the main menu—for accessing functions; however, you can use whatever method best suits your working needs.

- Page numbering is “online friendly.” Pages are numbered from 1 to x, *starting with the cover* and ending on the last page of the index.



Although numbering begins on the cover page, this number is not visible on the cover page or front matter pages. Page numbers are visible starting with the first page of the table of contents.

- This manual is intended for both print and online viewing.
 - If information appears in [blue](#), it is a hyperlink. Table of Contents and Index entries are also hyperlinks. Click the hyperlink to advance to the referenced information.

Assumptions for the manual

The *Mutation Surveyor User's Manual* assumes that:

- You are familiar with windows-based applications and basic Windows functions, navigational elements, and icons such as the Printer Setup icon on a Print Preview page.
- References to any third party standards or third party software functions were current during the development and release of this version of Mutation Surveyor.

Organization of the manual

In addition to this Preface, the *Mutation Surveyor User's Manual* contains the following chapters and appendices:

- [Chapter 1, “Getting Started with Mutation Surveyor,” on page 19](#) details the Mutation Surveyor installation requirements, and the procedures for installing and upgrading the application. It also describes how to launch Mutation Surveyor and it provides an overview of the major navigational elements in the application.
- [Chapter 2, “Project Setup,” on page 49](#) guides you through the necessary steps for setting up a project in Mutation Surveyor.
- [Chapter 3, “Graphical Analysis Display,” on page 75](#) details the layout and functions of the Graphical Analysis Display, which visually presents mutation project data.
- [Chapter 4, “Mutation Surveyor Graphical Reports,” on page 97](#) details the variety of graphic-based standard reports that are available in Mutation Surveyor for providing detailed information about a mutation project. The chapter contains a summary of descriptions, report field labels, report functions, and report caveats for all the various standard mutation reports.

- [Chapter 5, “Mutation Surveyor Standard Reports,” on page 125](#) details the variety of text-based standard reports that are available in Mutation Surveyor for providing detailed information about a mutation project. The chapter contains a summary of descriptions, report field labels, report functions, and report caveats for all the various standard mutation reports.
- [Chapter 6, “Mutation Surveyor Custom Report,” on page 163](#) details the Custom report, which contain options for building a tailored report that provides detailed information about your mutation projects in a format that best suits your specific working needs.
- [Chapter 7, “Mutation Surveyor Customer Reports,” on page 183](#) details the various customer report templates in Mutation Surveyor. Each report was initially developed as a template to provide custom content that met a specific customer’s needs. All these report templates are now available from the Reports > Customer Reports option on the Mutation Surveyor main menu.
- [Chapter 8, “GenBank File Editors,” on page 225](#) details the tools that are specifically designed for editing GenBank files and Sequence files.
- [Chapter 9, “Mutation Surveyor Tools,” on page 253](#) details the variety of tools that are available in Mutation Surveyor for optimizing input data and analyzing results.
- [Chapter 10, “Mutation Surveyor Specialized Applications,” on page 299](#) details the use of Mutation Surveyor for some specialized applications.
- [Chapter 11, “Multiple Projects,” on page 313](#) details Mutation Surveyor’s capabilities for modifying existing projects and for working with multiple projects.
- [Chapter 12, “Reviewing and Editing Mutation Projects,” on page 327](#) details the functions and the tools that are available in Mutation Surveyor for adding and editing mutation calls in a project.

Related document

In addition to this manual, you may find the following document helpful

- [Mutation Surveyor Reference and Track Manager Tool User’s Manual](#)

Preface

Chapter 1

Getting Started with Mutation Surveyor

The Mutation Surveyor software application uses patented anti-correlation technology to provide highly accurate discovery of DNA variants from Sanger sequencing traces. This software is ideal for the analysis of data from Applied Biosystems, Beckman-Coulter and MegaBACE systems. This chapter details the installation requirements, and the procedures for installing and upgrading the application. It also describes how to launch Mutation Surveyor and it provides an overview of the major navigational elements in the application. It also details how to configure user management for your Mutation Surveyor instance.

This chapter covers the following topics:

- [“Mutation Surveyor System Requirements” on page 21.](#)
- [“Installing Mutation Surveyor” on page 22.](#)
- [“Starting Mutation Surveyor” on page 23.](#)
- [“Mutation Surveyor Conventions” on page 24.](#)
- [“The Mutation Surveyor Main Window” on page 26.](#)
- [“Configuring User Management” on page 31.](#)
- [“Managing Groups in Mutation Surveyor” on page 38.](#)
- [“Managing Users in Mutation Surveyor” on page 43.](#)

Chapter 1

Getting Started with Mutation Surveyor

Mutation Surveyor System Requirements

Mutation Surveyor is supported only on Windows and Intel-powered Macintosh operating systems. It is not supported on Linux or UNIX-based operating systems. Mutation Surveyor recognizes only PC file formats.

System	Requirements
Windows PC	<ul style="list-style-type: none">• Operating System—Mutation Surveyor can function on Windows 32- or 64-bit operating systems that are within Microsoft Extended Support.• Processor—Dual core with 1.8 GHz• RAM: 2 GB• Available hard drive space: 60 GB <p>Note: Some optional features require a 64-bit operating system along with additional RAM and hard drive space.</p>
Intel Powered Macintosh	<ul style="list-style-type: none">• Operating System—Parallels desktop for Mac (Mac OS/virtual machine dependent) or Apple, Boot Camp or VMware Fusion (Mac OS/virtual machine dependent)• RAM: 2 GB• Available hard drive space: 60 GB

Installing Mutation Surveyor

Mutation Surveyor is licensed in three different ways, each of which follow slightly different installation procedures—Validation, Local, and Network:



You must have Local Administrator rights for the computer on which you are installing Mutation Surveyor.

- Validation license—The Validation license is a trial license that provides all the functionality of a purchased license. You can load data, create and save new files, analyze and visualize data, and so on. The Validation license does not require a registered license and you can install this version on an unlimited number of computers. The Validation license expires 35 calendar days from installation.



You must contact SoftGenetics to receive a disc that contains a fully functional, 35 day trial of the software.

- Local license—The Local license is for installation on a single computer.
- Network license—The Network license is for installation on multiple client computers that are connected to a license server computer.

To install Mutation Surveyor

The Mutation Surveyor Installation wizard guides you through the steps that are necessary to install a Mutation Surveyor license. The default installation location is

C:\Program Files (x86)\SoftGenetics\Mutation Surveyor

but you can always select a different installation directory.

When you are installing Mutation Surveyor, keep in mind the following:

Version	Comments
Validation	Accept the default Program Manager Group to ensure that program operability is not negatively affected.
Local	<ul style="list-style-type: none">• A dongle may be supplied with a Local license. Make sure that you follow the specific installation instructions that are included with the key.• Accept the default Program Manager Group to ensure that program operability is not negatively affected.• At the prompts to enter your registration information, enter the information exactly as SoftGenetics supplied it to you.
Network	<ul style="list-style-type: none">• At the prompts to enter your registration information on the server computer, enter the information exactly as SoftGenetics supplied it to you.

Starting Mutation Surveyor

After Mutation Surveyor has been installed on your computer, a shortcut icon for the application is placed on your desktop. An option for the application is also available from your Start menu. You can double-click the desktop icon to open Mutation Surveyor, or you can select the option from your Start menu.




If your Mutation Surveyor administrator has configured User Management for your Mutation Surveyor instance, then you are prompted to enter a user name and password before Mutation Surveyor opens. Contact your administrator if you need assistance.

Figure 1-1: Mutation Surveyor desktop icon



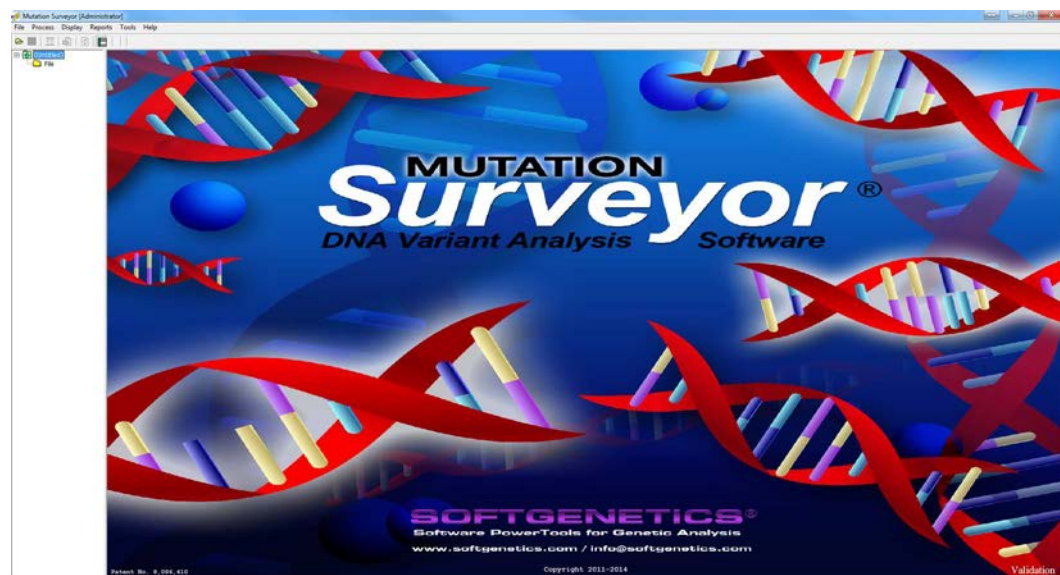Figure 1-2: Mutation Surveyor main window



Mutation Surveyor Conventions


Mutation Surveyor uses conventions in both its terminology ([genomic terms](#) and [mutation terms](#)) and its [output](#) to ensure consistency and accuracy when you are manipulating and analyzing mutation project data.

Genomic terms

In genomics, the same term or definition can have multiple meanings or interpretations. As a result, you must know how specific genomic terms are defined for use in Mutation Surveyor. [Figure 1-3](#) shows a sequence region which contains a gene model. Mutation Surveyor uses the following terms to define the different areas of a gene:

- (A)—The sequence, which is the complete order of nucleotides in a gene.
- (B)—mRNA, which is a region in a gene that contains part of the reading frame that codes for a specific portion of the complete protein. mRNA regions can include sequences that code for amino acids (red fill in [Figure 1-3](#)) and untranslated sequences (no fill in [Figure 1-3](#)).
- (C)—CDS, or the coding sequence, which is a DNA region in a gene that is translated into protein.
- (D)—Intron, which are the regions of the precursor mRNA that are spliced out to make the mature mRNA. These introns might be within the untranslated region of the mRNA or they might be between CDS regions of mRNA.

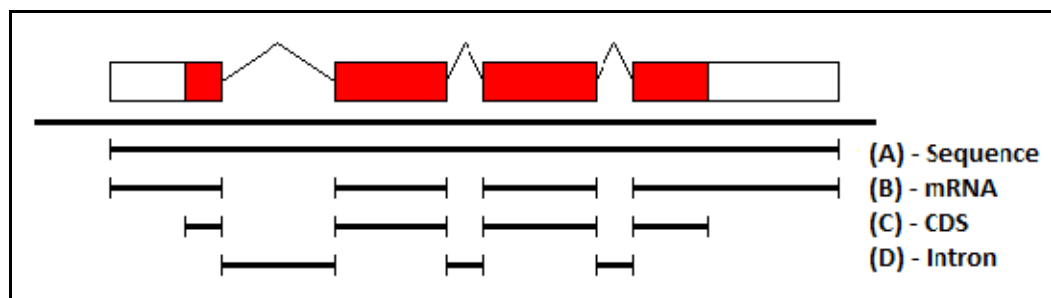Figure 1-3: Mutation Surveyor genomic terms



Mutation terms

The following list details how specific types of mutations are defined in Mutation Surveyor:

- Silent—Coding and non-coding mutations that do not affect the amino acid sequence.
- Synonymous—Mutations found only in the coding region that do not affect the amino acid sequence.

- Non-synonymous—Mutations found only in the coding region that affect the amino acid sequence.
- Nonsense—A non-synonymous mutation that results in a stop codon.
- Mosaic—Might be used in place of somatic mutations (2D small peak) or heteroplasmy.

Reports and graphical output

- When analyzing samples from a polyploid organism, Mutation Surveyor displays the results for only the two dominant alleles (the alleles with the two highest frequencies) in any report or graphical display.
- Mutation Surveyor supports both standard and non-standard genetic codes as defined in Version 4.4 of the NCBI Genetic Code table (see <ftp://ftp.ncbi.nih.gov/entrez/misc/data/gc.prt>). To analyze samples that require translation using a non-standard genetic code, the GenBank file must contain the correct Translation table number. If the GenBank file does not contain this information, Mutation Surveyor defaults to using standard genetic code.

The Mutation Surveyor Main Window

The Mutation Surveyor main window is your starting point for the Mutation Surveyor application. The window provides quick access to all the Mutation Surveyor functions and system tools. The Mutation Surveyor main window has five major components—the [title bar](#), the [main menu](#), the [toolbar](#), the [Browser pane](#), and the [Viewing and Reporting pane](#).


Figure 1-4: Mutation Surveyor main window



Title bar

The name “Mutation Surveyor” appears in the title bar at the top of the Mutation Surveyor main window. When a saved project is opened, the title bar also shows the version number in which the project was saved.

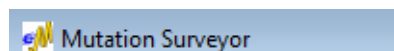
Figure 1-5: Title bar



Main menu

The main menu is set up in a standard Windows menu format with menu commands grouped into menus (File, Process, Display, Reports, Tools, and Help) across the menu bar. Some of these menu commands are available in other areas of the application.

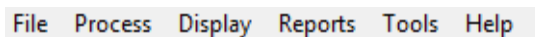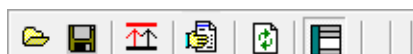
Figure 1-6: Main menu









Main window toolbar

The toolbar provides quick access to all the Mutation Surveyor functions. [Figure 1-7](#) below shows the Mutation Surveyor toolbar after the application first opens. The Mutation Surveyor toolbar is updated dynamically with other toolbar buttons based on the files that you load and/or the post-processing steps that you select. Place your mouse pointer over a toolbar button to open tooltip text for the button.

Figure 1-7: Mutation Surveyor Toolbar




Button	Description
	Open Files—Opens the Open Files dialog box which you use to load GenBank files, references files, and sample files for a project.
	Save Project—Saves all the project files to a single file mutation project file with an .sgp extension. You can open an entire project in a single step by opening this .sgp file.
	Raw Data to Processed Data—Enabled when AB1/ABI files without base call information are uploaded and Load Raw Data is selected for the Mutation Project settings. Converts the raw files so that base call information is displayed.
	Run—Runs the currently loaded project in Mutation Surveyor.
	Restart—Opens the Options dialog box which you use to modify the current project settings. You can then run the project again using these new settings.
	Show/Hide Browser—Shows/hides the Browser pane in the Mutation Surveyor main window.

Browser pane




The Browser (left) pane is arranged in a classic Windows Explorer-style folder structure. The top-level folders are the File folder and the Contig folder. The File folder contains up to three sub-level folders—GenBank File, Reference File, and Sample File—which contain the GenBank files, the reference files, and the sample files, respectively. The Contig folder contains subfolders for each of the contigs that were created from the analysis. After you run a project for the first time, an entry also exists for the Mutation Report. (See [Figure 1-8 on page 28.](#))

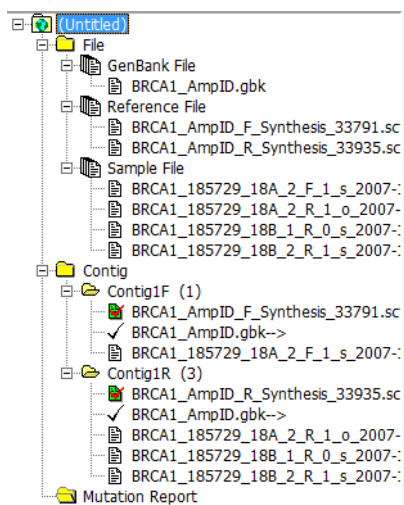Figure 1-8: Browser pane



For a detailed discussion about contigs and the Browser pane, see “[Contig directory for a Mutation Surveyor project](#)” on page 72.

To navigate the Browser pane

- To open a folder, click the Expand icon  next to the folder, or double-click the folder. To close a folder, click the Collapse icon  next to the folder.
- To view the raw trace for a file, double-click anywhere on the file entry. (The icon for the selected file turns green ) The trace opens in the Viewing and Reporting (right) pane. To clear the trace for a file, double-click anywhere on the file entry again.



To view corresponding traces at the same time, double-click individual samples.

Figure 1-9: Viewing a trace for a single file


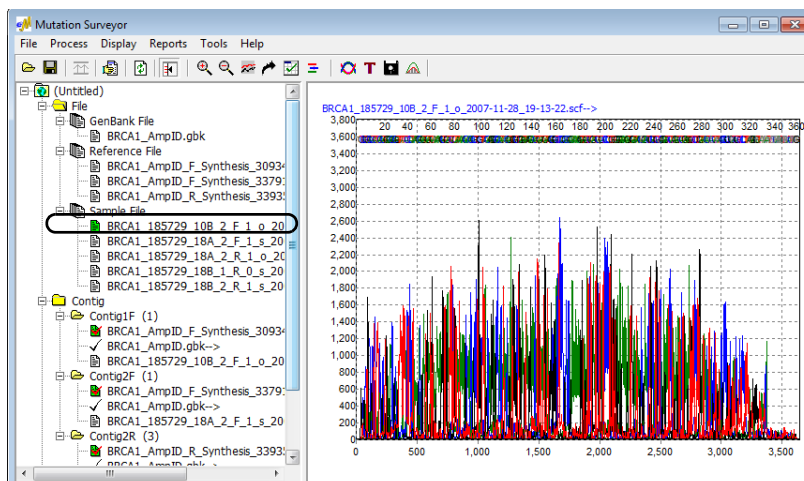


Figure 1-10: Viewing traces for multiple files

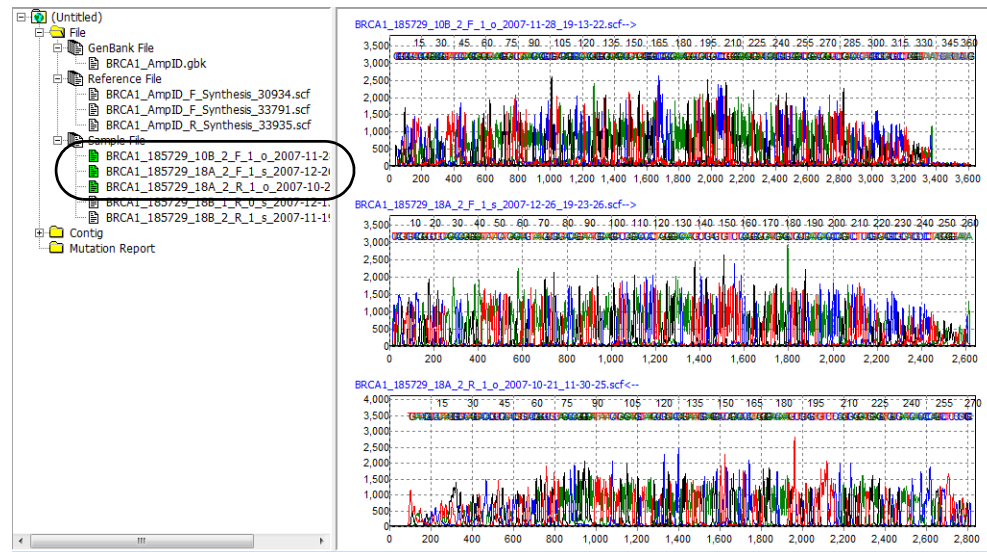

- To view the Mutation report, double-click the Mutation Report entry. The Mutation report opens in the Viewing and Reporting pane. (See “Mutation Report” on page 127.)


Figure 1-11: Mutation Report in the Viewing and Reporting pane

Sample File	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Qualit	Mut#	Mutation1
BRCA1_012345_10B_2_R_1_s_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30746	31027	282	11	1	30753A>AT,278T>T/S\$21
BRCA1_185729_10A_1_R_1_i_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30503	30852	350	26	0	
BRCA1_185729_10A_2_R_0_i_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30503	30852	350	26	0	
BRCA1_185729_10B_1_R_0_s_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30759	31027	269	11	1	30988A>AG,356Q>Q/R\$20
BRCA1_185729_10B_2_R_0_s_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30759	31027	269	11	1	30988A>AG,356Q>Q/R\$20
BRCA1_264381_10A_1_R_1_o_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30503	30852	350	25	0	
BRCA1_264381_10A_2_R_0_s_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30503	30852	350	25	0	
BRCA1_264381_10B_1_R_0_o_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30759	31041	283	10	1	30988A>AG,356Q>Q/R\$14
BRCA1_264381_10B_2_R_1_o_2007-1	BRCA1_AmpID_R_S1-R	BRCA1			2	30759	31041	283	10	1	30988A>AG,356Q>Q/R\$14
BRCA1_012345_11A_1_F_1_s_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	30977	31416	440	7	0	
BRCA1_012345_11A_2_F_1_o_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	30977	31416	440	7	0	
BRCA1_012345_11B_1_F_1_i_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	31143	31460	318	0	2	31389C>CG,490P>P/A\$22
BRCA1_012345_11B_2_F_1_i_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	31143	31460	318	0	2	31389C>CG,490P>P/A\$22
BRCA1_185729_11A_1_F_0_s_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	30975	31414	440	11	0	
BRCA1_185729_11A_2_F_0_o_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	30975	31414	440	11	0	
BRCA1_185729_11B_1_F_1_i_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	31138	31460	323	0	1	31407C>CT,496R>R/C\$32
BRCA1_185729_11B_2_F_1_i_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	31138	31460	323	0	1	31407C>CT,496R>R/C\$32
BRCA1_264381_11A_1_F_0_i_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	30975	31414	440	10	0	
BRCA1_264381_11A_2_F_0_s_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	30975	31414	440	10	0	
BRCA1_264381_11B_1_F_0_s_2007-1	BRCA1_AmpID_F_S2-F	BRCA1			2	31134	31460	327	0	1	31407C>CT,496R>R/C\$32

Viewing and Reporting pane

The Viewing and Reporting (right) pane displays graphics, reports, and other information depending on the post-processing options that you have selected.

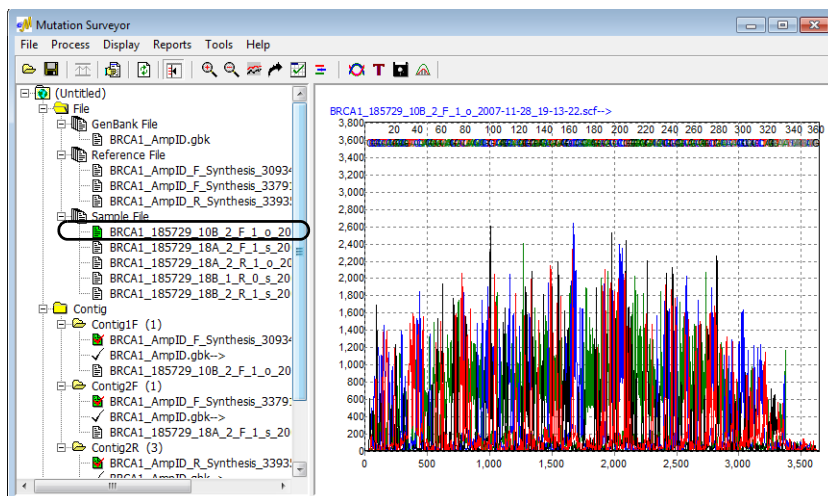Figure 1-12: Viewing a trace in the Viewing and Reporting pane



Reports and views are discussed in the appropriate chapters in this manual.

Configuring User Management

After Mutation Surveyor is installed, user management can be configured for it. User management requires that a user be authenticated before logging in and using the applications. You can [configure](#) user management independently for each computer (localhost) on which Mutation Surveyor is installed. In this configuration, the SoftGenetics Server service must be installed on each computer on which Mutation Surveyor is installed. Because the same user management configuration is part of the installation process for Geneticist Assistant and NextGENe, the steps that you must follow to install the SoftGenetics Server service depend on whether Geneticist Assistant and/or NextGENe has already been installed on the localhost. Alternatively, a single server can host the SoftGenetics Server service and you can configure each Mutation Surveyor host to connect to this single server to verify user credentials. When you configure user management, you must always configure the Administrator user account first. Only the Administrator user has all the necessary privileges for managing other users. All other users are standard users. After you configure user management, you must [turn on](#) user management. You can also always [turn off](#) user management at any time without deleting any of the user configuration information.



If you changed the directory for storing the MySQL information that Mutation Surveyor uses from the default directory (C:\ProgramData\MySQL\MySQL Server 5.1\Data), then before configuring user management, you must contact tech_support@softgenetics.com.

To configure user management



The following procedures details the configuration of user management independently for each computer (localhost) on which Mutation Surveyor is installed. To configure user management with a single server hosting the SoftGenetics Server service, contact tech_support@softgenetics.com.


1. If Geneticist Assistant or NextGENe is already installed on the computer on which you are configuring user management for Mutation Surveyor, go to [“To turn on user management” on page 34](#); otherwise, do the following:
 - Log on to the host computer as a Windows user that is a local Administrator.
 - To avoid issues with User Account Control settings, right-click on the Mutation Surveyor desktop shortcut and on the context menu that opens, select Run as administrator.
2. On the Mutation Surveyor main menu, click Help > User Management > Install Local Service.

The License page for the SoftGenetics Server Setup wizard opens. The page details the license agreement for installing the SoftGenetics Server service.



Be patient. It might take a few minutes for the SoftGenetics Server Setup wizard to open.

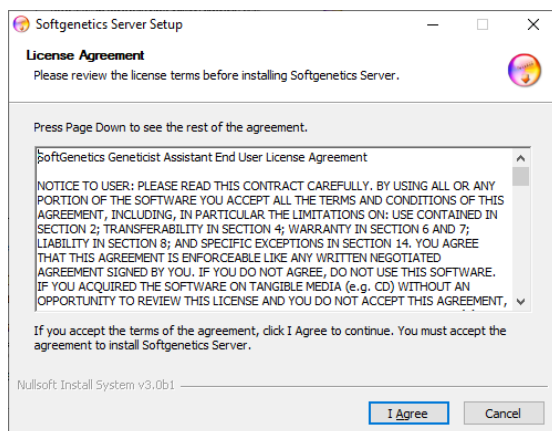
Figure 1-13: SoftGenetics Server Setup wizard, License Agreement page



3. Click I Agree to accept the license agreement.

The Settings page for the SoftGenetics Server Setup wizard opens. By default, the page is prepped for configuring the Administrator user.

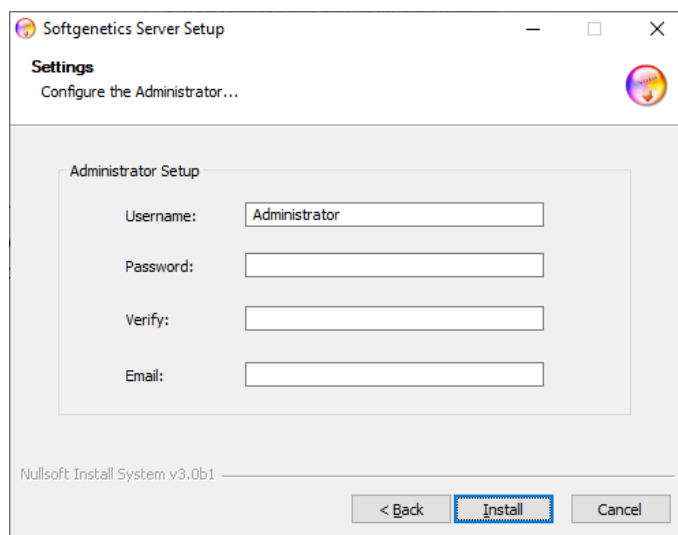Figure 1-14: SoftGenetics Server Setup wizard, Settings page



4. Do the following:
 - Leave the user name set to Administrator, or modify it as needed.
 - In the Password field, enter the password for the Administrator user.



The only invalid character for the password is a space. There are no other special requirements or restrictions for the Administrator password. It can adhere to your organization's standards and any other requirements as needed. If you forget or lose this password, it is not recoverable.

- In the Verify field, enter the Administrator password *exactly* as you entered it in the Password field.
- In the Email field, enter the email address for the Administrator user.




The current version of User Management does not support email notifications; however, an email address is still required.

5. Click Install.



After you select the server, the space requirements for installing the SoftGenetics Server service are displayed on the page. Make sure that you have sufficient space on the computer to install this service.

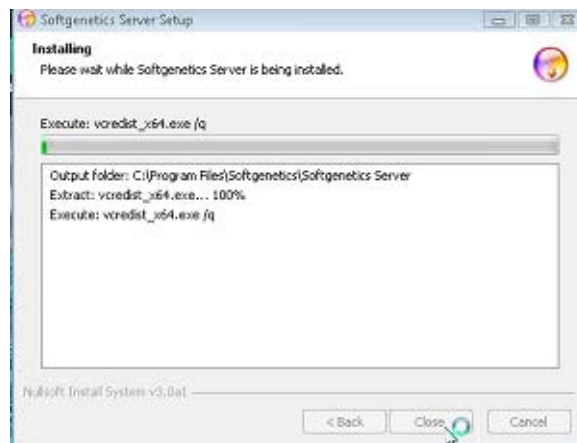
Figure 1-15: SoftGenetics Server Setup wizard, Installation page



Note the following about the installation:


- If MySQL has *not* already been installed on the localhost, then after installation of MySQL is complete, click Close at the prompt; otherwise, the installation begins with the installation of the other server components (Python, Django, and Apache).
- During the installation of the other server components, you might receive Security Alerts. The installation is set up to handle these alerts and with the exception of a Windows Security Alert for Apache (see below), no special action is required.
- After Apache is installed, a Windows Security Alert opens indicating that the Windows Firewall has blocked some features of the installation. Click Allow access to allow the Apache HTTP Server to operate correctly on the localhost.

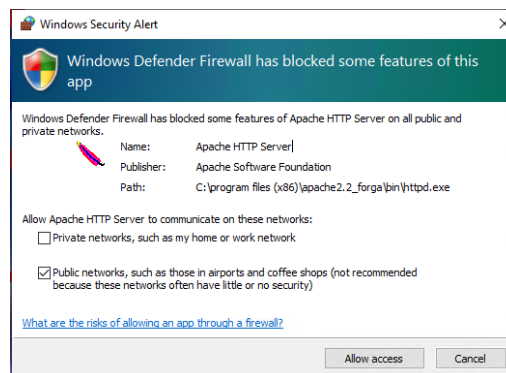Figure 1-16: Windows Security Alert for Apache



After installation is complete, Completed is displayed at the top of the Installation page.

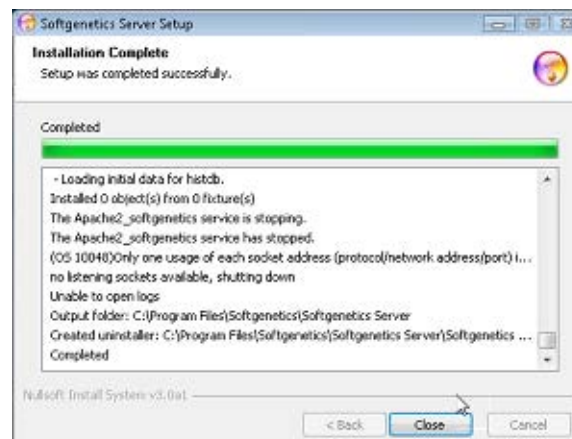Figure 1-17: SoftGenetics Server Setup wizard, Installation page for completed installation



6. Click Close.

The SoftGenetics Server Setup wizard closes. Mutation Surveyor remains open.


7. Continue to “To turn on user management” below.

To turn on user management

1. On the Mutation Surveyor main menu, click Help > User Management > Manage Settings.

The User Management Settings dialog box opens. The General tab is the open tab.

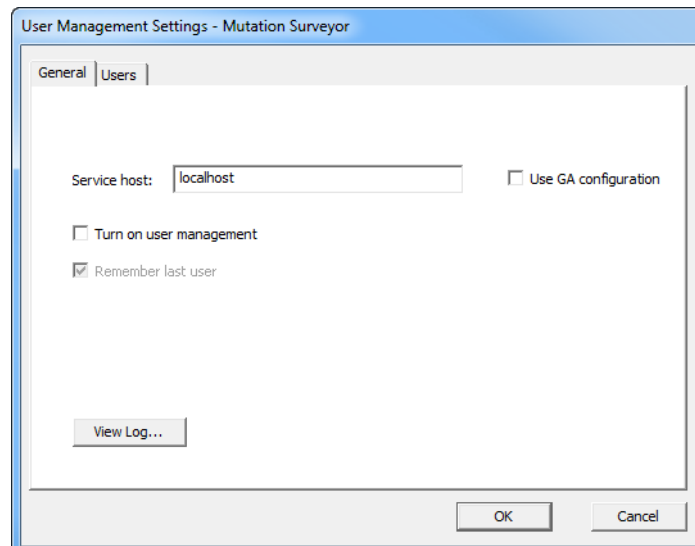Figure 1-18: User Management Settings dialog box, General tab



2. Leave Service host set to localhost.
3. Select Turn on user management.
Remember last user becomes available.
4. Leave Remember last user, or optionally, clear it.




If Remember last user is selected, then when a user logs into Mutation Surveyor, the Username field on the Login dialog box is automatically populated with the user name for the user who last logged into Mutation Surveyor.

5. Select Use GA configuration to integrate user accounts between multiple SoftGenetics products.
6. Click OK.

The Administrator Verification dialog box opens. The dialog box indicates that Administrator verification is required to apply the changes.

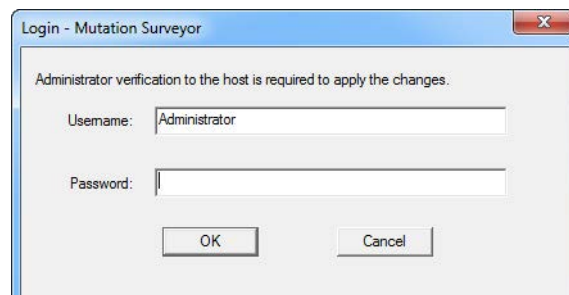Figure 1-19: Administrator Verification dialog box



7. In the Username field, leave the Administrator username as-is, or optionally, modify the name as needed.
8. In the Password field, enter the password for the Administrator user.
9. Click OK.

A message opens, indicating that to apply the changes that Mutation Surveyor must be closed and reopened and asking you if you want to close Mutation Surveyor now.


10. Click Yes.

The message closes.

11. Start Mutation Surveyor.

The Login dialog box opens.

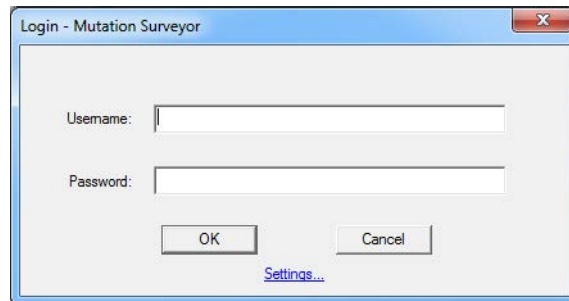Figure 1-20: Mutation Surveyor Login dialog box



12. Enter the Administrator username and password, and then click OK.

The Login dialog box closes and Mutation Surveyor opens. Now, every time a user opens Mutation Surveyor, they are prompted to enter a username and password before they can use the application.



If you are the Administrator user, you should now continue to setting up the needed groups and users for your Mutation Surveyor instance. See [“Managing Groups in Mutation Surveyor”](#) on page 38 and [“Managing Users in Mutation Surveyor”](#) on page 43.


To turn off user management

After configuring and turning on user management for your Mutation Surveyor instance, as the Administrator user, you always have the option of turning off user management. This does not delete any user configuration information. It simply means that users are not required to be authenticated before they log in to and use Mutation Surveyor. You can always turn user management back on.

1. Start Mutation Surveyor.

The Login dialog box opens.

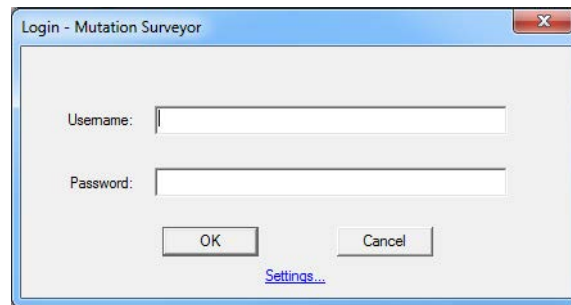Figure 1-21: Mutation Surveyor Login dialog box




2. Enter the Administrator username and password, and then click OK.

Mutation Surveyor opens.

3. On the Mutation Surveyor main menu, click Help > User Management > Manage Settings.

The User Management dialog box opens. The General tab is the open tab. Turn on user management is selected.

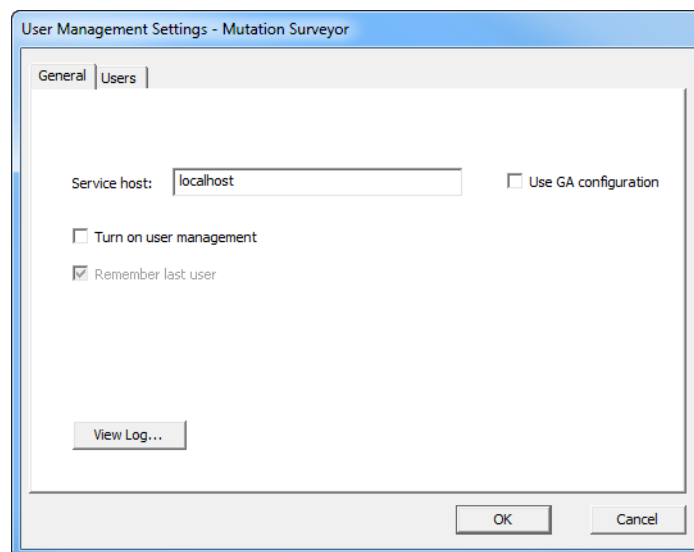Figure 1-22: User Management Settings dialog box, General tab



4. Clear Turn on user management.
5. Click OK.

A message opens, indicating that to apply the changes that Mutation Surveyor must be closed and reopened and asking you if you want to close Mutation Surveyor now.

6. Click Yes.

The message and Mutation Surveyor close. Now, any user can start Mutation Surveyor without any authentication. The user configuration information, however, is not deleted, so you can always turn user management back on if needed.

Managing Groups in Mutation Surveyor

Users are the people who log into Mutation Surveyor, whether they are adding and reviewing content, or just using the application in a read-only capacity. A *group* is a collection of users that have the same permissions in Mutation Surveyor. As the Administrator user for Mutation Surveyor, you are responsible for managing all the groups for your Mutation Surveyor instance and managing the users for these groups to ensure that your users have the appropriate permissions available to them in Mutation Surveyor. You can assign users to one of the four default groups that are installed with every instance of Mutation Surveyor, or you can create your own groups with the needed permissions, and then assign users to one of these groups.

	Mutation Surveyor Default User Group			
	Reporter	Technician	Analyst	Supervisor
Assigned Permissions				
View Project	Y	Y	Y	Y
Export Results	Y	Y	Y	Y
Create and Run Project	N	Y	Y	Y
Re-run Project	N	N	Y	Y
Edit Variants	N	N	Y	Y
Manage Analysis Settings	N	N	Y	Y
Manage Report Settings	N	N	Y	Y
Manage Knowledge Base	N	N	N	Y

[Managing](#) groups for Mutation Surveyor consists of adding new groups, editing existing groups, and deleting groups.

To manage groups in Mutation Surveyor


1. On the Mutation Surveyor main menu, click Help > User Management > Manage Settings.

The User Management Settings dialog box opens. The General tab is the open tab. See [Figure 1-22 on page 37](#).

2. Click the Groups tab to open it.

The tab lists the four default groups that are installed with every instance of Mutation Surveyor as well as any groups that have been configured for your Mutation Surveyor instance. If applicable, it also lists any groups that have been configured for your Geneticist Assistant and/or NextGENe instances. See [Figure 1-23 on page 39](#).

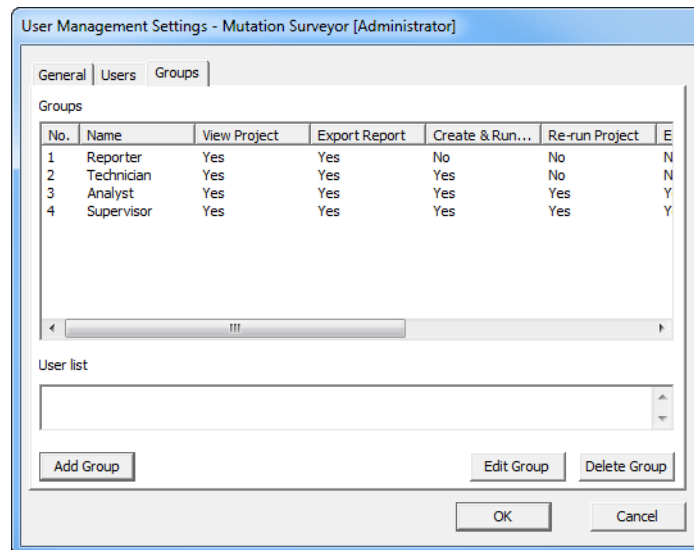Figure 1-23: User Management Settings dialog box, Groups tab



3. Optionally, to view a list of all users that are currently assigned to a group, select the group.

The users that are assigned to the selected group are displayed alphabetically by username in the User list pane.


4. Continue to one of the following:
 - [“To add a new group” on page 40.](#)
 - [“To edit a group” on page 40.](#)
 - [“To delete a group” on page 41.](#)

To add a new group

1. Click Add Group.

The Add Group dialog box opens.

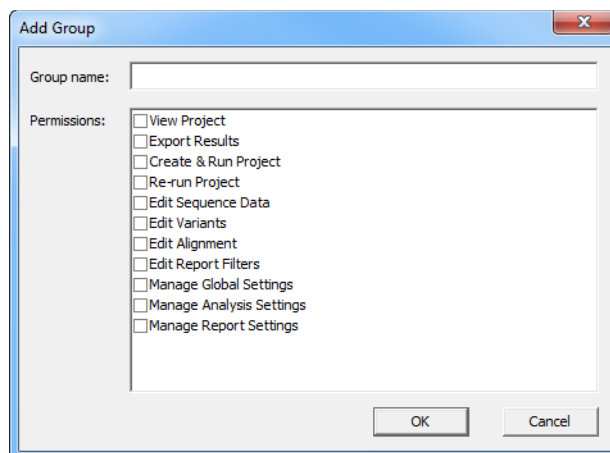Figure 1-24: Add Group dialog box



2. In the Group name field, enter the name for the new group.
3. On the Permissions list, select the permissions for the new group.
4. Click OK.

A message opens, indicating that the new group was successfully created.

5. Click OK.

The message closes. The Groups tab remains open with the newly added group displayed on the tab.

6. Click OK.

The User Management Settings dialog box closes.

To edit a group

Editing a group from the Groups tab consists of modifying the permissions for the group. If you want to edit a group by adding or deleting users, then you must do so from the Users tab. (See [“Managing Users in Mutation Surveyor” on page 43.](#)) Also, you cannot edit a group name. If you need to rename a group, then you must delete the current group, and create a new group with the new name, and then reassign all the users who were associated with the deleted group to this new group.




Although you can edit the permissions that are assigned to the Mutation Surveyor default groups, SoftGenetics strongly recommends that you not do so. Instead, you should create a new group with the appropriate permissions and then assign users to the new group.

1. Select the group for which you are modifying the permissions, and then click Edit Group.

The Edit Group dialog box opens. The group name is displayed in the Group name field, and you cannot edit it. The permissions that are currently assigned to the group are also displayed.

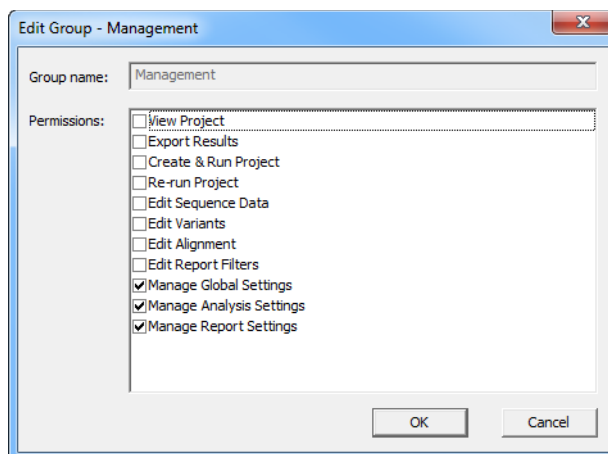Figure 1-25: Edit Group dialog box



2. Modify the permissions for the group as needed.
3. Click OK.

A message opens, indicating that the group was successfully edited.

4. Click OK.

The message closes. The Groups tab remain opens.

5. Click OK.

The User Management Settings dialog box closes.

To delete a group



Although you can delete any of the Mutation Surveyor default groups, SoftGenetics strongly recommends that you not do so. Instead, you should delete only those custom groups that you have added for your Mutation Surveyor installation.

1. Select the group that you are deleting, and then click Delete Group.

A message opens, indicating that you are deleting the selected group, and prompting you to click OK to confirm the deletion.

Chapter 1

Getting Started with Mutation Surveyor

2. Click OK.

The message closes, and a second message opens, indicating that you have successfully deleted the selected group.

3. Click OK.

The second message closes. The entry for the group is removed from the Groups tab. The Groups tab remains open.

4. Click OK.

The User Management Settings dialog box closes.

Managing Users in Mutation Surveyor

Users are the people who log into Mutation Surveyor, whether they are adding and reviewing content, or just using the application in a read-only capacity. If you are the Administrator user for Mutation Surveyor, then you are responsible for managing all the other users. [Managing](#) users for Mutation Surveyor consists of adding new users, editing existing users, and deleting users. You can also [view](#) the activity for your Mutation Surveyor users (logging in to or logging out of Mutation Surveyor) in a log file.

To manage users in Mutation Surveyor


1. On the Mutation Surveyor main menu, click Help > User Management > Manage Settings.

The User Management Settings dialog box opens. The General tab is the open tab. See [Figure 1-22 on page 37](#).

2. Optionally, to view the activity for your Mutation Surveyor users (logging in to or logging out of Mutation Surveyor) in a log file, click View Log.

The User Management Log file opens onscreen. The file lists all the activity for your Mutation Surveyor users, and if applicable, for your Geneticist Assistant and/or NextGENe users as well. You can click Save to File to save the log file with a name and a location of your choosing.

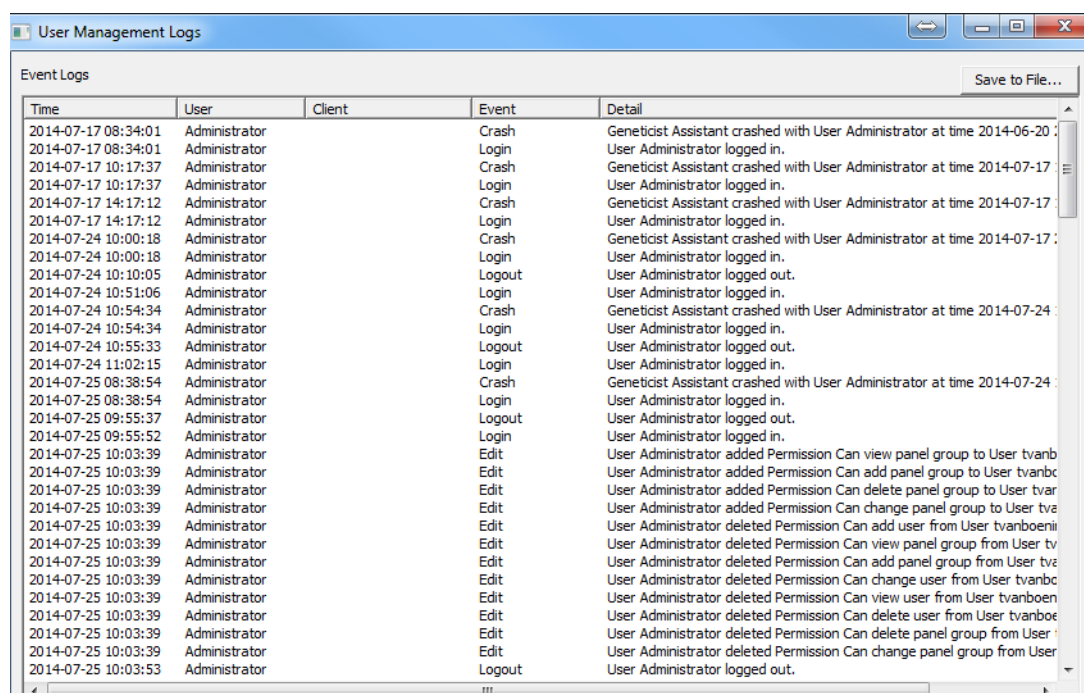Figure 1-26: User Management Log file




The screenshot shows a window titled "User Management Logs" with a "Save to File..." button. Below the title bar is a table with the following columns: Time, User, Client, Event, and Detail. The table contains a list of events, including logins, logouts, crashes, and permission changes, all performed by the Administrator user.

Time	User	Client	Event	Detail
2014-07-17 08:34:01	Administrator		Crash	Geneticist Assistant crashed with User Administrator at time 2014-06-20:
2014-07-17 08:34:01	Administrator		Login	User Administrator logged in.
2014-07-17 10:17:37	Administrator		Crash	Geneticist Assistant crashed with User Administrator at time 2014-07-17:
2014-07-17 10:17:37	Administrator		Login	User Administrator logged in.
2014-07-17 14:17:12	Administrator		Crash	Geneticist Assistant crashed with User Administrator at time 2014-07-17:
2014-07-17 14:17:12	Administrator		Login	User Administrator logged in.
2014-07-24 10:00:18	Administrator		Crash	Geneticist Assistant crashed with User Administrator at time 2014-07-17:
2014-07-24 10:00:18	Administrator		Login	User Administrator logged in.
2014-07-24 10:10:05	Administrator		Logout	User Administrator logged out.
2014-07-24 10:51:06	Administrator		Login	User Administrator logged in.
2014-07-24 10:54:34	Administrator		Crash	Geneticist Assistant crashed with User Administrator at time 2014-07-24:
2014-07-24 10:54:34	Administrator		Login	User Administrator logged in.
2014-07-24 10:55:33	Administrator		Logout	User Administrator logged out.
2014-07-24 11:02:15	Administrator		Login	User Administrator logged in.
2014-07-25 08:38:54	Administrator		Crash	Geneticist Assistant crashed with User Administrator at time 2014-07-24:
2014-07-25 08:38:54	Administrator		Login	User Administrator logged in.
2014-07-25 09:55:37	Administrator		Logout	User Administrator logged out.
2014-07-25 09:55:52	Administrator		Login	User Administrator logged in.
2014-07-25 10:03:39	Administrator		Edit	User Administrator added Permission Can view panel group to User tvanb
2014-07-25 10:03:39	Administrator		Edit	User Administrator added Permission Can add panel group to User tvanb
2014-07-25 10:03:39	Administrator		Edit	User Administrator added Permission Can delete panel group to User tvanb
2014-07-25 10:03:39	Administrator		Edit	User Administrator added Permission Can change panel group to User tvanb
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can add user from User tvanboeni
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can view panel group from User tv
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can add panel group from User tv
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can change user from User tvanb
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can view user from User tvanboen
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can delete user from User tvanboe
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can delete panel group from User
2014-07-25 10:03:39	Administrator		Edit	User Administrator deleted Permission Can change panel group from User
2014-07-25 10:03:53	Administrator		Logout	User Administrator logged out.

3. Click the Users tab to open it.

The tab lists all the user accounts that have been configured for your Mutation Surveyor instance, and if applicable, any user accounts that have been configured for your Geneticist Assistant and/or NextGENe instances.

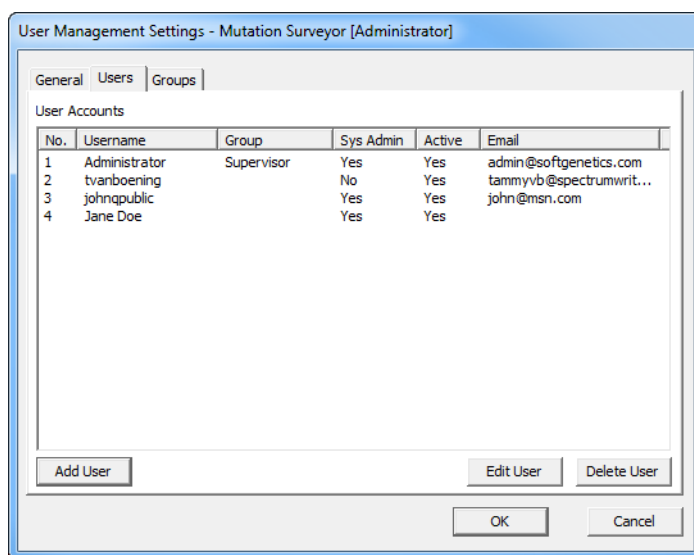Figure 1-27: User Management Settings dialog box, Users tab




4. Continue to one of the following:
 - “To add a user” on page 45.
 - “To edit a user” on page 46.
 - “To delete a user” on page 47.

To add a user

1. Click Add User.

The Add User dialog box opens.

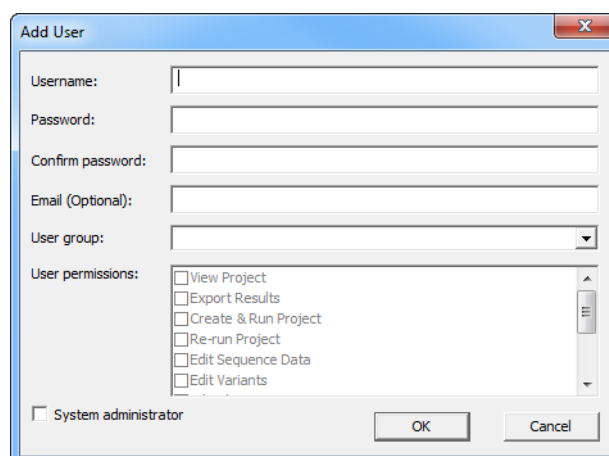Figure 1-28: Add User dialog box



2. Enter the information for the new user:

- In the Username field, enter the appropriate user name.
- In the Password field, enter the password for the user.



The only invalid character is a space. There are no other special requirements or restrictions for the user password. It can adhere to your organization's standards and any other requirements as needed. If you forget or lose this password, it is not recoverable.

- In the Verify field, enter the user password *exactly* as you entered it in the Password field.
- Optionally, in the Email field, enter the email address for the user.



The current version of User Management does not support email notifications; however, you can still enter an email address.

3. Assign the user to a selected group.



By assigning a user to a group assigns the user's permissions for Mutation Surveyor. If the appropriate group is not available, then you must add the group. See ["Managing Groups in Mutation Surveyor"](#) on page 38.

4. Optionally, if the user is to be responsible for User Management in Mutation Surveyor (managing groups and users), then select System administrator.

5. Click OK.

A message opens, indicating that the new user was created successfully.

6. Click OK.

The message closes. The entry for the new user is displayed on the Users tab. The Users tab remains open.

7. Click OK.

The User Management Settings dialog box closes.


To edit a user

You can edit the password, the email address, and the groups for a user. For any user other than the default Administrator user, you can edit the System administrator status. You cannot edit the username for any user. To edit the username, you must delete the user, and then create a new user with a different username. See [“To delete a user” on page 47](#).

1. Select the user that you are editing, and then click Edit User.

The Edit User dialog box opens.

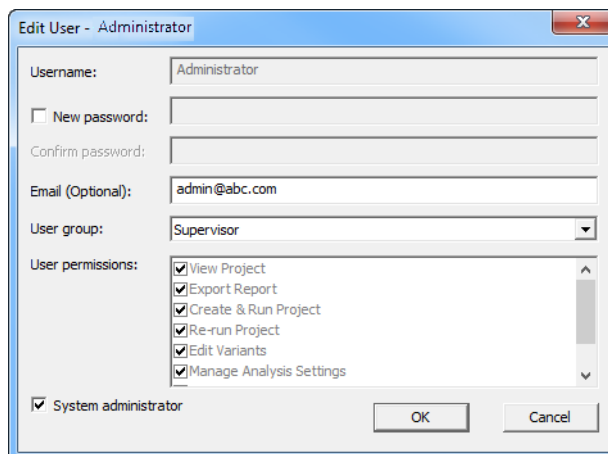Figure 1-29: Edit User dialog box

The screenshot shows the 'Edit User - Administrator' dialog box. It contains several fields: 'Username' is set to 'Administrator'; 'New password' and 'Confirm password' are empty, with a checkbox for 'New password' currently unchecked; 'Email (Optional)' is 'admin@abc.com'; 'User group' is set to 'Supervisor' in a dropdown menu; 'User permissions' are listed with checkboxes for 'View Project', 'Export Report', 'Create & Run Project', 'Re-run Project', 'Edit Variants', and 'Manage Analysis Settings', all of which are checked. At the bottom, there is a checkbox for 'System administrator' which is also checked. 'OK' and 'Cancel' buttons are at the bottom right.

2. Edit the information for the user as needed:

- To edit the password, select New password, and then do the following:
 - In the Password field, enter the password for the user.



The only invalid character is a space. There are no other special requirements or restrictions for the user password. It can adhere to your organization's standards and any other requirements as needed. If you forget or lose this password, it is not recoverable.

- In the Verify field, enter the user password *exactly* as you entered it in the Password field.

- Enter an email address for the user, or edit the existing address as needed.
 - Select a different group for the user.
 - Select or clear the System administrator status for the user.
3. Click OK.
A message opens, indicating that the new user was updated successfully.
 4. Click OK.
The message closes. The entry for the user is updated accordingly on the Users tab.
 5. Click OK.
The User Management Settings dialog box closes.

To delete a user

You cannot delete the default Administrator user. To edit the name for a user, you must delete the user, and then create a new user with a different user name. See [“To add a user” on page 45](#).

1. Select the user that you are deleting, and then click Delete User.
A message opens, indicating that you are deleting the user and asking you to click OK to continue.
2. Click OK.
The message closes, and a second message opens indicating that the selected user was successfully deleted.
3. Click OK.
The second message closes. The entry for the user is removed from the Users tab. The Users tab remains open.
4. Click OK.
The User Management Settings dialog box closes.

Chapter 1

Getting Started with Mutation Surveyor

Chapter 2

Project Setup

To set up a project in Mutation Surveyor, you must understand about the different file types, how to modify the default Mutation Surveyor project settings, and then using this information to set up a project that meets your working needs. You can set up single or multiple projects in Mutation Surveyor. This chapter guides you through the necessary steps for setting up a single project in Mutation Surveyor.



For information about working with multiple Mutation Surveyor projects, including processing and displaying results, see [Chapter 11, “Multiple Projects,”](#) on page 313.

This chapter covers the following topics:

- [“Project File Types” on page 50.](#)
- [“Modifying the Default Mutation Surveyor Project Settings” on page 52.](#)
- [“Setting up a Single Mutation Surveyor Project” on page 70.](#)

Project File Types

A project in Mutation Surveyor can have up to three different types of files—[sample files](#), [reference files](#), and [GenBank files](#). Each file serves a different purpose in the project, each file has a different source, and each file type has its own unique requirements.

Sample file

The sample file is always a required file for a Mutation Surveyor project. The Sample file is the file that contains an unknown base sequence that you want to determine. A sample file must be in one of the following formats—SCF, ABI/ABI, GZ, ESD, or RSD. A sample file can be a single one-directional file or it can be two separate files, with one file containing the forward read and the other file containing the overlapping reverse read.



If your sample files are human DNA, then you have the option of loading only sample files. See [“Automatic GenBank download” on page 51](#).

Reference file

The Reference file contains a known base sequence to which your sample files are compared. The Reference file is always a required file in a Mutation Surveyor project; however, the file source can vary. You can manually load a reference file, or you can have Mutation Surveyor create a synthetic reference file. The loaded file must be in one of the following formats—SCF, ABI/ABI, GZ, ESD, or RSD. If you do not load a reference file *and* your files do not meet the requirements for an Automatic GenBank download (see [“Automatic GenBank download” on page 51](#)) then you *must* have a GenBank file loaded so that Mutation Surveyor can automatically create a synthetic reference file for each contig based on the loaded GenBank file.

GenBank file

A GenBank file is a text-based file that shows the amino acid translation that corresponds to the base coding sequence (CDS). A GenBank file is always an optional file for a Mutation Surveyor project. The required file format is either GBK, GB, or SEQ. If you load a GenBank file, then Mutation Surveyor automatically creates a synthetic reference file for each contig based on the loaded GenBank file. Additionally, FASTA or FA files can be loaded as a GenBank file, but no CDS annotation will be provided.



You can load a GenBank file even you have loaded reference files for a project. This can be useful if you want to determine systematic errors in your amplicons such as PCR artifacts. See [“GenBank/Reference Comparison” on page 58](#).

Automatic GenBank download

If your sample files are human DNA, then you have the option of loading only sample files for a Mutation Surveyor project. When you run the project, by default, Mutation Surveyor automatically accesses a GenBank database on SoftGenetics's bwt server (this database contains Build 37 of the human genome) and downloads the correct GenBank reference file (gene) to which to match the sample files.



You can change the default location from which to download the GenBank reference file. See [“Automatically Download GenBank Files”](#) on the [“”](#) on page 66.

Modifying the Default Mutation Surveyor Project Settings

When you run a mutation project, you can use the default values for the mutation project settings, or you can [modify](#) any or all the settings as needed. You use the Mutation Project Settings dialog box to modify the settings for a mutation project. You can modify the value for two types of project settings—*analysis settings* and *output/display settings*.

- **Analysis settings**—Analysis settings determine how your data is analyzed and therefore, which positions in your sample traces are called as mutations and which are not. Any difference that is found between your sample traces and reference traces must meet *all* the project analysis settings to be called as a mutation. You must modify the analysis settings *before* running a project. The Input tab, the Contig tab, and the Database tab contain only analysis settings. Other analysis settings are found on the following tabs:

Tab	Project Analysis Settings
Mutation	<ul style="list-style-type: none">• Mutation Detection• Peak Fronting• Base Labeling• Data Processing
Display	<ul style="list-style-type: none">• Check 2D Small Peaks (Mosaic)
Others	<ul style="list-style-type: none">• Check Gene Direction• Methylation

- **Output/display parameters**—Output/display parameters do not affect the analysis of your data. They simply determine how your data is displayed to you in the different frames of the Graphical Analysis Display (GAD) for a project. For example, is an identified mutation displayed in red or blue in the Mutation table of the GAD? You can modify the output/display settings (found on the Mutation tab, the Output tab, the Display tab, the 2 Directions tab, the Others tab, and the Database tab) before *or* after running a project.



For detailed information about the Graphical Analysis Display for a project, see [Chapter 3, “Graphical Analysis Display,” on page 75.](#)

To modify the default Mutation Surveyor project settings

1. On the Mutation Surveyor main menu, click Process > Settings.
The Mutation Project Settings dialog box opens. The Contig tab is the active tab.
2. Do one of the following:
 - Click Load Settings to browse to and select a saved Settings file (.ini file) to analyze and display the project results according to the settings in the file.
 - Modify the project values as needed on any of the tabs. See:
 - “Input tab” on page 54.
 - “Contig tab” on page 55.
 - “Mutation tab” on page 59.
 - “Output tab” on page 62.
 - “Display tab” on page 64.
 - “2 Directions tab” on page 65.
 - “” on page 66.
 - “Database tab” on page 69.



To reset all values on all tabs to their default values in a single step, click Default.

3. Optionally, before you process the files, click Save Settings to save the settings that you have specified to a Settings.ini file.




You can always load this file at a later date and process other data files according to the saved settings in the file.

4. Click OK.
 - Unless you modify the values again, all *new* projects are carried out according to the options that you have loaded/specified.
 - If you load an existing project, the options revert to those set for the project. If you run a new project after loading an existing project, the new project is carried out according to the existing project settings.
5. Continue to “Setting up a Single Mutation Surveyor Project” on page 70.

Input tab

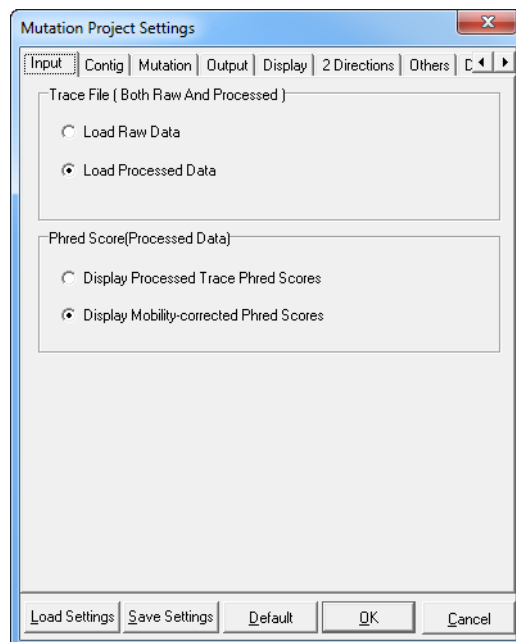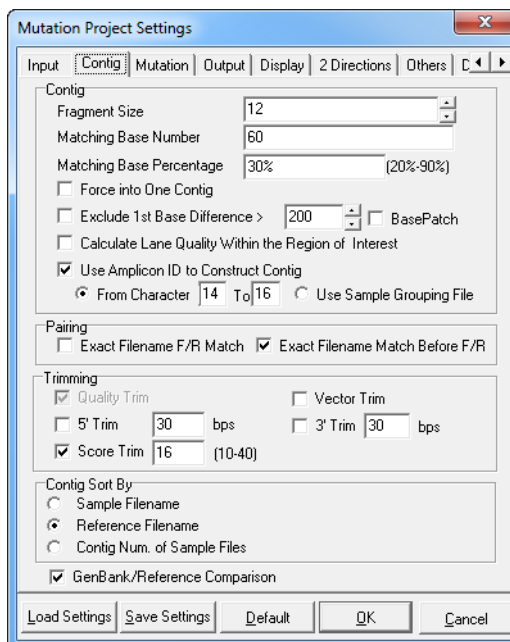
Figure 2-1: Mutation Project Settings dialog box, Input tab



Option	Description
Trace File (both Raw and Processed)	<p>These options are applicable only if you loaded AB1/ABI sample files.</p> <ul style="list-style-type: none"> Load Raw Data—Loads only the raw trace (no baseline subtraction, no smoothing, no mobility shift correction, or no base calling) in the AB1/ABI sample files for analysis. <p>Note: If the project contains only a few traces and some of the traces have heterozygous substitutions close to the start of the traces, then the mobility shift correction that Mutation Surveyor carries out might cause a single heterozygous peak to be corrected into two homozygous peaks, which results in a false positive of a single base insertion.</p> <ul style="list-style-type: none"> Load Processed Data—The default value. Loads only the processed data in the AB1/ABI sample files for analysis.
Phred Score (Processed Data)	<ul style="list-style-type: none"> Display Processed Trace Phred Scores—Load the Phred scores that are contained within the traces for each base. Display Mobility-corrected Phred Scores—Load the Phred scores that Mutation Surveyor has calculated for each base.

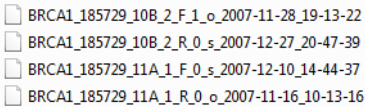
Contig tab

Figure 2-2: Mutation Project Settings dialog box, Contig tab



Option	Description
Contig	
Fragment Size	The number of sequential bases that are used for base matching. The default value is 12. The allowed range is 10 -16 bases. Mutation Surveyor evaluates a string of 12 bases in the sample trace and compares this string to the reference trace. If the string matches the reference trace, Mutation Surveyor moves to the next series of 12 bases and makes another comparison. Mutation Surveyor continues with these matching steps until it is able to determine whether the sample trace belongs to an existing contig or whether the sample trace belongs in its own contig.
Matching Base Number	The minimum number of bases that must match between two sample traces for the traces to be placed in the same contig. The default value is 60. The allowed range is 20 -100 bases.
Matching Base Percentage	The minimum percentage of bases that must match between two sample traces for the traces to be placed in the same contig. The default value is 30%. The allowed range is 20 - 90%.

Option	Description
Force Into One Contig	<p>Place sample traces that meet all the matching base criteria and that differ in the starting base positions by ≤ 300 bps into one contig.</p> <p>Note: Selected by default. If the quality of the sequences differs and sequences might be placed into either of two contigs, the samples might be placed into a different contig than expected. If samples that should be grouped into the same contig are not being grouped correctly using the Force Into One Contig option, then select Exclude 1st Base Difference $> []$ and accept the default value of 200 bps.</p> <p>Note: If you are using a Grouping File to correctly pair your sample files, make sure that this option is <i>not</i> selected. See "Pairing" on page 57.</p>
Exclude 1st Base Difference $> []$	<p>Place sample traces that meet all the matching base criteria but that differ in the starting base positions by more than "n" number of bases into two contigs. The default value is 200 bps.</p>
<p>Note: Force into One Contig and Exclude First Base Difference sort contigs using only base calling (peak quality is not evaluated) and therefore, are carried out before quality trimming. (See "Quality Trim" on page 58.) If neither Force into one Contig nor Exclude 1st Base Difference $> []$ is selected, then:</p> <ul style="list-style-type: none"> Place sample traces with > 500 bps that meet all the matching base criteria and that differ in the starting base positions by ≤ 250 bps into one contig. Place sample traces with > 300 bps and < 500 bps that meet all the matching base criteria and that differ in the starting base positions by $< \text{read length}/2$ bps into one contig. Place sample traces with < 300 bps that meet all the matching base criteria and that differ in the starting base positions by < 150 bps into one contig. 	
Base Patch	<p>If you are attempting to make mutation calls in the first 50 bases of sample traces, but the calls are being missed, then select Base Patch to correct the poor mobility shift at the starting of traces. With this option selected, Mutation Surveyor evaluates the spacing between each peak, and based on the results of this evaluation, Mutation Surveyor might make a different base call than what is included in the original trace.</p> <p>Note: If you are analyzing AB1/ABI files and these issues are still not corrected with Base Patch selected, select Load Raw Data on the Options tab. See "Trace File (both Raw and Processed)" on page 54.</p>
Calculate Lane Quality Within the Region of Interest	<p>Calculate the lane quality for the Region of Interest in the GenBank file, not the entire trace. If the ROI extends beyond the comparison region, then the ROI quality score is adjusted by the percent overlap of the aligned region and the ROI. For example, if the ROI length is 63 bp and the length of the overlap between the aligned region and the ROI is 8, then the percent overlap is $8/63 = 0.127$. The currently calculated quality score, which in this example is 88, is then adjusted by this percentage to give the new quality score or $(0.127)(88) = 11$.</p> <p>Note: To define a Region of Interest (ROI) in a GenBank file, see "Advanced GBK File Editor Tool" on page 231.</p>



Option	Description
Use Amplicon to Construct Contig	<p>Two options:</p> <ul style="list-style-type: none"> If all the filenames in a project contain an amplicon-specific designator, then select Select from Character X to Y and enter the starting and ending character positions of this designator. For example, Figure 2-3 shows a series of sample file names with the amplicon-specific designators of 10B and 11A located in positions 14 through 16 of the filename. <p><i>Figure 2-3: Amplicon-specific designators in sample file name</i></p>  <p>Note: If ROIs in the GenBank file are labeled with their Amplicon IDs, then selecting this option ensures that the appropriate ROI is evaluated for each amplicon.</p> <ul style="list-style-type: none"> If your sample file names do not follow any file-naming standards, and you used a Grouping File to load the sample files, then select Use Sample Grouping file to help organize the contigs correctly. See “2D Filename Match Editor tool” on page 261. <p>Note: If you do use a Grouping File, then when you are adding sample files to for your Mutation Surveyor project, make sure that Load Grouping File is selected on the Open Files dialog box. When you click Add in the Sample files pane, select the Grouping File. All the sample files that are defined in the Grouping File are added for the Mutation Surveyor project.</p>
Pairing	
Exact Filename F/R Match	<p>Select this option to pair your forward and reverse sample files if your filenames differ only by the directional identifier (f or r). For example:</p> <p>MMipOp260_f_018.ab1/MMipOp260_r_018.ab1</p>
Exact Filename Match Before F/R	<p>Select this option to pair your forward and reverse sample files if your filenames are identical before the directional identifier (f or r), but vary after the directional identifier. For example:</p> <p>MMipR260_f_017.ab1/MMipR260_r_018.ab1</p>
<p>Note: In some situations, neither of the above automatic file pairing options suits your current needs. For example, your files might require a complete naming overhaul before you can use either option. Additionally, you might want to set aside blocks of data for analysis at a later date. In such a situation, you should create a Grouping File to load your sample files. See “2D Filename Match Editor tool” on page 261.</p>	

Option	Description
Trimming	
Quality Trim	<p>Eliminates low quality data at the starting and the end of a sample trace. Quality trimming trims the sequence data based on the base quality. By default, when the S/N Ratio is greater than 15 for 7 out of 9 sequential bases, then mutation detection is carried out in this region.</p> <p>Note: Quality Trim is always implemented in Mutation Surveyor which is why the option is always selected and always unavailable. If you select other trimming options, then all the selected trimming methods are compared. The trimming method that results in the greatest number of bases being trimmed is the trim method that is applied. Quality Trimming is then applied after the selected trimming method.</p>
Vector Trim	Trims the sample sequence according to a starting vector sequence and an ending vector sequence that you define using the SEQ File Editor tool. See "SEQ File Editor Tool" on page 246.
5' Trim []	Trim the 5' end of the sample trace by the indicated number of bases.
3' Trim []	Trim the 3' end of the sample trace by the indicated number of bases.
Score Trim []	Reject regions for mutation detection in which the base quality for 7 out of 9 sequential bases falls below the specified value. For example, if the Score Trim is set to a value of 10, and the S/N Ratio is 8 for 7 out of 9 sequential bases, then mutation detection is not carried out in this region. The default value is 16.
Contig Sort By	
	<p>Determines the sort order of multiple contigs in a project.</p> <ul style="list-style-type: none"> • Sample Filename • Reference Filename • Contig Num. of Sample Files
GenBank/Reference Comparison	If you have both GenBank and reference files loaded, and this option is selected, then Mutation Surveyor creates a synthetic sample trace from the base sequence of the GenBank file and the reference trace to classify the variance between the GenBank file and your reference files. You can, in turn, use this variance to determine systematic errors in your amplicons such as PCR artifacts.

Mutation tab

Figure 2-4: Mutation Project Settings dialog box, Mutation tab

Option	Description												
Mutation													
Mutation Score	<p>Indicates the level of confidence that a called mutation is a true mutation. The score is based on the concept of Phred scores, where quality scores are logarithmically linked to error probabilities as shown in Figure 2-5.</p> <p><i>Figure 2-5: Phred scores and error probabilities</i></p> <table><tr><th>Phred quality score</th><th>Probability that the base is called wrong</th><th>Accuracy of the base call</th></tr><tr><td>10</td><td>1 in 10</td><td>90%</td></tr><tr><td>20</td><td>1 in 100</td><td>99%</td></tr><tr><td>30</td><td>1 in 1,000</td><td>99.9%</td></tr></table> <p>The values for “Overlapping Factor,” “Dropping Factor,” and “S/N Ratio” are used in the calculation of the Mutation Score. The lower the score, the greater number of mutation calls, but the confidence in the accuracy of the calls is reduced. The default value is 5.00.</p> <p>When calculating heterozygous and homozygous indel quality score values, note that:</p> <ul style="list-style-type: none">• Heterozygous Indel Score reflects the ratio of S/N before and after the start of the event.• Homozygous Indel Score evaluates the spacing between the flanking aligned bases as well as signal in the uncalled region.	Phred quality score	Probability that the base is called wrong	Accuracy of the base call	10	1 in 10	90%	20	1 in 100	99%	30	1 in 1,000	99.9%
Phred quality score	Probability that the base is called wrong	Accuracy of the base call											
10	1 in 10	90%											
20	1 in 100	99%											
30	1 in 1,000	99.9%											
Mutation Height	<p>The minimum height (in RFU) required for a peak in the mutation trace to be called as a mutation peak. The default value is 500 RFU.</p>												

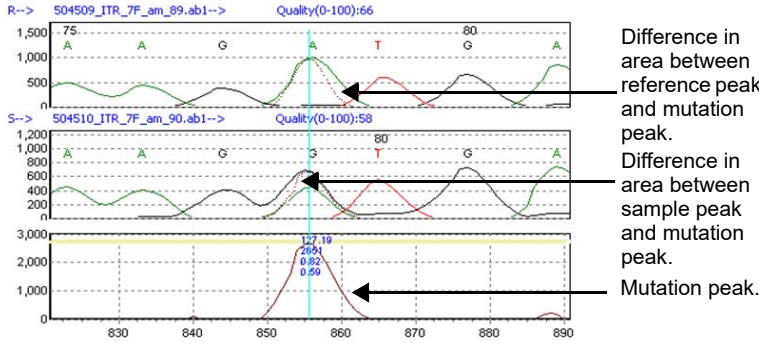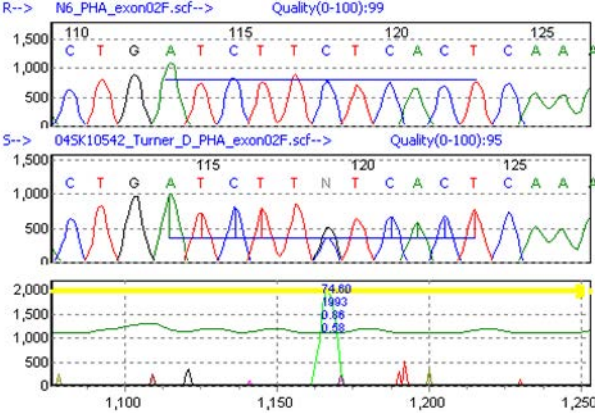
Option	Description
Overlapping Factor	<p>The <i>average</i> of the following two values:</p> <ul style="list-style-type: none"> The difference in the area between the reference peak and the mutation peak. The difference in the area between the sample peak and the mutation peak. <p>For an ideal mutation peak, the overlapping factor has a value of 1.0. In reality, this value is typically less than one. The overlapping factor must be greater than or equal to the specified value for the variant to be called a mutation. The default value is .20, or 20%.</p> <p>Figure 2-6: Calculating the Overlapping Factor</p>  <p>Difference in area between reference peak and mutation peak.</p> <p>Difference in area between sample peak and mutation peak.</p> <p>Mutation peak.</p>
Dropping Factor	<p>Measures the drop in peak intensity at the variant position by normalizing the peak intensity between the reference peak and the sample peak at positions two through five on both sides of the variant. The dropping factor for a homozygous mutation is 1, or 100%. The dropping factor for a heterozygous mutation is 0.5, or 50%. The dropping factor must be greater than or equal to the specified value for the variant to be called a mutation. The default value is .10, or 10%.</p> <p>Figure 2-7: Calculating the Dropping Factor</p> 

Option	Description
S/N Ratio	<p>Signal (S) is defined as the mutation peak intensity and noise (N) is defined as the median noise intensity for the mutation peak. The Signal to Noise Ratio (S/N Ratio) determines how large the signal must be for a mutation peak relative to the noise for the mutation to be registered.</p> <p>Note: The minimum noise at a given base position is 250 RFU.</p> <p>Note: S/N Ratio is output to some reports as SnRatio.</p>
Mutation Detection	
	<p>Determines the level of sensitivity that Mutation Surveyor uses to call mutations.</p> <ul style="list-style-type: none"> • High sensitivity • Medium sensitivity <p>Note: There are many functions that change the parameters between these two options. For example, if in 2D data, one trace indicates a mutation at a location, but the other trace does not, Mutation Surveyor checks for local base quality and decides whether to add or delete the mutation. In other situations, statistical analyses are carried out for the contigs to check for possible missed mutations, reported mutations are checked for missed mutations and so on. If you are concerned that your data might be showing a high number of false positives, then set the sensitivity to medium to reduce the number of false positives.</p>
Allow for Hypervariable Mutation Detection	<p>Select this option to have Mutation Surveyor automatically extend the starting and ending positions for the aligned comparison region if both of the following conditions are met:</p> <ul style="list-style-type: none"> • At least 4 of 10 consecutive bases outside of the aligned comparison region have a Phred-like score ≥ 15. • At least 8 of 12 consecutive bases outside of the aligned comparison region are identical to the reference bases. <p>Note: By extending this region, more potential mutations can be called, which is especially useful for samples that are known to have a high mutation frequency such as viruses and the human leukocyte antigen system (HLA).</p>
Delete Indels/het_indels Automatically	<p>Select this option to have Mutation Surveyor automatically delete all insertion and deletion calls from the project.</p> <p>Note: Selecting this option helps to eliminate false positives that are common to hypervariable sequences.</p>
Allow Software to Delete Mutations	<p>Select this option to have Mutation Surveyor automatically delete illogical mutation calls and mutation calls that have high background noise (> 250 RFU).</p>
Delete Indels Outside of ROI	<p>If selected, then Mutation Surveyor automatically deletes insertions and/or deletions that fall outside the ROI.</p>

Option	Description
Peak Fronting	
Correction	<p>Peak Fronting is often seen as a smaller peak that is in front of a major peak of <i>the same color</i>, which means that the peaks represent the same base in the sequence. This small peak is often the result of PCR reactions of a dye-labeled terminator when concentrated reagents are used. With Peak Fronting Correction, the minor peak is reduced in size relative to the major peak. The default value is 1%.</p> <p>Note: Peak Fronting Correction might reduce the mutation intensity of some peaks in the mutation trace.</p> <p>Note: Newer chemistries often have less peak fronting, so lowering this value can increase the sensitivity of mutation detection.</p>
Base Labeling - Determines the point on a peak that is used for determining peak height.	
<ul style="list-style-type: none"> Original Peak Center 	<ul style="list-style-type: none"> Use the base position that was defined by the original base calling software. Use the bases at the center of the peak to determine peak height.
Data Processing	
Dye blob baseline noise removal	<p>If the sequence traces exhibit broad peaks caused by a dye blob, then the baseline might introduce artifacts into the trace. Select this option to remove the baseline noise that is caused by a dye blob.</p> <p>Note: Baseline removal might reduce the mutation intensity of some peaks in the mutation trace.</p>

Output tab

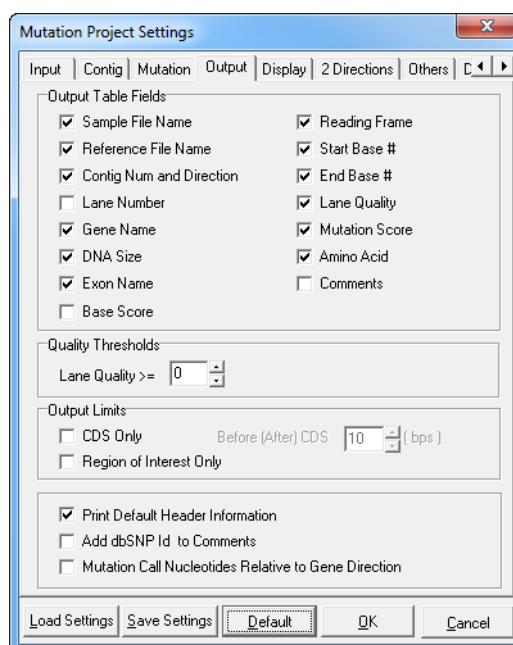
Figure 2-8: Mutation Project Settings dialog box, Output tab



Option	Description
Output Table Fields	<p>Select the fields that are to be displayed in the Viewing and Reporting pane for the different standard mutation reports.</p> <p>Note: Base Score - The average value of the Phred scores in the comparison region. The Phred score can be derived by either the Processed Trace Phred scores or the Mobility-corrected Phred scores.</p> <p>Note: The selections that you make here affect all standard mutation reports en masse. You cannot select different fields for different reports. See Chapter 5, "Mutation Surveyor Standard Reports," on page 125.</p>
Quality Thresholds	<p>Lane Quality >= []—Lane quality is a measure of the S/N Ratio within a sample trace prior to alignment to a reference trace. After the sample trace is aligned to the reference trace, peak spacing in the sample trace is adjusted, which often results in higher lane quality. Lane quality ranges from 0 to 100, with 100 being the highest quality. If multiple reference files are available, Mutation Surveyor chooses the reference file with the highest lane quality.</p> <p>Note: "Low Quality" is displayed in the Lane Quality column in reports for a trace with lane quality that falls below the Quality Threshold.</p>
Output Limits	
CDS Only • Before (After) CDS	<p>Limits the graphical output to the CDS.</p> <ul style="list-style-type: none"> Available only if CDS Only is selected. Specify the number of bases that are to precede/follow the coding region.
Region of Interest Only	Limits the graphical output to Regions of Interest.
Print Default Header Information	Selected by default. Prints the default Header information on a variety of standard mutation reports.
Add dbSNP Id to Comments	Displays the dbSNP identifier in the Comments field for a standard mutation report when Comments are selected for output.
Mutation Call Nucleotides Relative to Gene Direction	Make the mutation call based on the gene orientation. To make a mutation call for a gene on the reverse strand, a reverse complement is generated.

Display tab

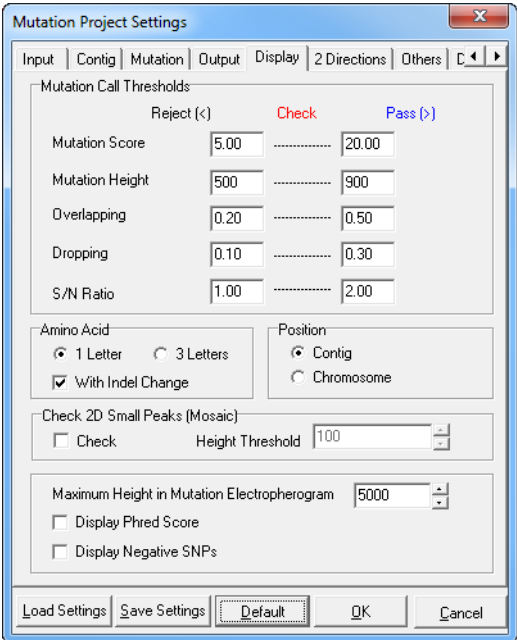Figure 2-9: Mutation Project Settings dialog box, Display tab




Option	Description
Mutation Threshold	<p>Determines how an identified mutation is displayed—red font (low confidence) or blue font (high confidence)—in the different mutation reports. The values in the Reject column indicate the upper rejection boundary. The values in the Pass column indicate the lower acceptance boundary.</p> <ul style="list-style-type: none">• If the value <i>for any</i> of the mutation settings is less than the rejection boundary, then the mutation is rejected and is not displayed in the different mutation reports.• If the values <i>for all</i> the mutation settings fall within the indicated ranges, the mutation is accepted and displayed in red in the different mutation reports.• If the values <i>for all</i> the mutation settings exceed their acceptance boundaries, then the mutation is accepted and displayed in blue in different mutation reports. <p>Note: For 2D data, the mutation scores (F and R) are summed and the sum score is used to determine if the mutation call is rejected, passed, or put into check.</p>
Amino Acid	<ul style="list-style-type: none">• 1 Letter—Display 1 Letter IUPAC amino acid code in the mutation reports and views.• 3 Letter—Display 3 Letter IUPAC amino acid code in the mutation reports and views.• With Indel Changes—Selected by default. Displays the amino acid output including indels. Clear this option to display the amino acid output without indels.

Option	Description
Position	<ul style="list-style-type: none"> Contig—Show the position within the GenBank file for the queried contig in the Nucleotide Sequence frame and Mutation table of the GAD, and in the Project Reviewer report. Chromosome—Show the chromosome position for the queried contig in the Nucleotide Sequence frame and Mutation table of the GAD, and in the Project Reviewer report. <p>Note: The Chromosome option works correctly only if the GenBank file contains chromosome information; otherwise, Mutation Surveyor defaults to showing the contig information.</p>
Check 2D Small Peaks (Mosaic)	<ul style="list-style-type: none"> Check—If selected, Mutation Surveyor searches for minor alleles that are present within the baseline of bi-directionally aligned peaks and indicates their presence with a short green bar. <p>Note: This option is useful for somatic mutation detection.</p> <ul style="list-style-type: none"> Height Threshold—Available only if Check 2D Small Peaks is selected. Specify the RFU value for the smallest peak that is to be detected.
Maximum Height in Mutation Electropherogram	If the height of a peak in the mutation trace exceeds this maximum value, then Mutation Surveyor normalizes the peak to this maximum height. The default value is 5000 RFU.
Display Phred Score	If selected, the Phred-like score for each peak in the mutation trace is displayed in the sample trace directly below the base call.
Display Negative SNPs	If selected, positions with reported variations that do not contain the SNP are indicated in the Graphical Analysis Display (GAD) by showing the possible alleles in green at the top of the mutation trace.

2 Directions tab

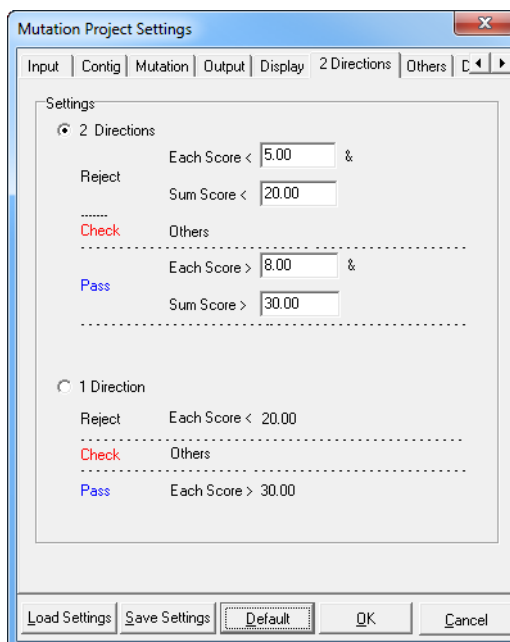Figure 2-10: Mutation Project Settings dialog box, 2 Directions tab



Option	Description
2 Directions	Select this option if 2D is being analyzed.
1 Direction	Select this option 1D data is being analyzed. Note: You can select the 2 Directions options for 1D data.

For 2D data, the mutation scores (F and R) are summed and the sum score is used to determine if the mutation call is rejected, passed, or put into check. [Figure 2-11](#) below illustrates how 2 Directions settings work based on the default values for a mutation project.

Figure 2-11: Examination of 2D mutation data

F	R	Sum	Result
3	3	6	Fail
6	10	16	Check
4	40	44	Check
12	18	30	Pass

For 1D data that is being analyzed with 2D settings, a mutation is called if it meets one of the following parameters; otherwise, it is rejected.

- Score > SumScoreReject
- Score > SumScoreReject*0.8 and Dropping Factor >= 0.3
- Score > SumScoreReject*0.6 and Dropping Factor >= 0.35
- Score > SumScoreReject*0.4 and Dropping Factor >= 0.4


[Figure 2-12](#) below illustrates these parameters based on the default value of 20 for the SumScoreReject.


Figure 2-12: Examination of 1D mutation data analyzed with 2D settings

Score	SumScore Reject Factor	SumScoreReject	Dropping Factor	Result
25		20	N/A	Pass
15	0.8	20	0.3	Fail
15	0.6	20	0.35	Pass
15	0.4	20	0.4	Pass

Others tab

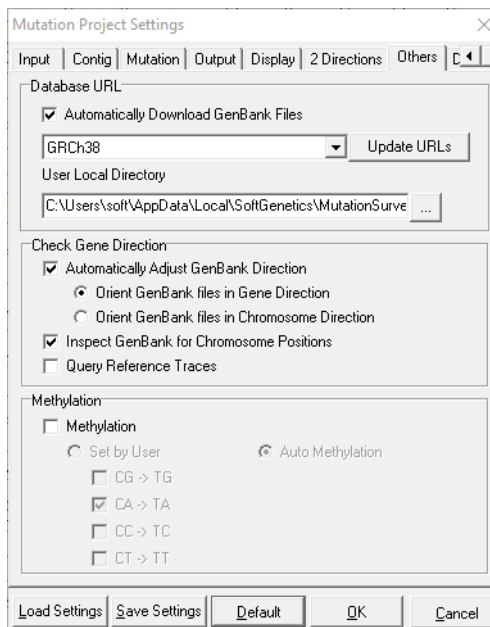Figure 2-13: Mutation Project Settings dialog box, Others tab



Option	Description
Database URL	
Automatically Download GenBank Files	<ul style="list-style-type: none"> If your sample files are human DNA, then you have the option of loading only sample files for a Mutation Surveyor project. When you run the project with this option selected, Mutation Surveyor automatically accesses the database that is specified by the URL and downloads the correct GenBank file (gene) to which to match the sample files. <p>Note: The default URL is GRCh38. Click on the drop-down arrow to change to GRCh37 or GRCh36. Click on Update URLs to manually add a new database URL.</p>
User Local Directory	<ul style="list-style-type: none"> If the GenBank file to which you want to match your sample files resides in a directory on your local computer, clear the Automatically Download GenBank Files option, and then click the Browse button  to browse to and select the local directory path.
Check Gene Direction	
Automatically Adjust GenBank Direction	<p>Selected by default. Determines whether the queried gene in the GenBank file is complemented based on its location (plus (+) or minus (-) strand).</p> <ul style="list-style-type: none"> Selected by default. Orient GenBank files in Gene Direction Orient GenBank files in Chromosome Direction

Option	Description
Inspect GenBank for Chromosome Positions	<p>Selected by default. Determines if the GenBank file header contains the genome build, the chromosome name, and the chromosome position. If this information is present, then chromosome information can be displayed in the Nucleotide Sequence frame and the Mutation table in the GAD, and the Project Reviewer report; otherwise, only the contig information can be displayed.</p> <p>Note: If you clear this selection, then Mutation Surveyor does not notify you when chromosome options will not work correctly; however, if chromosome information is contained in the GenBank file, then the chromosome options still work correctly.</p>
Query Reference Traces	<p>Not selected by default. If a GenBank file is not loaded for the project, but reference traces are, then the sample traces are aligned only to the reference traces. If selected, Mutation Surveyor downloads the appropriate GenBank file for the project, and aligns both the reference and sample traces to the GenBank file for annotation.</p>
Methylation	
Methylation <ul style="list-style-type: none"> • Set By User 	<p>Select this option if methylated samples are to be analyzed.</p> <ul style="list-style-type: none"> • Select this option to manually specify which forms of methylation Mutation Surveyor should detect.
<ul style="list-style-type: none"> • Automethylation 	<ul style="list-style-type: none"> • Selected by default when Methylation is selected. Mutation Surveyor automatically detects methylated sequences.
<p>Note: For information about methylation detection in Mutation Surveyor, see “Methylation Detection” on page 301.</p>	


Figure 2-14: Database URL, Update URLs button

The screenshot shows a window titled "Update Reference URLs". Inside the window, there are three labeled input fields. The first, "Reference Server", is a dropdown menu currently showing "GRCh38". The second, "Reference Server URL", is a text box containing "http://bwt38.softgenetics.com/query.php". The third, "Reference Server Name", is a text box containing "GRCh38". At the bottom right of the window are two buttons: "Delete" and "Update".

Database tab

Some GenBank files contain only NT-accession numbers and not chromosome information. You specify the database on the Database tab that is to be used to translate these contig positions into chromosome positions.

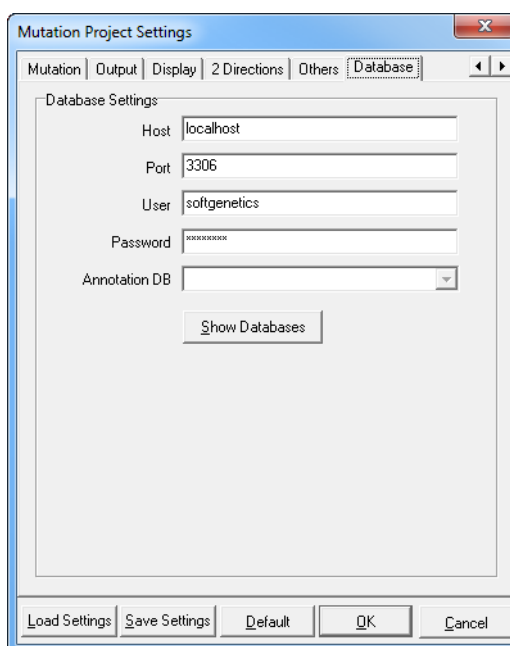Figure 2-15: Mutation Project Settings dialog box, Database tab



Option	Description
Host	The address for the Annotation Database server. The default value is set to localhost, which assumes that the server is installed on the same computer as Mutation Surveyor. If this is correct, then leave the default value as-is.; otherwise, modify the value accordingly.
Port	Leave this option set to the default value of 3306 only if no other application is already using MySQL on the port; otherwise, modify the value accordingly.
User	The user name for MySQL. Leave this option set to the default value of softgenetics unless your administrator has changed it.
Password	The password for MySQL. Leave this option set to the default value unless your administrator has changed it.
Show Database	To verify the MySQL connection and populate the list of available annotation databases, click this option. Note: If the MySQL connection cannot be verified, an error message opens indicating this, and that the list of annotation databases cannot be populated. Verify that you have imported the necessary annotation databases (see the Reference and Track Manager Tool User's Manual) and/or contact for your administrator for assistance.
Annotation DB	Select the appropriate annotation database.

Setting up a Single Mutation Surveyor Project

When you [set up](#) any Mutation Surveyor project, the following caveats apply:

- For one directional data, you can enter 48 or 2000 sample files depending on the version of Mutation Surveyor that you purchased. Only a single reference file is required.
- For two directional data, you can enter 24 or 1000 pairs of sample files depending on the version of Mutation Surveyor that you purchased. Both a forward reference file and a reverse reference file are required.
- Sample files and reference files are limited to a read length of 2500 bps. GenBank files are limited to a read length of 10 Mbp per file.
- If your amplicons do not overlap and your forward and reverse directional pairs have similar start and end locations, then the forward and reverse files with similar names that have been grouped in the same contig are automatically paired.

To set up a Mutation Surveyor project




The following procedure details how to set up a single new Mutation Surveyor project. For information about setting up and working with multiple Mutation Surveyor projects, see [Chapter 11, “Multiple Projects,”](#) on page 313.

1. Confirm that your Mutation Surveyor project options are correctly set, or modify them as needed. (See [“Modifying the Default Mutation Surveyor Project Settings”](#) on page 52.)
2. On the Mutation Surveyor main menu, click File > Open Files.

The Open Files dialog box opens.

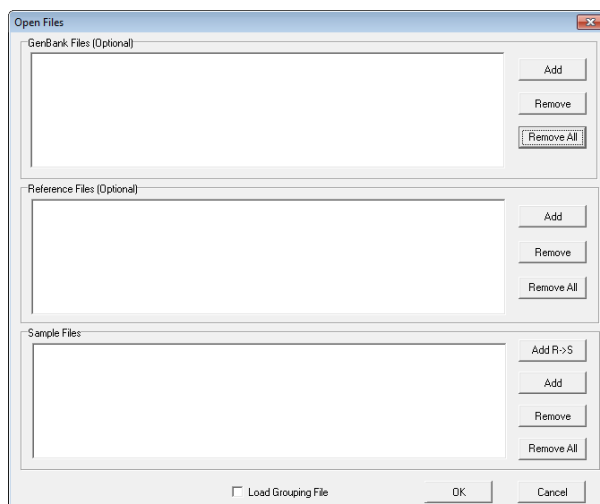Figure 2-16: Open Files dialog box




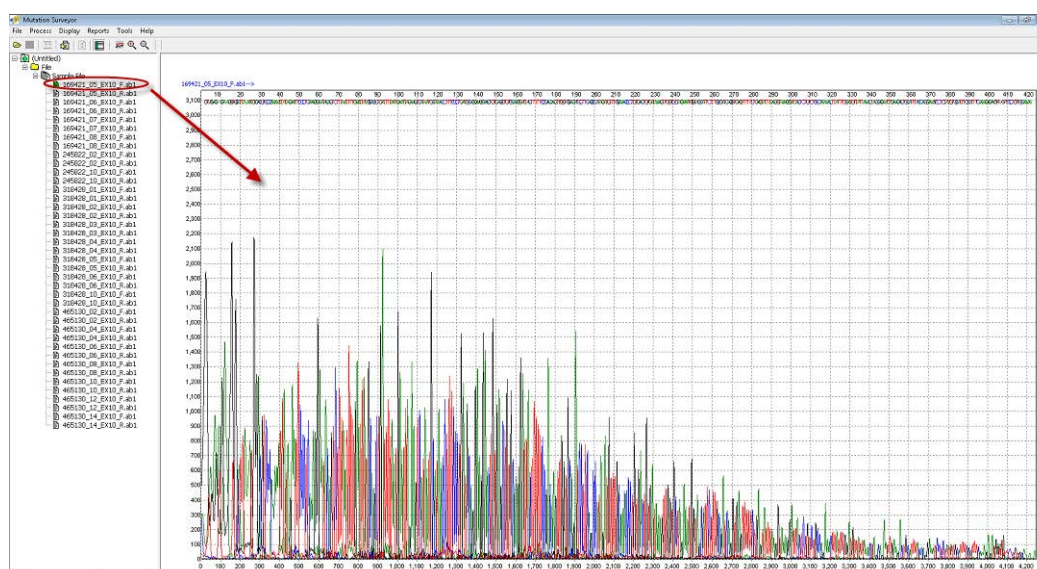
3. For each file type that you are adding to the project:
 - Click Add in the appropriate pane. (For example, to add a GenBank file, click Add in the GenBank Sequence File pane.)
 - In the Open <File> dialog box, browse to and select the files that you are adding to the project. (CTRL-click to select multiple files.)
 - Click Open to add the files to the project, and then click OK.
4. Optionally, to view the data before it is aligned, expand the Sample file folder in the Browser pane, and then double-click a sample file. You can do any of the following to change the view of the data:
 - Zoom in - Click and hold the left mouse button and draw a box from the upper left hand corner of the pane towards the lower right hand corner. A box is formed around the area that being reduced for viewing.
 - Zoom out - Click and hold the left mouse button and draw a box from the lower right hand corner of the pane towards the upper left hand corner. (The magnification for zooming out is always 100%.)
 - Move left/right - Click and hold the right mouse button, and then drag the display to the left or right.
 - If the data does not include basecalling, or if you selected the Mutation Project Setting “Load Raw Data”, click on the active “Raw Data to Processed Data” icon () to view the basecalling as processed by Mutation Surveyor. See “Input tab” on page 54.

Figure 2-17: Viewing unaligned data for a sample file



- On the Mutation Surveyor main menu, click Process > Run.

Mutation Surveyor aligns the nucleotide sequences of the sample traces to the reference traces and groups the sample traces into contigs based on their degree of overlap. (See “[Contig directory for a Mutation Surveyor project](#)” below.) The reference traces and sample traces are then compared at a trace by trace level to identify substitutions, insertions, and deletions. All the variants are aligned in the Mutation report, which is automatically displayed in the Viewing and Reporting pane after the project is run.

Figure 2-18: Mutation Report in Viewing and Reporting pane

No.	Sample File	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Qual	Mut#	Mutation1	Mutation2	Mutation3
1	BRCA1_185729_10	NT_010783_411956	1-F	BRCA1	10	2	31054	31392	339	25	0			
2	BRCA1_185729_10	NT_010783_411956	1-F	BRCA1	10	2	31054	31392	339	25	0			
3	BRCA1_264381_10	NT_010783_411956	1-F	BRCA1	10	2	31044	31392	349	25	0			
4	BRCA1_264381_10	NT_010783_411956	1-F	BRCA1	10	2	31044	31392	349	25	0			
5	BRCA1_185729_10	NT_010783_411956	1-R	BRCA1	10	2	31036	31383	348	26	0			
6	BRCA1_185729_10	NT_010783_411956	1-R	BRCA1	10	2	31036	31383	348	26	0			
7	BRCA1_264381_10	NT_010783_411956	1-R	BRCA1	10	2	31036	31384	349	25	0			
8	BRCA1_264381_10	NT_010783_411956	1-R	BRCA1	10	2	31036	31384	349	25	0			
9	BRCA1_185729_10	NT_010783_411956	2-F	BRCA1	10	2	31301	31575	275	6	1	31520A>AG_356		
10	BRCA1_185729_10	NT_010783_411956	2-F	BRCA1	10	2	31301	31575	275	6	1	31520A>AG_356		
11	BRCA1_264381_10	NT_010783_411956	2-F	BRCA1	10	2	31296	31575	280	6	1	31520A>AG_356		
12	BRCA1_264381_10	NT_010783_411956	2-F	BRCA1	10	2	31296	31575	280	6	1	31520A>AG_356		
13	BRCA1_185729_10	NT_010783_411956	2-R	BRCA1	10	2	31290	31559	270	11	1	31520A>AG_356		
14	BRCA1_185729_10	NT_010783_411956	2-R	BRCA1	10	2	31290	31559	270	11	1	31520A>AG_356		
15	BRCA1_264381_10	NT_010783_411956	2-R	BRCA1	10	2	31290	31573	284	10	1	31520A>AG_356		
16	BRCA1_264381_10	NT_010783_411956	2-R	BRCA1	10	2	31290	31573	284	10	1	31520A>AG_356		
17	BRCA1_185729_11	NT_010783_411956	3-F	BRCA1	10	2	31510	31948	439	11	0			
18	BRCA1_185729_11	NT_010783_411956	3-F	BRCA1	10	2	31510	31948	439	11	0			
19	BRCA1_264381_11	NT_010783_411956	3-F	BRCA1	10	2	31493	31948	456	10	0			
20	BRCA1_264381_11	NT_010783_411956	3-F	BRCA1	10	2	31493	31948	456	10	0			
21	BRCA1_185729_11	NT_010783_411956	3-R	BRCA1	10	2	31486	31934	449	9	0			
22	BRCA1_185729_11	NT_010783_411956	3-R	BRCA1	10	2	31486	31934	449	9	0			
23	BRCA1_264381_11	NT_010783_411956	3-R	BRCA1	10	2	31486	31934	449	9	0			
24	BRCA1_264381_11	NT_010783_411956	3-R	BRCA1	10	2	31486	31934	449	9	0			
25	BRCA1_185729_11	NT_010783_411956	4-F	BRCA1	10	2	31670	31995	326	1	1	31939C>CT_496		
26	BRCA1_185729_11	NT_010783_411956	4-F	BRCA1	10	2	31670	31995	326	1	1	31939C>CT_496		
27	BRCA1_264381_11	NT_010783_411956	4-F	BRCA1	10	2	31651	31995	345	2	1	31939C>CT_496		
28	BRCA1_264381_11	NT_010783_411956	4-F	BRCA1	10	2	31651	31995	345	2	1	31939C>CT_496		
29	BRCA1_185729_11	NT_010783_411956	4-R	BRCA1	10	2	31646	31983	338	16	1	31939C>CT_496		
30	BRCA1_185729_11	NT_010783_411956	4-R	BRCA1	10	2	31646	31983	338	16	1	31939C>CT_496		
31	BRCA1_264381_11	NT_010783_411956	4-R	BRCA1	10	2	31646	31974	329	15	1	31939C>CT_496		
32	BRCA1_264381_11	NT_010783_411956	4-R	BRCA1	10	2	31646	31974	329	15	1	31939C>CT_496		
33	BRCA1_185729_12	NT_010783_411956	5-F	BRCA1	10	2	31949	32130	182	0	0			
34	BRCA1_185729_12	NT_010783_411956	5-F	BRCA1	10	2	31949	32130	182	0	0			
35	BRCA1_264381_12	NT_010783_411956	5-F	BRCA1	10	2	31930	32130	201	0	0			
36	BRCA1_264381_12	NT_010783_411956	5-F	BRCA1	10	2	31930	32130	201	0	0			
37	BRCA1_185729_12	NT_010783_411956	5-R	BRCA1	10	2	31928	32126	199	28	0			
38	BRCA1_185729_12	NT_010783_411956	5-R	BRCA1	10	2	31928	32126	199	28	0			
39	BRCA1_264381_12	NT_010783_411956	5-R	BRCA1	10	2	31928	32118	191	29	0			
40	BRCA1_264381_12	NT_010783_411956	5-R	BRCA1	10	2	31928	32118	191	29	0			

- On the main menu, click File > Save Mutation Project.

The project is assigned a default name and has the type of Mutation Project Files (*.Sgp). You cannot change the file type, but you can change the project name and select any location in which to save the project.

Contig directory for a Mutation Surveyor project

Mutation Surveyor aligns the nucleotide sequences of the sample traces to the reference traces and groups the sample traces into contigs based on their degree of overlap. Because contigs are organized relative to the reference file, forward and reverse sequence pairs are grouped into two separate contig folders that are displayed in the Browser pane. The folders are labeled as ContignF and ContignR, where F stands for forward sequence, R stands for reverse sequence, and “n” stands for the contig file name, for example, Contig2F and Contig2R. The number in the parenthesis after the contig name indicates the number of

samples for the contig.




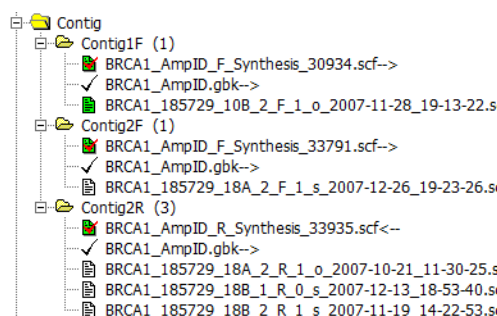
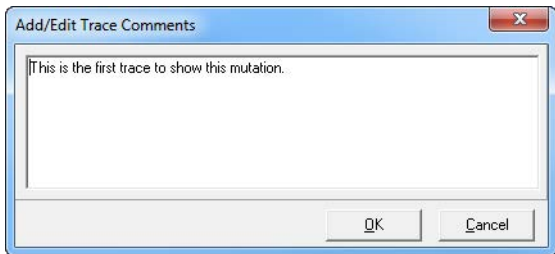
- A green File icon with a red checkmark  indicates the active reference file for the contig.
- If the reference file was derived from a GenBank file, a black checkmark  indicates this GenBank file.
- A green File icon  indicates the sample file that is currently selected. The corresponding data is displayed in the Viewing and Reporting pane of the GAD.

Figure 2-19: Contig directory in the Browser pane



You can right-click on the currently selected trace to open a context menu with the following options:

Option	Description
Copy Filename to Clipboard	Copies the complete file name for the selected trace to the clipboard. You can then paste the file name into a third-party application such as Word or Excel as needed.
Add/Edit Comment	<p>Opens the Add/Edit Comments dialog box in which you can add/edit the comments for the selected trace. The comments are displayed above the sample trace in the GAD.</p> <p><i>Figure 2-20: Add/Edit Trace Comments dialog box</i></p>  <p>Note: Trace comments can also be enabled in the Custom report. See “Display tab” on page 172.</p>
Delete Lane	<p>The Lane Quality for a deleted lane is set to “Low Quality” in the GAD. All the mutations for the deleted lane are removed from the Mutation table.</p> <p>Note: The Delete Lane option is available for a contig on both the Mutation table context menu and the Sample Trace name context menu. See “GAD analysis options” on page 91.</p>

Option	Description
Undo Lane Deletion	The Lane Quality is reset to its original value in the GAD. All the mutations for the deleted lane are again displayed in the Mutation table. Note: The Undo Lane Deletion option is available for a contig on both the Mutation table context menu and the Sample Trace name context menu. See “GAD analysis options” on page 91 .
Export Trace to Json	Saves the selected trace in JSON format.

Chapter 3

Graphical Analysis Display

The Graphical Analysis Display visually presents mutation project data. A variety of options are available for working with the display, including options for analyzing and manipulating mutation project data so that you identify the true mutations in a mutation project.

This chapter covers the following topics:

- [“Graphical Analysis Display Layout” on page 77.](#)
- [“Graphical Analysis Display Navigation and Analysis Options” on page 90.](#)
- [“Exporting the GAD” on page 94.](#)

Chapter 3


Graphical Analysis Display

Graphical Analysis Display Layout

The Graphical Analysis Display (GAD) visually presents mutation project data in the Viewing and Reporting pane. A variety of options are available for working with the display, including options for analyzing and manipulating project data to identify the true mutations in a project. To open the display for a mutation project, do one of the following:

- Double-click a sample file in the Mutation report.
- Double-click a sample file in the Browser pane.
- Double-click a specific mutation in the Mutation report.

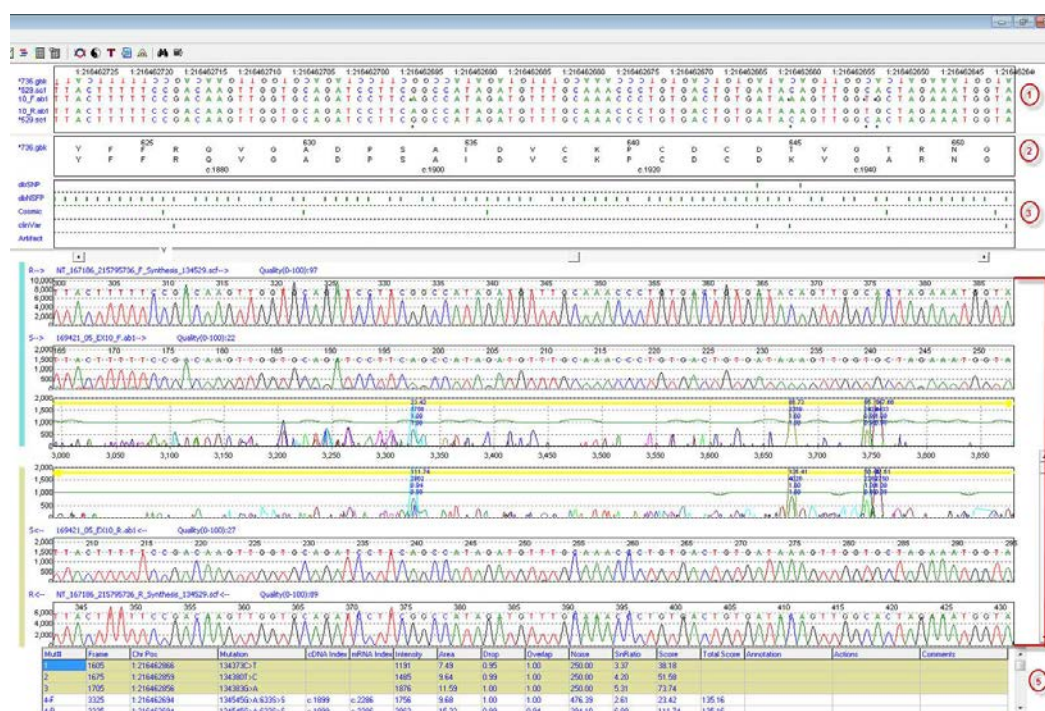Figure 3-1: Graphical Analysis Display in the Viewing and Reporting pane



- 1 - Nucleotide Sequence frame
- 2 - Amino Acid frame
- 3 - Tracks frame
- 4 - Sample Traces frame
- 5 - Mutation table

The Graphical Analysis Display for any mutation project has up to five major components. Four of the five components—the Nucleotide Sequence frame, the Amino Acid frame, the Sample Traces frame, and the Mutation table—are always displayed for a Mutation Surveyor project. The fifth component, the Tracks frame, is displayed only if information has been imported from an annotation database for the project, and/or false positives and/or artifacts have been manually added to the project. From top to bottom, these components are:

- The Nucleotide Sequence frame. See “[Nucleotide Sequence frame](#)” below.
- The Amino Acid frame. See “[Amino Acid frame](#)” on page 79.
- The Tracks frame. See “[Tracks frame](#)” on page 79.
- The Sample Traces frame. See “[Sample Traces frame](#)” on page 79.
- The Mutation table. See “[Mutation table](#)” on page 83.

The settings that you specify on the output/display tabs of the Mutation Project Settings dialog box determine some of the information that is displayed in the GAD. If you modify a setting on an output/display tab and then click OK, the GAD is dynamically updated to reflect this change. (See “[Modifying the Default Mutation Surveyor Project Settings](#)” on page 52.) In addition, the Mutation Surveyor toolbar is dynamically updated with buttons that are specific for navigating and working with the GAD. (See “[GAD Toolbar](#)” on page 84.)



The following sections describe the GAD components for a 2D mutation project in which a GenBank file was used to create a synthetic reference file. There are slight differences for 1D projects, or for projects that did not use a GenBank file.

Nucleotide Sequence frame

The Nucleotide Sequence frame displays the GenBank file, the reference traces, and the sample traces as called bases.

Figure 3-2: Nucleotide sequence frame



From top to bottom in the frame, the rows display the following information:

- Position Information—The option that is selected on the Display tab of the Mutation Project Settings dialog box determines the information that is displayed in the first row. The positions within the GenBank file for the queried contig are displayed if Contig is selected. Chromosome number:chromosome positions are displayed if Chromosome is selected. The location of the queried gene (plus (+) or minus strand (-)) and the option that is selected for Automatically Adjust GenBank Direction on the Others tab determines the order (ascending or descending) in which this position information is displayed.
- The nucleotide sequence for the GenBank file. Within a CDS, the GenBank nucleotides are displayed in uppercase. Outside the coding region, the GenBank nucleotides are displayed in lowercase. If the gene is on the minus (-) strand and Orient GenBank Files in Gene Direction is selected for Automatically Adjust GenBank Direction, then the nucleotide sequence is also displayed upside down.
- The nucleotide sequence for the forward reference trace.

- The nucleotide sequence for the forward sample trace.
- The nucleotide sequence for the reverse sample trace.
- The nucleotide sequence for the reverse reference trace.

An asterisk indicates inconsistencies in base calls between traces. A black asterisk (*) indicates an inconsistency in sequence between a reference trace and a sample trace. A red asterisk (*) indicates an inconsistency in sequence between a reference trace and the GenBank file.

Amino Acid frame

The Amino Acid frame displays the codon number and the IUPAC abbreviation (unless 3 letter abbreviation setting is selected) for the corresponding amino acid sequence for the called bases that fall within a coding region in the GenBank sequence (the first line in the frame) and in the sample file (the second line in the frame). Any differences between the amino acid translations as the result of mutations in the sample file are therefore visible in this frame. The frame also shows the coding DNA reference index. Just like the Nucleotide Sequence frame, the location of the queried gene (plus (+) or minus strand (-)) and the option that is selected for Automatically Adjust GenBank Direction on the Others tab determines the order (ascending or descending) in which the codon numbers and c.dna numbers are displayed.


Figure 3-3: Amino Acid frame



Tracks frame

The Tracks frame is displayed only if information has been imported from an annotation database for the project, and/or false positives and/or artifacts have been manually added to the project. A reported single base variant at a specific chromosome position is indicated by a green tick mark. A reported variant that spans multiple bases at a specific chromosome position is indicated by a green horizontal line. You can place your cursor over a green tick mark or horizontal line to open a popup that displays detailed information for the reported variant.

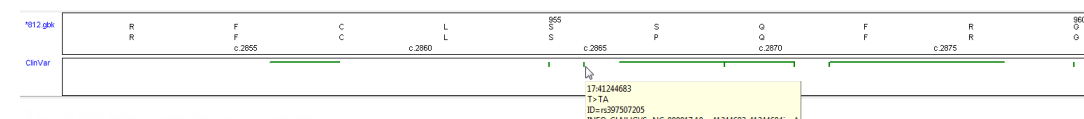Figure 3-4: Tracks frame with mutation information displayed for a reported variant



Sample Traces frame


The Sample Traces frame displays the traces for a selected sample grouped by direction.

From top to bottom, the forward reference trace, the forward sample trace, and the forward mutation trace are grouped together, and from bottom to top, the reverse reference trace, the reverse sample trace, and the reverse mutation trace are grouped together. The forward direction group is marked by a light blue vertical bar to the left of the group and the reverse direction group is marked by a light tan vertical bar to the left of the group. The trace name, the sequence direction, and the lane quality are displayed above each trace. If any comments have been added for a trace, they are also displayed above the trace. The location of the queried gene (plus (+) or minus strand (-)) and the option that is selected for Automatically Adjust GenBank Direction on the Others tab determines if the Reference trace is reverse complemented or not. The X-axis indicates the frame number and the Y-axis indicates peak intensity in Relative Fluorescence Units (RFU). The nucleotide numbering is from the 5' end to the 3' end on the forward trace and from the 3' end to the 5' end on the reverse trace. The following section identifies the different parts of the traces. In addition, the different types of peaks in the traces are identified with unique peak information. (See “[Mutation trace legend](#)” below and “[Mutation trace peak identification](#)” on page 82.)

Figure 3-5: Reference traces, sample traces and mutation traces



Mutation trace legend

A peak in the mutation trace is called as a *mutation peak* when its height (in RFU) is greater than or equal to the value specified for the Mutation Height in the Mutation Project settings. (See “[Mutation tab](#)” on page 59.) A mutation peak appears in frame with the reference trace. The peak color indicates the called base for the mutation. (To see the color that is assigned to a specific base or group of bases in the mutation trace, click the Show Lines button  on the GAD toolbar.)

Horizontal *location bars* run across the top of the mutation trace. The bar color indicates the location of the sequence that is being displayed in the trace.

- Yellow—Exon coding region
- Pink—Region of Interest (ROI)
- Light tan—Gene region
- Light blue—Primer region


- Green—Somatic (Mosaic) mutation



*Somatic mutations are marked in the mutation trace only if you have selected the **Check 2D Small Peaks (Mosaic)** option for the mutation project settings. See [“Display tab” on page 64](#).*

- Red—Homozygous deletion or duplication. See [“Mutation trace peak identification”](#) below.
- Red with green arrows—Homozygous insertion. See [“Mutation trace peak identification”](#) below.
- Brown—Heterozygous insertion or deletion. See [“Mutation trace peak identification”](#) below.

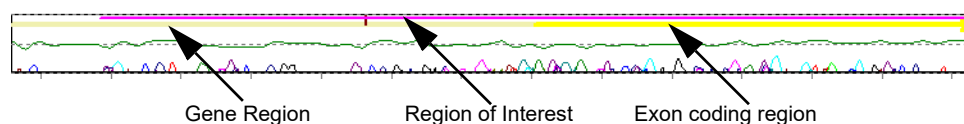Figure 3-6: Example of different colored Location bars in a mutation trace




Brown tick marks indicate reported variations in the GenBank file while tick marks of other colors indicate reported SNPs. Thick blue vertical bars define a *comparison region*, which is a region where the sample traces align to the reference traces, and all the mutation project *analysis* settings have been met, and therefore, mutations, with the exception of heterozygous indels, can be called in this region. (See [“Mutation trace peak identification”](#) below.)



*If **Display Negative SNPs** is selected for the mutation project, you can zoom in on the tick marks in the mutation trace to see the involved bases. See [“Display tab” on page 64](#).*

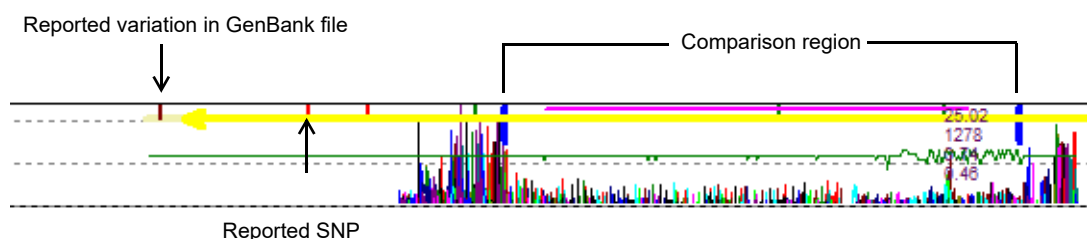
Figure 3-7: Comparison region in mutation trace



Mutation trace peak identification

If all the criteria are met for a substitution mutation event, the values for mutation score, peak height, overlapping factor, and dropping factor are displayed on the peak.

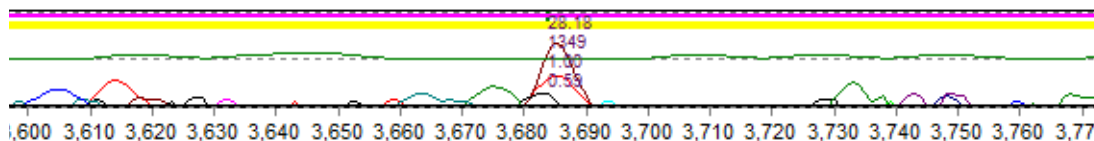Figure 3-8: Substitution mutation event



The *mobility line* is the horizontal line that runs through the center of the mutation trace. This line indicates the mobility of the sample trace relative to the reference trace—that is, by how much was the sample trace shifted so that the center of the peaks in the sample trace align to the center of the peaks in the reference trace. Typically, this line is green; however, the line changes color if an indel event is detected in the sample trace. If a *homozygous* deletion, duplication, or insertion event is detected in the sample trace, then the line is red. A positive peak in the mobility line spans a homozygous deletion event. A negative peak in the mobility line spans a homozygous duplication or insertion event.

Figure 3-9: Homozygous deletion event

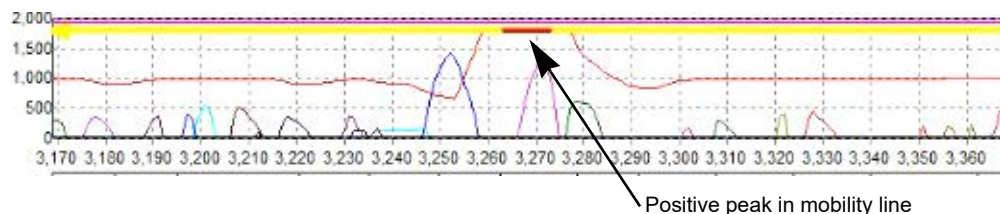


Figure 3-10: Homozygous duplication event



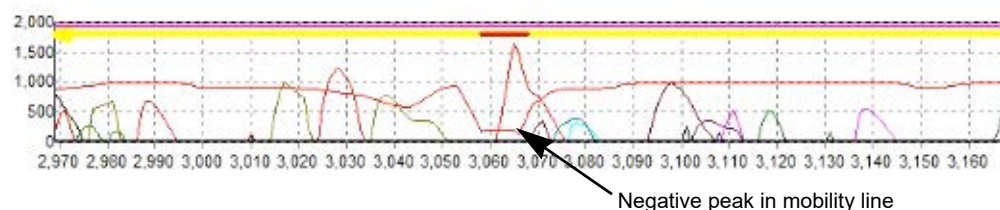


Figure 3-11: Homozygous insertion event

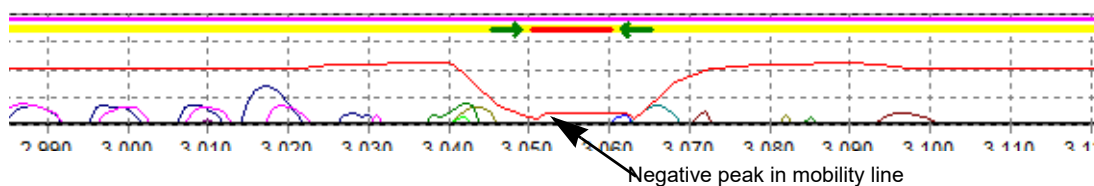

If a *heterozygous* insertion or deletion event is detected in the sample trace, then the line is brown and a blue vertical line indicates that automatic mutation detection is not possible past the point of the event.

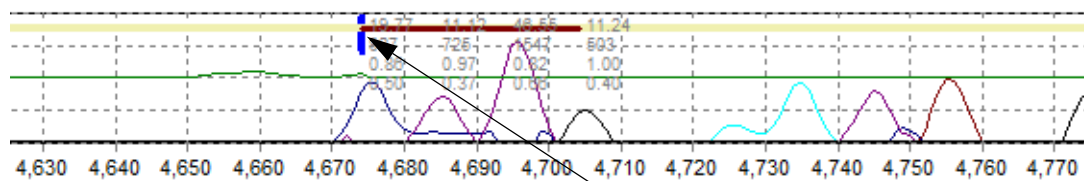Figure 3-12: Heterozygous mutation event



Mutation table

The Mutation table is the last component in the GAD. The Mutation table displays all the mutation data for the currently opened trace. See [“Modifying the Default Mutation Surveyor Project Settings”](#) on page 52.



For a detailed explanation about the code in the Mutation field, see [“Mutation Report”](#) on page 127.

Figure 3-13: Mutation table

Mut#	Frame	Chr Pos	Mutation	cDNA Index	mRNA Index	Intensity	Area	Drop	Overlap	Noise	SnRatio	Score
1	1605	1:216462866	134373C>T			1191	7.49	0.95	1.00	250.00	3.37	38.18
2	1675	1:216462859	134380T>C			1485	9.64	0.99	1.00	250.00	4.20	51.58
3	1705	1:216462856	134383G>A			1876	11.59	1.00	1.00	250.00	5.31	73.74
4-F	3325	1:216462694	134545G>A:6335>S	c.1899	c.2286	1756	9.68	1.00	1.00	476.39	2.61	23.42
4-R	3325	1:216462694	134545G>A:6335>S	c.1899	c.2286	2862	15.32	0.99	0.94	294.10	6.88	111.74











The color-coding for a substitution mutation event indicates a variety of information about the mutation.













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

- Text color—Blue text indicates that the mutation is called with high confidence. Red text indicates that the mutation is called with low confidence. (See [“Mutation Threshold”](#) on page 64.) Black text indicates that the mutation has been manually confirmed. (See [“Confirm”](#) on page 133.)
- Purple background with green text—A reported variation that was not found in the mutation project (Negative SNP).
- Pink background—A novel variant that results in a change to the amino acid sequence.
- Purple background—A reported variant that might or might not result in a change to the amino acid sequence.
- Blue background—A novel variant on the forward strand that does not result in a change to the amino acid sequence.
- Tan background—A novel variant on the reverse strand that does not result in a change to the amino acid sequence.
- No background color—A novel variant that is found in both directions that does not result




in a change to the amino acid sequence.

To view a mutation event in the mutation trace in the GAD, double-click the mutation in the table. If the mutation is a heterozygous indel, the Het-Indel tool also opens. See [“To add or edit a heterozygous indel \(The Het-Indel tool\)” on page 331.](#)

GAD Toolbar

Button	Description
	Open Files—Opens the Open Files dialog box which you use to load GenBank files, references files and sample files for a project.
	Save Project—Saves all the project files to a single file mutation project file with an .sgp extension. You can open an entire project in a single step by opening this SGP file.
	Raw Data to Processed Data—Enabled when AB1/ABI files without base call information are uploaded and Load Raw Data is selected for the Mutation Project settings. Converts the raw files so that base call information is displayed.
	Run—Runs the currently loaded project in Mutation Surveyor.
	Restart—Opens the Options dialog box which you use to modify the current project settings. You can then run the project again using these new settings.
	Show/Hide Browser—Shows/hides the Browser pane in the Mutation Surveyor main window. Also available on the Display menu (Display > Browser).
	Show/Hide Nucleotides—Shows/hides the Nucleotide Sequence frame in the GAD. Also available on the Display menu (Display > Nucleotides).
	<ul style="list-style-type: none"> Show/Hide Amino Acids—Shows/hides the Amino Acid frame in the GAD. Also available on the Display menu (Display > Amino Acids). Show/Hide Tracks—Shows/hides the Tracks frame in the GAD. Also available on the Display menu (Display > Tracks).
	Show/Hide Mutation table—Shows/hides the Mutation table in the GAD. Also available on the Display menu (Display > Peak table).
	<p>Show Lines—Provides options for:</p> <ul style="list-style-type: none"> Show All—Shows all nucleotides in the sample traces and mutation traces. Hide All—Hides all nucleotides in the sample traces and mutation traces. Showing/hiding selected nucleotides in the sample traces and mutation traces. Show/Hide Grid—Hides the background grid in all sample traces and mutation traces. Line Weight—Adjusts the weight of traces from 1 to 4 times. <p>Note: If you select Hide Grid, you can still select Show All.</p>

Button	Description
	Zoom In—Zooms in on the center of the GAD.
	Zoom Out—Zooms out from the center of the GAD.
Note: The GAD “remembers” <i>all</i> the Show/Hide settings that you last specified. Even if you close, and then reopen Mutation Surveyor, the Show/Hide settings remain the same for any project that you open.	
	Select Reference File—Opens the Select Reference file dialog box in which you can select a different reference file to which to align samples traces in a contig.
	Custom Report—Opens the Custom Report Settings dialog box which you use to specify the settings for a custom report. See “Custom Report” on page 164 .
	Graphic Display—Opens the Graphic Display Report settings dialog box in which you specify the settings for generating a Graphic Display report. See “Graphic Display Report” on page 108 .
	GAD report—Prints the currently opened GAD in a report format. See “GAD Report” on page 114 .
	Clinical Report—Generates a Clinical report. See “Clinical Report” on page 116 .
 	Display Mutations with 1D/2D Settings—Toggles the GAD between the 1D and 2D views. Note: With 1D setting selected, more low confidence mutation calls might be included in reports.
	Detect Heterozygous Indels—Opens the Heterozygous Indel Detection window. See “To add or edit a heterozygous indel (The Het-Indel tool)” on page 331 .
	Nucleotide Text Display Report—Generates a Nucleotide Text Display report. See “Nucleotide Text Display Report” on page 120 .
	Save Base-Modified Samples—Opens the Browser for Folder dialog box in which you can specify the directory in which to save modified traces. This button appears only after you edit a base call in a sample file or reference file. See “Edit Base” on page 93 .
	Re-process Modified Files—Displayed only after you edit a base call in a sample file or reference file. Reprocesses the mutation project based on the modified files. See “Edit Base” on page 93 .
	Export GAD—Exports selected GAD regions to a Word document, to JPG files, or directly to a printer. See “Exporting the GAD” on page 94 .

Button	Description
	Quantify—Opens the Mutation Quantification function. See “Mutation Quantifier Function” on page 303 .
	Search the GenBank Sequence—Opens the Search dialog box. You use the options on this dialog box to search for a specific sequence in the loaded GenBank file for a project. See “Search GenBank sequence” below.
	Variation Tracks Settings icon - Opens the Variation Tracks Settings dialog box. The Variation Tracks Settings dialog box displays the available tracks settings for the Mutation table based on the variation databases that were imported for the project. See “Variation Tracks Settings dialog box” on page 87 . Note: After being imported into Mutation Surveyor, a variation database is referred to as a <i>track</i> .


Search GenBank sequence

You use the Search function to search for a specific sequence or position in the GenBank file for the displayed project.

1. On the GAD toolbar, click the Search the GenBank Sequence icon .

The Search dialog box opens.

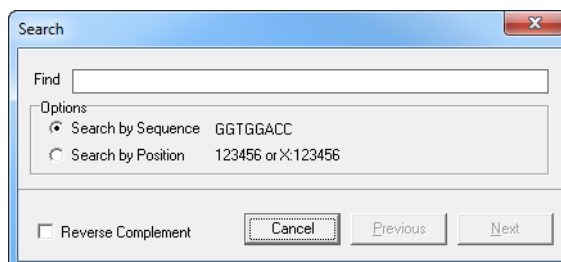Figure 3-14: Search dialog box



2. In the Find field, enter the appropriate search string.
3. Under Options, indicate whether to search by sequence (the default value), or by position.
4. Optionally, if you select Search by Sequence, then to search by a reverse complement sequence, click Reverse Complement.
5. Click Previous or Next to search through the GenBank sequence in the Nucleotide Sequence frame.

Matching sequences or positions are highlighted in yellow.


Variation Tracks Settings dialog box

The Variation Tracks Settings dialog box contains the tracks settings for the Mutation table based on the variation databases that were imported for the project. (After being imported into Mutation Surveyor, a variation database is referred to as a *track*. See the [Reference and Track Manager Tool User's Manual](#). You can select what information to display for the tracks in the Annotation field, and you can filter the data that is displayed in the Mutation table based on the tracks, or you can choose not to filter the data based on any of the tracks.

1. On the GAD toolbar, click the Variation Tracks Settings icon .

The Variation Tracks Settings dialog box opens. The Tracks pane is the left pane of the dialog box. The pane displays all the variation databases, or *tracks*, that were imported for the selected project.

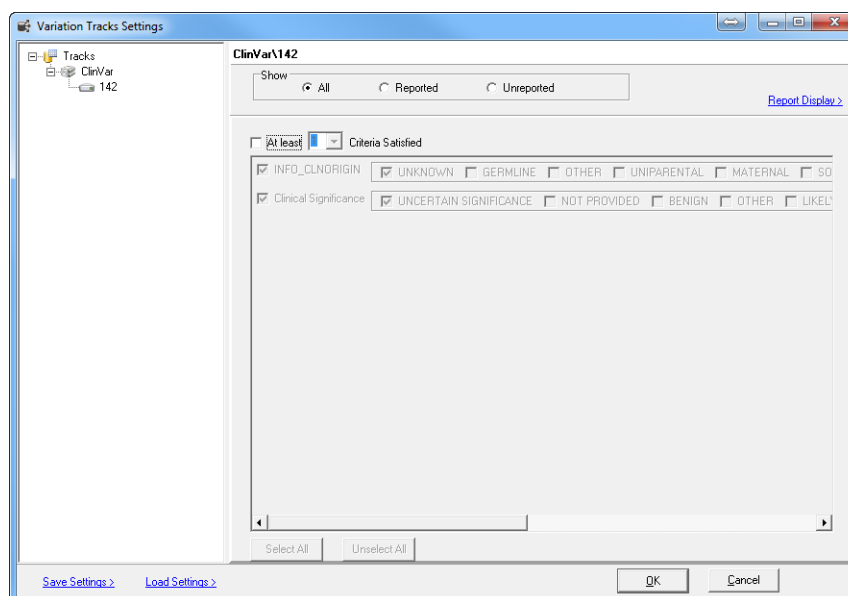Figure 3-15: Variation Tracks Settings dialog box, Filter Settings pane



2. If you do not want to filter the data for the project based on any of the tracks, then click Load Settings > Clear all track filters, and then click OK; otherwise, go to [Step 3](#).
3. In the Tracks pane, select a track, and then do the following:
 - a. Indicate the types of variants that are to be included in the Mutation report.
 - All - By default, all variants that meet all the filtering criteria are displayed in the Mutation report, whether they are included in the selected track.
 - Reported - Select Reported to display only those variants that meet all the filtering criteria and that are included in the selected track.
 - Unreported - Select Unreported to display only those variants that are not included in the selected track.

- b. Specify the filter settings for the track.




The available settings vary for each track.

- Click Report Display to open the Report Display Settings dialog box.

The Report Display Settings pane displays all the columns that can be included in the Annotation field of the Mutation table. By default, no columns are selected.

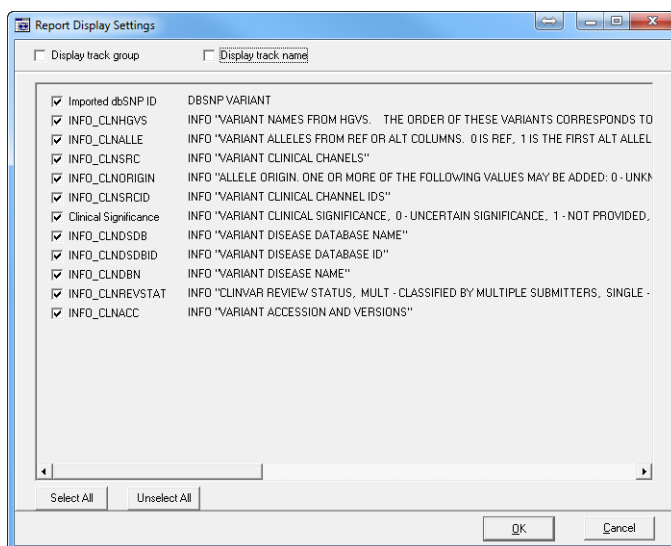Figure 3-16: Mutation Tracks Settings dialog box, Report Display Settings pane



4. Select the columns that are to be included in the Annotation field of Mutation table, and then optionally, select one or both of the following:
 - Display track group—Displays the track group in the Annotation field, for example, ClinVar for the ClinVar database.
 - Display track name—Displays the track name in the Annotation field. (This is the name that you assigned to the track during its import with the Track Manager tool. See the [Reference and Track Manager Tool User's Manual](#).
5. Click OK to close the Report Display Settings dialog box.
6. Repeat [Step 3](#) through [Step 5](#) for each track that is to be used to filter the Mutation table.
7. Do one of the following to save your settings and close the Variation Tracks Settings dialog box:
 - Click OK.

Going forward, the Mutation table is generated according to these saved settings until you change them.

- Save Settings > Save User Defaults, and then click OK.

The settings that you have specified for all the tracks are saved as your (the logged in user's) default settings. Going forward, any new project that *you* run in Mutation Surveyor uses these settings by default. If you change the settings for a project, and want to generate the Mutation table display based on your default settings, then you can click Load Settings > Load User Defaults to restore your default settings.

- Save Settings > Save To File, and then click OK.

The settings that you have specified for all the tracks are saved to a Settings (.ini) file. Going forward, you can click Load Settings > Load From File to load this saved Settings file and generate the Mutation table display according to the settings in the file.

Graphical Analysis Display Navigation and Analysis Options

You can easily change the [view](#) in the GAD or [navigate](#) the GAD viewer using some of the toolbar icons (see “[GAD Toolbar](#)” on page 84) or your mouse and some keyboard hotkeys. Options are also available for [analyzing](#) mutation project data so that you identify the true mutations in a mutation project. Miscellaneous options are also available for [manipulating](#) the mutation project data so that you can work with the data in outside programs if needed.

GAD viewing and navigation options


Action	Description
Zoom in	Click and hold the left mouse button and draw a box from the upper left hand corner of the pane towards the lower right hand corner. A box is formed around the area that being reduced for viewing.
Zoom out	Click and hold the left mouse button and draw a box from the lower right hand corner of the pane towards the upper left hand corner. Note: The magnification for zooming out is always 100%.
Moving an image in the GAD	To move any image (frame or trace) in the GAD, right-click and hold the mouse button on the image, and then move the mouse.
Note: The frames and traces are linked in the GAD. If you zoom in on a region in one of the frames or traces or zoom out from a region in one of the frames or traces, then this action is automatically duplicated in the same region in all the other frames and traces. If you move a frame or a trace, then all other frames and traces are moved to the same position as well.	
Moving to a mutation peak	Press and hold the [Ctrl] key and then press: <ul style="list-style-type: none">• F to move forward to the next mutation peak.• B to move back to the previous mutation peak. Note: You move to each mutation peak on a trace by trace basis. For example, if Sample Trace #1 has five mutation peaks, Sample Trace #2 has no mutation peaks, and Sample Trace #3 has one mutation peak, you must press CTRL+F five times to move to each mutation peak in the Sample Trace #1. When you press CTRL+F the 6th time, you move to the mutation peak in Sample Trace #3.
Moving to a trace	Use the Browse function on the right side of the Traces component in the GAD to move to the next or previous trace.

GAD analysis options

Analysis options are available from the context menu for the Mutation table and from various context menus in the GAD. Any edit action (addition, substitution, or deletion) that you carry out for a mutation is displayed in the Actions column for the mutation in the Mutation table.

Action	Description
Mutation table context menu analysis options	
Edit Comments	<p>You can manually add or edit comments for a mutation call.</p> <ul style="list-style-type: none"> Right-click the entry for the mutation call on the Mutation table, and on the context menu that opens, click Edit Comments. You can add original comments in the Comments field, or you can select from a list of up to 200 customizable Comment templates. Right-click the entry for the mutation call on the Mutation table, and on the context menu that opens, click Edit. The Edit Substitution dialog box or the Edit Indel dialog box opens. Add comments in the Comments field in this dialog box. <p>Note: For detailed information about editing a mutation call, see Chapter 12, "Reviewing and Editing Mutation Projects," on page 327.</p>
Edit	<p>Right-click the entry for the mutation in the Mutation table, and on the context menu that opens, click Edit. The Edit Substitution dialog box or the Edit Indel dialog box opens. Edit the mutation call as appropriate.</p> <p>Note: For detailed information about editing a mutation call, see Chapter 12, "Reviewing and Editing Mutation Projects," on page 327.</p>
Delete Mutation	<ul style="list-style-type: none"> Right-click the entry for the mutation in the Mutation table, and on the context menu that opens, click Delete Mutation. Click the entry for the mutation in the Mutation table, and then click [Ctrl]-D. Press and hold the [Ctrl] key, and then right-click the mutation peak in the trace to open a Warning message about deleting the mutation call. Click OK in the message to delete the mutation call. <p>The text for the mutation code turns gray in the Mutation table. The comment <<Deleted>> is displayed in the Actions field in the Mutation table.</p>
Delete Mutation and Add to Database	<p>This action is available only when project meets all requirements for UKDB. Either click OK to delete the mutation without adding it to the User Knowledge database, or if the GenBank file that was used for the project contains chromosome information, then you can add the mutation to the User Knowledge database. (See "Creating a User Knowledge Database for Known Deletions" on page 342.)</p>
Undo Mutation Deletion	<ul style="list-style-type: none"> Right-click the entry for the deleted mutation in the Mutation table, and on the context menu that opens, click Undo Mutation Deletion. Press and hold the [Ctrl] key, and then right-click the deleted mutation in the trace to open a Warning message about undoing the deletion of the mutation call. Click OK in the message to undo the deletion of the mutation call.

Action	Description
Delete Lane	<p>Make sure that the appropriate trace is open in the GAD, and then right-click any entry in the Mutation table and on the context menu that opens, click Delete Lane.</p> <p>Note: The Lane Quality for a deleted lane is set to “Low Quality” in the GAD. All the mutations for the deleted lane are removed from the Mutation table.</p> <p>Note: The context menu for a contig in the Browser pane also has a Delete Lane option. See “Contig directory for a Mutation Surveyor project” on page 72.</p>
Undo Lane Deletion	<p>Make sure that the appropriate trace is open in the GAD, and then right-click on the trace's title and on the context menu that opens, and click Undo Lane Deletion.</p> <p>Note: The Lane Quality for the lane is reset to its original value. All the mutations that were deleted for the lane are again displayed in the Mutation table.</p> <p>Note: The context menu for a contig in the Browser pane also has a UnDo Lane Deletion option. See “Contig directory for a Mutation Surveyor project” on page 72.</p>
Confirm/Undo Confirmation	Right-click the entry for the mutation in the in the Mutation table, and on the context menu that opens, click the appropriate option.
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See “Viewing the Edit History (Audit Trail) for a Mutation” on page 344.
Hyperlink to NCBI's dbSNP	Opens the NCBI's database entry (via the web) for the variant.
GAD context menu analysis options	
Delete Lane	<p>Make sure that the appropriate trace is open in the GAD, and then right-click the Sample Trace name, and on the context menu that opens, click Delete Lane.</p> <p>Note: The Lane Quality for a deleted lane is set to “Low Quality” in the GAD. All the mutations for the deleted lane are removed from the Mutation table.</p> <p>Note: The context menu for a contig in the Browser pane also has a Delete Lane option. See “Contig directory for a Mutation Surveyor project” on page 72.</p>
Undo Lane Deletion	<p>Make sure that the appropriate trace is open in the GAD, and then right-click the Sample Trace name, and on the context menu that opens, click Undo Lane Deletion.</p> <p>Note: The Lane Quality for the lane is reset to its original value. All the mutations that were deleted for the lane are again displayed in the Mutation table.</p> <p>Note: The context menu for a contig in the Browser pane also has a Undo Lane Deletion option. See “Contig directory for a Mutation Surveyor project” on page 72.</p>

Action	Description
Add Substitution Call	<p>Right-click in the mutation trace at the position at which you want to add the substitution, and on the context menu that opens, click Add Substitution. The Add Substitution dialog box opens. Add the substitution as appropriate.</p> <p>Note: For detailed information about modifying a mutation call, see Chapter 12, “Reviewing and Editing Mutation Projects,” on page 327.</p> <p>Note: If you click a peak for a substitution mutation event, a message opens indicating that substitution already exists at the selected position and asking you if you want to replace it. Click OK to continue with the substitution.</p>
Add Insertion/ Deletion Call	<p>Press and hold the [Ctrl] key, and then click the mobility line at the position of the indel. The Add Indel dialog box opens. Add the Indel call as appropriate.</p> <p>Note: For detailed information about modifying a mutation call, see Chapter 12, “Reviewing and Editing Mutation Projects,” on page 327.</p>
Edit Base	<ul style="list-style-type: none"> Right-click a reference or sample trace and click Edit Base. A dialog box opens that displays the complete base sequence for the trace. Click the base that you are editing and enter the text for the new base (A,C, G, or T.) Click the Close button (“X”) to close the dialog box, and then click “Yes” at the prompt to save the base call changes. <p>Note: If you do not want to override the original traces, you can save the modified traces to a different location. See Save Base-Modified Samples on page 85.</p> <ul style="list-style-type: none"> Click the Re-process Modified Files button  to run the project with the modified sample and/or reference traces.
Copy Sequence	<p>Right-click a reference or sample trace and click Copy Sequence. The base calls for the portion of the trace that is in view will be copied to the clipboard.</p>

Miscellaneous GAD options

Action	Description
Copy Image	<p>Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Copy Text	<p>Right-click a table cell in the Mutation table and click Copy to copy the text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p>

Exporting the GAD

You use the Export GAD function to export mutation traces and other information for selected mutations to a Word document, to separate JPG files, or directly to a printer.


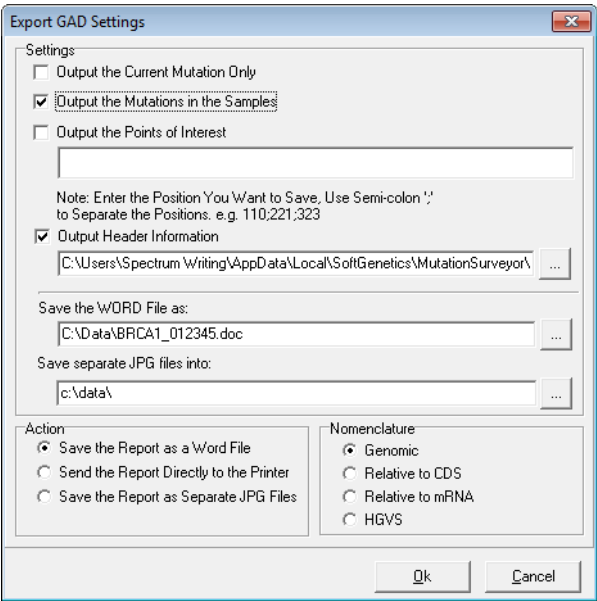



 A variety of graphic-based reports that present part or all the GAD in a report format are also available in Mutation Surveyor. See [Chapter 4, “Mutation Surveyor Graphical Reports,”](#) on page 97.

Figure 3-17: Export GAD dialog box



Option	Description
Settings	
Output the Current Mutation Only	Exports the GAD for only those mutations (+-16 bp) that are currently displayed in the GAD for the selected sample file.
Output the Mutations in the Samples	Exports the GADs for all the mutations in all the sample files for every contig.
Output Points of Interest	<p>Exports the GAD only for the specified base positions. (The Base position must be the contig position where at least one sample contains a mutation call at the point of interest.)</p> <ul style="list-style-type: none">• If Output the Mutations in the Samples is also selected, then images are captured for not only all mutation calls, but also for positive and negative mutations at all Points of Interest.• If Output the Mutations in the Samples is not selected, then images are captured for only positive and negative mutations at all Points of Interest.

Option	Description
Output Header Information	<p>Automatically populated with the default Mutation Surveyor header information file. You can click the Browse button  to open the Open File dialog box, and browse to and select a custom header to add to the report.</p> <p>Note: A header is a .inf file. For a discussion of Custom headers and how to create one, see “Print Header Editor Tool” on page 267.</p>
Save the Word File As	<p>Available only if Save the Report as a Word file is the selected action. Automatically populated with the default location (c:\data) and filename for the Word file to which the mutations traces for the selected mutations are being exported. You can click the Browse button  to open the Save As dialog box, and specify a different location, name, or both for the Word files.</p> <p>Note: This function works only when Microsoft Word is installed on the client computer. If the client's copy of Word has additional add-ons, such as End Note, installed, this function might not work properly.</p>
Save separate JPG files into	<p>Available only if Saving Report as Separate JPG Files is the selected action. Automatically populated with the default location (c:\data) to which the mutation traces for the selected mutations are being exported.</p> <p>You can click the Browse button  to open the Browse for Folder dialog box, and specify a different location for the JPG files.</p>
Action	
<ul style="list-style-type: none"> Save the Report as a Word File Send the Report Directly to the Printer Save the Report as Separate JPG Files 	<ul style="list-style-type: none"> Exports all the mutation traces to a single Word file. Exports the files to directly to the system's default printer Export each mutation trace as a separate JPG file
Nomenclature	
<ul style="list-style-type: none"> Genomic Relative to CDS Relative to mRNA 	<ul style="list-style-type: none"> Displays the mutation call positions based on the genomic position referenced in the GenBank file. Displays the mutation call positions relative to the CDS numbering in the GenBank reference file. A mutation that is called in the CDS is prefixed with a "c." A mutation that is called in a non-coding region is prefixed with "IVS." Its position is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide Displays the mutation call positions relative to the mRNA numbering in the GenBank reference file. A mutation that is called in the mRNA region is prefixed with a "c." A mutation that is called outside of an mRNA region is identified by the closest nucleotide in the mRNA region plus or minus the number of bps by which it resides away from the closest nucleotide.

Chapter 3
Graphical Analysis Display

Option	Description
Nomenclature (cont'd)	
<ul style="list-style-type: none"> • HGVS 	<ul style="list-style-type: none"> • Displays the mutation call positions relative to its position in the coding sequence. A mutation that is called in a non-coding region is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide. Both alleles are shown in separate bracketed statements
<p>Note: For any of the four standard Nomenclature options, if a GenBank file was not loaded as the reference file, then the Base Number is enclosed in parenthesis and it indicates the position of the base in the reference file. If both a GenBank file and a reference file were loaded for the mutation project and the mutation was called at a position that does not align to the GenBank file, then the Base Number is enclosed in parenthesis and is displayed as +/- the number of base pairs by which the mutation was shifted in comparison to the GenBank file.</p>	

Chapter 4

Mutation Surveyor Graphical Reports

A variety of graphic-based reports that provide detailed information about a mutation project are available in Mutation Surveyor. All these reports, which are available from the Display option on the Mutation Surveyor main menu, present part or all the GAD in a report format. All these reports rely heavily on the use of colored legends to indicate critical information about the mutation project data. You can print the information in many of these reports and you can save and store the report information for use at a later date.

This chapter covers the following topics:

- [“Project Reviewer Report” on page 99.](#)
- [“Graphic Display Report” on page 108.](#)
- [“GAD Report” on page 114.](#)
- [“Clinical Report” on page 116.](#)
- [“Nucleotide Text Display Report” on page 120.](#)


Chapter 4

Mutation Surveyor Graphical Reports

Project Reviewer Report

You use the Project Reviewer report to view mutation project data for one contig at a time, or for all contigs at the same time. The data that you can view includes the reference nucleotide sequence, the consensus sequence for all the sample traces, the sequence for each sample trace, and all the mutation calls for all the samples. You can use the Project Reviewer functions to customize the viewing of this information for your sample files. For example, you can use the Sample Grouping function to group all your sample files by an identifier and view any overlapping contigs that result. You can also use the Project Reviewer to edit mutation calls.

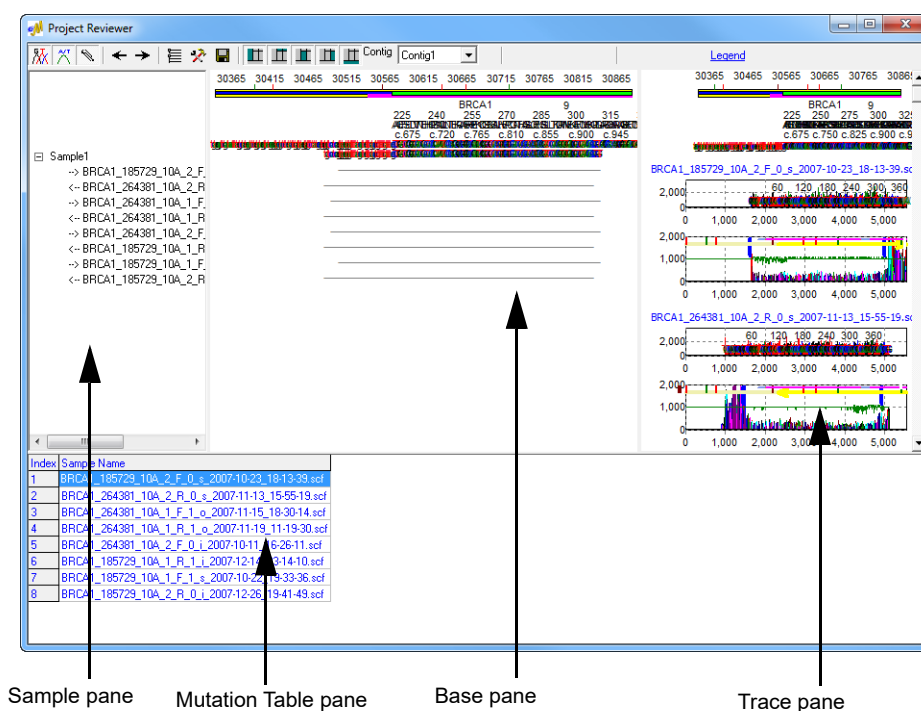Figure 4-1: Project Reviewer report



Project Reviewer report layout

To view the color coding legends that are used in the report, click Legend at the top of the report. The report displays mutation project data according to these legends in four panes.

- **Sample pane**—The Sample pane is the top left pane. This pane lists all the sample files that are contained in the currently selected contig.
- **Base pane**—The Base pane is the top center pane. The top row of numbers indicates the genomic base position in the GenBank file. The color of the horizontal *location bars* that run below the genomic position numbers indicates the location of the sequence that is being displayed in the trace. Colored tick marks indicate reported variations. The first row of nucleotides in the pane is the GenBank nucleotide sequence. The second row of

nucleotides in the pane is the consensus sequence for the sample traces. The remaining rows show the sequence for each sample trace. A black dash (-) indicates a match between the reference file and the sample file. If a different base is at the same position in the sample file and reference file, then the standard IUB/IUPAC code for the base is shown in the sample file at the position. Color-coding is used to indicate the mutation type. No information is displayed for areas that fall outside the comparison region. A gray square indicates a mutation that has been deleted either by Mutation Surveyor or by a user.

- Trace pane—The Trace pane is the top right pane. This pane displays the sample traces, the mutation traces, or both for all the selected samples in the sample and base panes. Blue highlighted bases indicate positions of called mutations.
- Mutation Table pane—The Mutation Table pane is the lower pane. This pane displays all the mutation calls for all the samples in the selected contig. The mutation call codes and color-coding that are used in this table are identical to the mutation call codes and color-coding that are used in the Mutation standard report. See [“Mutation Report” on page 127](#).

The report is interactive:

- To zoom in on a selected region for the sample files, in the Base pane, click and hold the left mouse button and draw a box from the upper left hand corner of the region towards the lower right hand corner of the region. A box is formed around the area that is being reduced for viewing.
- To zoom in on a selected region for a trace, in the Trace pane, click and hold the left mouse button and draw a box from the upper left hand corner of the region towards the lower right hand corner of the region. A box is formed around the area that is being reduced for viewing.
- To zoom out from the sample files or trace, click and hold the left mouse button and draw a box from the lower right hand corner of the file sequence or trace towards the upper left hand corner of the sequence or trace. The magnification for zooming out is always 100%.




By default, the Base pane and Trace pane are linked in the Project Reviewer. If you zoom in on a region in one of the panes or zoom out from a region in one of the panes, then this zooming action is automatically duplicated in the same region in the other pane. You can turn off this automatic linking. See [“Project Reviewer report toolbar” on page 102](#).

- To change the view in the reviewer to a selected mutation, double-click the mutation in the Mutation Table pane.
- To change the view to that for a different contig, select the contig on the Contig dropdown list.




If you select All, then any overlap among the contigs is displayed in the Base pane and in the Trace pane

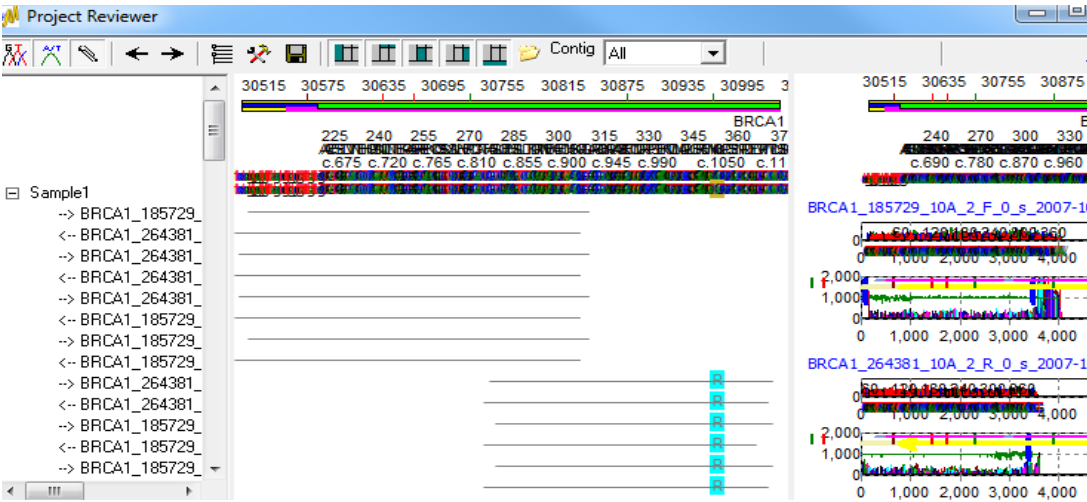Figure 4-2: Overlapping contigs displayed in the Base pane



- To manipulate and analyze mutation project data, right-click a mutation call in the Mutation Table pane and on the context menu that opens, click the appropriate option.

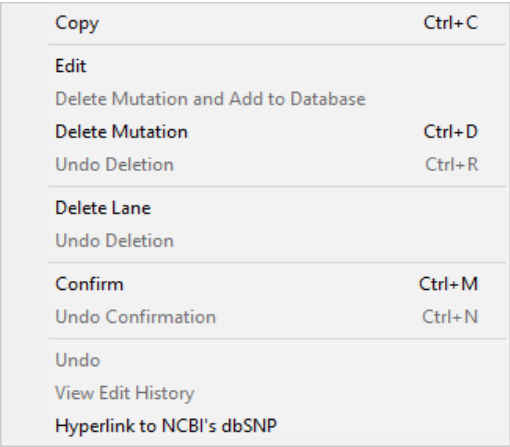Figure 4-3: Mutation Table pane context menu



Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Project Reviewer report as an image. Press and hold the [Shift] key and the [Ctrl] key, and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>

Option	Description
Edit	Opens the Edit Substitution or Edit Indel dialog box in which you can edit the selected mutation call. Note: For detailed information about modifying a mutation call, see Chapter 12, "Reviewing and Editing Mutation Projects," on page 327.
Delete	Deletes the selected mutations. Note: SHIFT-click to select multiple contiguous mutations. The text for the mutation code turns gray in both the Project Reviewer report and the Mutation Table pane.
Undo Deletion	Reverses the deletion of the selected mutation. Note: SHIFT-click to select multiple contiguous mutations.
Delete Mutation and Add to Database	Deletes the selected mutation and adds it to the database.
Delete Lane	Note: Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to "Low Quality." Note: SHIFT-click to select multiple contiguous samples.
Undo Deletion	Reverses the deletion of the selected lane. Note: SHIFT-click to select multiple contiguous samples.
Confirm	Confirms the selected mutations. The text for the mutation code turns black in both the Project Reviewer report and the Mutation Table pane. Note: SHIFT-click to select multiple contiguous mutations.
Undo Confirmation	Removes the "Confirmed" status for the mutation. Note: SHIFT-click to select multiple contiguous mutations.
Hyperlink to NCBI's dbSNP	Opens the NCBI's database entry (via the web) for the variant.

Project Reviewer report toolbar

The Project Reviewer report toolbar contains buttons that are specific for navigating and working with the Project Reviewer report. The buttons that are displayed on the toolbar depend on whether a single contig is selected or all contigs are selected.

Figure 4-4: Project Reviewer report toolbar for a selected contig

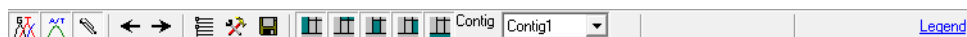



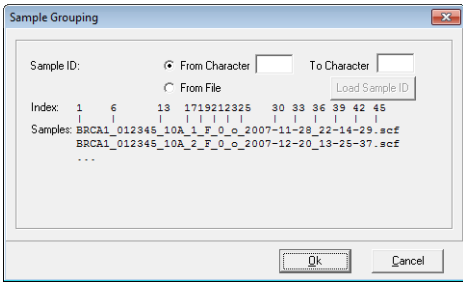
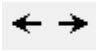










Figure 4-5: Project Reviewer report toolbar for all contigs



Button	Description
	Show Sample Trace—Selected by default. Shows the sample traces in the Trace pane.


Button	Description
	Show Mutation Trace—Selected by default. Shows the mutation traces in the Trace pane.
	Link—Selected by default. Links the Base pane and the Trace pane. If you zoom in on a region in one of the panes or zoom out from a region in one of the panes, then this zooming action is automatically duplicated in the same region in the other pane. You can turn off this automatic linking.
	<p>Sample Grouping—Opens the Sample Grouping dialog box in which you can specify how sample traces are to be grouped in the report.</p> <p><i>Figure 4-6: Sample Grouping dialog box</i></p>  <ul style="list-style-type: none"> From Character [] To Character []—Select this option if all the filenames in your project contain the same character string/designator in the same positions. From File—Select this option if traces that are to be grouped together contain a common Sample ID in their filenames. (The Sample IDs do not need to be in the same location in the filename.) After you click From File, click Load Sample ID to open the Load Sample ID dialog box and specify the different Sample IDs, such as Patient #, by which to group the traces. You can also click Open in the Load Sample ID dialog box to load an existing Sample ID text file. <p>Note: By default, the Load Sample ID dialog box contains rows for up to ten different Sample IDs. You can right-click and click Insert or Delete as needed.</p>
	Move Left/Move Right—Moves the display in the Base pane and Trace pane in the indicated direction.
	Report Settings—Opens the Project Reviewer Report Settings dialog box, in which you can specify what information is to be displayed in the report, what color-coding is to be used in the report, and what nomenclature is to be used in the report. See “Project Reviewer Report Settings dialog box” on page 104.
	Saves current view as a JPG file.
	Show/Hide Sample—Shows/hides the Sample pane.
	Show/Hide Annotation—Shows/hides the gene/CDS annotation in the Base pane and the Trace pane.

Button	Description
	Show/Hide Base—Shows/hides the Base pane.
	Show/Hide Trace—Shows/hides the Trace pane.
	Show/Hide Mutation Table—Shows/hides the Mutation Table pane.
	Add GenBank File—Available only if All contigs are selected. If you used SEQ files as reference files, click the this button to open the Add GenBank file dialog box and select a GenBank to add to the project. The Project Reviewer then aligns the SEQ files to the GenBank and as a result, the SEQ files are displayed in genomic order.

Project Reviewer Report Settings dialog box

The Project Reviewer Report Settings dialog box contains four tabs—[Mutation](#), [Color](#), [Nomenclature](#), and [Display](#)—on which you can specify what information is to be displayed in the report, what color-coding is to be used in the report, what nomenclature is to be used in the report, and how many traces to view together, respectively.

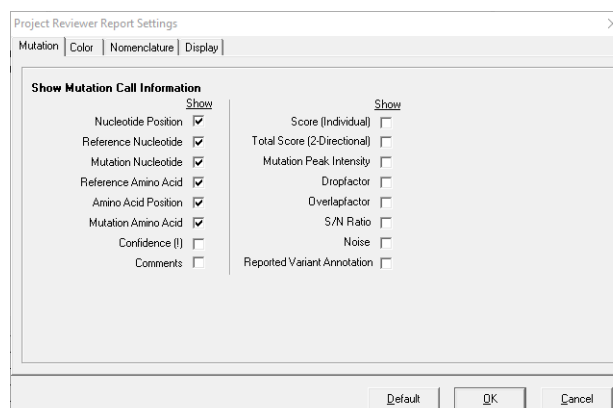Figure 4-7: Project Reviewer Report Settings dialog box, Mutation tab



Mutation tab



The majority of options on this tab are self-explanatory, and therefore, not all these options are discussed in this section.

- Nucleotide Position—Nucleotide positions are displayed in the nomenclature that you select on the Nomenclature tab. (See “[Nomenclature tab](#)” on page 106.)
- Amino Acid Position—By default, amino acid information is organized with the reference amino acid displayed first, the amino acid position displayed next, and the mutation amino acid displayed last. If you do not select Amino Acid Position, then a slash (/) is displayed between the reference amino acid and the mutation amino acids.

- **Confidence(!)**—Displays a Confidence mark (!) for mutation calls that exceed certain thresholds as defined in the MutationSurveyor.ini file.




Although you can modify these thresholds, SoftGenetics does not recommend that you do so on your own. Contact SoftGenetics for assistance.

- **Comments**—Displays all comments that were added to a mutation call, either automatically by Mutation Surveyor or manually by a user.

Color tab

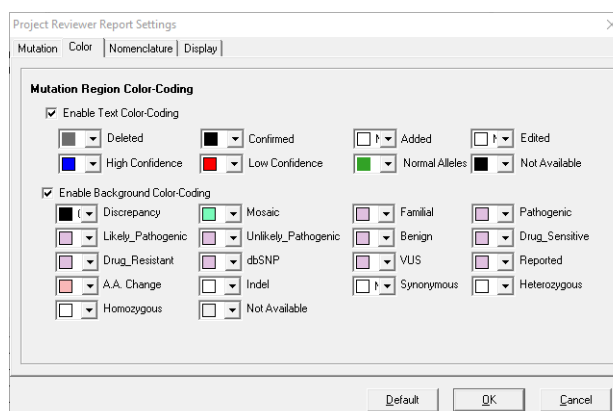Figure 4-8: Project Reviewer Report Settings dialog box, Color tab




The Color tab provides a variety of color-coding options to assist you with identifying different types of mutation calls and potential analysis problems. Color-coding is applied according to the following:

- If a mutation call meets several color-coding criteria, the call is displayed in the report in the color of highest priority.
- If None is selected for a color-coding option, then the color of the next highest priority is used.
- For Mutation Region Text color-coding, the priorities from highest to lowest are as follows—Deleted, Confirmed, Added, Edited, High Confidence, Low Confidence, Normal Alleles, and Not Available.
- For Mutation Region Background color coding, priorities from highest to lowest are as follows—Mosaic, Reported, A.A. Change, Indel, Synonymous, Heterozygous, Homozygous, and Not Available, where *Synonymous* identifies the coding and non-coding substitutions that do not affect the amino acid sequence.

Nomenclature tab

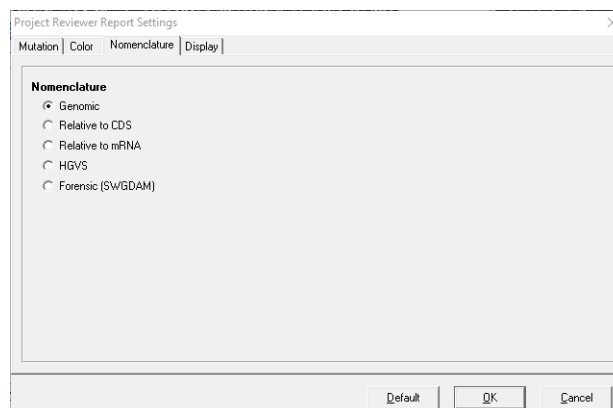Figure 4-9: Project Reviewer Report Settings dialog box, Nomenclature tab



If you select Nucleotide positions on the Mutation tab, then the positions are displayed in the nomenclature that you select on the Nomenclature tab.

Option	Description
Genomic	Displays the mutation call positions based on the genomic position referenced in the GenBank file.
Relative to CDS	Displays the mutation call positions relative to the CDS numbering in the GenBank reference file. A mutation that is called in the CDS is prefixed with a "c." A mutation that is called in a non-coding region is prefixed with "IVS." Its position is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide.
Relative to mRNA	Displays the mutation call positions relative to the mRNA numbering in the GenBank reference file. A mutation that is called in the mRNA region is prefixed with a "c." A mutation that is called outside of an mRNA region is identified by the closest nucleotide in the mRNA region plus or minus the number of bps by which it resides away from the closest nucleotide.
HGVS	Displays the mutation call positions relative to its position in the coding sequence. A mutation that is called in a non-coding region is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide. Both alleles are shown in separate bracketed statements.
Forensic (SWGAM)	Displays the mutation call using SWGDAM Guidelines.
Note: If any of the five standard Nomenclature options is selected and a GenBank file was not loaded as the reference file, then the Base Number is enclosed in parenthesis and it indicates the position of the base in the reference file. If both a GenBank file and a reference file were loaded for the mutation project and the mutation was called at a position that does not align to the GenBank file, then the Base Number is enclosed in parenthesis and is displayed as +/- the number of base pairs by which the mutation was shifted in comparison to the GenBank file.	


Figure 4-10: Nomenclature Examples

			Genomic	Relative to CD5	Relative to mRNA	HGV5	Forensic (SWGDAM)
Substitution	Homozygous	non-coding	873C>T	IV541-14C>T	c.28-14C>T	c.[41-14C>T]+[41-14C>T]	C873T
		coding	77A>T	c.31A>T	c.18A>T	c.[31A>T]+[31A>T]	A77T
	Heterozygous	non-coding	821C>CT	IV541-66C>CT	c.[=,28-66C>T]	c.[41-66C>T]+[=]	C821Y
		coding	80C>CA	c.34C>CA	c.[=,21C>A]	c.[34C>A]+[=]	C80M
Indel	Homozygous	non-coding	211delT	IV571+4delT	c.58+4delT	c.71+4delT	211d
		coding	18_20delTTT	c.72_74delTTT	c.59_61delTTT	c.72_74delTTT	18d,19d,20d
	Heterozygous	non-coding	133_134het_delTA	IV571+16_71+17het_delTA	c.58+16_58+17het_delTA	c.71+16_71+17het_delTA	133d,134d
		coding	14_15het_insT	c.68_69het_insT	c.55_56het_insT	c.68_69het_insT	15.1T

Display tab

Select the number of traces to display in the Trace Pane. Default is 4. A maximum of 8 can be selected.

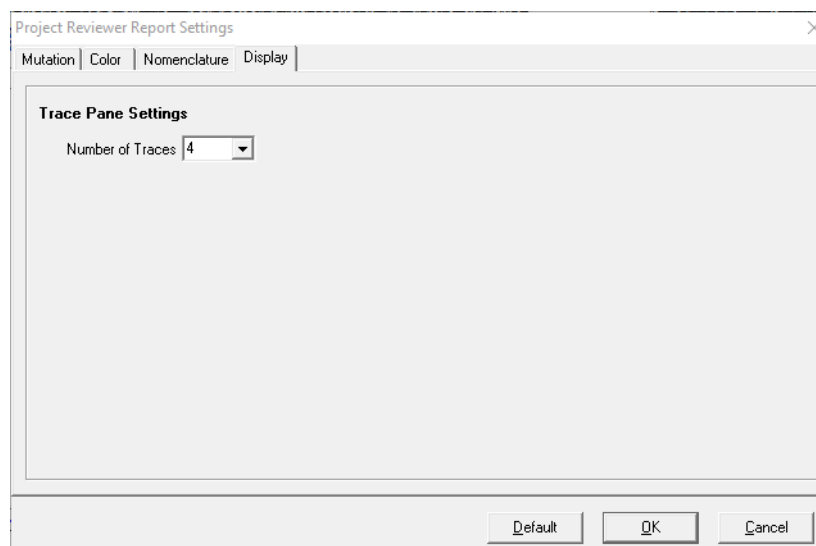Figure 4-11: Project Reviewer Report Settings dialog box, Display tab



Graphic Display Report


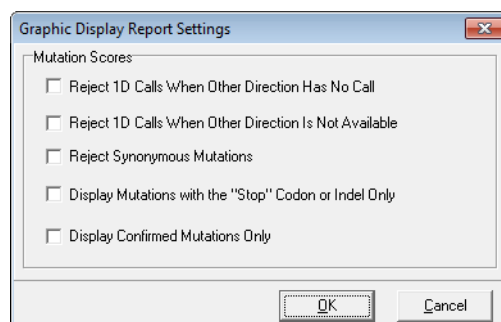
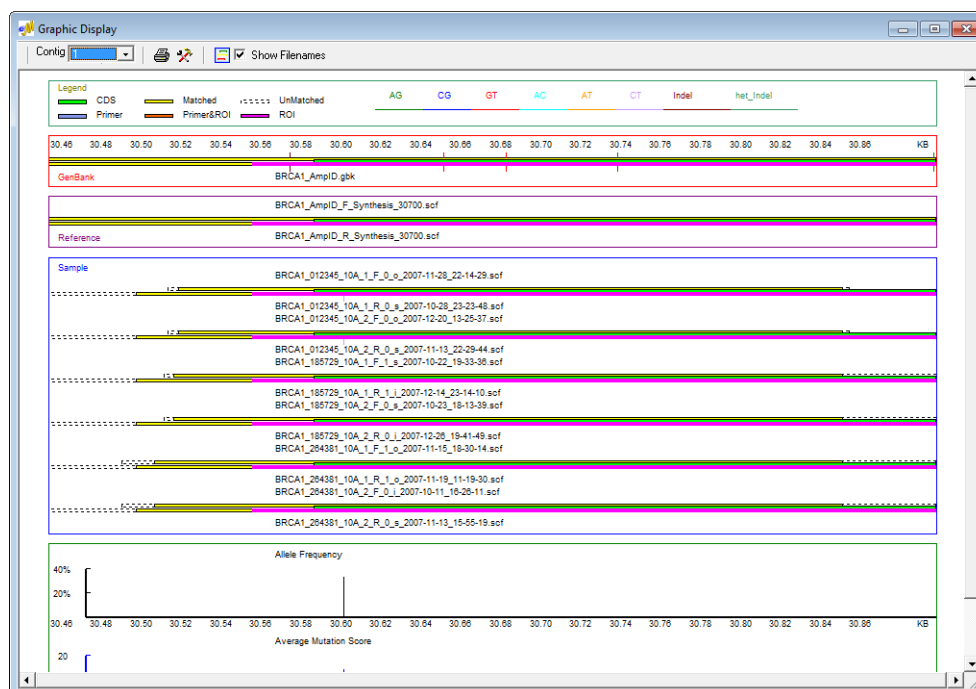
You use the Graphic Display report to examine sample alignment and mutation calls for a single contig. The Graphic Display report is available from the Display option on the Mutation Surveyor main menu or you can click the Graphic Display icon  on the GAD. After you select either option, the Graphic Display Report Settings dialog box opens. You specify the options in this dialog box, and then click OK to generate the report. (See [Figure 4-13 on page 109.](#))

Figure 4-12: Graphic Display Report Settings dialog box



Option	Description
Reject 1D Calls When Other Direction has No Call	Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction.
Reject 1D Calls When Other Direction is Not Available	Reject mutation calls that occur outside the region where the sample traces in both directions are aligned within the comparison region.
Reject Synonymous Mutations	Coding region substitutions that do not affect amino acid calls are not displayed in the report.
Display Mutations with the "Stop" Codon or Indel Only	Display only nonsense mutations and indels in the report.
Display Confirmed Mutations Only	Display only those mutations that you have manually confirmed in the report.

Figure 4-13: Graphic Display report



Graphic Display report layout

The top row of the Graphic Display report details the color coding legends that are used in the report. The report displays mutation project data below these legends in four panes. From top to bottom, these panes are as follows:

- **GenBank pane**—The GenBank pane displays information about the loaded GenBank file. Contig positions are indicated in kilobases (kb). The color of the horizontal *location bars* that run below the contig position numbers indicates the location of the sequence that is being displayed in the trace. Colored tick marks indicate reported variations.
- **Reference pane**—The Reference pane displays information about the reference trace that was used for the selected contig. The trace can be a reference file that you manually added to the project, or it can be a synthetic reference file that Mutation Surveyor created from the loaded GenBank file.
- **Sample pane**—The Sample pane displays all the sample files that are contained within the selected contig. Tick marks represent called mutations. Color-coding is used to indicate the mutation type.
- **Allele Frequency pane**—The Allele Frequency pane shows the percentage of samples in the contig that have the mutation call at the indicated position.
- **Mutation Score pane**—The Mutation Score pane shows the average mutation score for all the calls at the indicated position for the contig.

The report is interactive:

- To zoom in on a selected region, in any pane, click and hold the left mouse button and draw a box from the upper left hand corner of the region towards the lower right hand corner of the region. A box is formed around the area that is being reduced for viewing.
- To zoom out from any region, click and hold the left mouse button and draw a box from the lower right hand corner of the region towards the upper left hand corner of the region. The magnification for zooming out is always 100%.



All the panes in the Graphic Display are linked. If you zoom in on a region in any pane, or zoom out from a region in any pane, then this zooming action is automatically duplicated in the same regions in all the other panes.




- To change the view to that for a different contig, select the contig on the Contig dropdown list.
- By default, the filenames for all the files (GenBank, reference files, and sample files) are displayed on the report. To remove the filenames from the report display, clear the Show Filenames option.

Graphic Display report toolbar

The Graphic Display report toolbar contains buttons that are specific for navigating and working with the Graphic Display report.

Figure 4-14: Graphic Display report toolbar




Button	Description
	Print Preview—Opens the Graphic Display Print Preview window, which displays a preview of what is to be printed for the report. Note: Before you print the Graphic Display report, make sure to adjust the zoom level so that the area of the report that you want to print is displayed in the report window.
	Report Settings—Opens the Graphic Display Report Settings dialog box, in which you can specify what information is to be displayed in the report. See Figure 4-12 on page 108 .
	Project Display Report—Opens the Project Display Report Settings dialog box, in which you specify the information that is to be displayed in the Project Display subreport, as well as how the sample files are to be grouped in the subreport. See “Project Display subreport” on page 111 .
Show Filenames	Selected by default. Shows the filenames on the report.

Project Display subreport

You use the Project Display subreport as another way of graphically viewing mutation project data. Options are available for filtering mutation calls that are displayed in the report and for grouping sample files based on the character string/designator that is used in your sample filenames. This subreport, which has a layout that is almost identical to that of the Graphic Display report, is ideal for comparing overlapping sample traces that were analyzed in separate contigs. The subreport opens after you click OK in the Project Display Report Settings dialog box.

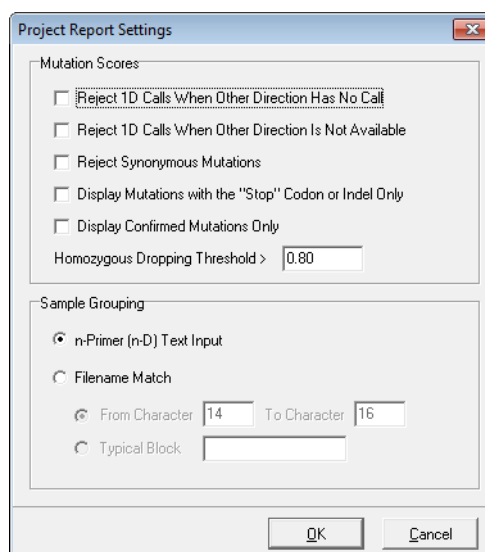Figure 4-15: Project Display Report Settings dialog box




Option	Description
Mutation Scores	
Reject 1D Calls When Other Direction has No Call	Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction.
Reject 1D Calls When Other Direction is Not Available	Reject mutation calls that occur outside the region where the sample traces in both directions are aligned within the comparison region.
Reject Synonymous Mutations	Coding region substitutions that do not affect amino acid calls are not displayed in the report.
Note: In cases where the entire column was not omitted from the report because of rejection by one of these criteria, the rejected mutations have no background color and an “n.a.” is displayed in the Mutation cell instead of the mutation code.	
Display Mutations with the “Stop” Codon or Indel Only	Display only nonsense mutations and indels in the report.
Display Confirmed Mutations Only	Display only those mutations that you have manually confirmed in the report.

Chapter 4

Mutation Surveyor Graphical Reports

Option	Description
Homozygous Dropping Threshold	<p>Indicates the value for the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions.</p> <p>You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.</p>
Sample Grouping	
	<p>Specify the filename substring by which Mutation Surveyor groups all the mutation project samples files in the subreport.</p> <ul style="list-style-type: none"> n-Primer (n-D) Text Input—Use the ND Grouping file to group the sample files. See "ND Filename Match Editor Tool" on page 264. Filename Match (based on one of the following): <ul style="list-style-type: none"> From Character [] To Character []—If all the filenames in your project contain the same characters/designator in the same positions, then select this option and enter the starting and ending character positions. <p>Note: Typical Block—If all the filenames in your project contain the same characters/designator but in different positions, then select this option and enter the value for the character/designator block.</p>

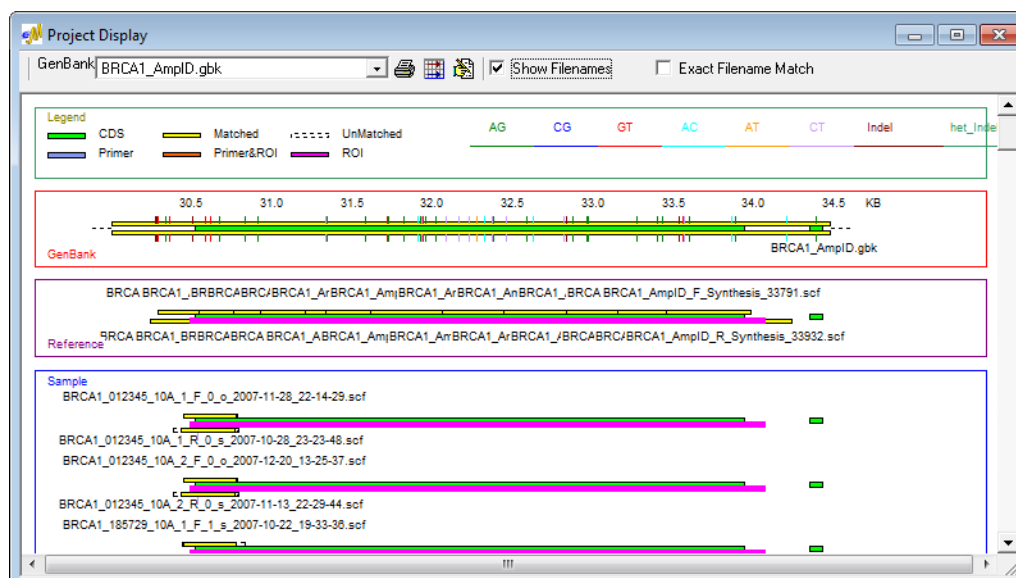Figure 4-16: Project Display subreport






Project Display subreport toolbar

The Project Display subreport toolbar contains options and buttons that are specific for navigating and working with the subreport.

Figure 4-17: Project Display subreport toolbar



Option/Button	Description
Select GenBank File	If your mutation project used multiple GenBank files, then select one of these files from this list to view the file and the sample files that align to it.
	Print Preview—Opens the Preview window, which displays a preview of what it to be printed for the report. Note: Before you click the Print Preview button, make sure to adjust the zoom level so that the area of the report that you want to print will be displayed in the Preview window.
	Project Report—Generates a Project report for the sample files. See “Project Report” on page 146 .
	Modify Mutation Projects—See “Modifying Mutation Surveyor Projects” on page 315 .
Show Filenames	By default, the filenames for all the files (GenBank, reference files, and sample files) are displayed on the subreport. To remove the filenames from the report display, clear the Show Filenames option.
Exact Filename Match	Select this option to pair sample files in the report display when the filenames for the forward and reverse sample files are very similar. Note: This sorting option is used <i>in conjunction</i> with the sorting options that you specified on the Project Display Report Settings dialog box to further group sample files in the report display.

GAD Report

The GAD report displays the Graphical Analysis Display (GAD) for the sample files that are contained in a single contig in a report format. You can specify whether the report is to display all the traces for all the sample files in the contig, or whether the report is to display only those traces for the sample files that contained mutation calls (the default setting for the report). After you click GAD Report on the Display menu, the GAD Report Settings dialog box opens. You can leave the default option of Print Samples with Mutations Only selected, or you can clear this option to display all the traces for all the sample files for a selected contig in the report. After you click OK, the GAD report opens.

Figure 4-18: GAD Report Settings dialog box


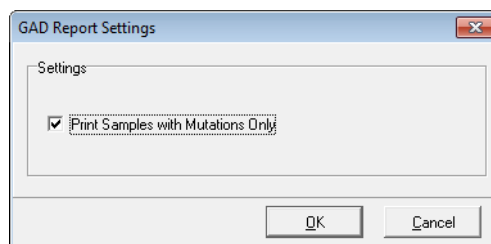


Figure 4-19: GAD report

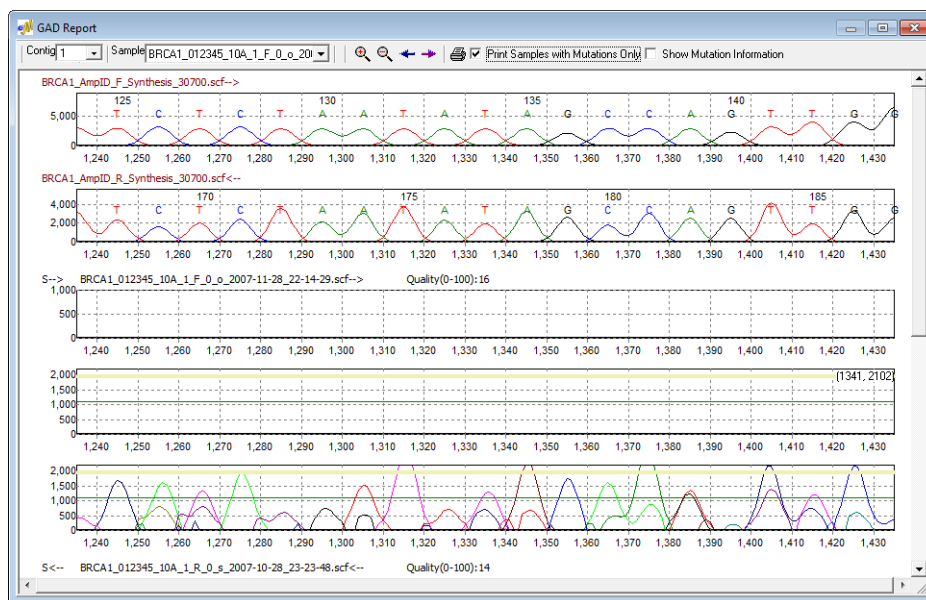






The first two traces that are displayed in the report are the forward and reverse reference traces used for the selected contig. (The names of the traces are displayed in brown font to indicate that they are reference traces.) The remaining traces are the traces for the sample files in the selected contig. The traces are either the traces for all the sample files that are contained in the selected contig, or they are those traces for display the trace regions for only those samples in the selected contig that contain mutation calls. To view the traces for a different contig, select the contig on the Contig dropdown list.

GAD Report toolbar

The GAD report toolbar contains options and buttons that are specific for navigating and working with the report.

Figure 4-20: GAD report toolbar




Option/Button	Description
Sample	Indicates the sample for which the sample traces are currently in view for the report. You can select a different sample from the list, or you can use the Scroll feature on the right side of the report window to change the view.
	Zoom In—Zooms in on a region of the report for viewing. Note: You can also manually zoom in on a region for viewing. Click and hold the left mouse button and draw a box from the upper left hand corner of the pane towards the lower right hand corner. A box is formed around the area that being reduced for viewing.
	Zoom Out—Zooms out from a region of the report for viewing. The magnification for zooming out is always 100%. Note: You can also manually zoom out from a region for viewing. Click and hold the left mouse button and draw a box from the lower right hand corner of the pane towards the upper left hand corner.
	Move Left—Moves the currently displayed report view to the left. Note: You can also manually move the report view to the left. Right-click and hold the mouse button on the report view, and then move the mouse.
	Move Right—Moves the currently displayed report view to the right. Note: You can also manually move the report view to the left. Right-click and hold the mouse button on the report view, and then move the mouse.
	Print Preview—Opens the Preview window, which displays a preview of what it to be printed for the GAD report. You can use the options on the Preview window's toolbar to adjust variables such as the color scale and the page layout before you print the report. Note: Before you click the Print Preview button, make sure to adjust the zoom level so that the area of the report that you want to print will be displayed in the Preview window.
Print Samples with Mutations Only	Select this option to display only those traces for the sample files that contained mutation calls.
Show Mutation Information	Select this option to display the code for each mutation call on the report. The code for each call is displayed directly above the mutation peak in the sample trace.

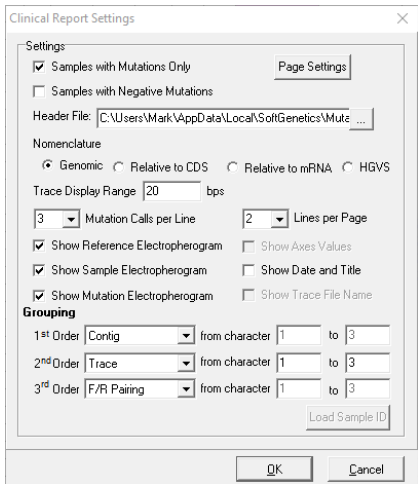
Clinical Report

The Clinical report is a custom-formatted report that contains options for displaying only those sample files with mutation calls or for displaying all positions with reported annotation information, including negative SNPs. The report shows one pair of sample files in a contig per page, and the report includes the mutation code for each called mutation. To generate a Clinical report, modify any of the settings as needed and click OK. The report opens in the Print Preview window. You can use the options on the Preview window’s toolbar to adjust variables such as the color scale and the page layout. See [Figure 4-23 on page 119](#).

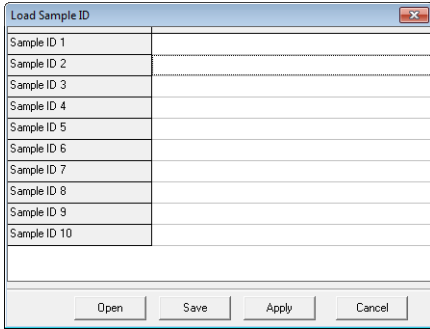
The Clinical Report Settings dialog box opens.

Figure 4-21: Clinical Report Settings dialog box




Option	Description
Print Samples with mutations only	Show only those samples that contain called mutations.
Print Samples with negative mutations	Show those samples that contain negative mutation calls, where a negative SNP is defined as position in the GenBank file at which a variation has been reported, but no variation was detected at the same position in the sample trace. Sample traces that have negative SNPs might still contain mutation calls.
Header File	Click the Browse button  to open the Open File dialog box, and browse to and select a custom header to add to the report. Note: A header is a .inf file. For a discussion of Custom headers and how to create one, see “Print Header Editor Tool” on page 267 .

Option	Description
Nomenclature	<ul style="list-style-type: none"> • Genomic—Displays the mutation call positions based on the genomic position referenced in the GenBank file. • Relative to CDS—Displays the mutation call positions relative to the CDS numbering in the GenBank reference file. A mutation that is called in the CDS is prefixed with a “c.” A mutation that is called in a non-coding region is prefixed with “IVS.” Its position is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide. • Relative to mRNA—Displays the mutation call positions relative to the mRNA numbering in the GenBank reference file. A mutation that is called in the mRNA region is prefixed with a “c.” A mutation that is called outside of an mRNA region is identified by the closest nucleotide in the mRNA region plus or minus the number of bps by which it resides away from the closest nucleotide. • HGVS—Displays the mutation call positions relative to its position in the coding sequence. A mutation that is called in a non-coding region is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide. Both alleles are shown in separate bracketed statements.
Trace Display Range <> bp	Total number of bases in range of mutation call, with call centered in this range. The value can be an integer from 1 to 255. The default value is 20.
Mutation Calls per Line	<p>Maximum number of mutation calls that would be displayed in a single row of the report. Options are:</p> <ul style="list-style-type: none"> • 1 • 2 • 3 (Default value)
Lines per Page	<p>Maximum number of lines of mutation calls that would be displayed on a single page of the report. If a trace contains more than would fit on a single page, the report would include additional pages. Options are:</p> <ul style="list-style-type: none"> • 1 • 2 (Default value)
Show Reference Electropherogram	Option to display the bases of reference trace at the location of the mutation call. This option is selected by default.
Show Sample Electropherogram	Option to display the bases of the sample trace at the location of the mutation call. This option is selected by default.
Show Mutation Electropherogram	Option to display the peaks of the mutation trace at the location of the mutation call. This option is selected by default.
Show Axes Values	Option to display or hide the y-axis labels and gridlines. This option is unavailable for 3 mutation calls per line. This option is deselected by default.
Show Date and Title	Option to display or hide the current date/time and the first order grouping type at the top of report's heading section. This option is deselected by default.
Show Trace File Name	Option to display each trace's file name of its image. This option may not be available for 2 lines per page, depending on electropherograms to be displayed. This option is deselected by default.

Option	Description
Grouping —Traces are initially grouped based on the 1st Order Grouping option. Further sorting is then carried out based on the 2nd Order Grouping option, and then on the 3rd Order Grouping option.	
<ul style="list-style-type: none"> Contig Trace Grouping File Sample ID 	<ul style="list-style-type: none"> The default value for 1st Order. Groups are sorted based on the Contigs that Mutation Surveyor created for determining mutation calls. When Trace is selected, the From Character X to Y option is enabled. Traces are sorted into groups based on character matching in the specified positions in the trace filenames. Traces are sorted into groups based on the grouping method that is specified in the Load Grouping text file. See “2D Filename Match Editor tool” on page 261. When Sample ID is selected, both the Load Sample ID button and the From Character X to Y option are enabled. Click Load Sample ID to open the Load Sample ID dialog box and specify the different Sample IDs, such as Patient #, by which to group the traces. You can also click Open to load an existing Sample ID file. <p>Note: By default, the Load Sample ID dialog box contains rows for up to ten different Sample IDs. You can right-click and click Insert or Delete as needed.</p> <p><i>Figure 4-22: Load Sample ID dialog box</i></p>  <p>The traces that are to be grouped together must contain at least one of these Sample IDs in their filenames but they do not need to be in the same location in the filename. When grouping by Sample ID, the trace filenames are first checked based on the Sample IDs defined in the Load Sample ID dialog box. The values specified in the From Character X to Y fields are ignored if all the trace filenames contain at least one of these Sample IDs. Traces that do not contain any of these Sample IDs in their filenames are sorted into groups based on the values specified in the From Character X to Y fields. Traces that do not contain any of these Sample IDs and that do not contain the characters specified in the From Character X to Y fields are sorted into their own group.</p>
<ul style="list-style-type: none"> F/R pairing 	<ul style="list-style-type: none"> The default value for the 2nd Order. Places two-directional traces in adjacent rows.

Option	Description
<ul style="list-style-type: none"> Trace Position 	<ul style="list-style-type: none"> The default value for the 3rd Order. Traces are sorted into groups based on where they match to the GenBank reference file. Each trace is matched at its starting genomic base number. Groups can be sorted so that traces that match to the earliest reference base are displayed first.
<ul style="list-style-type: none"> Gene 	<ul style="list-style-type: none"> Traces are sorted into groups based on the gene names defined in the GenBank reference file.
<ul style="list-style-type: none"> Exon 	<ul style="list-style-type: none"> Traces are sorted into groups based on the exons defined in the GenBank reference file.
<ul style="list-style-type: none"> Trace Quality 	<ul style="list-style-type: none"> Traces are sorted into groups based on the lane quality for the trace.
<ul style="list-style-type: none"> None 	<ul style="list-style-type: none"> Ignore the sorting level.
<ul style="list-style-type: none"> In One Group 	<ul style="list-style-type: none"> Traces are placed in a single group. Only available for the 1st Order.

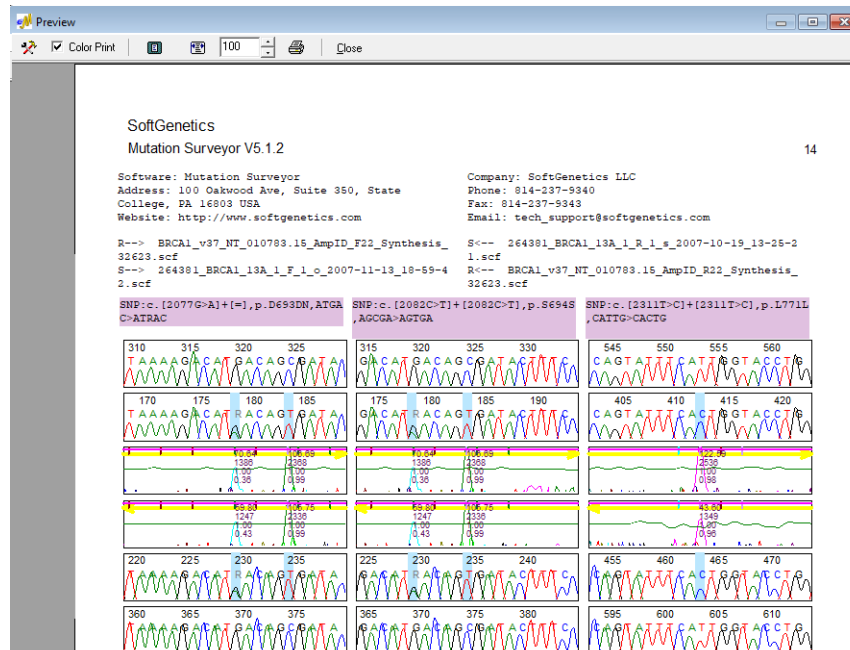Figure 4-23: Clinical report




Nucleotide Text Display Report

The Nucleotide Text Display report is a color-coded report of the text base calls of all the sample file sequences for a selected contig. The report has options for displaying every nucleotide position for every sample trace in a selected contig or for displaying only those nucleotide positions at which a mutation was called. To generate a Nucleotide Text Display report, do the following:

- 1. On the Display menu, click Nucleotide Text Display.
The Nucleotide Text Display Report Settings dialog box opens.

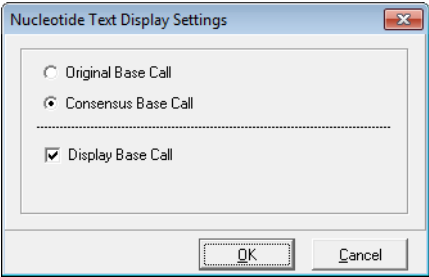Figure 4-24: Nucleotide Text Display Report Settings dialog box




- 2. Select the options as appropriate for your report:

Option	Description
Original Base Call	Display the bases as called in the original sample trace at locations where variations exist between the sample and the reference.
Consensus Base Call	Display the base relative to the mutation call at locations where mutation calls were identified.
Display Base Call	Selected by default. Display every nucleotide position for every sample trace in a selected contig. If you clear this option, then only the variations relative to the Original Base Call or the Consensus Base Call are displayed for the sample trace in the selected contig.

- 3. Click OK.

The Nucleotide Text Display report opens. See [Figure 4-25 on page 121](#).

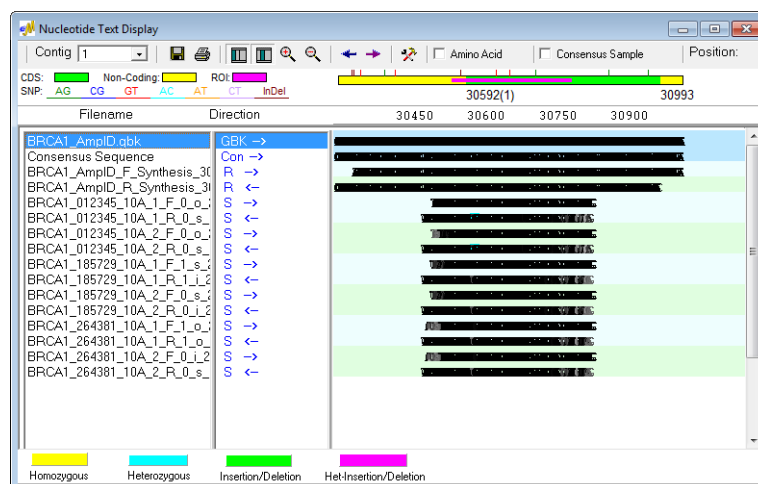Figure 4-25: Nucleotide Text Display report



Nucleotide Text Display report layout

The top row of the Nucleotide Text Display report details the color coding legends that are used in the report. The color of the horizontal *location bars* indicates the location of the sequence that is being displayed in the trace. Colored tick marks indicate reported variations. The report displays mutation project data below these legends in three panes. From left to right these panes are:


- **Filenames pane**—The first row in the Filenames pane is the name of the reference nucleotide sequence that was used to align the sample files. The second row in the Filenames pane is the consensus sequence for the sample traces. The remaining rows in the pane shows the name for each reference and sample trace contained in the selected contig.
- **Direction pane**—The Direction pane shows the trace direction for the reference sequence, for the consensus sequence, and for each sample trace for the selected contig.
- **Nucleotide Text pane**—The Nucleotide Text pane shows the textual representation of the reference sequence, the consensus sequence, and for each sample trace contained in the selected contig.

The report is interactive:

- To view the nucleotide text display for a different contig, select the contig on the Contig dropdown list.
- To view the mutation electropherograms for sample files at a given mutation position, double-click the mutation in the report. The Contig Trace Figure window opens. (See [Figure 4-26 on page 122.](#)) You can use the Scroll function on this window to scroll through all the sample traces at the selected mutation position. Each mutation call is highlighted with either a blue or purple vertical line, for novel or reported variants

respectively.

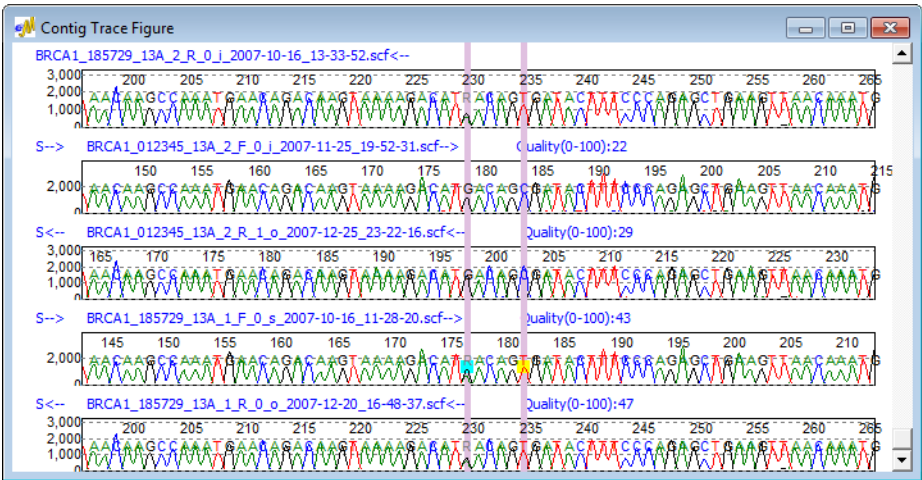Figure 4-26: Contig Trace Figure window



- To highlight the information for a sequence in the report, click the sequence in the Filenames pane.
- To view the Position (Frame #, Nucleotide #, and Row #) for a base, point the cursor on the base in the Nucleotide Text pane. The Position information is updated on the report toolbar.




These last two options work best after you have zoomed in on a region in the Nucleotide Text pane for viewing. See the Zoom In function on the “Nucleotide Text Display report toolbar.”









Nucleotide Text Display report toolbar

The Nucleotide Text Display report toolbar contains options and buttons that are specific for navigating and working with the report.

Figure 4-27: Nucleotide Text Display report toolbar



Option/Button	Description
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as a tab-delimited text file.

Option/Button	Description
	<p>Print Preview—Opens the Preview window, which displays a preview of what it to be printed for the Nucleotide Text Display report. You can use the options on the Preview window's toolbar to adjust variables such as the color scale and the page layout before you print the report.</p> <p>Note: Before you click the Print Preview button, make sure to adjust the zoom level so that the area of the report that you want to print will be displayed in the Preview window.</p>
	Show/Hide Filenames—Shows/hides the Filenames pane.
	Show/Hide Direction—Shows/hides the Direction pane.
	<p>Zoom In—Zooms in on a region of the report for viewing.</p> <p>Note: You can also manually zoom in on a region for viewing. Click and hold the left mouse button and draw a box from the upper left hand corner of the pane towards the lower right hand corner. A box is formed around the area that being reduced for viewing.</p>
	<p>Zoom Out—Zooms out from a region of the report for viewing. The magnification for zooming out is always 100%.</p> <p>Note: You can also manually zoom out from a region for viewing. Click and hold the left mouse button and draw a box from the lower right hand corner of the pane towards the upper left hand corner.</p>
	<p>Move Left—Moves the currently displayed report view to the left.</p> <p>Note: You can also manually move the report view to the left. Right-click and hold the mouse button on the report view, and then move the mouse.</p>
	<p>Move Right—Moves the currently displayed report view to the right.</p> <p>Note: You can also manually move the report view to the left. Right-click and hold the mouse button on the report view, and then move the mouse.</p>
	Report Settings—Opens the Nucleotide Text Display Report Settings dialog box in which you can specify the information that is to be displayed in the report. See Figure 4-24 on page 120 .
Amino Acid	Select this option to display the amino acid sequence for which the consensus sequence codes. The code is displayed in the top row of the Nucleotide Text pane as the 1 Letter IUPAC code.
Consensus Sample	Displays the consensus sequence for each of the sample pairs. An extra line marked as "Consensus Sample" is displayed below the sample pairs in all three of the report panes.

Chapter 4

Mutation Surveyor Graphical Reports

Chapter 5

Mutation Surveyor Standard Reports

A variety of text-based standard reports that provide detailed information about a mutation project are available in Mutation Surveyor. All these reports, which are available from the Reports option on the Mutation Surveyor main menu, display mutation project data according to a predefined format. You can print the report information and you can save and store the report output for use at a later date.

This chapter covers the following topics:

- [“Mutation Report” on page 127.](#)
- [“Two Direction Report” on page 134.](#)
- [“Advanced Two Direction Report” on page 140.](#)
- [“Project Report” on page 146.](#)
- [“HGVS Report” on page 157.](#)

Chapter 5

Mutation Surveyor Standard Reports

Mutation Report

Mutation Surveyor aligns the nucleotide sequences of the sample traces to the reference traces and groups the sample traces into contigs based on their degree of overlap. The reference traces and sample traces are then compared at a trace by trace level to identify substitutions, insertions, and deletions. All the variants are aligned in the Mutation report, which is automatically displayed in the Viewing and Reporting pane after the project is run. In addition, the Mutation Surveyor toolbar is updated with buttons that are specific for working with the Mutation report.








You can also open the Mutation report from the Mutation Surveyor main menu (Reports> Mutation Report), or you can double-click the Mutation Report entry in the Browser pane.

Mutation Surveyor toolbar

Figure 5-1: Mutation Surveyor toolbar



Button	Description
	Open Files—Opens the Open Files dialog box which you use to load GenBank files, references files and sample files for a project.
	Save Project—Saves all the project files to a single file mutation project file with an .sgp extension. You can open an entire project in a single step by opening this SGP file.
	Raw Data to Processed Data—Enabled when AB1/ABI files without base call information are uploaded and Load Raw Data is selected for the Mutation Project settings. Converts the raw files so that base call information is displayed.
	Run—Runs the currently loaded project in Mutation Surveyor.
	Restart —Opens the Options dialog box which you use to modify the current project settings. You can then run the project again using these new settings.
	Show/Hide Browser—Shows/hides the Browser pane in the Mutation Surveyor main window.
	Save Report—Opens the Save Report As dialog box in which you can specify the location for saving the Mutation report as a .txt file.
	Custom Report—Opens the Custom Report Settings dialog box which you use to specify the settings for a custom report. See Chapter 6, "Mutation Surveyor Custom Report," on page 163.

Button	Description
	Graphic Display—Opens the Graphic Display Report settings dialog box in which you specify the settings for generating a Graphic Display report. See “Graphic Display Report” on page 108 .
	Display Mutations with 1D/2D Settings—Toggles the Mutation report between the 1D and 2D views.
	Show Nucleotide Text—Generates a Nucleotide Text Display report. See “Nucleotide Text Display Report” on page 120 .
	Quantify—Opens the Mutation Quantification function. See “Mutation Quantifier Function” on page 303 .
	Variation Tracks Settings icon - Opens the Variation Tracks Settings dialog box. The Variation Tracks Settings dialog box displays the available tracks settings for the Mutation report based on the variation databases that were imported for the project. See “Variation Tracks Settings dialog box” on page 87 . Note: After being imported into Mutation Surveyor, a variation database is referred to as a <i>track</i> .

Mutation report fields

Figure 5-2 shows the default Mutation report in the Viewing and Reporting pane.

Figure 5-2: Mutation report

No.	Sample File	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Qualit	Mut#	Mutation1
16	BRCA1_012345_10B_2_R_1_s_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30746	31027	282	11	1		30753A>AT,278T>T/S\$21
17	BRCA1_185729_10A_1_R_1_i_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30503	30852	350	26	0		
18	BRCA1_185729_10A_2_R_0_i_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30503	30852	350	26	0		
19	BRCA1_185729_10B_1_R_0_s_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30759	31027	269	11	1		30988A>AG,356Q>Q/R\$20
20	BRCA1_185729_10B_2_R_0_s_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30759	31027	269	11	1		30988A>AG,356Q>Q/R\$20
21	BRCA1_264381_10A_1_R_1_o_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30503	30852	350	25	0		
22	BRCA1_264381_10A_2_R_0_s_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30503	30852	350	25	0		
23	BRCA1_264381_10B_1_R_0_o_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30759	31041	283	10	1		30988A>AG,356Q>Q/R\$14
24	BRCA1_264381_10B_2_R_1_o_2007-1	BRCA1_AmplD_R_S1-R	BRCA1		2	30759	31041	283	10	1		30988A>AG,356Q>Q/R\$14
25	BRCA1_012345_11A_1_F_1_s_2007-1	BRCA1_AmplD_F_S2-F	BRCA1		2	30977	31416	440	7	0		
26	BRCA1_012345_11A_2_F_1_o_2007-1	BRCA1_AmplD_F_S2-F	BRCA1		2	30977	31416	440	7	0		
27	BRCA1_012345_11B_1_F_1_i_2007-1	BRCA1_AmplD_F_S2-F	BRCA1		2	31143	31460	318	0	2		31389C>CG,490P>P/A\$22
28	BRCA1_012345_11B_2_F_1_s_2007-1	BRCA1_AmplD_F_S2-F	BRCA1		2	31143	31460	318	0	2		31389C>CG,490P>P/A\$22

The table below summarizes the default fields that are displayed in the Mutation report.

Field	Description
Sample File	A list of all the sample traces in the mutation project, including any unmatched traces.
Reference File	The reference trace to which the sample trace was matched.
Direction	The contig number and the direction of the trace (Forward (F) or Reverse (R)).
Gene	The gene name as noted in the GenBank reference file.

Field	Description
Exon	The exon as noted in the GenBank reference file.
RF	The reading frame of the exon.
Start/End	The starting and ending bases for the comparison region.
Size	The number of bps in the comparison region.
Quality	Lane quality for the trace. See “Quality Thresholds” on page 63.
Mut #	The number of mutations found in the sample trace.
Mutation	Displays the mutation code for the called mutation. See “Mutation report codes.” Note: If Comments is selected on the Output tab on the Mutation Project Options dialog box, then Comments are displayed in this field after the mutation code.

To change the fields that are displayed in the Mutation report, select or clear the appropriate options for the Output Table Fields on the Output tab on the Mutation Project Options dialog box. (See [“Output tab” on page 62.](#))


Mutation report codes

The *mutation code* is the unique code that identifies each mutation in the sample trace. The code format indicates whether the mutation is a [substitution](#) event or whether it is an [insertion/duplication](#) or [deletion](#) event.

Substitution mutation code


The mutation code for a substitution event has four components—the base number, the mutation, the amino acid sequence that is affected by the mutation, and the mutation score.

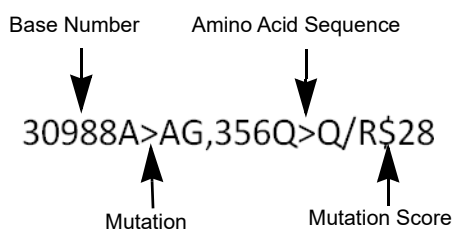Figure 5-3: Example of a substitution mutation code



- **Base Number**—If a GenBank file was loaded, the Base Number is the contig position in the GenBank file. If a GenBank file was not loaded, then the Base Number is enclosed in parenthesis and it indicates the position of the base in the reference file. If both a GenBank file and a reference file were loaded for the mutation project and the mutation was called at a position that does not align to the GenBank file, then the Base Number is enclosed in parenthesis and is displayed as +/- the number of base pairs by which the mutation was shifted in comparison to the GenBank file.

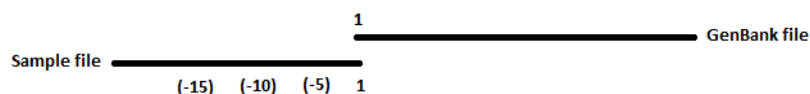Figure 5-4: Mutation called at a position that does not align to GenBank file



- **Mutation**—The mutation is written as reference base > variant base, where the *reference base* represents the base called at the position in the reference trace and the *variant base* represents the base called at the position in the sample trace. The mutation can be *homozygous*, such as A > C, or the mutation can be *heterozygous*, such as A > AC.
- **Amino Acid**—If the mutation falls within the coding region of the GenBank file, the amino acid portion of the mutation code shows the codon number, the normal amino acid that is affected by the mutation, and the resulting amino acid.
- **Mutation Score**—The mutation score is the 1D score for the mutation. See “[Mutation Score](#)” on page 59.

Figure 5-5: Example of a mutation code for a homozygous mutation


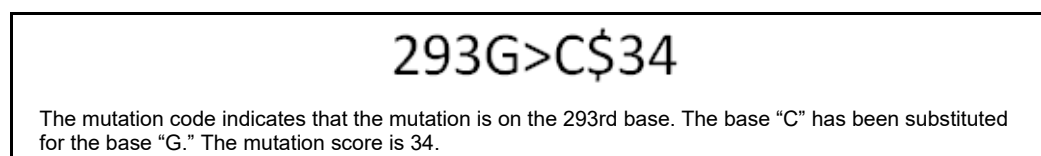


Figure 5-6: Example of a mutation code for a heterozygous mutation

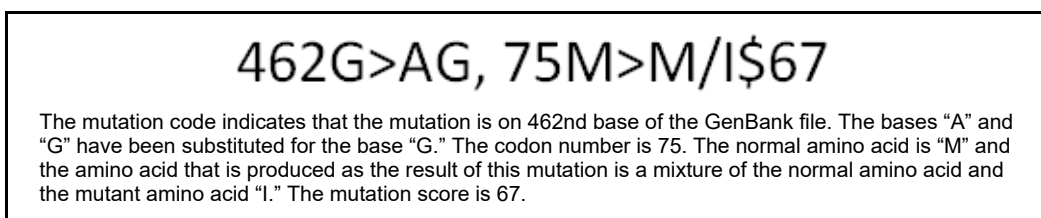

Color codes

The color of a substitution mutation code indicates a variety of information about the mutation.

- Text color—Blue text indicates that the mutation is called with high confidence. Red text indicates that the mutation is called with low confidence. (See “**Mutation Threshold**” on [page 64](#).) Black text indicates that the mutation has been manually confirmed. (See “**Confirm**” on [page 133](#).)
- Purple background with green text—A reported variation that was not found in the mutation project.
- Pink background—The mutation is a novel variant that results in a change to the amino acid sequence.
- Purple background—The mutation is a reported variant that might or might not result in

a change to the amino acid sequence.

- No background color—The mutation is a novel variant that is found on either the forward or reverse trace and that does not result in a change in the amino acid sequence.

Insertion, duplication, and deletion mutation codes

Mutation Surveyor monitors the mobility of a sample fragment relative to the reference. When the sequences no longer align, Mutation Surveyor continues to attempt alignment until it again occurs. If a homozygous insertion or duplication event is detected, Mutation Surveyor artificially gaps the reference trace. If a homozygous deletion event is detected, Mutation Surveyor artificially gaps the sample trace. If a heterozygous insertion or deletion event is detected, Mutation Surveyor automatically deconvolutes the mixed trace into two clean traces.

Figure 5-7: Example of a mutation code for a homozygous insertion event

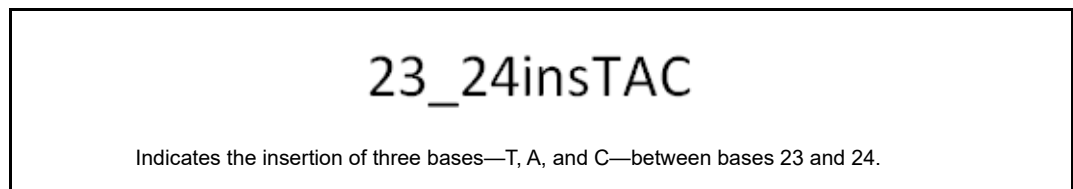


Figure 5-8: Example of a mutation code for a homozygous duplication event

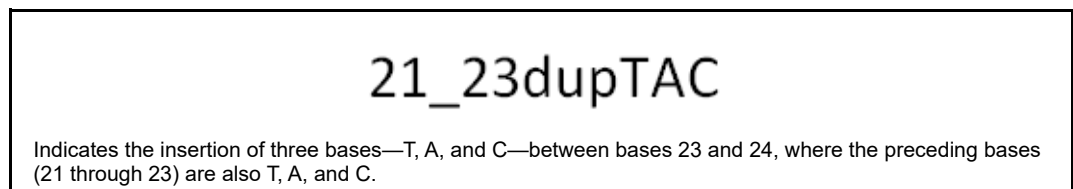


Figure 5-9: Example of a mutation code for a homozygous deletion event

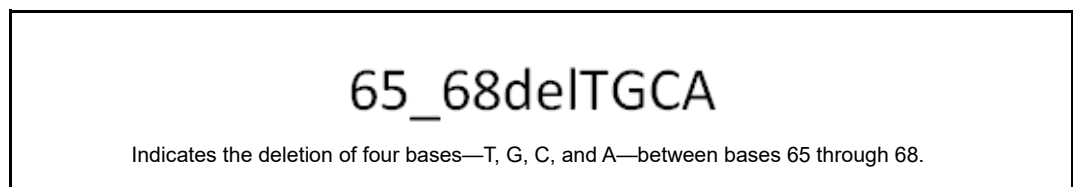


Figure 5-10: Example of a mutation code for a heterozygous insertion event

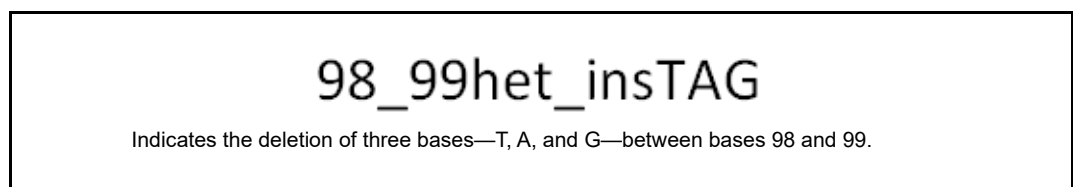


Figure 5-11: Example of a mutation code for a heterozygous duplication event

1316het_dupT

Indicates the insertion of one base - T - between bases 1316 and 1317, where the preceding base (1316) is also T.

Figure 5-12: Example of a mutation code for a heterozygous deletion event

123_127het_delCCTGA

Indicates the heterozygous deletion of five bases—C, C, T, G, and A—between bases 123 through 127.

Mutation report viewing and analysis options

A context menu is available for the Mutation report. You use the options on this menu to change the view of the Mutation report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the Mutation report.

Figure 5-13: Mutation report context menu

Shrink	Ctrl+S
Expand	Ctrl+U
Copy	Ctrl+C
Sort	▶
Confirm	Ctrl+M
Undo Confirmation	Ctrl+N
Delete Lane	
Undo Lane Deletion	
Undo	
View Edit History	
Hyperlink to NCBI's dbSNP	

Option	Description
Shrink	Fit the complete Mutation report in the Viewing and Reporting pane.
Expand	Expand all columns in the Mutation report to their default widths. Note: You might need to Browse the Viewing and Reporting pane to view the entire report.

Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Mutation report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Sort	Click in the column by which you want to sort the report, and then click Sort > Descending or Sort > Ascending.
Confirm	Confirms the selected mutation. The text for the mutation code turns black in both the Mutation report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.
Undo Confirmation	Removes the “Confirmed” status for the mutation.
Delete Lane	Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to “Low Quality.”
Undo Lane Deletion	The Lane Quality for the lane is reset to its original value.
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See “Viewing the Edit History (Audit Trail) for a Mutation” on page 344.
Hyperlink to NCBI's dbSNP	Opens the NCBI's database entry (via the web) for the variant.

You can also view the information for a mutation in the Mutation report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.



If you click a trace in the GAD with the Mutation report open, the GAD is brought to the front. To bring the Mutation report to the front again, press [CTRL]-T.

Two Direction Report

The Two Direction report is automatically generated after you click Reports > Two Direction Report on the Mutation Surveyor main menu. The Two Direction report shows mutation information organized by paired sample traces. The fields that are displayed by default in the report and the color coding of the mutation calls in the report are identical to the default fields and color coding in the Mutation report. (See “Mutation report fields” on page 128 and “Color codes” on page 130.) Also, just like the Mutation report, you can change the fields that are displayed in the report by selecting or clearing the appropriate options on the Output tab on the Mutation Project Options dialog box. (See “Output tab” on page 62.) Figure 5-14 shows the default Two Direction report. The report opens in a window that is separate from the Viewing and Reporting pane.

Figure 5-14: Two Direction report

No.	Sample File	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Qual	Mult	Mutation1	Mutation2	Mutation
1	BRCA1_185729_10/BRCA1_AmpD_R_5.1-R	BRCA1				2	30503	30852	350	26	0		n.a.	n.a.
2	BRCA1_185729_10/BRCA1_AmpD_F_5.1-F	BRCA1				2	30522	30857	336	25	0		n.a.	n.a.
3	BRCA1_185729_10/BRCA1_AmpD_R_5.1-R	BRCA1				2	30503	30852	350	26	0		n.a.	n.a.
4	BRCA1_185729_10/BRCA1_AmpD_F_5.2-F	BRCA1				2	30768	31041	274	6	1	30988A>AG.356	n.a.	n.a.
5	BRCA1_185729_10/BRCA1_AmpD_R_5.2-R	BRCA1				2	30759	31027	268	11	1	30988A>AG.356	n.a.	n.a.
6	BRCA1_185729_10/BRCA1_AmpD_F_5.3-F	BRCA1				2	30975	31414	440	11	0		n.a.	n.a.
7	BRCA1_185729_10/BRCA1_AmpD_R_5.3-R	BRCA1				2	30954	31403	450	9	0		n.a.	n.a.
8	BRCA1_185729_11/BRCA1_AmpD_F_5.4-F	BRCA1				2	31136	31460	323	0	1	31407C>CT.496	n.a.	n.a.
9	BRCA1_185729_11/BRCA1_AmpD_R_5.4-R	BRCA1				2	31114	31451	338	16	1	31407C>CT.496	n.a.	n.a.
10	BRCA1_185729_11/BRCA1_AmpD_F_5.4-F	BRCA1				2	31136	31460	323	0	1	31407C>CT.496	n.a.	n.a.
11	BRCA1_185729_11/BRCA1_AmpD_R_5.4-R	BRCA1				2	31114	31451	338	16	1	31407C>CT.496	n.a.	n.a.
12	BRCA1_185729_12/BRCA1_AmpD_F_5.5-F	BRCA1				2	31417	31595	178	0	0		n.a.	n.a.
13	BRCA1_185729_12/BRCA1_AmpD_R_5.5-R	BRCA1				2	31397	31591	195	23	0		n.a.	n.a.
14	BRCA1_185729_12/BRCA1_AmpD_F_5.5-F	BRCA1				2	31417	31595	178	0	0		n.a.	n.a.
15	BRCA1_185729_12/BRCA1_AmpD_R_5.5-R	BRCA1				2	31397	31591	195	23	0		n.a.	n.a.
16	BRCA1_185729_12/BRCA1_AmpD_F_5.6-F	BRCA1				2	31541	31902	362	11	1	31608C>CT.563	n.a.	n.a.
17	BRCA1_185729_12/BRCA1_AmpD_R_5.6-R	BRCA1				2	31529	31902	374	9	1	31608C>CT.563	n.a.	n.a.
18	BRCA1_185729_12/BRCA1_AmpD_F_5.6-F	BRCA1				2	31541	31902	362	11	1	31608C>CT.563	n.a.	n.a.
19	BRCA1_185729_12/BRCA1_AmpD_R_5.6-R	BRCA1				2	31529	31902	374	9	1	31608C>CT.563	n.a.	n.a.
20	BRCA1_185729_13/BRCA1_AmpD_F_5.7-F	BRCA1				2	31834	32312	479	43	4	31882delA	31988G>GA.693.32003C	31988G>GA.693.32003C
21	BRCA1_185729_13/BRCA1_AmpD_R_5.7-R	BRCA1				2	31820	32306	487	47	3	31882delA	31988G>GA.693.32003C	31988G>GA.693.32003C
22	BRCA1_185729_13/BRCA1_AmpD_F_5.7-F	BRCA1				2	31834	32312	479	43	4	31882delA	31988G>GA.693.32003C	31988G>GA.693.32003C
23	BRCA1_185729_13/BRCA1_AmpD_R_5.7-R	BRCA1				2	31820	32306	487	47	3	31882delA	31988G>GA.693.32003C	31988G>GA.693.32003C
24	BRCA1_185729_14/BRCA1_AmpD_F_5.8-F	BRCA1				2	32270	32711	442	48	1	32533C>CT.871	n.a.	n.a.
25	BRCA1_185729_14/BRCA1_AmpD_R_5.8-R	BRCA1				2	32253	32705	453	48	1	32533C>CT.871	n.a.	n.a.
26	BRCA1_185729_14/BRCA1_AmpD_F_5.8-F	BRCA1				2	32270	32711	442	48	1	32533C>CT.871	n.a.	n.a.
27	BRCA1_185729_14/BRCA1_AmpD_R_5.8-R	BRCA1				2	32253	32705	453	48	1	32533C>CT.871	n.a.	n.a.
28	BRCA1_185729_15/BRCA1_AmpD_F_5.9-F	BRCA1				2	32602	33090	488	41	2	32787T>C.956S.33040G>GA.104	n.a.	n.a.
29	BRCA1_185729_15/BRCA1_AmpD_R_5.9-R	BRCA1				2	32631	33082	452	42	2	32787T>C.956S.33040G>GA.104	n.a.	n.a.

Two Direction report layout

The Two Direction report provides the same general information as the Mutation report with the following exceptions:

- Forward and reverse samples traces are paired next to each other in the report based on the contig in which they were found. Alternating colors of light blue and light tan are used to differentiate the sample pairs.
- Mutations are grouped in the report based on their position in the contig. For a position that shows a mutation call for some of the sample traces:
 - The Mutation # cell is blank for other sample traces that do not show a mutation at this position in the aligned region.






- “n.a.” is displayed in the Mutation # cell for other sample traces that show the mutation at this position, but is filtered due to report settings, or the position is outside the aligned region.
- A row below each group of paired sample files shows the sum of all sizes in the contig, the average lane quality, and the allele frequency (as a percentage).

Two Direction report toolbar

The Two Direction report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 5-15: Two Direction report toolbar



Button	Description
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
	Report Settings—Opens the Two Direction Report Settings dialog box which contains options for filtering the information that is to be displayed in the report. See “Two Direction report settings” on page 136 .
	Advanced Two Direction Report—Opens the Advanced Two Direction Report Settings dialog box which contains options for filtering the information that is to be displayed in the report. It also contains options for specifying the formatting of the mutation code in the report. See “Advanced Two Direction Report” on page 140 .
	Undo Latest Deletions—Restores the deletion that you last carried out in the Two Direction report. Allows for multiple undos.
Lock Sample File in View	Freezes the Sample Filename column while scrolling.

Two Direction report mutation codes

The mutation codes that are used in the Two Direction report are identical to the codes that are used in the Mutation report with one exception—by default, the substitution mutation code also indicates the Dropping Factor and Overlapping Factor for the mutation call.

Figure 5-16: Substitution mutation code in Two Direction report

30988A>AG,356Q>Q/R\$22,0.54,1.00

Dropping factor Overlapping factor

To prevent one or both of these values from being displayed in the mutation code, open the Two Direction Report Settings dialog box (see [“Two Direction report settings” on page 136](#)) and clear the appropriate options.

Two Direction report viewing and analysis options

The Two Direction report has multiple options available for changing the view of the report and for manipulating and analyzing mutation project data. These options are:

- The Two Direction report settings. See [“Two Direction report settings”](#) below.
- The Two Direction report context menu. See [“Two Direction report context menu” on page 137](#).

Two Direction report settings


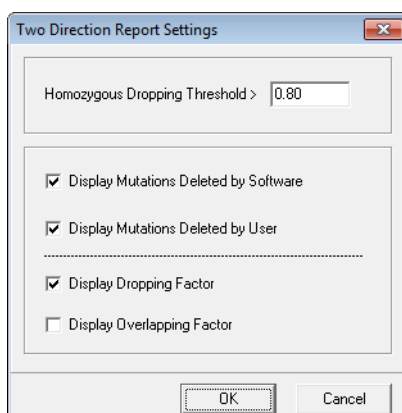
To open the Two Direction Report Settings dialog box, click the Report Settings button  on the Two Direction report toolbar.

Figure 5-17: Two Direction Report Settings dialog box



Option	Description
Homozygous Dropping Threshold	Indicates the value for the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions. Note: You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.
Display the Mutations Deleted by the Software	By default, mutations that are deleted automatically by Mutation Surveyor are displayed in gray text in the Two Direction report. Clear this option to remove these mutation calls from the report display.
Display Mutations Deleted by the User	By default, mutations that are deleted by a user are displayed in gray text in the Two Direction report. Clear this option to remove these mutation calls from the report display.
Display Dropping Factor	By default, the substitution mutation code in the Two Direction report indicates the Dropping factor for the mutation call. Clear this option to remove this value from the mutation code.
Display Overlapping Factor	By default, the substitution mutation code in the Two Direction report indicates the Overlapping factor for the mutation call. Clear this option to remove this value from the mutation code.


Two Direction report context menu

A context menu is available for the Two Direction report. You use the options on this menu to change the view of the Two Direction report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the Two Direction report.

Figure 5-18: Two Direction report context menu

Copy	Ctrl+C
Confirm	▶
Undo Confirmation	▶
Sort By Score(Descending)	
Sort By Score(Ascending)	
Delete Mutations	▶
Delete Lane	
Undo Deletion	
Undo	
View Edit History	
Hyperlink to NCBI's dbSNP	

Chapter 5
Mutation Surveyor Standard Reports

Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Two Direction report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the Two Direction report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Undo Confirmation	<p>Removes the “Confirmed” status for the mutation.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p>
Sort By Score (Descending) Sort by Score (Ascending)	<p>Click the appropriate option to sort the all the data by descending mutation score or ascending mutation score.</p>
Delete Mutations	<p>Deletes the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations, to select all mutations in a column, or to select multiple rows.</p> <p>The text for the mutation code turns gray in both the Two Direction report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p> <p>Note: To undo the last deletion that you carried out, click the Undo Latest Deletions button  on the report toolbar.</p>
Delete Lane	<p>Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to “Low Quality.”</p>
Undo Lane Deletion	<p>The Lane Quality for the lane is reset to its original value.</p>
Undo	<p>Undo the <i>last</i> edit action that was carried out for the mutation.</p>
View Edit History	<p>See “Viewing the Edit History (Audit Trail) for a Mutation” on page 344.</p>
Hyperlink to NCBI's dbSNP	<p>Opens the NCBI's database entry (via the web) for the variant.</p>

You can also view the information for a mutation in the Two Direction report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.




If you click a trace in the GAD with the Two Direction report open, the GAD is brought to the front. To bring the Two Direction report to the front again, press [CTRL]-T.

Advanced Two Direction Report


Like the Two Direction report, the Advanced Two Direction report shows mutation information organized by paired sample traces. Also, the layout of the Advanced Two Direction report is identical to the layout of the Two Direction report. (See [“Two Direction report layout” on page 134.](#)) The Advanced Two Direction report, however, has different settings for changing the mutation information that is displayed in the report.

The fields that are displayed by in the report, the mutation codes, and the color coding of the mutation calls in the report are identical to the fields, mutation codes, and color coding in the Mutation report (See [“Mutation report fields” on page 128](#), [“Mutation report codes” on page 129](#), and [“Color codes” on page 130.](#)) Just like the Mutation report and the Two Direction report, you can change the fields that are displayed in the report by selecting or clearing the appropriate options on the Output tab on the Mutation Project Options dialog box. (See [“Output tab” on page 62.](#))

1. To open the Advanced Two Direction report, do one of the following:
 - Click the Advanced Two Direction Report button  on the Two Direction report toolbar.
 - On the Mutation Surveyor main menu, click Reports > Adv. Two Direction Report.

The Advanced Two Direction Report Settings dialog box opens.

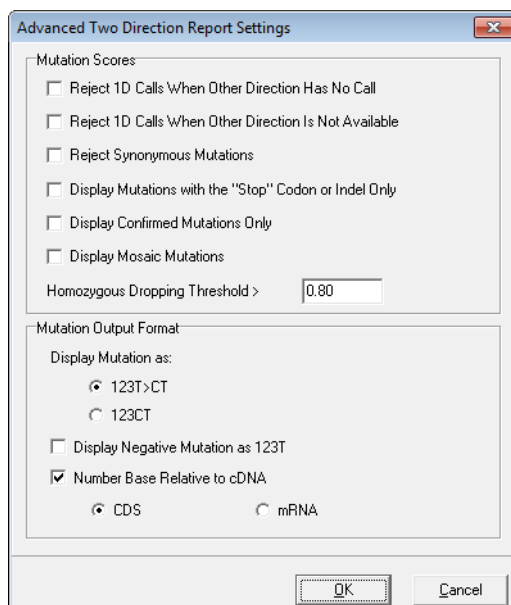Figure 5-19: Advanced Two Direction Report Settings dialog box



2. Accept the default settings for the report, or make any changes as needed.

Option	Description
Mutation Scores	
Reject 1D Calls When Other Direction has No Call	Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction.
Reject 1D Calls When Other Direction is Not Available	Reject mutation calls that occur outside the region where the sample traces in both directions are aligned within the comparison region.
Reject Synonymous Mutations	Coding region substitutions that do not affect amino acid calls are not displayed in the report.
Note: In cases where the entire column was not omitted from the report because of rejection by one of these criteria, the rejected mutations have no background color and an "n.a." is displayed in the Mutation cell instead of the mutation code.	
Display Mutations with the "Stop" Codon or Indel Only	Display only nonsense mutations and indels in the report.
Display Confirmed Mutations Only	Display only those mutations that you have manually confirmed in the report.
Display Mosaic Mutations	Display mosaic mutations in the report.
Homozygous Dropping Threshold	<p>This value is the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions.</p> <p>Note: You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.</p>
Mutation Output Format	
Display Mutation As	<p>Select one option to set the format of the mutation code in the report.</p> <ul style="list-style-type: none"> 123T>CT (Default value) 123CT
Display Negative Mutation as 123T	Display the position and the normal allele for a negative mutation in the report.
Number Bases Relative to cDNA	<ul style="list-style-type: none"> If this option is cleared, the numbering is the genomic position within the GenBank file. If this option is selected: <ul style="list-style-type: none"> Select CDS to number the bases relative to the CDS numbering in the GenBank reference file. Select mRNA to number the bases relative the mRNA numbering in the GenBank reference file.

3. Click OK.

The Advanced Two Direction report is generated and displayed in a window that is separate from the Viewing and Reporting pane.

Figure 5-20: Advanced Two Direction report




No.	Sample File	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Quali	Mut#	Mutation1	Mutation2	Muta
1	BRCA1_012345_10	BRCA1_AmplD_F_S	1-F	BRCA1		2	30524	30857	334	16	0			n.a.
2	BRCA1_012345_10	BRCA1_AmplD_R_S1-R	BRCA1			2	30503	30851	349	14	1	c.686C>CT.p.S2		n.a.
3	BRCA1_012345_10	BRCA1_AmplD_F_S1-F	BRCA1			2	30524	30857	334	16	0			n.a.
4	BRCA1_012345_10	BRCA1_AmplD_R_S1-R	BRCA1			2	30503	30851	349	14	1	c.686C>CT.p.S2		n.a.
5	BRCA1_185729_10	BRCA1_AmplD_F_S1-F	BRCA1			2	30522	30857	336	25	0			n.a.
6	BRCA1_185729_10	BRCA1_AmplD_R_S1-R	BRCA1			2	30503	30852	350	26	0			n.a.
7	BRCA1_185729_10	BRCA1_AmplD_F_S1-F	BRCA1			2	30522	30857	336	25	0			n.a.
8	BRCA1_185729_10	BRCA1_AmplD_R_S1-R	BRCA1			2	30503	30852	350	26	0			n.a.
9	BRCA1_264381_10	BRCA1_AmplD_F_S1-F	BRCA1			2	30512	30857	346	25	0			n.a.
10	BRCA1_264381_10	BRCA1_AmplD_R_S1-R	BRCA1			2	30503	30852	350	25	0			n.a.
11	BRCA1_264381_10	BRCA1_AmplD_F_S1-F	BRCA1			2	30512	30857	346	25	0			n.a.
12	BRCA1_264381_10	BRCA1_AmplD_R_S1-R	BRCA1			2	30503	30852	350	25	0			n.a.
12									4130	22	0.17	33.3%		
13	BRCA1_012345_10	BRCA1_AmplD_F_S2-F	BRCA1			2	30768	31039	272	8	0	n.a.	n.a.	
14	BRCA1_012345_10	BRCA1_AmplD_R_S2-R	BRCA1			2	30746	31027	282	11	2	c.826A>AC.p.T2 c.832A>AT.p.T2		
15	BRCA1_012345_10	BRCA1_AmplD_F_S2-F	BRCA1			2	30768	31039	272	8	0	n.a.	n.a.	
16	BRCA1_012345_10	BRCA1_AmplD_R_S2-R	BRCA1			2	30746	31027	282	11	2	c.826A>AC.p.T2 c.832A>AT.p.T2		
17	BRCA1_185729_10	BRCA1_AmplD_F_S2-F	BRCA1			2	30768	31041	274	6	1	n.a.	n.a.	c.10f
18	BRCA1_185729_10	BRCA1_AmplD_R_S2-R	BRCA1			2	30759	31027	269	11	1	n.a.	n.a.	c.10f
19	BRCA1_185729_10	BRCA1_AmplD_F_S2-F	BRCA1			2	30768	31041	274	6	1	n.a.	n.a.	c.10f
20	BRCA1_185729_10	BRCA1_AmplD_R_S2-R	BRCA1			2	30759	31027	269	11	1	n.a.	n.a.	c.10f


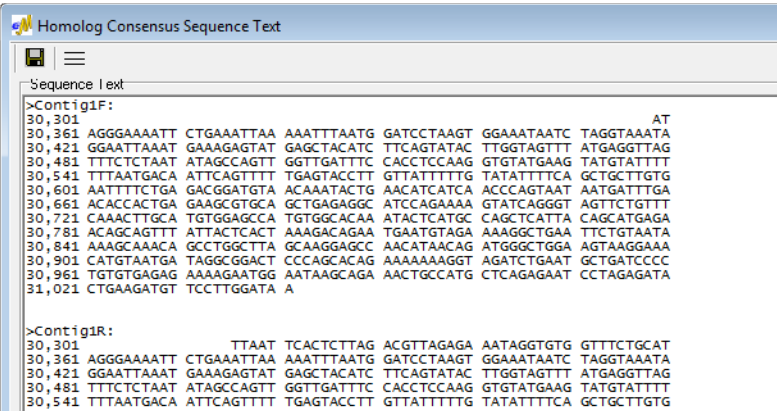

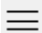
Advanced Two Direction report toolbar

The Advanced Two Direction report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 5-21: Advanced Two Direction report toolbar




Button	Description
	Copy—Copies the selected the text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
	Report Settings—Opens the settings dialog box to allow adjustment of the Homozygous Dropping Factor.

Button	Description
	<p>Homolog Consensus Sequence Text—Generates an onscreen text file of the consensus base calls for the forward samples in a contig, for the reverse samples in a contig, and for the forward and reverse samples together.</p> <p><i>Figure 5-22: Homolog Consensus Sequence Text window</i></p>  <p>The screenshot shows a window titled "Homolog Consensus Sequence Text". Inside, there are two sections: ">Contig1F:" and ">Contig1R:". Each section displays a list of sequence reads with their positions and base calls. For example, under ">Contig1F:", the first line is "30,301 AGGGAAAATT CTGAAATTAA AAATTTAATG GATCCTAAGT GGAAATAATC TAGGTAATAA". The second section, ">Contig1R:", shows reverse sequence reads.</p> <ul style="list-style-type: none"> • Save button  —Opens the Save As dialog box in which you can specify a name and location for saving the sequence text as a .txt file. • Display Sequence in One Line button  —Displays the consensus sequences in fasta format.
Lock Sample File in View	Freezes the Sample Filename column while scrolling.

Advanced Two Direction report viewing and analysis options

A context menu is available for the Advanced Two Direction report. You use the options on this menu to change the view of the Advanced Two Direction report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the Advanced Two Direction report.

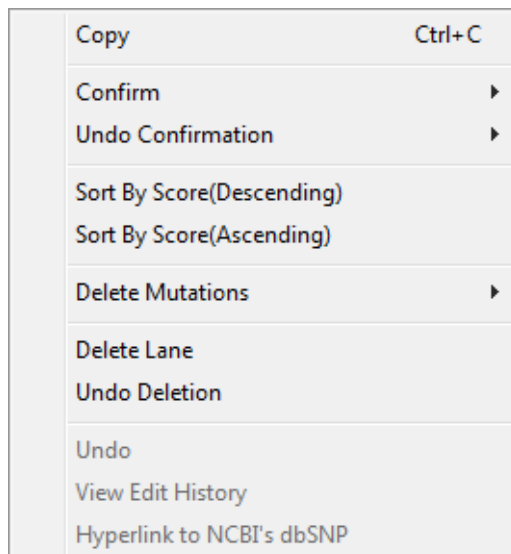Figure 5-23: Advanced Two Direction report context menu



Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Advanced Two Direction report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the Advanced Two Direction report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Undo Confirmation	<p>Removes the "Confirmed" status for the mutation.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p>

Option	Description
Sort By Score (Descending) Sort by Score (Ascending)	Click the appropriate option to sort the all the data by descending mutation score or ascending mutation score.
Delete Mutations	Deletes the selected mutations. Note: SHIFT-click to select multiple contiguous mutations, to select all mutations in a column, or to select multiple rows. The text for the mutation code turns gray in both the Advanced Two Direction report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.
Delete Lane	Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to "Low Quality."
Undo Deletion	The Lane Quality for the lane is reset to its original value.
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See "Viewing the Edit History (Audit Trail) for a Mutation" on page 344.
Hyperlink to NCBI's dbSNP	Opens the NCBI's database entry (via the web) for the variant.

You can also view the information for a mutation in the Advanced Two Direction report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.



If you click a trace in the GAD with the Advanced Two Direction report open, the GAD is brought to the front. To bring the Advanced Two Direction report to the front again, press [CTRL]-T.


Project Report

Like the directional reports, the Project report shows mutation information organized by paired sample traces. Although the layout of the Project report is very similar to the layout of the directional reports, the report has different settings for working with the report and for changing the information that is displayed in the report, and reports samples aligning to a single GenBank file. The fields that are displayed in the report, the mutation codes, and the color coding of the mutation calls in the report are identical to the fields, mutation codes, and color coding in the Mutation report (See [“Mutation report fields” on page 128](#), [“Mutation report codes” on page 129](#), and [“Color codes” on page 130](#).) Just like the other standard mutation reports, you can change the fields that are displayed in the report by selecting or clearing the appropriate options on the Output tab on the Mutation Project Options dialog box. (See [“Output tab” on page 62](#).)

1. To open the Project report, on the Mutation Surveyor main menu, click Reports > Project report.

The Project Report Settings dialog box opens.

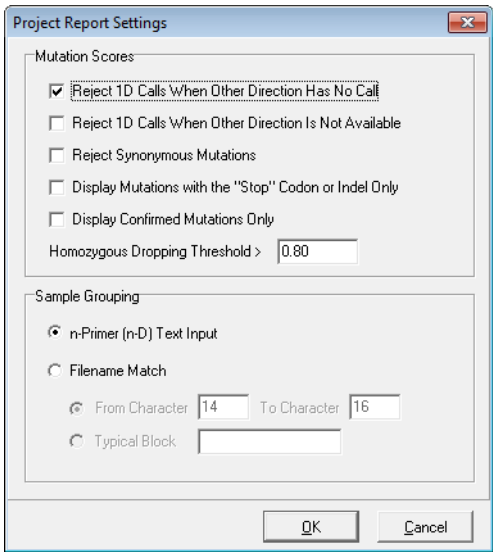Figure 5-24: Project Report Settings dialog box



2. Accept the default settings for the report, or make any changes as needed.


Option	Description
Mutation Scores	
Reject 1D Calls When Other Direction has No Call	Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction.
Reject 1D Calls When Other Direction is Not Available	Reject mutation calls that occur outside the region where the sample traces in both directions are aligned within the comparison region.

Option	Description
Reject Synonymous Mutations	Coding region substitutions that do not affect amino acid calls are not displayed in the report.
Note: In cases where the entire column was not omitted from the report because of rejection by one of these criteria, the rejected mutations have no background color and an “n.a.” is displayed in the Mutation cell instead of the mutation code.	
Display Mutations with the “Stop” Codon or Indel Only	Display only nonsense mutations and indels in the report.
Display Confirmed Mutations Only	Display only those mutations that you have manually confirmed in the report.
Display Mosaic Mutations	Display mosaic mutations in the report. This option requires using Check 2D Small Peaks in the Mutation Project Settings “Display tab” on page 64 .
Homozygous Dropping Threshold	This value is the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions. You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.
Sample Grouping	
	<p>Specify the filename substring by which Mutation Surveyor groups all the mutation project samples files in the report. (These groupings are also used in the Project Sample Assembly subreport and in the Comparison functions. See “Project report toolbar” on page 148.)</p> <ul style="list-style-type: none"> • n-Primer (n-D) Text Input—Use the ND Grouping file to group the sample files. See “ND Filename Match Editor Tool” on page 264. • Filename Match (based on one of the following): <ul style="list-style-type: none"> • From Character [] To Character []—If all the filenames in your project contain the same characters/designator in the same positions, then select this option and enter the starting and ending character positions. <p>Note: Typical Block—If all the filenames in your project contain the same characters/designator but in different positions, then select this option and enter the value for the character/designator block.</p>

3. Click OK.

The Project report is generated and displayed in a window that is separate from the Viewing and Reporting pane.

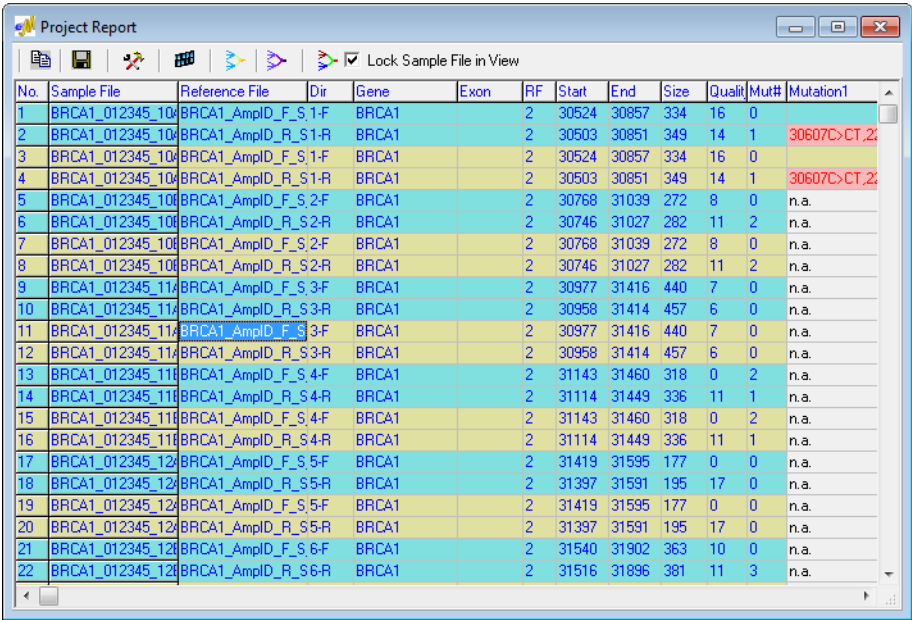Figure 5-25: Project report










Project report toolbar

The Project report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report. It also contains buttons for generating subreports in which you can view mutations in a matrix and view the alleles for each mutation.

Figure 5-26: Project report toolbar



Button	Description
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).

Button	Description
	Report Settings—Opens the Project Report Settings dialog box. See Step 2 of “Project Report” on page 146 .
	Project Sample Assembly—Opens the Project Sample Assembly subreport. The Project Sample Assembly subreport creates a consensus of all the samples traces as defined by the Sample Grouping parameters set in the Project Report Settings dialog box. This subreport also lists the alleles for each mutation. See “Project Sample Assembly subreport” on page 150 .
	Consensus Sequence Text—Opens the Consensus Sequence Text dialog box. The Consensus Sequence Text function searches sample files that have been grouped based on a specific filename for consensus sequence strings. See “Consensus Sequence Text function” on page 151 .
	Sequence Text Comparison—Opens the Sequence Text Comparison dialog box. The Sequence Text Comparison function compares the sequence text for sample files that have been grouped based on a specific filename substring to the sequence text in the reference file. See “Sequence Text Comparison function” on page 152 .
	Amino Acid Text Comparison—Opens the Amino Acid Text Comparison dialog box. The Amino Acid Text Comparison function compares the amino acid sequence for sample files that have been grouped based on a specific filename substring to the amino acid sequence in the reference file. See “Amino Acid Text Comparison function” on page 154 .
Lock Sample File in View	Freezes the Sample Filename column while scrolling.

Project Sample Assembly subreport

The Project Sample Assembly subreport displays a separate row for each substring, the total number of mutations found in all the files that contain the specified substring, and the code for each called mutation. The subreport can be saved as a text file.

Figure 5-27: Project Sample Assembly subreport

No.	Sample	Mut#	Mutation1(33.3%)	Mutation2(25.0%)	Mutation3(25.0%)	Mutation4(33.3%)	Mutation5(16.7%)	Mutation6(33.3%)	Mutation7
1	BRCA1_012345_1041	30607C>CT,229S>S				n.a.	n.a.	n.a.	n.a.
2	BRCA1_012345_1041	30607C>CT,229S>S				n.a.	n.a.	n.a.	n.a.
3	BRCA1_012345_10E2	n.a.	30747A>AC,276T>T	30753A>AT,278T>T			n.a.	n.a.	n.a.
4	BRCA1_012345_10E2	n.a.	30747A>AC,276T>T	30753A>AT,278T>T			n.a.	n.a.	n.a.
5	BRCA1_012345_1140	n.a.	n.a.	n.a.	n.a.				
6	BRCA1_012345_1140	n.a.	n.a.	n.a.	n.a.				
7	BRCA1_012345_11E3	n.a.	n.a.	n.a.	n.a.	n.a.	31389C>CG,490P>P	31407C>CT,496R>R	31407C>C
8	BRCA1_012345_11E3	n.a.	n.a.	n.a.	n.a.	n.a.	31389C>CG,490P>P	31407C>CT,496R>R	31407C>C
9	BRCA1_012345_1240	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
10	BRCA1_012345_1240	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.

- To view the group consensus results within a matrix, click Display Results on the subreport toolbar.

Figure 5-28: Results matrix view of mutations

No.	Mutations	BRCA1_012345_104	BRCA1_012345_10E	BRCA1_012345_10E	BRCA1_012345_10E	BRCA1_012345_114	BRCA1_012345_11E	BRCA1_012345_124
1	30607C>CT,p.S229S	YES?	YES?	n.a.	n.a.	n.a.	n.a.	n.a.
2	30747A>AC,p.T276T	NO	NO	YES	YES	n.a.	n.a.	n.a.
3	30753A>AT,p.T278T	NO	NO	YES	YES	n.a.	n.a.	n.a.
4	30988A>AG,p.Q356L	n.a.	n.a.	NO	NO	NO	NO	n.a.
5	31389C>CG,p.P490P	n.a.	n.a.	n.a.	n.a.	NO	NO	YES?
6	31407C>CT,p.R496R	n.a.	n.a.	n.a.	n.a.	NO	NO	YES?
7	31407C>CG,p.R496R	n.a.	n.a.	n.a.	n.a.	NO	NO	YES?
8	31518A>AC,p.N533I	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
9	31526A>AG,p.G535V	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
10	31608C>CT,p.Q563L	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.

- To view a matrix of the alleles for each group consensus for each mutation, click Display Alleles on the subreport toolbar.

Figure 5-29: Allele matrix view of mutations

No.	Mutations	BRCA1_012345_104	BRCA1_012345_10E	BRCA1_012345_10E	BRCA1_012345_10E	BRCA1_012345_114	BRCA1_012345_11E	BRCA1_012345_124
1	30607C>CT,p.S229S	YES?	YES?	n.a.	n.a.	n.a.	n.a.	n.a.
2	30747A>AC,p.T276T	AA	AA	AC	AC	n.a.	n.a.	n.a.
3	30753A>AT,p.T278T	AA	AA	AT	AT	n.a.	n.a.	n.a.
4	30988A>AG,p.Q356L	n.a.	n.a.	AA	AA	AA	AA	n.a.
5	31389C>CG,p.P490P	n.a.	n.a.	n.a.	n.a.	CC	CC	YES?
6	31407C>CT,p.R496R	n.a.	n.a.	n.a.	n.a.	CC	CC	YES?
7	31407C>CG,p.R496R	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	YES?
8	31518A>AC,p.N533I	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
9	31526A>AG,p.G535V	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
10	31608C>CT,p.Q563L	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.

Consensus Sequence Text function

The Consensus Sequence Text function searches sample files that have been grouped based on a specific filename for consensus sequence strings. You can also specify a minimum lane quality to exclude low quality samples

Figure 5-30: Consensus Sequence Text Settings dialog box


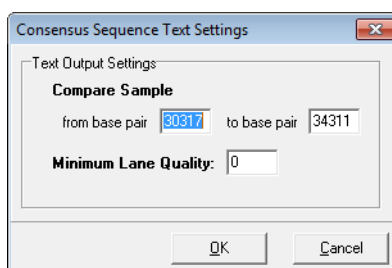



Figure 5-31 is an example of the Consensus Sequence Text report. The reference information is displayed at the top of the report. The consensus sequence for the base pair region is displayed below the reference information and is grouped by filename substring.

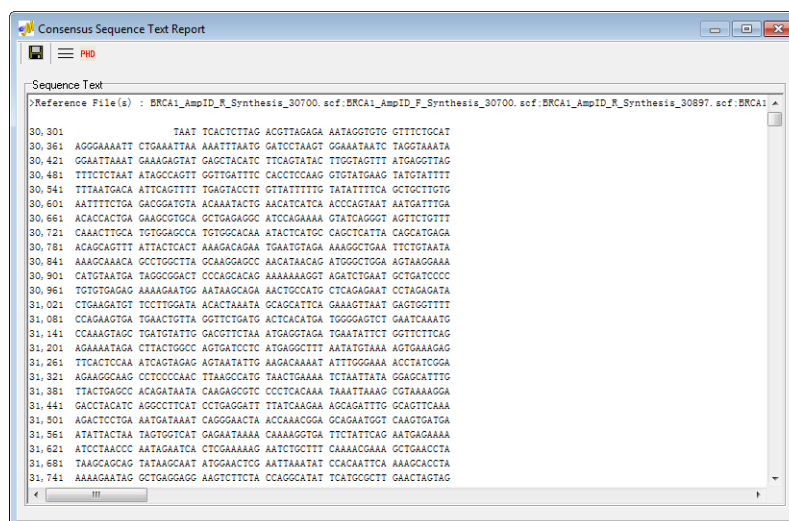

Figure 5-31: Consensus Sequence Text report



Consensus Sequence Text window toolbar

The Consensus Sequence Text window toolbar contains buttons that are specific for working with the output and for changing the output display.


Button	Description
	Save Sequence Text—Opens the Save As dialog box in which you can specify a name and location for saving the output as a .txt file.

Button	Description
	Display Sequence in One Line—Displays the consensus sequences in fasta format.
	Save As PHD file—Opens the Browse for Folder dialog box in which you can specify the location for saving the output as a .phd file. A .phd file contains three columns: <ul style="list-style-type: none"> • Column 1-Base call • Column 2-Quality score for the called base • Column 3-Raw data frame for called base You can open a .phd file in Notepad.

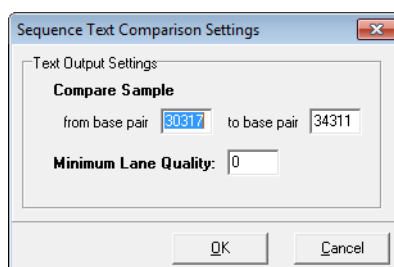Sequence Text Comparison function

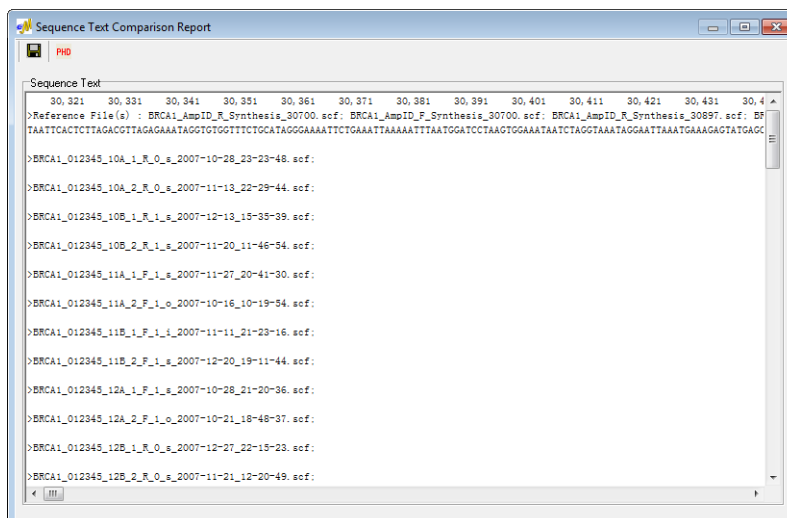
The Sequence Text Comparison function compares the sequence text for sample files that have been grouped based on a specific filename substring to the sequence text in the reference file. You can also specify a minimum lane quality to exclude low quality samples.

Figure 5-32: Sequence Text Comparison Settings dialog box





[Figure 5-33 on page 153](#) is an example of the Sequence Text Comparison report. The reference sequence is displayed at the top of the report as a single continuous string of characters without numbering or spacing. Sample files are grouped by substring and are displayed below the reference sequence. In a sample file sequence, a black dot (.) indicates a match between the reference file and the sample file. A black dash (-) indicates an artificial gap added to a sample file so that all the sample files remain aligned. This gap can be the result of a deletion in the gapped sample or of an insertion in another sample. If a different base is at the same position in the sample file and reference file, then the standard IUB/IUPAC code for the base is shown in the sample file at the position.

Figure 5-33: Sequence Text Comparison report



Sequence Text Comparison window toolbar

The Sequence Text window toolbar contains buttons that are specific for working with the output.

Button	Description
	Save Sequence Text—Opens the Save As dialog box in which you can specify a name and location for saving the output as a .txt file.
	Save As PHD file—Opens the Browse for Folder dialog box in which you can specify the location for saving the output as a .phd file. A .phd file contains three columns: <ul style="list-style-type: none"> • Column 1-Base call • Column 2-Quality score for the called base • Column 3-Raw data frame for called base You can open a PHD file in Notepad.

Amino Acid Text Comparison function

The Amino Acid Text Comparison function compares the amino acid sequence for sample files that have been grouped based on a specific filename substring to the amino acid sequence in the reference file. You can also specify a minimum lane quality to exclude low quality samples.

Figure 5-34: Amino Acid Text Comparison Settings dialog box


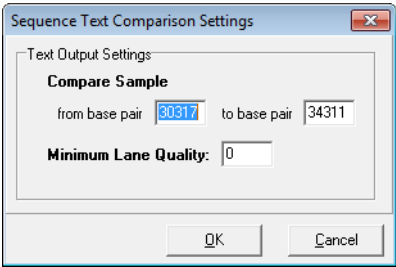



Figure 5-35 is an example of the Amino Acid Text Comparison report. The amino acid sequence for the reference file displayed at the top of the report using standard IUB/IUPAC amino acid codes. Sample files are grouped by substring and are displayed below the reference sequence. In a sample file sequence, a black dot (.) indicates a match between the reference file and the sample file. If a different amino acid is at the same position in the sample file and reference file, then the standard IUB/IUPAC code for the amino acid is shown in the sample file at the position.

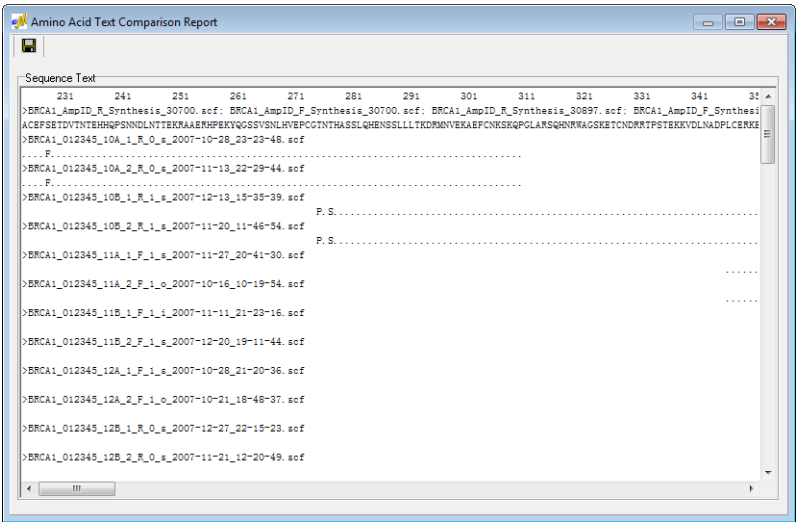Figure 5-35: Amino Acid Text Comparison report



Amino Acid Text Comparison window toolbar

The Amino Acid Text Comparison window toolbar contains a single button that is specific for working with the output.

Button	Description
	Save Sequence Text—Opens the Save As dialog box in which you can specify a name and location for saving the output as a .txt file.

Project report viewing and analysis options

A context menu is available for the Project report. You use the options on this menu to change the view of the Project report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the Project report.

Figure 5-36: Project report context menu

Copy	Ctrl+C
Confirm	Ctrl+M
Undo Confirmation	Ctrl+N
Sort By Score (Descending)	
Sort By Score (Ascending)	

Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Project report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the Project report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>

Chapter 5

Mutation Surveyor Standard Reports

Option	Description
Undo Confirmation	Removes the “Confirmed” status for the mutation. Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.
Sort By Score (Descending) Sort by Score (Ascending)	Click the appropriate option to sort the all the data by descending mutation score or ascending mutation score.

You can also view the information for a mutation in the Project report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.



If you click a trace in the GAD with the Project report open, the GAD is brought to the front. To bring the Project report to the front again, press [CTRL]-T.

HGVS Report

Like the directional reports, the HGVS report shows mutation information organized by paired sample traces. Also, the layout of the HGVS report is very similar to the layout of the directional reports. (See [“Two Direction report layout” on page 134.](#)) The HGVS report, however, displays mutation calls according to the guidelines established by the Human Genome Variations Society. (The color coding of these calls is still the same as the color coding that is used in the other standard Mutation Surveyor reports. See [“Color codes” on page 130.](#)) In addition to the standard report mutation fields specified for the project (see [“Output tab” on page 62](#)), the HGVS report shows many quality values for the project, including quality values for the entire trace, the comparison region, and the ROIs.

1. To open the HGVS report, on the Mutation Surveyor main menu, click Reports > HGVS report.

The HGVS Report Settings dialog box opens.

Figure 5-37: HGVS Report Settings dialog box

HGVS Report Settings

Mutation Scores

- ☐ Reject 1D Calls When Other Direction Has No Call
- ☐ Reject 1D Calls When Other Direction Is Not Available
- ☐ Reject Synonymous Mutations
- ☐ Display Mutations with the "Stop" Codon or Indel Only
- ☐ Display Confirmed Mutations Only
- ☐ Display Mosaic Mutations

Homozygous Dropping Threshold >

Phred Score Threshold

Trace Quality

- ☒ Entire Trace ☐ Comparison Region ☐ ROI (CDS)
- ☒ Phred Score ☐ S/N Ratio

Quality Threshold

Low Quality Threshold is: 0

Amino Acid

- ☒ 1 Letter ☐ 3 Letters

Display Start and End

- ☒ Genomic ☐ HGVS Nomenclature

2. Accept the default settings for the report, or make any changes as needed.


Option	Description
Mutation Scores	
Reject 1D Calls When Other Direction has No Call	Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction.
Reject 1D Calls When Other Direction is Not Available	Reject mutation calls that occur outside the region where the sample traces in both directions are aligned within the comparison region.
Reject Synonymous Mutations	Coding region substitutions that do not affect amino acid calls are not displayed in the report.
Note: In cases where the entire column was not omitted from the report because of rejection by one of these criteria, the rejected mutations do not have a background color and an "n.a." is displayed in the Mutation cell instead of the mutation code.	
Display Mutations with the "Stop" Codon or Indel Only	Display only nonsense mutations and indels in the report.
Display Confirmed Mutations Only	Display only those mutations that you have manually confirmed in the report.
Display Mosaic Mutations	Display mosaic mutations in the report. This option requires using Check 2D Small Peaks in the Mutation Project Settings. See "Display tab" on page 64 .
Homozygous Dropping Threshold	<p>Indicates the value for the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions. You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.</p> <p>Note: You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.</p>
Phred Score Threshold	A measure of base quality, all bases that are in the ROI and that have a Phred score <= to the specified value are displayed in the Bases Below Threshold (ROI) report column. Reducing this value reduces the number of bases that are displayed. Conversely, increasing this value increased the number of based that are displayed.
Trace Quality	

Option	Description
	Display the trace quality for one or more or all the following: <ul style="list-style-type: none"> Entire Trace Comparison Region ROI (CDS) and set the trace quality relative to <i>one</i> of the following: <ul style="list-style-type: none"> Phred Score S/N Ratio
Quality Threshold	A filtering option to visually indicate all the quality values that fall below this threshold. Quality values that fall below this threshold are highlighted in red in the report.
Low Quality Threshold is: []	Any lanes with quality that is less than this threshold are not displayed in the report. Note: The value shown is the value that you specified for the Quality Threshold on the Output tab of the Mutation Project Settings dialog box. To change this value, see "Output tab" on page 62 .
Display Start and End	
	Display start and end base numbering as <i>one</i> of the following: <ul style="list-style-type: none"> Genomic HGVS nomenclature
Amino Acid	
<ul style="list-style-type: none"> 1 Letter 3 Letter 	<ul style="list-style-type: none"> Display 1 Letter IUPAC amino acid code in the report. Display 3 Letter IUPAC amino acid code in the report.

3. Click OK.

The HGVS report is generated and displayed in a window that is separate from the Viewing and Reporting pane.

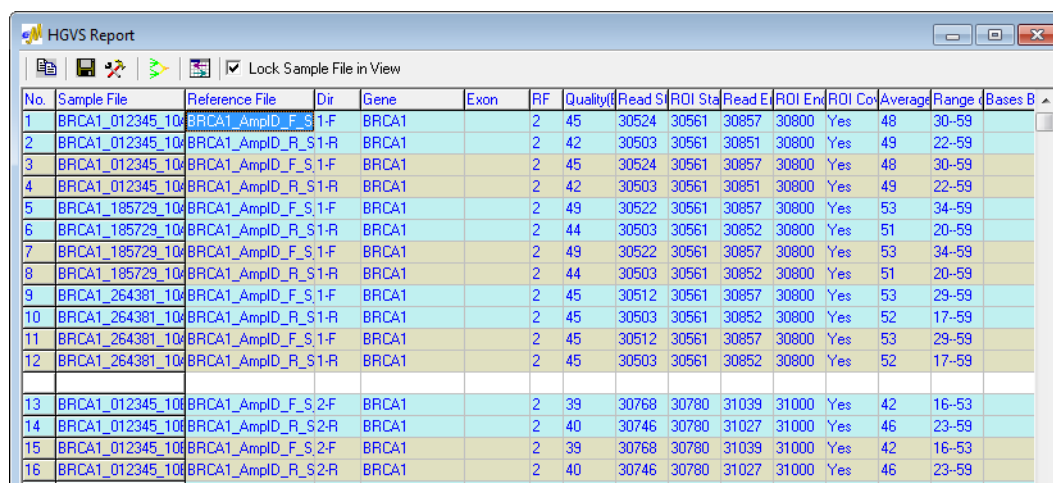Figure 5-38: HGVS report

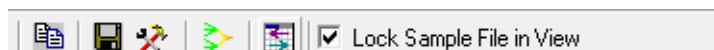






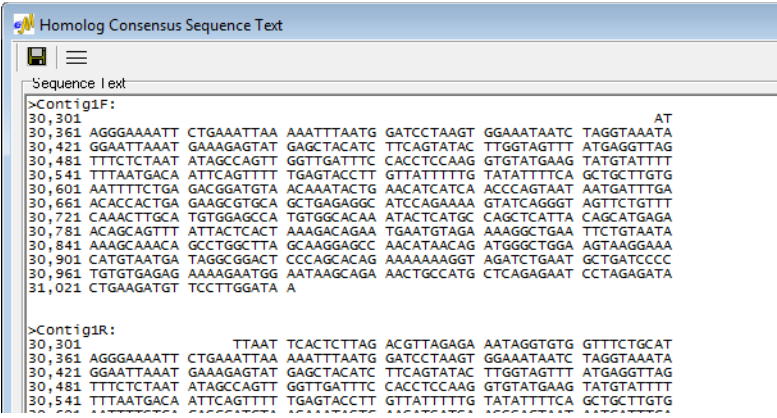

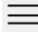
No.	Sample File	Reference File	Dir	Gene	Exon	RF	Quality	Read S	ROI Sta	Read E	ROI End	ROI Co	Average	Range	Bases B
1	BRCA1_012345_10	BRCA1_AmplD_F_S1-F	1-F	BRCA1		2	45	30524	30561	30857	30800	Yes	48	30-59	
2	BRCA1_012345_10	BRCA1_AmplD_R_S1-R	1-R	BRCA1		2	42	30503	30561	30851	30800	Yes	49	22-59	
3	BRCA1_012345_10	BRCA1_AmplD_F_S1-F	1-F	BRCA1		2	45	30524	30561	30857	30800	Yes	48	30-59	
4	BRCA1_012345_10	BRCA1_AmplD_R_S1-R	1-R	BRCA1		2	42	30503	30561	30851	30800	Yes	49	22-59	
5	BRCA1_185729_10	BRCA1_AmplD_F_S1-F	1-F	BRCA1		2	49	30522	30561	30857	30800	Yes	53	34-59	
6	BRCA1_185729_10	BRCA1_AmplD_R_S1-R	1-R	BRCA1		2	44	30503	30561	30852	30800	Yes	51	20-59	
7	BRCA1_185729_10	BRCA1_AmplD_F_S1-F	1-F	BRCA1		2	49	30522	30561	30857	30800	Yes	53	34-59	
8	BRCA1_185729_10	BRCA1_AmplD_R_S1-R	1-R	BRCA1		2	44	30503	30561	30852	30800	Yes	51	20-59	
9	BRCA1_264381_10	BRCA1_AmplD_F_S1-F	1-F	BRCA1		2	45	30512	30561	30857	30800	Yes	53	29-59	
10	BRCA1_264381_10	BRCA1_AmplD_R_S1-R	1-R	BRCA1		2	45	30503	30561	30852	30800	Yes	52	17-59	
11	BRCA1_264381_10	BRCA1_AmplD_F_S1-F	1-F	BRCA1		2	45	30512	30561	30857	30800	Yes	53	29-59	
12	BRCA1_264381_10	BRCA1_AmplD_R_S1-R	1-R	BRCA1		2	45	30503	30561	30852	30800	Yes	52	17-59	
13	BRCA1_012345_10	BRCA1_AmplD_F_S2-F	2-F	BRCA1		2	39	30768	30780	31039	31000	Yes	42	16-53	
14	BRCA1_012345_10	BRCA1_AmplD_R_S2-R	2-R	BRCA1		2	40	30746	30780	31027	31000	Yes	46	23-59	
15	BRCA1_012345_10	BRCA1_AmplD_F_S2-F	2-F	BRCA1		2	39	30768	30780	31039	31000	Yes	42	16-53	
16	BRCA1_012345_10	BRCA1_AmplD_R_S2-R	2-R	BRCA1		2	40	30746	30780	31027	31000	Yes	46	23-59	


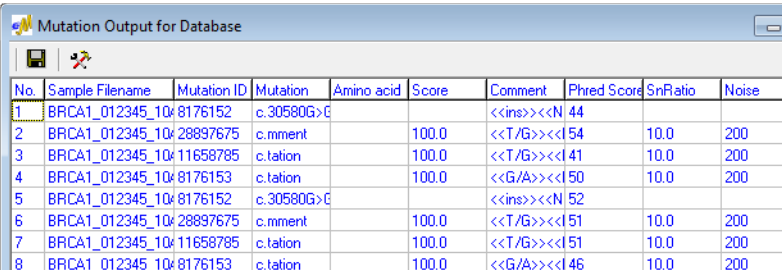


HGVS report toolbar

The HGVS report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 5-39: HGVS report toolbar



Button	Description
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
	Report Settings—Opens the HGVS Report Settings dialog box. See Step 2 of "HGVS Report" on page 157 .
	<p>Homolog Consensus Sequence Text—Generates an onscreen text file of the consensus base calls for the forward samples in a contig, for the reverse samples in a contig, and for the forward and reverse samples together.</p> <p>Figure 5-40: Homolog Consensus Sequence Text window</p>  <p>The screenshot shows a window titled "Homolog Consensus Sequence Text". It contains two sections: ">Contig1F:" and ">Contig1R:". Each section displays a list of sequence positions (30,301 to 31,021) and the corresponding consensus base calls for multiple samples. The sequences are aligned, showing variations between samples.</p> <ul style="list-style-type: none"> • Save button  —Opens the Save As dialog box in which you can specify a name and location for saving the .txt file. • Display Sequence in One Line button  —Displays the consensus sequences in fasta format.

Button	Description
	<p>Mutation Output for Database—Outputs the HGVS report to a format that can be exported to a database.</p> <p><i>Figure 5-41: Mutation Output for Database</i></p>  <p>• Save button  —Opens the Save As dialog box in which you can specify a name and location for saving the mutation output as a .txt file.</p> <p>• Report Settings button  —Opens the Mutation Output dialog box. You use the options on this dialog box to select the fields that are to be included in the Mutation output. By default, when the dialog box first opens, all standard mutation report fields are selected.</p>
Lock Sample File in View	Freezes the Sample Filename column while scrolling.

HGVS report viewing and analysis options

A context menu is available for the HGVS report. You use the options on this menu to change the view of the HGVS report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the HGVS report.

Figure 5-42: HGVS report context menu

Copy	Ctrl+C
Confirm	▶
Undo Confirmation	▶
Sort By Score (Descending)	
Sort By Score (Ascending)	
Delete Mutations	▶
Delete Lane	
Undo Deletion	
Undo	
View Edit History	

Option	Description
Copy	<p>Copies the selected the text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the HGVS report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the HGVS report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Undo Confirmation	<p>Removes the “Confirmed” status for the mutation.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p>
Sort By Score (Descending) Sort by Score (Ascending)	Click the appropriate option to sort the all the data by descending mutation score or ascending mutation score.
Delete Mutations	<p>Deletes the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations, to select all mutations in a column, or to select multiple rows.</p> <p>The text for the mutation code turns gray in both the HGVS report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Delete Lane	Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to “Low Quality.”
Undo Deletion	The Lane Quality for the lane is reset to its original value.
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See “Viewing the Edit History (Audit Trail) for a Mutation” on page 344.

You can also view the information for a mutation in the HGVS report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.



If you click a trace in the GAD with the HGVS report open, the GAD is brought to the front. To bring the HGVS report to the front again, press [CTRL]-T.

Chapter 6

Mutation Surveyor Custom Report

A text-based custom report is available in Mutation Surveyor. This reports contain options for building a tailored report that provides detailed information about your mutation projects in a format that best suits your specific working needs. In addition, you can print the report information and you can save and store the report output for use at a later date.

This chapter covers the following topics:

- [“Custom Report” on page 164.](#)

Custom Report

The Custom report offers a variety of report formats in which to display your mutation project data. A variety of options are also available for tailoring the report to best suit your working needs, including sample grouping options, report heading options, options for filtering the variations that are displayed in the report, display options for samples and mutation calls, color-coding options for mutation calls, and nomenclature options for mutation calls. You can [create](#) a custom report template from scratch, or you can load an existing custom report template, modify the settings as needed, and then save the modified template.

To create a Custom Report template

1. On the Mutation Surveyor main menu, click Reports > Custom Report.

The Custom Report Settings dialog box opens. The Format tab is the active tab.

Figure 6-1: Custom Report Settings dialog box, Format tab

The screenshot shows the 'Custom Report Settings' dialog box with the 'Format' tab selected. The dialog has several sections: 'Table Format', 'Headings', 'Grouping', and 'Group Summary'. The 'Table Format' section includes options for 'One Trace Per Row', 'Generate Forward/Reverse Consensus', 'Generate Group Consensus', 'Mutation Position Specific for Column' (with sub-options 'By Group' and 'By Table'), 'One Mutation Per Row', 'Transpose', and 'Report One Nucleotide Position Per Column'. The 'Headings' section includes checkboxes for 'Show Display Region Heading', 'Show Mutation Number', 'Show Position Information', and 'Show Default Header Information', along with a text field for 'Custom Header File' and a 'Create Header' button. The 'Grouping' section includes three rows for '1st Order', '2nd Order', and '3rd Order', each with a dropdown menu and 'from character' and 'to' fields. There is also a checkbox for 'Add Blank Row Between Groups' and a 'Load Sample ID' button. The 'Group Summary' section includes options for 'Do Not Show Group Summary Information', 'Show Allele Information' (with sub-options 'Show Allele Frequency' and 'Show Advanced Allele Information'), and 'Show Mutation Summary'. At the bottom are buttons for 'Default', 'Load', 'Save', 'Ok', and 'Cancel'.

Section	Option	Value/Status
Table Format	One Trace Per Row	Selected
	Generate Forward/Reverse Consensus	Not Selected
	Generate Group Consensus	Not Selected
	Mutation Position Specific for Column	Selected
	By Group	Selected
	By Table	Not Selected
	One Mutation Per Row	Not Selected
	Transpose	Not Selected
Table Format	Report One Nucleotide Position Per Column	Not Selected
	Headings	
Headings	Show Display Region Heading	Selected
Headings	Show Mutation Number	Selected
Headings	Show Position Information	Not Selected
Headings	Show Default Header Information	Selected
Headings	Custom Header File	[Empty]
Headings	Create Header	Button
Grouping	1st Order	Contig
	from character	1
	to	3
	2nd Order	F/R Pairing
	from character	1
	to	3
3rd Order	Trace Position	
from character	1	
to	3	
Grouping	Add Blank Row Between Groups	Selected
Grouping	Load Sample ID	Button
Group Summary	Do Not Show Group Summary Information	Not Selected
	Show Allele Information	Selected
	Show Allele Frequency	Selected
	Show Advanced Allele Information	Not Selected
	Show Mutation Summary	Not Selected
Default		
Load		
Save		
Ok		
Cancel		

2. Do one of the following:

- To set up a new custom report from scratch, select your options and specify the appropriate values on the different dialog box tabs.
- To generate the report based on existing settings, click Load to browse to and select the MS Report Template.ini file that you want to use. Optionally, before you generate the report, modify the settings for the loaded report as needed.

See:

- “Format tab” on page 166.
- “Filters tab” on page 170.
- “Display tab” on page 172.
- “Mutation tab” on page 173.
- “Color tab” on page 176.
- “Nomenclature tab” on page 177.
- “Other tab” on page 179.



Click Default at any time to reset all the values on all the tabs to the system default values.

3. Do one of the following:

- To generate an onscreen report according to the settings that you have specified, click OK.

The report opens in the Custom Report window.

Figure 6-2: Custom Report window

Trace #	Sample Name	Reference N	Direction	Gene	Exon	Reading Frame	Lane Quality	# of
1	BRCA1_012345_10A_1_F_0_o_2007-11-28_2	BRCA1_Amp 1-F		BRCA1	2	16	0	
2	BRCA1_012345_10A_1_R_0_s_2007-10-28_2	BRCA1_Amp 1-R		BRCA1	2	14	0	
3	BRCA1_012345_10A_2_F_0_o_2007-12-20_1	BRCA1_Amp 1-F		BRCA1	2	16	0	
4	BRCA1_012345_10A_2_R_0_s_2007-11-13_2	BRCA1_Amp 1-R		BRCA1	2	14	0	
5	BRCA1_185729_10A_1_F_1_s_2007-10-22_1	BRCA1_Amp 1-F		BRCA1	2	25	0	
6	BRCA1_185729_10A_1_R_1_i_2007-12-14_2	BRCA1_Amp 1-R		BRCA1	2	26	0	
7	BRCA1_185729_10A_2_F_0_s_2007-10-23_1	BRCA1_Amp 1-F		BRCA1	2	25	0	
8	BRCA1_185729_10A_2_R_0_i_2007-12-26_1	BRCA1_Amp 1-R		BRCA1	2	26	0	
9	BRCA1_264381_10A_1_F_1_o_2007-11-15_1	BRCA1_Amp 1-F		BRCA1	2	25	0	
10	BRCA1_264381_10A_1_R_1_o_2007-11-19_1	BRCA1_Amp 1-R		BRCA1	2	25	0	
11	BRCA1_264381_10A_2_F_0_i_2007-10-11_16	BRCA1_Amp 1-F		BRCA1	2	25	0	
12	BRCA1_264381_10A_2_R_0_s_2007-11-13_1	BRCA1_Amp 1-R		BRCA1	2	25	0	

Chapter 6

Mutation Surveyor Custom Report

- To save the settings for loading and generating a report at a later date, click Save.



The Save Mutation Surveyor Report Template dialog box opens. By default, a Custom report is saved as MutationSurveyor_CRSettings.ini file in the %localappdata%\SoftGenetics\MutationSurveyor\Report folder. If you do not want to override the default custom report file supplied with Mutation Surveyor, change the name of the file, the location of the file, or both.


Format tab

The Format tab contains options for the formatting and layout of the custom report.

Figure 6-3: Custom Report Settings dialog box, Format tab

Option	Description
Table Format	
One Trace Per Row	Display all mutation information for each trace in its own row.
Generate Forward/Reverse Consensus	Create a single consensus by merging the forward consensus and reverse consensus.
Generate Group Consensus	Create a single consensus by merging all the consensus information for the sample group.
Mutation Position Specific for Column	Aligns mutations by column within each group (By Group) or aligns mutations by column across all groups (By Table).
One Mutation per Row	Display information for each mutation call in its own row. Note: Select if mutation information is to be exported to a database.
Transpose	Generates a table that has base positions in rows and sample traces in columns.

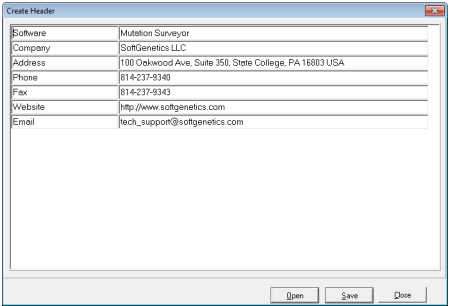Option	Description
Headings	
Show Display Region Heading	Selected by default. If you clear this option, then column headings are not displayed in the report.
Show Mutation Number	Labels mutation call columns with Variant number.
Show Position Information	Labels mutation call columns by the genomic position referenced in the GenBank file. Note: A particularly useful option if you display mutation calls by a nomenclature other than Genomic.
Show Default Header Information	Displays the Mutation Surveyor version number, the date and time that the report was generated, the project name (if loading a saved project) and the report settings file name. "Default" is displayed as the report settings filename when the report is generated using the system default settings. The actual report filename is displayed when the report is generated using a saved report settings file. "Customized" is displayed in all other cases. (The settings were not loaded using a saved report settings file and the settings are not the system default settings.)
Create Header/Custom Header File	<p>Click Create Header to open the Create Header dialog box which contains a template for creating a custom header that can be displayed at the top of a custom report.</p> <p>Figure 6-4: Create Header dialog box</p>  <p>Do one of the following:</p> <ul style="list-style-type: none"> • Modify any of the default template information in either column. • Click Open to select an existing custom header file to modify. <p>You can add or delete rows of information as needed.</p> <ul style="list-style-type: none"> • To delete a row from the header, right-click the row, and then click Delete a Row. • To insert a row into the header, right-click the row that is to be located <i>below</i> the inserted row, and then click Insert a Row. • To add a row as the last row in the header, right-click any row, and then click Add a Row. <p>Click Save to save the custom header as an .inf file, and then click the Browse button  to browse to and load the saved header file for the custom report.</p>

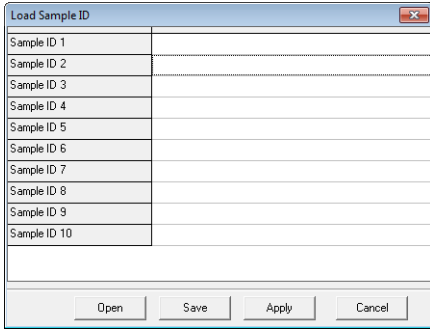Option	Description
Grouping —Traces are initially grouped based on the 1st Order Grouping option. Further sorting is then carried out based on the 2nd Order Grouping option, and then on the 3rd Order Grouping option.	
<ul style="list-style-type: none"> Contig Trace Grouping File Sample ID 	<ul style="list-style-type: none"> The default value for 1st Order. Groups are sorted based on the Contigs that Mutation Surveyor created for determining mutation calls. When Trace is selected, the From Character X to Y option is enabled. Traces are sorted into groups based on character matching in the specified positions in the trace filenames. Traces are sorted into groups based on the grouping method that is specified in the Load Grouping text file. See “2D Filename Match Editor tool” on page 261. When Sample ID is selected, both the Load Sample ID button and the From Character X to Y option are enabled. Click Load Sample ID to open the Load Sample ID dialog box and specify the different Sample IDs, such as Patient #, by which to group the traces. You can also click Open to load an existing Sample ID file. <p>Note: By default, the Load Sample ID dialog box contains rows for up to ten different Sample IDs. You can right-click and click Insert or Delete as needed.</p> <p><i>Figure 6-5: Load Sample ID dialog box</i></p>  <p>The traces that are to be grouped together must contain at least one of these Sample IDs in their filenames but they do not need to be in the same location in the filename. When grouping by Sample ID, the trace filenames are first checked based on the Sample IDs defined in the Load Sample ID dialog box. The values specified in the From Character X to Y fields are ignored if all the trace filenames contain at least one of these Sample IDs. Traces that do not contain any of these Sample IDs in their filenames are sorted into groups based on the values specified in the From Character X to Y fields. Traces that do not contain any of these Sample IDs and that do not contain the characters specified in the From Character X to Y fields are sorted into their own group.</p>
<ul style="list-style-type: none"> F/R pairing 	<ul style="list-style-type: none"> The default value for the 2nd Order. Places two-directional traces in adjacent rows.

Option	Description
<ul style="list-style-type: none"> Trace Position Gene Exon Trace Quality None All in One Group 	<ul style="list-style-type: none"> The default value for the 3rd Order. Traces are sorted into groups based on where they match to the GenBank reference file. Each trace is matched at its starting genomic base number. Groups can be sorted so that traces that match to the earliest reference base are displayed first. Traces are sorted into groups based on the gene names defined in the GenBank reference file. Traces are sorted into groups based on the exons defined in the GenBank reference file. Traces are sorted into groups based on the lane quality for the trace. Ignore the sorting level. Traces are placed in a single group. Only available for the 1st Order.
Add Blank Rows between Groups	Selected by default. Clear this option if you do not want a blank row to be displayed between groups of traces in the custom report.
Group Summary	
<ul style="list-style-type: none"> Do Not Show Group Summary Information Show Allele Information Show Mutation Summary 	<ul style="list-style-type: none"> Do not display a row below each group that shows such information as the average frequency of the call among the traces. Show Allele Frequency—Adds a row below each group that displays the average frequency of the call among all the traces. Traces for which data is not available for the indicated position are not used in the frequency calculation. Show Advanced Allele Information—Applicable only if you are analyzing diploid samples. Displays the frequency of each allele by traces and as a grand total. <ul style="list-style-type: none"> Total—Shows the frequency of the two alleles within each sample trace. Grand Total (2D data) <ul style="list-style-type: none"> If both the forward and reverse traces show a mutation at the position, then the alleles in the Forward direction mutation call are used in the calculation of the allele frequency. If data is not available for the position in one of the traces, then the alleles in the mutation call in the other direction are used in the calculation of the allele frequency. If data is available for the position in both traces, but only one of the traces has a mutation call, then Mutation Surveyor ignores this pair and the alleles in the sample trace with the mutation call are not used in the calculation of the allele frequency. Grand Total (1D data)—The alleles in each trace are used in the calculation of the allele frequency. Applicable only if you have loaded a GenBank file with reported variations. Displays a listing of variants within each group, with variants placed in one of two columns—Unreported and Reported. The Unreported column displays novel variants that were found in the sample files, but were not reported in the GenBank file. The Reported column displays the reported variants that were found in the sample files.

Filters tab

The Filters tab contains options for specifying which traces are to be displayed in the custom report.



Some of the options on this tab are self-explanatory, and therefore, not all these options are discussed in this section.

Figure 6-6: Custom Report Settings dialog box, Filters tab

Option	Description
Traces	
Show Selected Traces	<p>Filtering options that limit the traces that are displayed in the report to only those that meet the selected criteria.</p> <ul style="list-style-type: none">Containing No Mutation Calls<ul style="list-style-type: none">Containing only normal alleles—Traces without mutations that can have mosaics, negative SNPs, and/or negative mutations.Containing Negative SNPs—A <i>negative SNP</i> is defined as position in the GenBank file at which a variation has been reported, but no variation was detected at the same position in the sample trace. Sample traces that have negative SNPs might still contain mutation calls.Containing Negative Mutations—A <i>negative mutation</i> is defined as a position in a sample trace at which a variation was detected, but no variation was detected at the same position for other sample traces in the group. Sample traces that have negative mutations might still contain mutation calls.Unmatched Traces—Clear this option to remove sample traces from the Custom report that do not match to the reference trace.

Option	Description
Nucleotide Quality	
<ul style="list-style-type: none"> Phred Score Intensity 	Define nucleotide quality thresholds based on the nucleotide Phred Score and the nucleotide Intensity. Used to mitigate the calling of false positives when the data quality is poor.
Mutation Calls	
Show All Mutations	Show all mutations of all types in all regions.
Show Selected Mutations	<p>Filtering options that limit the mutations that are displayed in the report to only those that meet the selected criteria of type and region. (Show only mutations of the selected type in the selected regions.)</p> <ul style="list-style-type: none"> Regions—Limit the mutations that are displayed to only those that are called within regions of interest (ROI) that are defined in the GenBank reference file. Mutation Types <ul style="list-style-type: none"> Selected Types—You must choose at least one type from each group. <p>Note: When choosing at least one type from each group:</p> <ul style="list-style-type: none"> Homozygous Dropping Threshold—Indicates the value for the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions. You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions. Reject 1D Calls When Other Direction has No Call—Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction. Reject 1D Calls When Other Direction is Not Available—Reject a mutation that is called in only one direction when the position falls outside the comparison region of the other direction.

Display tab

The Display tab contains options for specifying the columns to display in the custom report.



The majority of the options on this tab are self-explanatory, and therefore, not all these options are discussed in this section.

Figure 6-7: Custom Report Settings dialog box, Display tab

The screenshot shows the 'Custom Report Settings' dialog box with the 'Display' tab selected. The 'Display Region Information' section has 'Show Selected Information' checked. Under 'General Information', 'Sample Name' is checked. Under 'Comparison Region', 'Read Start' and 'Read End' are checked. Under 'Region of Interest', 'ROI Start', 'ROI End', 'ROI Phred String', and 'ROI Coverage' are checked. Under 'Quality Information', 'Entire Trace' is checked. Under 'Nucleotide Quality Information', 'Phred Score' is checked. The 'Quality Threshold' is set to 15 and the 'Nucleotide Quality Threshold' is set to 20.00. The 'Default', 'Load', 'Save', 'OK', and 'Cancel' buttons are at the bottom.

Option	Description
Region of Interest	<ul style="list-style-type: none"> ROI Coverage—Identifies when an ROI has not aligned entirely within the Comparison Region. ROI Phred String—Display a single ASCII character representation of the Phred score for each nucleotide position within an ROI. The ASCII character that is displayed for a Phred Score is the Phred Score + 32. For example, if the Phred Score = 1, then 1 + 32 = 33, or the 33rd character in the ASCII character set, which is an exclamation point (!). <p>Note: For an ASCII character map, see http://www.asciitable.com/.</p>
Quality Information	<ul style="list-style-type: none"> Trace Quality Information—Options in this section are used to identify traces of low quality. The Low Quality Threshold is the value that is specified for Lane Quality in the Mutation project settings. (See “Quality Trim” on page 58.) The trace is also identified with the background color that has been selected for Quality Failure on the Colors tab. (See “Color tab” on page 176.) Nucleotide Quality Information—Options in this section are used to identify nucleotide positions of poor quality. (See “Score Trim []” on page 58.)

Mutation tab

The Mutation tab contains options for displaying information for the various mutation calls as well as options for displaying information about positions where mutations were not called.



The majority of options on this tab are self-explanatory, and therefore, not all these options are discussed in this section.

Figure 6-8: Custom Report Settings, Mutation tab

Option	Description
Show Mutation Call Information	
Nucleotide Position	Nucleotide positions are displayed in the nomenclature that you select on the Nomenclature tab. (See “Nomenclature tab” on page 177.)
Amino Acid Position	By default, amino acid information is organized with the reference amino acid appearing first, the amino acid position appearing next, and the mutation amino acid appearing last. If you do not select Amino Acid Position, then a slash (/) is displayed between the reference amino acid and the mutation amino acids.
Confidence(!)	Displays a Confidence mark (!) for mutation calls that exceed certain thresholds that are defined in the MutationSurveyor.ini file. Caution: Although you can modify these thresholds, SoftGenetics does not recommend that you do so on your own. Contact SoftGenetics for assistance.
Comments	Displays all comments that were added to a mutation call, either automatically by Mutation Surveyor or manually by a user.


Option	Description
Show Non-Mutation Call Information	
Show Normal Alleles	<ul style="list-style-type: none"> • Show Negative SNP—Display selected Show Normal Allele Information for nucleotides that do not show the variations reported in the GenBank or SEQ file. For indels, “no_ins” or “no_del” is displayed. • Show Negative Mutations—Display selected Show Normal Allele Information for a position in a sample trace at which a variation was <i>not</i> detected, but the variation was detected at the same position for other sample traces in the group. • Show Normal Allele Information—Specify the information that is to be displayed for positions that do not have an identified mutation call. <p>Note: When Normal Allele Information is displayed and One Mutation per Row is selected (see “Table Format” on page 166), all options selected in the Show Mutation Call Information section are exported for both mutation calls and normal alleles.</p>
General Information	
Reported Variant Annotation	Displays all the variations’ annotations that are included in the GenBank file such as the dbSNP identifier as well as the annotation fields from the Variation Tracks selected to report.
Genotype	With this option selected, Mutation Surveyor assumes that the analysis is for a diploid organism and displays the two alleles at the nucleotide position. For indels, a homozygous insertion of “A,” for example, is displayed as “insA insA,” and a heterozygous insertion of “A,” is displayed as “normal ins A.” If there is no insertion, then “Normal Normal” is displayed.
NM_Score	Assigns a score to each sample nucleotide that indicates the contribution of the normal allele at the position compared to the contribution of the possible mutant alleles at the same position. The score is based on the height ratio of the four different dye colors and the shape of the four peaks at each nucleotide position. The range is 0-100. A high score indicates a high degree of similarity between the sample trace and the reference trace, which, in turn, means that the sample trace more than likely contains a normal allele at the position. A low score indicates a high degree of difference between the sample trace and the reference trace, which, in turn, means that the sample trace more than likely contains a homozygous substitution. Scores in the middle indicate high levels of noise and error or heterozygous substitutions in the sample trace.

Option	Description
Display Information —Specifies how the mutation call information is to be distributed in table cells for display.	
<ul style="list-style-type: none"> One Field Multiple Fields 	<ul style="list-style-type: none"> All the information displayed in a single field. Report the Deletion and Insertion as Delins—If an insertion and deletion occur at the same position, then they are merged into a single call at the position. Mutation Call Nucleotides Relative to Gene Direction—Make the mutation call based on the gene orientation. To make a mutation call for a gene on the reverse strand, a reverse complement is generated. Multiple Fields (Group Nts, A.A., and Score into One Field or Group Nts and A.A. into Relevant Fields)—Specific pieces of information grouped into fields. All Information in Unique Fields—Each piece of information displayed in its own column.
Show Advanced Display	<ul style="list-style-type: none"> [] # of nucleotides Displayed Before/After Mutation—Displays the specified normal number of alleles on either side of the variant. The position is displayed with a slash (/) between the normal allele and the mutation alleles. Display Position—Displays the reference nucleotide and its position, the consensus nucleotides for the group at this position, the amino acid information (if applicable), and the reported variants' annotations. This information can be displayed in One Field or in Multiple Fields.

Color tab

The Color tab provides a variety of color-coding options to assist you with identifying different types of mutation calls and potential analysis problems.

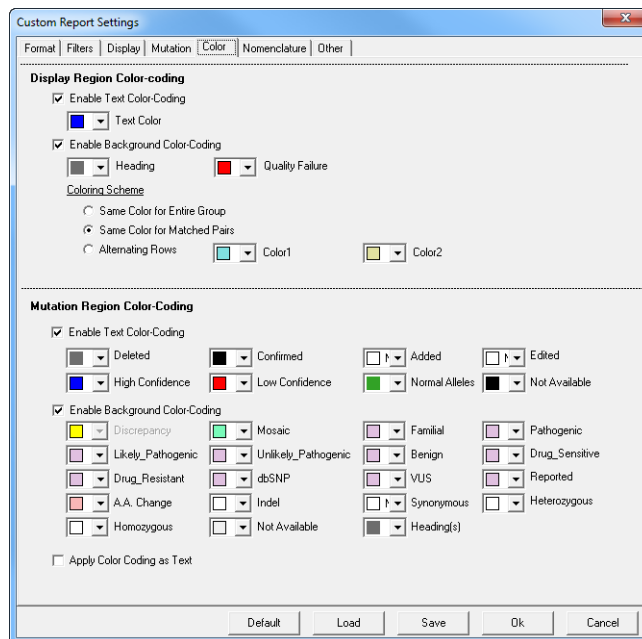Figure 6-9: Custom Report Settings dialog box, Color tab



Color-coding is applied according to the following:

- If a mutation call meets several color-coding criteria, the call is displayed in the report in the color of highest priority.
- If None is selected for a color-coding option, then the color of the next highest priority is used.
- For Mutation Region color-coding, Quality Failure has the highest priority.
- For Mutation Region Text color-coding, the priorities from highest to lowest are as follows—Deleted, Confirmed, Added, Edited, High Confidence, Low Confidence, Normal Alleles, and Not Available.
- For Mutation Region Background color coding, priorities from highest to lowest are as follows—Discrepancy, Mosaic, Familial, Pathogenic, Likely_Pathogenic, Unlikely_Pathogenic, Benign, Drug_Sensitive, Drug_Resistant, dbSNP, VUS, Reported, A.A. Change, Indel, Synonymous, Heterozygous, Homozygous, Not Available, and Heading(s) where:
 - *Discrepancy* identifies the variants within a consensus of traces that have at least one member of the consensus that differs from the others. The Discrepancy indicator is available for the Generate Forward/ Reverse Consensus table format and the Generate Group Consensus table format. (See “Table Format” on page 166.)

- *Synonymous* identifies the coding region substitutions that do not affect the amino acid sequence.




When a consensus is formed, discrepancies are identified only for positions that have data available. For example, if Generate Forward/Reverse Consensus is the selected table format, and the Forward trace contains a mutation call at position #320 but the Reverse trace does not at this position because data is not available, then a discrepancy is not detected.

If you select Apply Color Coding as text, then when you save the Custom report as a .txt file, the color-coding is shown as a comment in the file.

Nomenclature tab

The Nomenclature tab contains options for specifying the nomenclature that is to be used for the mutation calls.

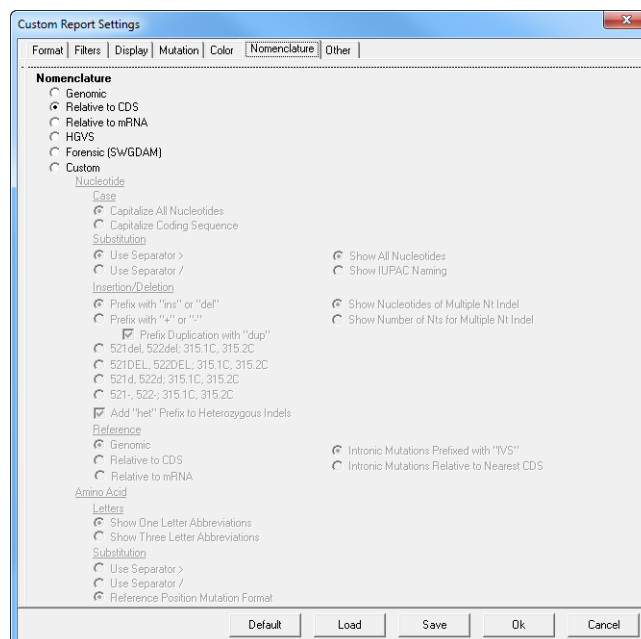Figure 6-10: Custom Report Settings dialog box, Nomenclature tab



Option	Description
Genomic	Displays the mutation call positions based on the genomic position referenced in the GenBank file.
Relative to CDS	Displays the mutation call positions relative to the CDS numbering in the GenBank reference file. A mutation that is called in the CDS is prefixed with a "c." A mutation that is called in a non-coding region is prefixed with "IVS." Its position is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide.

Option	Description
Relative to mRNA	Displays the mutation call positions relative to the mRNA numbering in the GenBank reference file. A mutation that is called in the mRNA region is prefixed with a "c." A mutation that is called outside of an mRNA region is identified by the closest nucleotide in the mRNA region plus or minus the number of bps by which it resides away from the closest nucleotide.
HGVS	Displays the mutation call positions relative to its position in the coding sequence. A mutation that is called in a non-coding region is identified by the closest nucleotide in the coding sequence plus or minus the number of bps by which it resides away from the closest nucleotide. Both alleles are shown in separate bracketed statements.
Forensic (SWGDAM)	Displays the mutation call using SWGDAM Guidelines.
Note: If any one of the four standard Nomenclature options is selected and a GenBank file was not loaded as the reference file, then the Base Number is enclosed in parenthesis and it indicates the position of the base in the reference file. If both a GenBank file and a reference file were loaded for the mutation project and the mutation was called at a position that does not align to the GenBank file, then the Base Number is enclosed in parenthesis and is displayed as +/- the number of base pairs by which the mutation was shifted in comparison to the GenBank file.	
Custom	Allows you to tailor the nomenclature for mutation calls to best suit your working needs.


Figure 6-11: Nomenclature examples

			Genomic	Relative to CD5	Relative to mRNA	HGVS	Forensic (SWGDAM)
Substitution	Homozygous	non-coding	873C>T	IV541-14C>T	c.28-14C>T	c.[41-14C>T]+[41-14C>T]	C873T
		coding	77A>T	c.31A>T	c.18A>T	c.[31A>T]+[31A>T]	A77T
	Heterozygous	non-coding	821C>CT	IV541-66C>CT	c.[=,28-66C>T]	c.[41-66C>T]+[=]	C821Y
		coding	80C>CA	c.34C>CA	c.[=,21C>A]	c.[34C>A]+[=]	C80M
Indel	Homozygous	non-coding	211delT	IV571+4delT	c.58+4delT	c.71+4delT	211d
		coding	18_20delTTT	c.72_74delTTT	c.59_61delTTT	c.72_74delTTT	18d,19d,20d
	Heterozygous	non-coding	133_134het_delTA	IV571+16_71+17het_delTA	c.58+16_58+17het_delTA	c.71+16_71+17het_delTA	133d,134d
		coding	14_15het_insT	c.68_69het_insT	c.55_56het_insT	c.68_69het_insT	15.1T

Other tab

The Other tab contains options that affect the report display as well as processing options for the report.

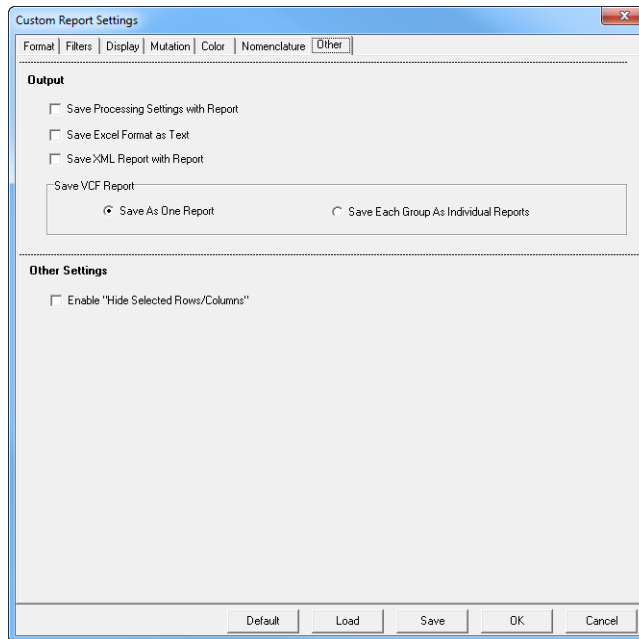Figure 6-12: Custom Report Settings dialog box, Other tab






Option	Description
Output	
Save Processing Settings with Report	Saves all the settings that you selected for the report as part of the report output.
Save Excel Format as Text	Saves the Custom report as an Excel file with all the data columns formatted as Text.
Save XML Report with Report	Saves the Custom report as an XML file.
Save VCF Report	<ul style="list-style-type: none"> Save as One Report—Selected by default. Generate a single VCF report that includes the information for all the sample groupings. Save Each Group as Individual Reports—Generate a VCF report for each sample grouping. <p>Note: Remember, you specify the sample grouping on the Format tab. See “Format tab” on page 166.</p>
Other Settings	
Enable “Hide Selected Rows/Columns”	Enables on the Hide Columns and Hide Rows function on the Custom report context menu. See “Custom Report viewing and analysis options” on page 181 .

Custom Report toolbar

The Custom report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

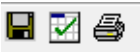Figure 6-13: Custom report toolbar








Button	Description
	Save—Opens the Save Report As dialog box in which you can specify a name for the saved report as well as its format (.vcf, .xls/.xlsx, or .txt) The default location is %localappdata%\SoftGenetics\MutationSurveyor\Report, but you can select a different location.
	Settings—Opens the Custom Report Settings dialog box in which you can adjust your report settings.

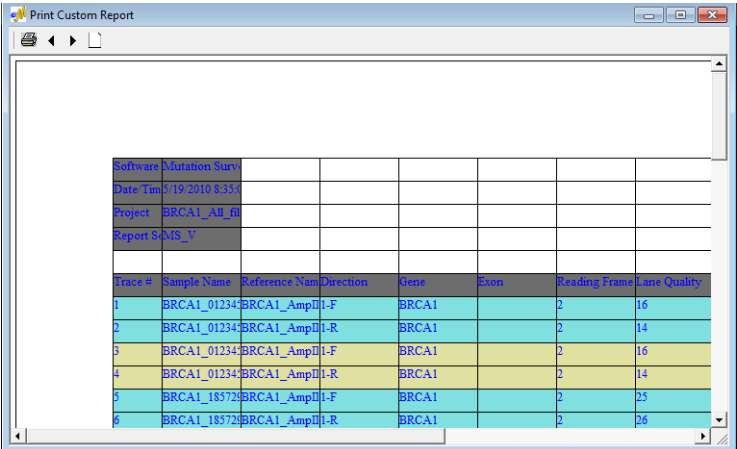

Print—Opens the custom report in the Print Custom Report window. The window toolbar contains buttons for working with the report and for printing the report.

Figure 6-14: Print Custom Report window




- Print button  —Opens the Print dialog box in which you can specify the options for printing the Custom report. By default, only the page that is currently displayed in the Print Custom Report window is selected for printing. You can leave the default as is, or you can select the entire report or a range of pages for printing.
- Display Previous  and Display Next  buttons—Use these buttons to Browse forward and backward through the Custom report one page at a time.
- Page Setup button  —Opens the Page Setup dialog box in which you can specify the page properties of the report pages (such as page size and orientation) for printing.

Custom Report viewing and analysis options

A context menu is available for the Custom report. You use the options on this menu to change the view of the Custom report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the Custom report.

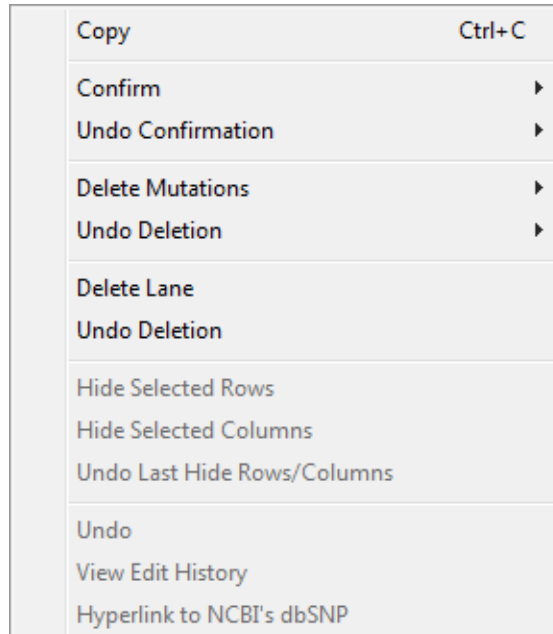Figure 6-15: Custom report context menu



Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Custom report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations for confirmation or to select all mutations in a column for confirmation.</p> <p>The text for the mutation code turns black in both the Custom report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>

Option	Description
Undo Confirmation	Removes the “Confirmed” status for a selected mutation or a selected column. Note: SHIFT-click to select multiple contiguous mutations for confirmation or to select all mutations in a column for confirmation.
Delete Mutations	Deletes the selected mutations. Note: SHIFT-click to select multiple contiguous mutations for deletion, to select all mutations in a column for deletion, or to select multiple rows for deletion. The text for the mutation code turns gray in both the Custom report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.
Undo Deletion	Removes the “Deleted” status for a selected mutation, a selected column, or a selected row.
Delete Lane	Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to “Low Quality.”
Undo Lane Deletion	The Lane Quality for the lane is reset to its original value.
Hide Selected Rows Hide Selected Columns	Enabled if “Enable Hide Selected Rows/Columns” is selected on the Other tab. (See “Other tab” on page 179.)
Undo Last Hide Rows/ Columns	Enabled if selected rows and/or columns are hidden in the Custom report.
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See “Viewing the Edit History (Audit Trail) for a Mutation” on page 344.
Hyperlink to NCBI's dbSNP	Opens the NCBI's database entry (via the web) for the variant.

Other options include the following:

- To open the GAD for the trace that contains a mutation, double-click the mutation in the Custom report.



If you click a trace in the GAD with the Custom report open, the GAD is brought to the front. To bring the Custom report to the front again, press [CTRL]-T.

- To select all cells in the Custom report in a single step, press [CTRL]-A.
- To auto-scale the width of a column to its largest entry, double-click the first cell in the column.



If you auto-scale the report after selecting all the cells in the report, you can resize the entire report easily in a single step.

Chapter 7

Mutation Surveyor Customer Reports

Customer reports are another type of custom report in Mutation Surveyor. Each report was initially developed as a template to provide custom content that met a specific customer's needs. All these report templates are now available from the Reports > Customer Reports option on the Mutation Surveyor main menu

This chapter covers the following topics:

- [“JHU report” on page 185.](#)
- [“Tassel report” on page 190.](#)
- [“Emory report” on page 193.](#)
- [“Bentley report and MSM_Song report” on page 196.](#)
- [“PrettyBase report” on page 200.](#)
- [“Methylation report” on page 203.](#)
- [“Vertex report” on page 207.](#)
- [“Vertex2 report” on page 209.](#)
- [“Vertex3 report” on page 212.](#)
- [“Full Gene report” on page 215.](#)
- [“FullBase report” on page 222.](#)


Chapter 7

Mutation Surveyor Customer Reports

JHU report

The JHU report displays various quality scores for a defined ROI organized by paired sample traces. The layout of the JHU Output report is identical to the layout of the Two Direction report. (See [“Two Direction report layout” on page 134.](#)) The mutation codes, and the color coding of the mutation calls in the report are identical to the mutation codes and color coding in the Mutation report (See [“Mutation report fields” on page 128,](#) [“Mutation report codes” on page 129,](#) and [“Color codes” on page 130.](#)) The JHU report contains fields that are specific to the report as well as all the fields that are displayed in any of the standard mutation reports.

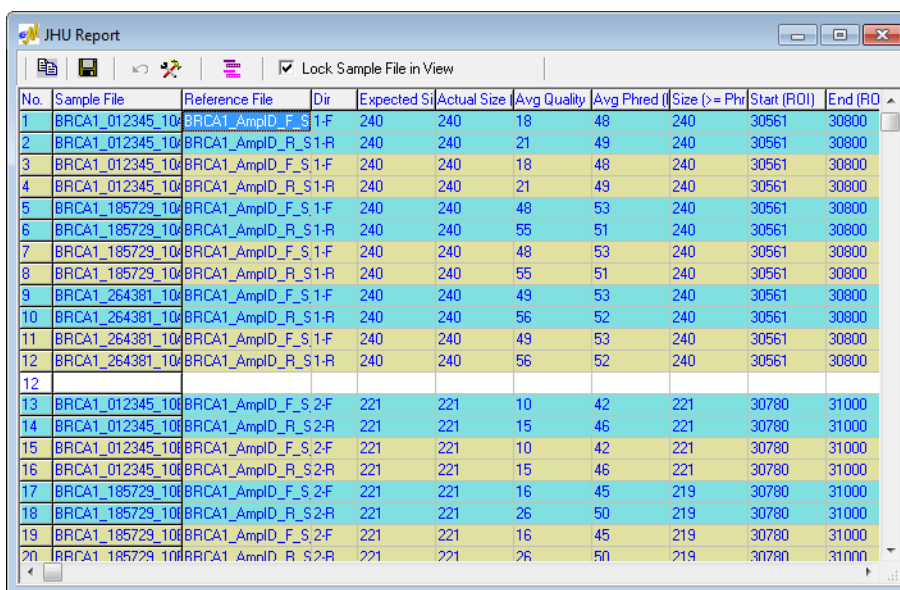Figure 7-1: JHU report



The screenshot shows a window titled "JHU Report" with a toolbar and a table of data. The table has columns for No., Sample File, Reference File, Dir, Expected Size, Actual Size, Avg Quality, Avg Phred, Size (>= Phr), Start (ROI), and End (ROI). The data is organized into two groups of paired sample traces (1-F and 1-R) for different sample files and reference files.






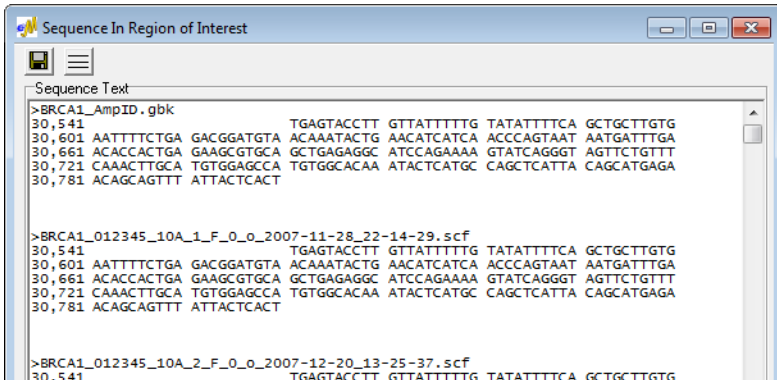


No.	Sample File	Reference File	Dir	Expected Size	Actual Size	Avg Quality	Avg Phred	Size (>= Phr)	Start (ROI)	End (ROI)
1	BRCA1_012345_10x	BRCA1_AmplD_F_S	1-F	240	240	18	48	240	30561	30800
2	BRCA1_012345_10x	BRCA1_AmplD_R_S	1-R	240	240	21	49	240	30561	30800
3	BRCA1_012345_10x	BRCA1_AmplD_F_S	1-F	240	240	18	48	240	30561	30800
4	BRCA1_012345_10x	BRCA1_AmplD_R_S	1-R	240	240	21	49	240	30561	30800
5	BRCA1_185729_10x	BRCA1_AmplD_F_S	1-F	240	240	48	53	240	30561	30800
6	BRCA1_185729_10x	BRCA1_AmplD_R_S	1-R	240	240	55	51	240	30561	30800
7	BRCA1_185729_10x	BRCA1_AmplD_F_S	1-F	240	240	48	53	240	30561	30800
8	BRCA1_185729_10x	BRCA1_AmplD_R_S	1-R	240	240	55	51	240	30561	30800
9	BRCA1_264381_10x	BRCA1_AmplD_F_S	1-F	240	240	49	53	240	30561	30800
10	BRCA1_264381_10x	BRCA1_AmplD_R_S	1-R	240	240	56	52	240	30561	30800
11	BRCA1_264381_10x	BRCA1_AmplD_F_S	1-F	240	240	49	53	240	30561	30800
12	BRCA1_264381_10x	BRCA1_AmplD_R_S	1-R	240	240	56	52	240	30561	30800
13	BRCA1_012345_10x	BRCA1_AmplD_F_S	2-F	221	221	10	42	221	30780	31000
14	BRCA1_012345_10x	BRCA1_AmplD_R_S	2-R	221	221	15	46	221	30780	31000
15	BRCA1_012345_10x	BRCA1_AmplD_F_S	2-F	221	221	10	42	221	30780	31000
16	BRCA1_012345_10x	BRCA1_AmplD_R_S	2-R	221	221	15	46	221	30780	31000
17	BRCA1_185729_10x	BRCA1_AmplD_F_S	2-F	221	221	16	45	219	30780	31000
18	BRCA1_185729_10x	BRCA1_AmplD_R_S	2-R	221	221	26	50	219	30780	31000
19	BRCA1_185729_10x	BRCA1_AmplD_F_S	2-F	221	221	16	45	219	30780	31000
20	BRCA1_185729_10x	BRCA1_AmplD_R_S	2-R	221	221	26	50	219	30780	31000

JHU Report toolbar

The JHU report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 7-2: JHU report toolbar



Button	Description
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
	Undo Latest Deletions—Restores the deletion that you last carried out in the JHU report. Allows for multiple undos.
	Report Settings—Opens the JHU Report Settings dialog box which contains options for filtering the information that is to be displayed in the report. See “JHU report settings” on page 187 .
	<p>Display the Sequence of ROI—Generates an onscreen text file of the sequences in the GenBank ROIs.</p> <p><i>Figure 7-3: Sequence in Region of Interest window</i></p>  <ul style="list-style-type: none"> Save button  —Opens the Save Sequence Text As dialog box in which you can specify a name and location for saving the .txt file. Display Sequence in One Line button  —Displays the sequences in fasta format.
Lock Sample File in View	Freezes the Sample Filename column while scrolling.

JHU report viewing and analysis options

The JHU report has multiple options available for changing the view of the report and for manipulating and analyzing mutation project data. These options are:

- The JHU report settings. See [“JHU report settings”](#) below.
- The JHU report context menu. See [“JHU report context menu”](#) on page 188.

JHU report settings


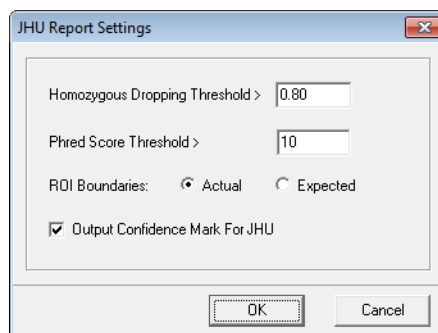
To open the JHU Report Settings dialog box, click the Report Settings button  on the JHU report toolbar.

Figure 7-4: JHU Report Settings dialog box




Option	Description
Homozygous Dropping Threshold	Indicates the value for the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions. Note: You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.
Phred Score Threshold	Call a mutation only if the Phred score for the variant base or bases is greater than the indicated threshold.
ROI Boundaries	Affects the values that are output to the Start (ROI) and End (ROI) column for traces that do not have complete coverage of the Region of Interest. For example, an ROI is set for positions 50 to 150 (Expected value). The sample trace's comparison region is positions 100 to 200. Only positions 100 to 150 of the ROI are within the comparison region (Actual value) whereas positions 50-99 are not covered.
Output Confidence Mark for JHU	Displays a Confidence mark (!) for mutation calls that exceed certain thresholds that are defined in the MutationSurveyor.ini file. Note: Although you can modify these thresholds, SoftGenetics does not recommend that you do so on your own. Contact SoftGenetics for assistance.


JHU report context menu

A context menu is available for the JHU report. You use the options on this menu to change the view of the JHU report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the JHU report.

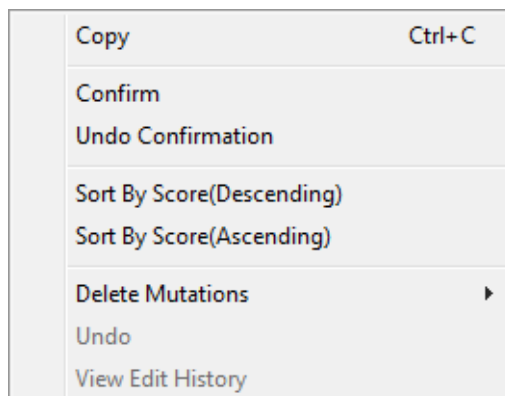Figure 7-5: JHU report context menu



Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the JHU report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the JHU report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Undo Confirmation	<p>Removes the “Confirmed” status for the mutation.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p>
Sort By Score (Descending) Sort by Score (Ascending)	<p>Click the appropriate option to sort all the data by descending mutation score or ascending mutation score.</p>


Option	Description
Delete Mutations	<p>Deletes the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations, to select all mutations in a column, or to select multiple rows.</p> <p>The text for the mutation code turns gray in both the JHU report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p> <p>Note: To undo the last deletion that you carried out, click the Undo Latest Deletions button  on the report toolbar.</p>
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See “Viewing the Edit History (Audit Trail) for a Mutation” on page 344.

You can also view the information for a mutation in the JHU report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

Tassel report

The Tassel report is ideal for identifying mutations in your sample traces if your traces are grouped by Sample ID (see “[Sample ID](#)” on page 168) or if all the filenames in your project contain the same characters/designator but in different positions. After you click Tassel Report on the Customer Reports menu, the Tassel Report Settings dialog box opens.

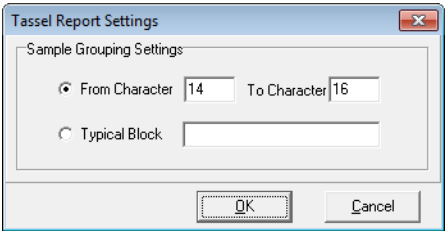Figure 7-6: Tassel Report Settings dialog box



1. Do one of the following:
 - If all the filenames in your project contain the same character string/designator in the same positions, then select From Character [] To Character [] and enter the starting and ending character positions of the characters/designator.
 - If all the filenames in your project contain the same character string /designator but in different positions, then select Typical Block and enter the character string/designator.
2. Click OK.

The Tassel report is generated.

Figure 7-7: Tassel report

No.	Sample Name	Sample ID	30607C		
1	BRCA1_012345_10A_1_10A	y		n.a.	n.a.
2	BRCA1_012345_10A_2_10A	y		n.a.	n.a.
3	BRCA1_185729_10A_1_10A	C		n.a.	n.a.
4	BRCA1_185729_10A_2_10A	C		n.a.	n.a.
5	BRCA1_264381_10A_1_10A	C		n.a.	n.a.
6	BRCA1_264381_10A_2_10A	C		n.a.	n.a.
			30747A	30753A	30988A
7	BRCA1_012345_10B_1_10B	m	w	A	
8	BRCA1_012345_10B_2_10B	m	w	A	
9	BRCA1_185729_10B_1_10B	n.a.	n.a.	R	
10	BRCA1_185729_10B_2_10B	n.a.	n.a.	R	
11	BRCA1_264381_10B_1_10B	n.a.	n.a.	R	
12	BRCA1_264381_10B_2_10B	n.a.	n.a.	R	
13	BRCA1_012345_11A_1_11A		n.a.	n.a.	n.a.
14	BRCA1_012345_11A_2_11A		n.a.	n.a.	n.a.
15	BRCA1_185729_11A_1_11A		n.a.	n.a.	n.a.
16	BRCA1_185729_11A_2_11A		n.a.	n.a.	n.a.
17	BRCA1_264381_11A_1_11A		n.a.	n.a.	n.a.
18	BRCA1_264381_11A_2_11A		n.a.	n.a.	n.a.

The layout of the Tassel report is somewhat similar to the layout of the Two Direction Output report. (See “[Two Direction report layout](#)” on page 134.) The similarities between the Tassel report and any of the other standard reports, however, ends there. In the Tassel report:


- Cells showing a novel variant that results in a change to the amino acid sequence are displayed in pink.
- Cells showing a reported variant that results in a change to the amino acid sequence are displayed in lime green.
- A lowercase letter indicates a 1D mutation, while an uppercase letter indicates a 2D mutation.
- If a mutation is a substitution, then the standard IUPAC code for the variant base or bases is displayed for the mutation.
- If a mutation is a deletion, then a zero (0) is displayed for the mutation.



You can also view the information for a mutation in the Tassel report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

Tassel report toolbar

The Tassel report toolbar contains just two buttons for working with the report.

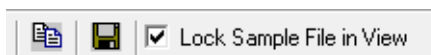Figure 7-8: Tassel report toolbar



Button	Description
	<p>Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Tassel report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
	<p>Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).</p>

Tassel report viewing and analysis options

Although the Tassel report does not have a context menu, you can still copy the report as an


image. You can also view the information for a mutation in the report in the GAD.

- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

Emory report

The layout of the Emory report is very similar to the layout of the Two Direction report. (See [“Two Direction report layout” on page 134.](#)) The mutation codes and the color coding of the mutation calls in the report are identical to the mutation codes and color coding in the Mutation report (See [“Mutation report codes” on page 129](#) and [“Color codes” on page 130.](#)) The Emory report, however, displays only the quality score and the number of mutations for each sample.

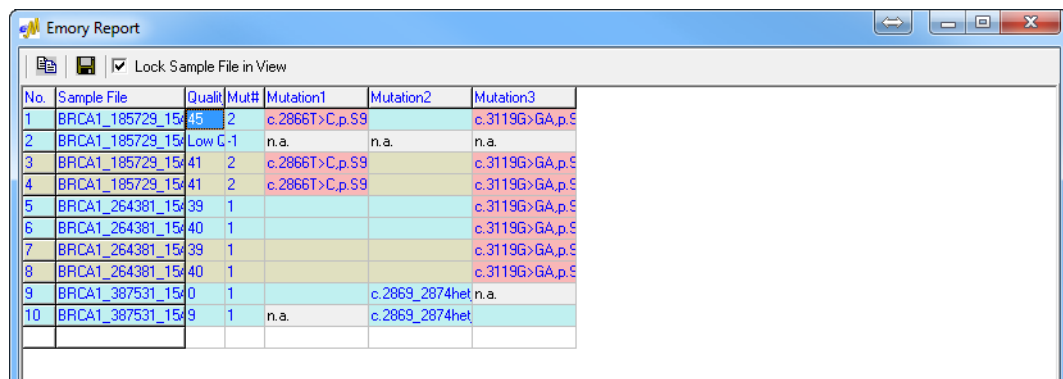Figure 7-9: Emory report






No.	Sample File	Quality	Mut#	Mutation1	Mutation2	Mutation3
1	BRCA1_185729_15	45	2	c.2866T>C.p.S9		c.3119G>GA.p.S9
2	BRCA1_185729_15	Low C-1		n.a.	n.a.	n.a.
3	BRCA1_185729_15	41	2	c.2866T>C.p.S9		c.3119G>GA.p.S9
4	BRCA1_185729_15	41	2	c.2866T>C.p.S9		c.3119G>GA.p.S9
5	BRCA1_264381_15	39	1			c.3119G>GA.p.S9
6	BRCA1_264381_15	40	1			c.3119G>GA.p.S9
7	BRCA1_264381_15	39	1			c.3119G>GA.p.S9
8	BRCA1_264381_15	40	1			c.3119G>GA.p.S9
9	BRCA1_387531_15	0	1		c.2869_2874het	n.a.
10	BRCA1_387531_15	9	1	n.a.	c.2869_2874het	

Emory report toolbar

The Emory report toolbar contains just two buttons for working with the report and an option to control the display while scrolling the report.

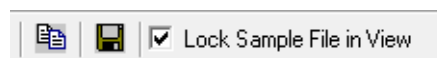Figure 7-10: Emory report toolbar




Button	Description
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
Lock Sample File in View	Freezes the Sample Filename column while scrolling.

Emory report viewing and analysis options

A context menu is available for the Emory report. You use the options on this menu to change the view of the Emory report and to manipulate and analyze mutation project data so that you identify the true mutations in a mutation project. To open the context menu, right-click anywhere in the Emory report.

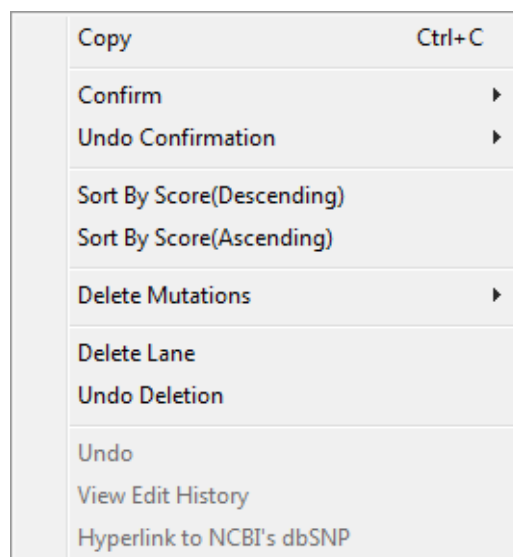Figure 7-11: Emory report context menu



Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Emory report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the Emory report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Undo Confirmation	<p>Removes the "Confirmed" status for the mutation.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p>

Option	Description
Sort By Score (Descending) Sort by Score (Ascending)	Click the appropriate option to sort all the data by descending mutation score or ascending mutation score.
Delete Mutations	Deletes the selected mutations. Note: SHIFT-click to select multiple contiguous mutations, to select all mutations in a column, or to select multiple rows. The text for the mutation code turns gray in both the Emory report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.
Delete Lane	Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to "Low Quality."
Undo Lane Deletion	The Lane Quality for the lane is reset to its original value.
Undo	Undo the <i>last</i> edit action that was carried out for the mutation.
View Edit History	See "Viewing the Edit History (Audit Trail) for a Mutation" on page 344.
Hyperlink to NCBI's dbSNP	Opens the NCBI's database entry (via the web) for the variant.

You can also view the information for a mutation in the Emory report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

Bentley report and MSM_Song report

The Bentley and MSM_Song reports are useful when carrying out analyses of diploid samples. Both reports show the genomic position of the mutation, the genotype, and the frequency of each allele by traces and as a grand total.



For a detailed explanation of allele frequency by trace and grand total allele frequency, see “[Show Allele Information](#)” on page 169.

The layout of the reports is very similar to the layout of the Two Direction Output report. (See “[Two Direction report layout](#)” on page 134.) Both of the reports contain the same fields that are displayed in any of the standard mutation reports. There are few minor difference between the Bentley report and the MSM_Song report:

- By default, the MSM_Song report displays the entire sequence of the GenBank file in the top row of the report. The Bentley does not.
- In both reports, the mutation call is split between two columns, with the first column displaying a summary for a specific position and the second column displaying the genotype for each sample file. The format, however, for these two columns differs slightly between the reports.

Figure 7-12: Bentley report

No.	Sample ID	Well Pos	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Quality	Mult	BRCA1	(Mutation 1) GENOTYPE
1	BRCA1_012345_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30524	30857	334	16	0	30607	CC
2	BRCA1_012345_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30851	349	14	1	ttcagCTGCTTG1CT	
3	BRCA1_012345_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30524	30857	334	16	0		CC
4	BRCA1_012345_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30851	349	14	1		CT
5	BRCA1_185729_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30522	30857	336	25	0		CC
6	BRCA1_185729_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	26	0		CC
7	BRCA1_185729_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30522	30857	336	25	0		CC
8	BRCA1_185729_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	26	0		CC
9	BRCA1_264381_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30512	30857	346	25	0		CC
10	BRCA1_264381_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	25	0		CC
11	BRCA1_264381_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30512	30857	346	25	0		CC
12	BRCA1_264381_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	25	0		CC
													Total by Traces:	24
													C	91.7%
													T	8.3%
													Grand Total:	12
													C	100.0%

Figure 7-13: MSM_Song report




No.	Sample ID	Well Pos	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Quality	Mult	BRCA1	(Mutation 1) GENOTYPE
1	BRCA1_012345_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30524	30857	334	16	0	30607	CC
2	BRCA1_012345_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30851	349	14	1	ttcagCTGCTTG1CT	
3	BRCA1_012345_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30524	30857	334	16	0	C/T	CC
4	BRCA1_012345_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30851	349	14	1	TGAGACGGATCCT	
5	BRCA1_185729_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30522	30857	336	25	0		CC
6	BRCA1_185729_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	26	0		CC
7	BRCA1_185729_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30522	30857	336	25	0		CC
8	BRCA1_185729_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	26	0		CC
9	BRCA1_264381_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30512	30857	346	25	0		CC
10	BRCA1_264381_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	25	0		CC
11	BRCA1_264381_10/10A		BRCA1_AmplD_F_S1-F		BRCA1		2	30512	30857	346	25	0		CC
12	BRCA1_264381_10/10A		BRCA1_AmplD_R_S1-R		BRCA1		2	30503	30852	350	25	0		CC
													Total by Traces:	24
													C	91.7%
													T	8.3%
													Grand Total:	12
													C	100.0%

Bentley/MSM_Song report toolbar

The Bentley toolbar and the MSM_Song toolbar contain the same buttons that are specific for working with the reports and for changing the information that is displayed in the reports.

Figure 7-14: Bentley/MSM_Song report toolbar



Button	Description
	<p>Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also right-click and select Copy to copy the selected text to your clipboard.</p> <p>Note: You can also copy both output reports as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
	<p>Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).</p>
	<p>Report Settings—Opens the Report Settings dialog box which contains options for filtering the information that is to be displayed in the report. See “Bentley/MSM_Song report settings” on page 198.</p>


Bentley/MSM_Song reports viewing and analysis options

The Bentley and MSM_Song reports have several options available for changing the view of the reports and for manipulating and analyzing mutation project data. These options are:

- The Bentley/MSM_Song report settings. See [“Bentley/MSM_Song report settings” on page 198](#).
- The Bentley/MSM_Song report context menu. See [“Bentley/MSM_Song reports context menu” on page 198](#).

Bentley/MSM_Song report settings

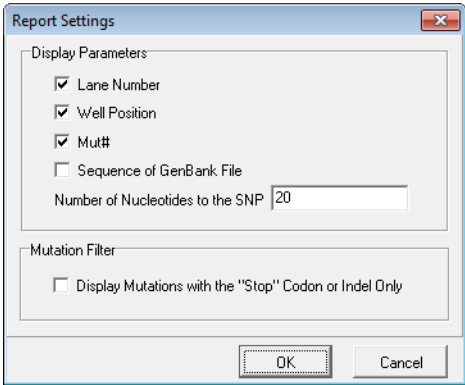Figure 7-15: Bentley/MSM_Song Report Settings dialog box



Option	Description
Display Parameters	
	<ul style="list-style-type: none">• Lane Number—Selected by default. Clear this option if you do not want the Lane Number to be displayed in the report.• Well Position—Selected by default. Clear this option if you do not want the Well Position to be displayed in the report.• Mut#—Selected by default. Clear this option if you do not want the mutation number to be displayed in the report.• Sequence of GenBank File<ul style="list-style-type: none">• Not selected by default for the Bentley Output report—Select this option if you want the entire sequence of the GenBank file to be displayed in the top row of the report.• Selected by default for the MSM_Song Output report—Clear this option if you do not want the entire sequence of the GenBank file to be displayed in the top row of the report.
Mutation Filter	
Display Mutations with the “Stop” Codon or Indel Only	Display only nonsense mutations and indels in the report.

Bentley/MSM_Song reports context menu

Although the Bentley and MSM_Song reports do not have a full context menu, you can still copy text from each report, or copy each report as an image. You can also view the information for a mutation in the reports in the GAD.

- Select the text that you want to copy or to copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Right-click and select Copy to copy the selected text to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then


click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.

- To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

PrettyBase report

The PrettyBase format was developed by the SeattleSNPs project for easily representing SNP information for groups of subjects over a given span of reference sequence. The report displays the following information (in order from left to right)—SNP Position, Sample Name, Normal Allele, Allele 1, Allele 2, IUPAC code for the mutation call, Mutation Call, Comments, Phred Score, and NM_Score.

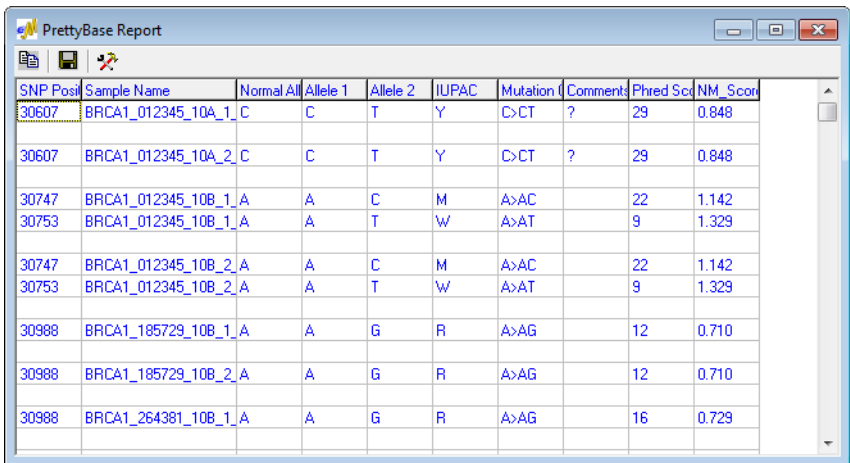Figure 7-16: PrettyBase report






SNP Pos	Sample Name	Normal Al	Allele 1	Allele 2	IUPAC	Mutation C	Comments	Phred Sc	NM_Score
30607	BRCA1_012345_10A_1_C	C	C	T	Y	C>CT	?	29	0.848
30607	BRCA1_012345_10A_2_C	C	C	T	Y	C>CT	?	29	0.848
30747	BRCA1_012345_10B_1_A	A	A	C	M	A>AC		22	1.142
30753	BRCA1_012345_10B_1_A	A	A	T	W	A>AT		9	1.329
30747	BRCA1_012345_10B_2_A	A	A	C	M	A>AC		22	1.142
30753	BRCA1_012345_10B_2_A	A	A	T	W	A>AT		9	1.329
30988	BRCA1_185729_10B_1_A	A	A	G	R	A>AG		12	0.710
30988	BRCA1_185729_10B_2_A	A	A	G	R	A>AG		12	0.710
30988	BRCA1_264381_10B_1_A	A	A	G	R	A>AG		16	0.729

PrettyBase report toolbar

The PrettyBase report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 7-17: PrettyBase report toolbar



Button	Description
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
	Report Settings—Opens the PrettyBase Report Settings dialog box which contains options for filtering the information that is to be displayed in the report. See “PrettyBase report settings” on page 201 .

PrettyBase report viewing and analysis options

The PrettyBase report has multiple options available for changing the view of the report and for manipulating and analyzing mutation project data. These options are:

- The PrettyBase report settings. See [“PrettyBase report settings”](#) below.
- The PrettyBase report context menu. See [“PrettyBase report context menu”](#) on page 202.

PrettyBase report settings


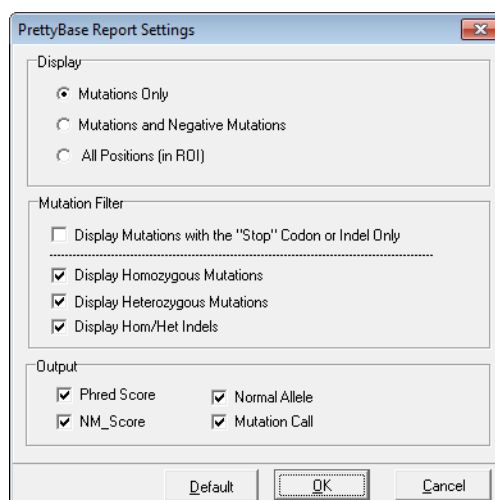
To open the PrettyBase Report Settings dialog box, click the Report Settings button  on the PrettyBase report toolbar.

Figure 7-18: PrettyBase Report Settings dialog box



Option	Description
Display	
Mutations Only	<ul style="list-style-type: none"> • Mutations Only—Selected by default. Show only all called mutations • Mutations and Negative Mutations—Select this option to show all called mutations and negative mutations. • All Positions (in ROI)—Select this option to show all nucleotide positions within defined ROIs.
Mutation Filter	
Display Mutations with the “Stop” Codon or Indel Only	Display only nonsense mutations and indels in the report.

Option	Description
	<ul style="list-style-type: none"> • Display Homozygous Mutations—Selected by default. Clear this option if you do not want to show homozygous mutations in the report. • Display Heterozygous Mutations—Selected by default. Clear this option if you do not want to show heterozygous mutations in the report. • Display Hom/Het Indels—Selected by default. Clear this option if you do not want to show homozygous and heterozygous indels in the report.
Output	
	<ul style="list-style-type: none"> • Phred Score • NM_Score • Normal Allele • Mutation Call <p>All four of these columns are displayed by default in the PrettyBase Output report. Clear any or all options that you do want to be displayed in the report.</p> <p>Note: For information about the NM_Score, see “NM_Score” on page 174.</p>

PrettyBase report context menu

Although the PrettyBase report does not have a full context menu, you can still copy text from the report, or copy the report as an image. You can also view the information for a mutation in the report in the GAD.

- Select the text that you want to copy or to copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Right-click and select Copy to copy the selected text to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

Methylation report

How Mutation Surveyor calls mutations for methylated DNA assumes 100% conversion of the non-methylated cytosines to uracil during bisulfite treatment. The evaluation of the methylation level in CpG islands as well as the level of incomplete conversion of unmethylated cytosines, however, is critical for ensuring that your methylation project has the necessary high quality data that is required for accurate mutation calls. The Methylation report displays information for methylation detection projects. The report distinguishes between cytosines that are contained in CpG islands and cytosines that are contained in other sequences. The report also displays a Conversion ratio, which is the ratio of the number of samples that were successfully converted at a given position to the total number of samples with the position. A Conversion Ratio of 100% indicates that the conversion was successful for all samples at the given position.

Figure 7-19: Methylation report


			5505_V578	7373_V578	7708_V578	5505_V584	7373_V584	7708_V584
58040	C	t	S					
58041	T	t	-					
58042	C	c	M					
58043	G	g	-					
58044	G	g	-					
58045	C	t	S	S	S			
58046	C	t	S	S	S			
58047	A	a	-	-	-			
58048	G	g	-	-	-			
58049	C	t	S	S	S			
58050	A	a	-	-	-			
58051	G	g	-	-	-			
58052	G	g	-	-	-			
58053	A	a	-	-	-			
58054	T	t	-	-	-			
58055	C	t	S	S	S			
58056	T	t	-	-	-			
58057	T	t	-	-	-			
58058	A	a	-	-	-			
58059	T	t	-	-	-			
58060	T	t	-	-	-			
58061	C	t	S	S	S			
58062	T	t	-	-	-			
58063	G	g	-	-	-			
58064	C	t	S	S	S			
58065	T	t	-	-	-			
58066	G	g	-	-	-			



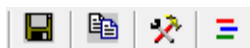
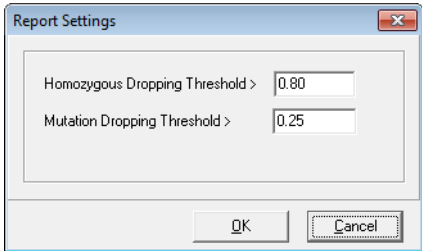

Code	Description
Cytosines contained in CpG islands	<ul style="list-style-type: none"> M—Methylated U—Non-methylated
Cytosines contained in other sequences	<ul style="list-style-type: none"> S—Successful conversion of cytosine (C) to thymine (T) I—Incomplete conversion

Methylation report toolbar

The Methylation report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 7-20: Methylation report toolbar



Button	Description
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as a tab-delimited .txt file.
	Copy—Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
	<p>Report Settings—Opens the Methylation Report Settings dialog box which contains options for filtering the information that is to be displayed in the report.</p> <p><i>Figure 7-21: Methylation Report Settings dialog box</i></p>  <ul style="list-style-type: none"> • Homozygous Dropping Threshold—Indicates the value for the dropping factor at which a mutation is identified as either a homozygous substitution or a heterozygous substitution. For example, is a mutation A>C or A>AC? Because Mutation Surveyor evaluates the traces rather than the base calls, noise below a major peak can sometimes be interpreted as a heterozygous substitution. This value assists in your analysis by converting these calls to homozygous substitutions. <p>Note: You can increase this value to view the exact call that Mutation Surveyor made, but an increased value might result in some calls remaining as heterozygous substitutions.</p> <ul style="list-style-type: none"> • Mutation Dropping Threshold—The dropping threshold must be greater than or equal to the specified value for the variant to be called a mutation. See “Dropping Factor” on page 60.
	Graphic Display—Opens the Graphic Display of Methylation report. See “Graphic Display of Methylation report” on page 205 .

Graphic Display of Methylation report

The Graphic Display of Methylation report graphically shows the pattern of methylated and non-methylated cytosines along the DNA fragments used for analysis. By default, when the report first opens, only cytosines that are part of a CpG island are indicated in the report. Red tick marks indicate methylated sites. Dark blue tick marks indicate non-methylated sites. The Methylation Ratio graph represents the percentage of samples that are methylated at the indicated position in a CpG island. The Conversion Ratio graph represents the percentage of samples that were converted at the indicated position outside of a CpG island.

Figure 7-22: Graphic Display of Methylation report



The report is interactive:

- To view the cytosines that are contained in other sequences, click Display S/I. Light blue tick marks indicate successful conversion of cytosine (C) to uracil (U). Gray tick marks indicate incomplete conversion.
- To zoom in on a selected region, click and hold the left mouse button and draw a box from the upper left hand corner of the region towards the lower right hand corner of the region. A box is formed around the area that is being reduced for viewing.
- To zoom out from any region, click and hold the left mouse button and draw a box from the lower right hand corner of the region towards the upper left hand corner of the region. The magnification for zooming out is always 100%.
- If your mutation project used multiple GenBank files, then select one of these files from the Select GenBank file to view the display for the file and the sample files that align to it.
- By default, the filenames for all the files (GenBank, reference files, and sample files) are

displayed on the report. To remove the filenames from the report display, clear the Show Filenames option.

Methylation report context menu


Although the Methylation report does not have a full context menu, you can still copy text from the report, or copy the report as an image. You can also view the information for a mutation in the report in the GAD.

- Select the text that you want to copy or to copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Right-click and select Copy to copy the selected text to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To open the GAD for the trace that contains a mutation, double-click the mutation in the report.


Vertex report

The Vertex report displays mutations in all the sample sequences for a selected contig. By default, the mutations are shown as the 1 Letter IUPAC representation of the normal amino acid that is affected by the mutation and the resulting amino acid. Also by default, the entire sequence of the GenBank file is displayed at the top of the report.

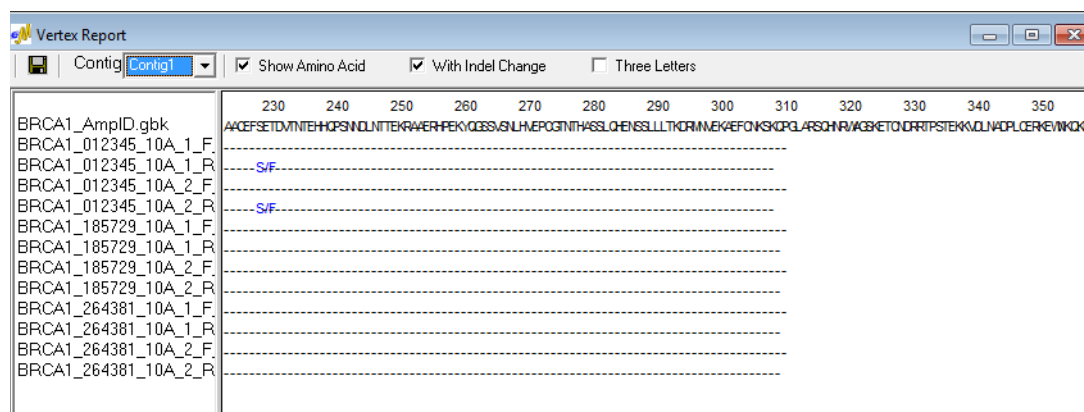Figure 7-23: Vertex report



The report is interactive:

- To view the amino acid sequences for the sample traces in a different contig, select the contig from the Select Contig list.
- To display the 3 Letter IUPAC representation for the amino acids, select Three Letter.
- To toggle between the amino acids and the called bases for the mutation, clear/select Show Amino Acid. (Like the amino acid codes, the called bases are represented by their IUPAC codes.)
- To zoom in on a selected region of the GenBank file, click and hold the left mouse button and draw a box from the upper left hand corner of the region towards the lower right hand corner of the region. A box is formed around the area that is being reduced for viewing.
- To zoom out from the GenBank file, click and hold the left mouse button and draw a box from the lower right hand corner of the sequence towards the upper left hand corner of the sequence. The magnification for zooming out is always 100%.
- To save the report as an .xls/.xlsx or .txt file, click the Save Report icon  on the report toolbar. A Save Report As dialog box opens in which you can specify the name, format and location for the saved report. (The .txt format is available only when saving the nucleotide sequence for the report. The .xls/.xlsx format is available only when saving the amino acid sequence for the report.)
- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to

Chapter 7


Mutation Surveyor Customer Reports

copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.

Vertex2 report

You use the Vertex2 report to assemble variations from multiple sequences or samples within one or multiple projects. The Vertex2 report is especially useful for analyzing hypervariable sequences. After you click Vertex2 Report on the Customer Reports menu, the Vertex Report Settings dialog box opens. You can leave the default values as is, or you can modify any of the values as needed, and then click OK to generate the report.

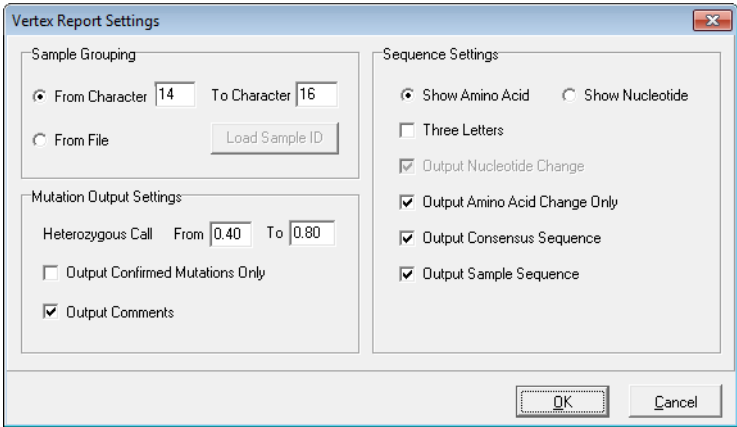Figure 7-24: Vertex Report Settings dialog box




Option	Description
Sample Grouping —Determines how traces are to be grouped in the report.	
	<ul style="list-style-type: none">From Character [] To Character []—Select this option if all the filenames in your project contain the same character string/designator in the same positions.From File—Select this option if traces that are to be grouped together contain a common Sample ID in their filenames. (The Sample IDs do not need to be in the same location in the filename.) After you click From File, click Load Sample ID to open the Load Sample ID dialog box and specify the different Sample IDs, such as Patient #, by which to group the traces. You can also click Open in the Load Sample ID dialog box to load an existing Sample ID text file. <p>Note: By default, the Load Sample ID dialog box contains rows for up to ten different Sample IDs. You can right-click and click Insert or Delete as needed.</p> <p>Figure 7-25: Load Sample ID dialog box</p>

Chapter 7
Mutation Surveyor Customer Reports

Option	Description
Mutation Output Settings	
Heterozygous Call From [] To []	A mutation call with a dropping factor that falls within the specified range is classified as a heterozygous mutation call. It is classified as a homozygous normal allele if its dropping factor falls below the lower threshold and as a homozygous mutation if its dropping factor exceeds the upper threshold.
Output Confirmed Mutations Only	Display only manually confirmed mutations in the report.
Output Comments	Display automatically added and manually added comments in the report.
Sequence Settings	
	<ul style="list-style-type: none"> • Show Amino Acid—Show the 1 Letter IUPAC representation of the normal amino acid that is affected by the mutation and the resulting amino acid. • Show Nucleotide—Show the IUPAC codes for the called bases for the mutation. • Three Letters—Available only if Show Amino Acid is selected. Show the 3 Letter representation of the normal amino acid that is affected by the mutation and the resulting amino acid. • Output Nucleotide Change—Display only those traces with called mutations in the report. • Output Amino Acid Change Only—Display only those traces with called mutations in the report, and for these traces, display the 1 Letter or 3 Letter IUPAC representation of the normal amino acid that is affected by the mutation and the resulting amino acid. • Output Consensus Sequence—Display the consensus sequence for all sample traces in the report. • Output Sample Sequences—Display the sequence for each individual sample trace in the report.

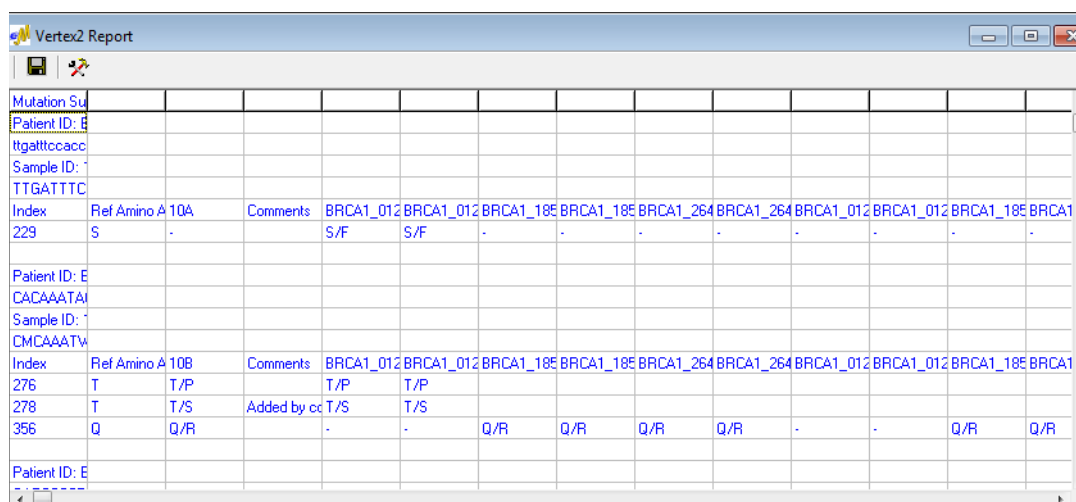Figure 7-26: Vertex2 report



The screenshot shows a window titled "Vertex2 Report" with a table of mutation data. The table has columns for Index, Ref Amino A, Comments, and multiple columns for BRCA1 variants (BRCA1_012, BRCA1_012, BRCA1_185, BRCA1_185, BRCA1_264, BRCA1_264, BRCA1_012, BRCA1_012, BRCA1_185, BRCA1_185). The data is organized into sections for different patient IDs and sample IDs.



Index	Ref Amino A	Comments	BRCA1_012	BRCA1_012	BRCA1_185	BRCA1_185	BRCA1_264	BRCA1_264	BRCA1_012	BRCA1_012	BRCA1_185	BRCA1_185
229	S	-	S/F	S/F	-	-	-	-	-	-	-	-
276	T	T/P	T/P	T/P	-	-	-	-	-	-	-	-
278	T	T/S	T/S	T/S	-	-	-	-	-	-	-	-
356	Q	Q/R	-	-	Q/R	Q/R	Q/R	Q/R	-	-	Q/R	Q/R

Vertex2 report toolbar

The Vertex2 report toolbar contains just two buttons for working with the report.

Figure 7-27: Vertex2 report toolbar



Button	Description
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.xls/.xlsx or .txt).
	Report Settings—Opens the Vertex2 Report Settings dialog box which contains options for grouping and filtering the information that is to be displayed in the report. See "Vertex2 report" on page 209 .

Vertex2 report viewing and analysis options


Although the Vertex2 report does not have a context menu, you can copy the report as an image. You can also view the information for a mutation in the report in the GAD.

- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

Vertex3 report

Like the Vertex2 report, you can use the Vertex3 report to assemble variations from multiple sequences or samples within one or multiple projects. After you click Vertex3 Report on the Customer Reports menu, the Vertex3 Report Settings dialog box opens. You can leave the default values as they are, or you can modify any of the values as needed, and then click OK to generate the report.

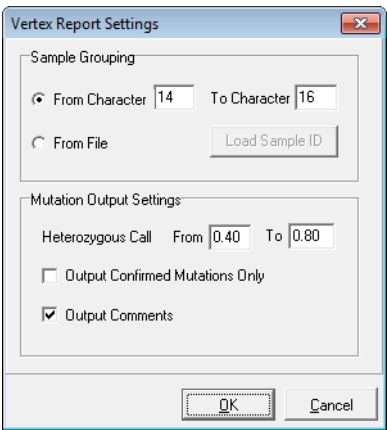Figure 7-28: Vertex Report Settings dialog box



Option	Description
Sample Grouping —Determines how traces are to be grouped in the report.	
	<ul style="list-style-type: none">From Character [] To Character []—Select this option if all the filenames in your project contain the same character string/designator in the same positions.From File—Select this option if traces that are to be grouped together contain a common Sample ID in their filenames. (The Sample IDs do not need to be in the same location in the filename.) After you click From File, click Load Sample ID to open the Load Sample ID dialog box and specify the different Sample IDs, such as Patient #, by which to group the traces. You can also click Open in the Load Sample ID dialog box to load an existing Sample ID text file. <p>Note: By default, the Load Sample ID dialog box contains rows for up to ten different Sample IDs. You can right-click and click Insert or Delete as needed.</p> <p>Figure 7-29: Load Sample ID dialog box</p>

Option	Description
Mutation Output Settings	
Heterozygous Call From [] To []	A mutation call with a dropping factor that falls within the specified range is classified as a heterozygous mutation call. It is classified as a homozygous normal allele if its dropping factor falls below the lower threshold and as a homozygous mutation if its dropping factor exceeds the upper threshold.
Output Confirmed Mutations Only	Display only manually confirmed mutations in the report.

Vertex3 report layout


The Vertex3 report is organized by Clone (Sample) ID and contains three sections for each Sample ID.

Section	Description
Nucleotide Sequences	Displays the reference nucleotide sequence for the samples and the consensus nucleotide sequence for the samples.
Nucleotide Changes	Displays a list of each location that contains a variation between the reference nucleotide and the sample nucleotide, the reference nucleotide at the location, the sample nucleotide at the location and a column denoting whether the variant is synonymous (S) or non-synonymous (N).
Amino Acid	Displays the codon number, the reference amino acid and the consensus amino acid for each non-synonymous variation.

Figure 7-30: Vertex3 report

[illegible]

Vertex3 report toolbar

The Vertex3 Report toolbar contains a single button on the toolbar—the Save Report button . Click this button to open a Save Report As dialog box in which you can specify the name and location for saving the report as a .txt file.

Vertex3 report viewing and analysis options

Although the Vertex3 report does not have a full context menu, you can still copy text from the report, or copy the report as an image.

- Select the text that you want to copy or to copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Right-click and select Copy to copy the selected text to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.

Full Gene report



Of all the Customer reports, the Full Gene report is the most customizable report. It has options for:

- Organizing sample traces into groups based on matching filenames and then further sort the sample traces within the groups based on specific character positions in the filenames.
- Custom color-coding to assist you with identifying different types of mutation calls and potential analysis problems.
- Creating and displaying a custom header above each group of sample traces.

After you click Full Gene Report on the Customer Reports menu, the Full Gene Report Settings dialog box opens. You can leave the default values as they are, or you can modify any of the values as needed, and then click OK to generate the report. See:


- “Full Gene Report Settings dialog box, Format tab.”
- “Full Gene Report Settings, Header tab” on page 217.
- “Full Gene Report Settings, Output tab” on page 217.

Figure 7-31: Full Gene report

Full Gene Report																																																																					
<div><div></div><div><table><tr><td>Software</td><td>Mutation Surveyor</td><td>Date</td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td></tr><tr><td>Date/Time</td><td>5/24/2010 2:01</td><td>Batch/Run ID</td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td></tr><tr><td>Technologist</td><td></td><td>Result</td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td></tr><tr><td>Sample ID</td><td>012345</td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td><td></td></tr></table></div></div>														Software	Mutation Surveyor	Date												Date/Time	5/24/2010 2:01	Batch/Run ID												Technologist		Result												Sample ID	012345												
Software	Mutation Surveyor	Date																																																																			
Date/Time	5/24/2010 2:01	Batch/Run ID																																																																			
Technologist		Result																																																																			
Sample ID	012345																																																																				
No.	Sample File	Reference File	Dir	Gene	Exon	RF	Start	End	Size	Quality	Variant1	Variant2	Variant3	Variant4																																																							
1	BRCA1_012345	BRCA1_AmpD_1-F		BRCA1		2	30524	30957	334	16		n.a.	n.a.	n.a.																																																							
2	BRCA1_012345	BRCA1_AmpD_1-R		BRCA1		2	30503	30951	349	14		n.a.	n.a.	n.a.																																																							
3	BRCA1_012345	BRCA1_AmpD_1-F		BRCA1		2	30524	30957	334	16		n.a.	n.a.	n.a.																																																							
4	BRCA1_012345	BRCA1_AmpD_1-R		BRCA1		2	30503	30951	349	14		n.a.	n.a.	n.a.																																																							
5	BRCA1_012345	BRCA1_AmpD_2-F		BRCA1		2	30768	31039	272	8	n.a.	n.a.	n.a.	n.a.																																																							
6	BRCA1_012345	BRCA1_AmpD_2-R		BRCA1		2	30746	31027	282	11	c.1496C>T	n.a.	n.a.	n.a.																																																							
7	BRCA1_012345	BRCA1_AmpD_2-F		BRCA1		2	30768	31039	272	8	n.a.	n.a.	n.a.	n.a.																																																							
8	BRCA1_012345	BRCA1_AmpD_2-R		BRCA1		2	30746	31027	282	11	c.1496C>T	n.a.	n.a.	n.a.																																																							
9	BRCA1_012345	BRCA1_AmpD_3-F		BRCA1		2	30977	31416	440	7	n.a.																																																										
10	BRCA1_012345	BRCA1_AmpD_3-R		BRCA1		2	30958	31414	457	6	n.a.																																																										
11	BRCA1_012345	BRCA1_AmpD_3-F		BRCA1		2	30977	31416	440	7	n.a.																																																										
12	BRCA1_012345	BRCA1_AmpD_3-R		BRCA1		2	30958	31414	457	6	n.a.																																																										
13	BRCA1_012345	BRCA1_AmpD_4-F		BRCA1		2	31143	31450	318	0	n.a.		c.1498C>CT.p.R																																																								
14	BRCA1_012345	BRCA1_AmpD_4-R		BRCA1		2	31114	31449	336	11	n.a.			c.1498C>CT.p.R																																																							
15	BRCA1_012345	BRCA1_AmpD_4-F		BRCA1		2	31143	31450	318	0	n.a.	c.1498C>CT.p.R	c.1498C>CT.p.R																																																								
16	BRCA1_012345	BRCA1_AmpD_4-R		BRCA1		2	31114	31449	336	11	n.a.			c.1498C>CT.p.R																																																							
17	BRCA1_012345	BRCA1_AmpD_5-F		BRCA1		2	31419	31595	177	0	n.a.	n.a.	n.a.	n.a.																																																							
18	BRCA1_012345	BRCA1_AmpD_5-R		BRCA1		2	31397	31591	195	17	n.a.	n.a.																																																									
19	BRCA1_012345	BRCA1_AmpD_5-F		BRCA1		2	31419	31595	177	0	n.a.	n.a.	n.a.	n.a.																																																							
20	BRCA1_012345	BRCA1_AmpD_5-R		BRCA1		2	31397	31591	195	17	n.a.	n.a.																																																									
21	BRCA1_012345	BRCA1_AmpD_6-F		BRCA1		2	31540	31902	363	10	n.a.	n.a.	n.a.	n.a.																																																							
22	BRCA1_012345	BRCA1_AmpD_6-R		BRCA1		2	31516	31896	381	11	n.a.	n.a.	n.a.	n.a.																																																							
23	BRCA1_012345	BRCA1_AmpD_6-F		BRCA1		2	31540	31902	363	10	n.a.	n.a.	n.a.	n.a.																																																							

Full Gene Report Settings dialog box, Format tab

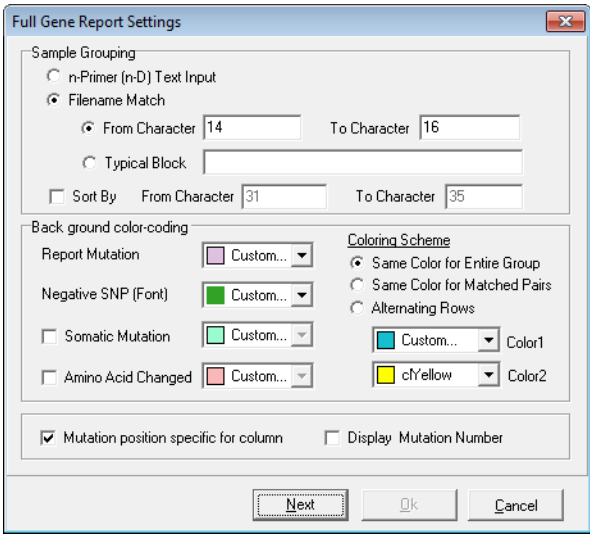Figure 7-32: Full Gene Report Settings dialog box, Format tab




Option	Description
Sample Grouping	
n-Primer (n-D) Text input	Select this option if you have defined primers when using a GenBank reference file to group sample files in forward and reverse pairs.
Filename match	<ul style="list-style-type: none">From Character [] To Character []—Select this option if all the filenames in your project contain the same character string/designator in the same positions.Typical Block—If all the filenames in your project contain the same character string/designators but in different positions, then select this option and enter the character string/designator.
Sort By	From Character [] To Character []—Select this option to further sort sample traces within a group based on specific character positions in the filenames.
Background Color-Coding	
Accept the default colors for the different types of mutations, or modify them as you want.	
	<ul style="list-style-type: none">Mutation position specific for column—Aligns mutations by column within each group.Display Mutation Number—Display the number of mutations found in the sample trace.

Full Gene Report Settings, Header tab

The Header tab opens after you click Next on the Format tab.

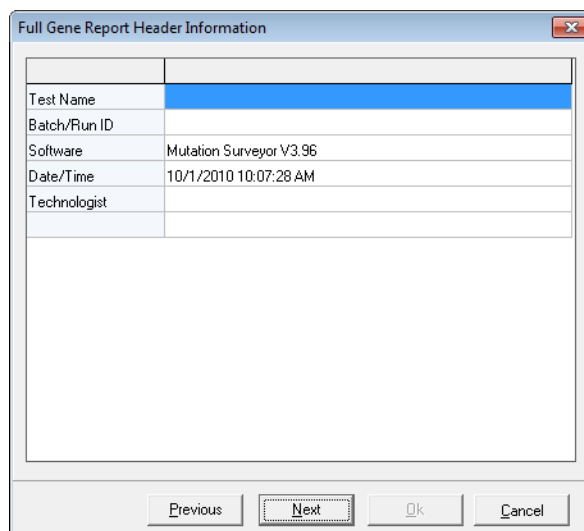Figure 7-33: Full Gene Settings dialog box, Header information



The dialog box titled "Full Gene Report Header Information" contains a table with the following fields:


Test Name	
Batch/Run ID	
Software	Mutation Surveyor V3.96
Date/Time	10/1/2010 10:07:28 AM
Technologist	

At the bottom of the dialog box are four buttons: Previous, Next, Ok, and Cancel. The "Next" button is highlighted with a blue border.

Full Gene Report Settings, Output tab

The Output tab opens after you click Next on the Header tab.

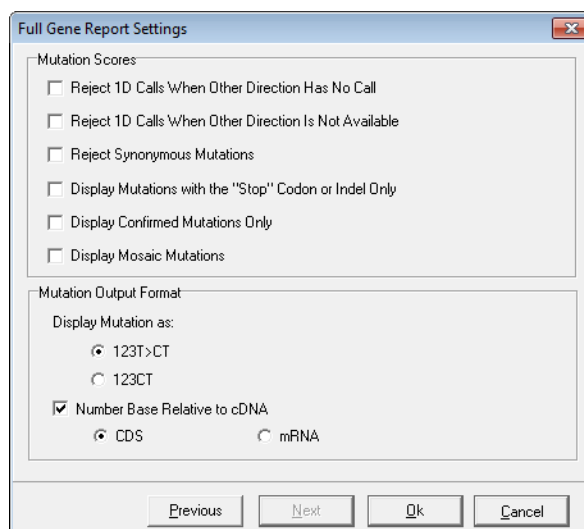Figure 7-34: Full Gene Report Settings dialog box, Mutation Score and Output options



The dialog box titled "Full Gene Report Settings" contains two sections:

Mutation Scores

- ☐ Reject 1D Calls When Other Direction Has No Call
- ☐ Reject 1D Calls When Other Direction Is Not Available
- ☐ Reject Synonymous Mutations
- ☐ Display Mutations with the "Stop" Codon or Indel Only
- ☐ Display Confirmed Mutations Only
- ☐ Display Mosaic Mutations

Mutation Output Format

Display Mutation as:

- ☒ 123T>CT
- ☐ 123CT

☒ Number Base Relative to cDNA

- ☒ CDS
- ☐ mRNA

At the bottom of the dialog box are four buttons: Previous, Next, Ok, and Cancel. The "Next" button is highlighted with a blue border.

Chapter 7
Mutation Surveyor Customer Reports

Option	Description
Mutation Scores	
Reject 1D Calls When Other Direction has No Call	Reject a mutation that is called in only one direction when the position falls within the comparison region of the other direction.
Reject 1D Calls When Other Direction is Not Available	Reject a mutation that is called in only one direction when the position falls outside the comparison region of the other direction.
Reject Synonymous Mutations	Coding region substitutions that do not affect amino acid calls are not displayed in the report.
Note: An "n.a." is displayed in the Mutation cell for rejected mutations instead of the mutation code.	
Display Mutations with the "Stop" Codon or Indel Only	Display only nonsense mutations and indels in the report.
Display Confirmed Mutations Only	Display only those mutations that you have manually confirmed in the report.
Display Mosaic Mutations	Display mosaic mutations in the report.
Mutation Output Format	
Display Mutation As	Select one option to set the format of the mutation code in the report. <ul style="list-style-type: none"> • 123T>CT (Default value) • 123CT
Number Bases Relative to cDNA	<ul style="list-style-type: none"> • If this option is cleared, the numbering is the genomic position within the GenBank file. • If this option is selected: <ul style="list-style-type: none"> • Select CDS to number the bases relative to the CDS numbering in the GenBank reference file. • Select mRNA to number the bases relative the mRNA numbering in the GenBank reference file.

Figure 7-35: Full Gene report



Full Gene Report														
	Software	Mutation Surveyor		Date										
	Date/Time	5/24/2010 2:01		Batch/Run ID										
	Technologist			Result										
	Sample ID	012345												
No.	Sample File	Reference File	Dr	Gene	Exon	RF	Start	End	Size	Quality	Variant1	Variant2	Variant3	Variant4
1	BRCA1_012345	BRCA1_AmpID_1F		BRCA1		2	30524	30557	334	15		n.a.	n.a.	n.a.
2	BRCA1_012345	BRCA1_AmpID_1R		BRCA1		2	30503	30551	349	14		n.a.	n.a.	n.a.
3	BRCA1_012345	BRCA1_AmpID_1F		BRCA1		2	30524	30557	334	15			n.a.	n.a.
4	BRCA1_012345	BRCA1_AmpID_1R		BRCA1		2	30503	30551	349	14			n.a.	n.a.
5	BRCA1_012345	BRCA1_AmpID_2F		BRCA1		2	30769	31039	272	8		n.a.	n.a.	n.a.
6	BRCA1_012345	BRCA1_AmpID_2R		BRCA1		2	30746	31027	282	11		n.a.	n.a.	n.a.
7	BRCA1_012345	BRCA1_AmpID_2F		BRCA1		2	30769	31039	272	8		n.a.	n.a.	n.a.
8	BRCA1_012345	BRCA1_AmpID_2R		BRCA1		2	30746	31027	282	11		n.a.	n.a.	n.a.
9	BRCA1_012345	BRCA1_AmpID_3F		BRCA1		2	30977	31416	440	7		n.a.		
10	BRCA1_012345	BRCA1_AmpID_3R		BRCA1		2	30958	31414	457	6		n.a.		
11	BRCA1_012345	BRCA1_AmpID_3F		BRCA1		2	30977	31416	440	7		n.a.		
12	BRCA1_012345	BRCA1_AmpID_3R		BRCA1		2	30958	31414	457	6		n.a.		
13	BRCA1_012345	BRCA1_AmpID_4F		BRCA1		2	31143	31450	318	0		n.a.		c.1498C>CT p.R
14	BRCA1_012345	BRCA1_AmpID_4R		BRCA1		2	31114	31449	336	11		n.a.		c.1498C>CT p.R
15	BRCA1_012345	BRCA1_AmpID_4F		BRCA1		2	31143	31450	318	0		n.a.		c.1498C>CT p.R
16	BRCA1_012345	BRCA1_AmpID_4R		BRCA1		2	31114	31449	336	11		n.a.		c.1498C>CT p.R
17	BRCA1_012345	BRCA1_AmpID_5F		BRCA1		2	31419	31595	177	0		n.a.	n.a.	n.a.
18	BRCA1_012345	BRCA1_AmpID_5R		BRCA1		2	31397	31591	195	17		n.a.	n.a.	n.a.
19	BRCA1_012345	BRCA1_AmpID_5F		BRCA1		2	31419	31595	177	0		n.a.	n.a.	n.a.
20	BRCA1_012345	BRCA1_AmpID_5R		BRCA1		2	31397	31591	195	17		n.a.	n.a.	n.a.
21	BRCA1_012345	BRCA1_AmpID_6F		BRCA1		2	31540	31902	363	10		n.a.	n.a.	n.a.
22	BRCA1_012345	BRCA1_AmpID_6R		BRCA1		2	31516	31936	381	11		n.a.	n.a.	n.a.
23	BRCA1_012345	BRCA1_AmpID_6F		BRCA1		2	31540	31902	363	10		n.a.	n.a.	n.a.

Full Gene report toolbar

The Full Gene report toolbar contains just two buttons for working with the report.

Figure 7-36: Full Gene report toolbar




Button	Description
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as well as its format (.xls/.xlsx, .xls/.xlsx (Excel sample in different sheet) or .txt).
	Report Settings—Opens the Full Gene Report Settings dialog box which contains options for grouping and filtering the information that is to be displayed in the report. Note: The Sample Grouping options are not available on the Format tab when you open the Full Gene Report Settings dialog box from the report toolbar. This prevents you from changing the report format after you have added custom text. To change the report format, close and then re-open the report. See “Full Gene Report Settings dialog box, Format tab,” “Full Gene Report Settings, Header tab,” and “Full Gene Report Settings, Output tab.”

Full Gene viewing and analysis options

A context menu that contains options for changing the view of the Full Gene report and for manipulating and analyzing mutation project data. To open the context menu, right-click anywhere in the Full Gene report.

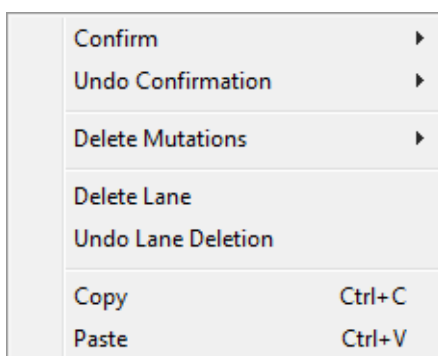Figure 7-37: Full Gene report context menu



Option	Description
Confirm	<p>Confirms the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p> <p>The text for the mutation code turns black in both the Full Gene report and the Mutation table in the GAD. The comment <<Checked>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Undo Confirmation	<p>Removes the “Confirmed” status for the mutation.</p> <p>Note: SHIFT-click to select multiple contiguous mutations or to select all mutations in a column.</p>
Delete Mutations	<p>Deletes the selected mutations.</p> <p>Note: SHIFT-click to select multiple contiguous mutations, to select all mutations in a column, or to select multiple rows.</p> <p>The text for the mutation code turns gray in both the Full Gene report and the Mutation table in the GAD. The comment <<Deleted>> is displayed in the Actions field for the mutation call in the Mutation table in the GAD.</p>
Delete Lane	<p>Deletes all the mutations for a selected sample in a single step. The Lane Quality for a deleted lane is set to “Low Quality.”</p>
Undo Lane Deletion	<p>The Lane Quality for the lane is reset to its original value.</p>

Option	Description
Copy	<p>Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.</p> <p>Note: You can also copy the Full Gene report as an image. Press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.</p>
Paste	Pastes copied text into an editable field (Date, Batch Run/ID or/or Result fields).

The Full Gene report is interactive.

- You can double click in the Date, Batch Run/ID or/or Result fields and enter any text as needed.
- You can also view information for a mutation in the report in the GAD. To open the GAD for the trace that contains a mutation, double-click the mutation in the report.

FullBase report

The FullBase report shows the forward and reverse consensus sequences for each base within specified ROIs. Amplicon IDs must be defined in GenBank or SEQ reference files and these IDs must be present in all the respective sample filenames. The Full Base report groups replicated positions into a single cell. After you click FullBase Report on the Customer Reports menu, the FullBase Report Settings dialog box opens. You can leave the default values as they are, or you can modify any of the values as needed, and then click OK to generate the report.



You use GenBank File Editors to define Amplicon IDs. See [Chapter 8, “GenBank File Editors,”](#) on page 225 for information about defining Amplicon IDs.

Figure 7-38: FullBase Report Settings dialog box

Option	Description
Amplicon ID from Character [] To Character []	Displays the starting and ending positions of the amplicon-specific starting designator in your reference filenames.
Sample ID from Character [] To Character []	Enter the starting and ending positions for the range in your sample filenames that contain the amplicon-specific designator.
Expected Number of Reads per Group	The expected total number of traces (forward and reverse) to pass the discrepancy check for a group of sample files. Note: More traces can be in the consensus without being highlighted with a yellow background.
Analyst 1	Optional information to further clarify your analysis setup.
Analyst 2	Optional information to further clarify your analysis setup.
Reviewer	Optional information to further clarify your analysis setup.

FullBase report layout


Mutation calls are displayed in cells where a variation has been identified with the font color indicating the call confidence. Forward and reverse samples traces are paired next to each other with alternating colors of light tan and light blue differentiating the sample pairs. Positions that do not have enough samples for determining mutation calls are highlighted in yellow. *Discrepancies* (identifies the variants within a consensus of traces that have at least one member of the consensus that differs from the others) are also highlighted in yellow. The Comments field is a free text field for entering additional information about your project.

Figure 7-39: Full Base report

The screenshot shows the 'FullBase Report' window. It contains a metadata section at the top with fields for Project, Software, Report Date/Time, Processing Settings, and Expected number of samples. Below this is a section for Analysts with fields for Printed Name, Signature, and Signature Date. The main part of the report is a large table with columns for Nucleotide Position, Reference Sequence, Amplicon ID, and multiple columns for different sample pairs (e.g., BRCA1_012345_10, BRCA1_012345_11). The table displays mutation calls with varying font colors (black, red, green, blue) and background colors (light tan, light blue, yellow) to indicate confidence and discrepancies. A toolbar at the bottom left includes a 'Save Report' button.

Project: BRCA1_All_files.SGF		Comments:					
Software: Mutation Surveyor V							
Report Date/Time: 09/03/2010 14:17:5							
Processing Settings: MutationSurveyor.ini							
Expected number of	3	Analyst 1	Analyst 2				
Analyst 1:		Printed Name					
Analyst 2:		Signature					
		Signature Date					
Nucleotide Position	Reference Sequence	Amplicon ID	BRCA1_012345_10	BRCA1_012345_10	BRCA1_012345_10	BRCA1_012345_10	BRCA1_012345_11
30561	t	10A	t:34;...;t:52;...	t:34;...;t:52;...
30562	g	10A	g:34;...;g:52;...	g:34;...;g:52;...
30563	a	10A	a:34;...;a:52;...	a:34;...;a:52;...
30564	g	10A	g:34;...;g:52;...	g:34;...;g:52;...
30565	t	10A	t:38;...;t:50;...	t:38;...;t:50;...
30566	a	10A	a:44;...;a:50;...	a:44;...;a:50;...
30567	c	10A	c:44;...;c:50;...	c:44;...;c:50;...
30568	c	10A	c:44;...;c:52;...	c:44;...;c:52;...
30569	t	10A	t:44;...;t:52;...	t:44;...;t:52;...
30570	t	10A	t:53;...;t:52;...	t:53;...;t:52;...
30571	g	10A	g:38;...;g:53;...	g:38;...;g:53;...
30572	t	10A	t:44;...;t:53;...	t:44;...;t:53;...
30573	t	10A	t:44;...;t:53;...	t:44;...;t:53;...

FullBase report toolbar

The FullBase Report toolbar contains a single button on the toolbar—the Save Report button . Click this button to open a Save As dialog box in which you can specify the name and location for saving the report as an .xls/.xlsx or .txt file.

FullBase viewing and analysis options

Although the FullBase report does not have a full context menu, you can still copy text from the report, or copy the report as an image, and you can paste copied text into the editable Comments field. You can also view the information for a mutation in the Full Base report in the GAD.

- Select the text that you want to copy or to copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Right-click and select Copy to copy the selected text to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application, or use the Paste command on the Context menu to paste copied text into the editable Comments field.
- To copy the report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.
- Because the FullBase report shows every position and not just mutations, double-click a cell in the report to open the relevant position in the first trace of the consensus.

Chapter 8

GenBank File Editors

Mutation Surveyor provides many tools for optimizing input data and analyzing results. This chapter details the three tools that are specifically designed for editing GenBank files and Sequence files—the GBK File Editor, the Advanced GenBank Editor, and the SEQ File Editor. You use all these tools to view, add, and edit CDS information in multi-exon GenBank files, which are essential functions for creating and using GenBank-derived reference traces.



Although Mutation Surveyor provides many tools for optimizing input data and analyzing results, the GenBank Editor tools are the most commonly used of all the tools and that is why they are afforded their own chapter. Other Mutation Surveyor tools are discussed in appropriate chapters, including [Chapter 9](#), “Mutation Surveyor Tools,” on page 253.

This chapter covers the following topics:

- “Overview of GenBank Files” on page 227.
- “Advanced GBK File Editor Tool” on page 231.
- “SEQ File Editor Tool” on page 246.

Chapter 8

GenBank File Editors

Overview of GenBank Files

The *GenBank sequence database* is an open access, annotated collection of all publicly available nucleotide sequences and their protein translations. This database is produced at the National Center for Biotechnology Information (NCBI) as part of the International Nucleotide Sequence Database Collaboration, or INSDC. A *GenBank file* is a text-based file that shows the amino acid translation that corresponds to the base coding sequence (CDS). A GenBank file is always an optional file for a Mutation Surveyor project. If you do use a GenBank file as a reference file for Mutation Surveyor project, then the file must be in one of two formats—.gbk or .gb, which represents the sequence text of multiple CDSs or .seq, which represents the sequence text of a single CDS. All three file formats contain the same information and have similar conventions.

- A comment line is preceded with the greater than (>) symbol or a forward slash (/) and it contains the following information—the gene name, the CDS number, the reading frame number of the starting CDS base, the starting base number, the coding sequence (CDS), the starting amino acid number, the starting and ending vector sequences, the translation protein sequence from the CDS, and the DNA variations. Any text in a comment line is not read into the DNA sequence.
- The coding sequence is in uppercase. All other sequences are in lowercase.



Mutation Surveyor is designed to work with GenBank files that are obtained solely from NCBI. You should not use GenBank files that are obtained from other sources. See “To download a GenBank file.”

To download a GenBank file

1. Open the NCBI’s website (www.ncbi.nlm.nih.gov).
2. On the Search dropdown menu, select Gene.
3. In the search field, enter the official name or symbol for your gene of interest, and then click Search.

A list of all organisms that contain the specified gene is displayed.

Figure 8-1: Searching for a gene (NCBI)



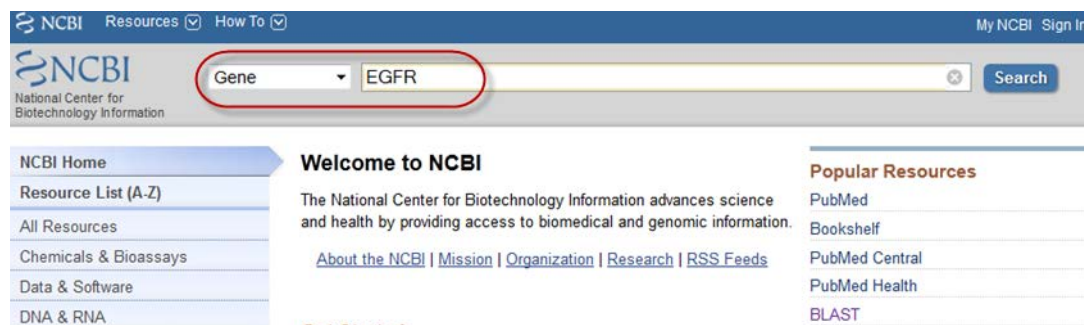


Figure 8-2: Gene search results

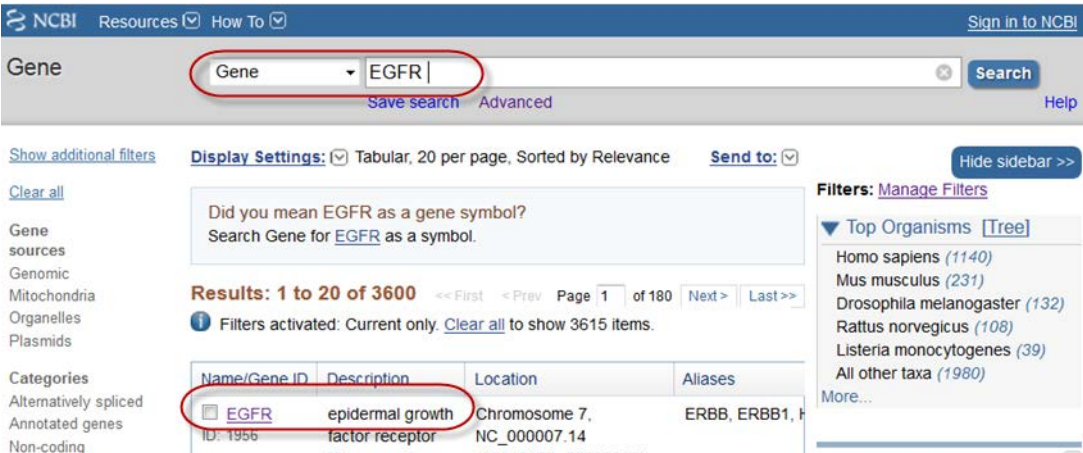

4. Scroll to the Genomics region section in the search results and select the correct assembly for your project.

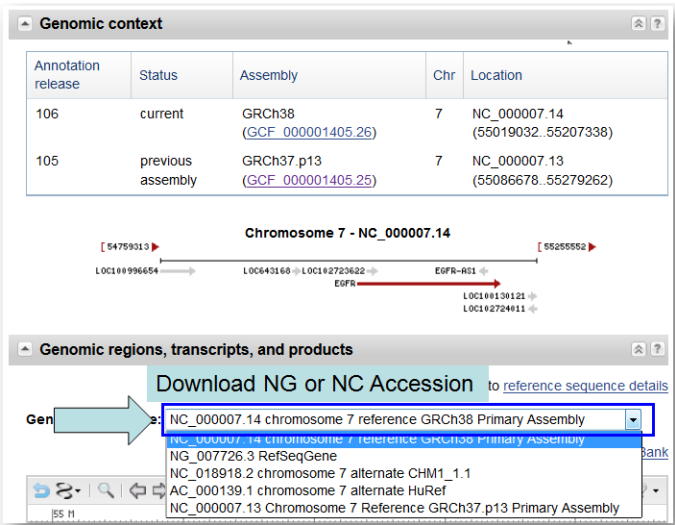Figure 8-3: Assembly options in the Gene report



5. Download the selected assembly. Note the following about the different assemblies:

Assembly	Description
NG Accession	<ul style="list-style-type: none"> The RefSeq. Most RefSeq GenBank files contain only a single transcript. NG Accession can show dbSNP. Nucleotide positions do not relate to chromosome coordinates. Click the GenBank link to download your selected NG Accession. See Figure 8-4 below.
NC Accession	<ul style="list-style-type: none"> NC accessions often contain several transcripts. NC accessions do not have dbSNP. Nucleotide positions are chromosome coordinates. Click the GenBank link to download your selected NC Accession. See Figure 8-5 below.
NT Accession	<ul style="list-style-type: none"> NT Accessions often contain several transcripts Nucleotide positions can be translated to chromosome coordinates. Available for download through Mutation Surveyor

Figure 8-4: NG Accession selected



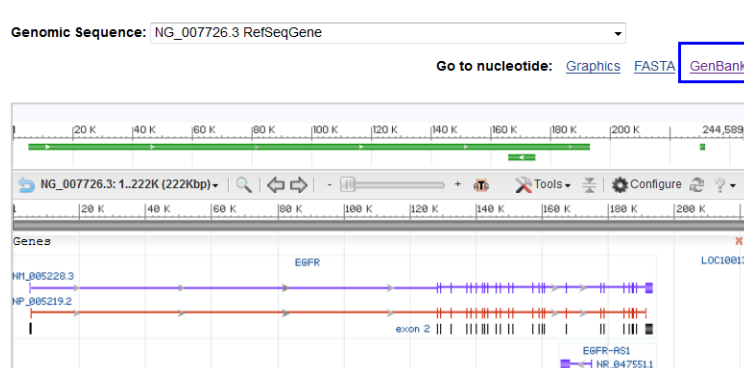


Figure 8-5: NC Accession selected

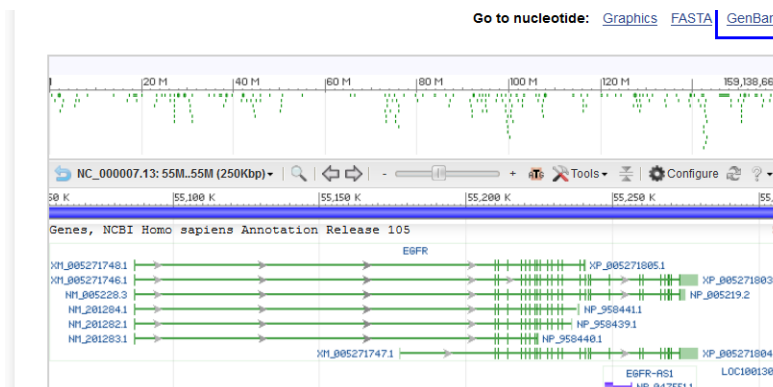

6. Optionally, before you download and save the GenBank file, do one or more of the following:
 - To include additional upstream and/or downstream sequences (for example, to include promoter regions), under Change Region Shown, click Selected Region, enter the new starting and ending base numbers, and then click Update View.
 - To include SNPs for an NG Accession, open the Customize View menu, and under Features added by NCBI, select the [] SNPs option, and then click Update View.
 - To change the view of the GenBank file for a gene that is on the minus strand, open the Customize View menu, and under Display Options, deselect Show Reverse Complement, and then click Update View.

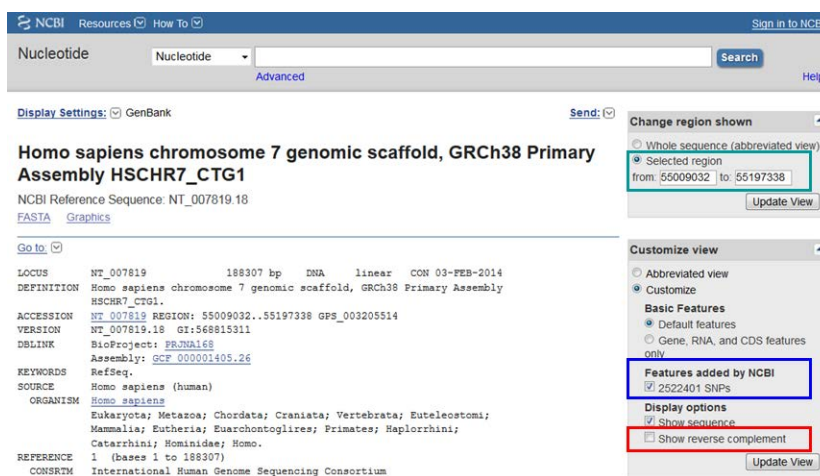Figure 8-6: Optional actions for a GenBank file



7. To save the GenBank file, open the Send menu, and then do the following.
 - a. Select Complete Record.
 - b. For Choose Destination, select File.
 - c. For Format, select GenBank.
 - d. Click Create File.

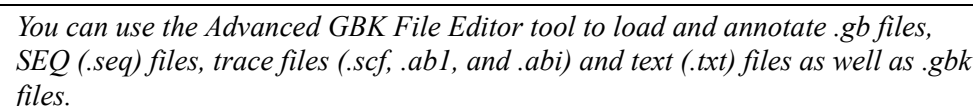
A File Download dialog box opens with options to Open or Save the selected GenBank file.

8. Save the file in a location of your choosing.

You can now use the Advanced GBK Editor to view, edit, and annotate this GenBank file and you can also use the Advanced GBK Editor to save the file as a SEQ file. See “Advanced GBK File Editor Tool” on page 231.

Advanced GBK File Editor Tool

You [use](#) the Advanced GBK Editor tool to view, edit, or annotate a GenBank file. If the GenBank file contains multiple protein and mRNA isoforms, the Advanced GBK File Editor tool displays all the proteins in the file and it displays the corresponding mRNA information for each isoform. You can also use the Advanced GenBank File Editor tool to add, view, or edit custom tags and identifiers to the reported variations in the file.




To load a GenBank file into the Advanced GBK File Editor tool

1. On the Mutation Surveyor main menu, click Tools > Advanced GBK File Editor.
2. On the Advanced GBK File Editor window main menu, click File > Open and in the Open File dialog box, select the appropriate file, and then click Open.

The file is loaded into the Advanced GBK File Editor. The Advanced GBK File Editor window has a toolbar with buttons that are specific for buttons for working with the loaded file and it has two panes—the GenBank Tree File pane and the Tabs pane. When the window first opens, the Sequence tab is the active tab in the Tabs pane.



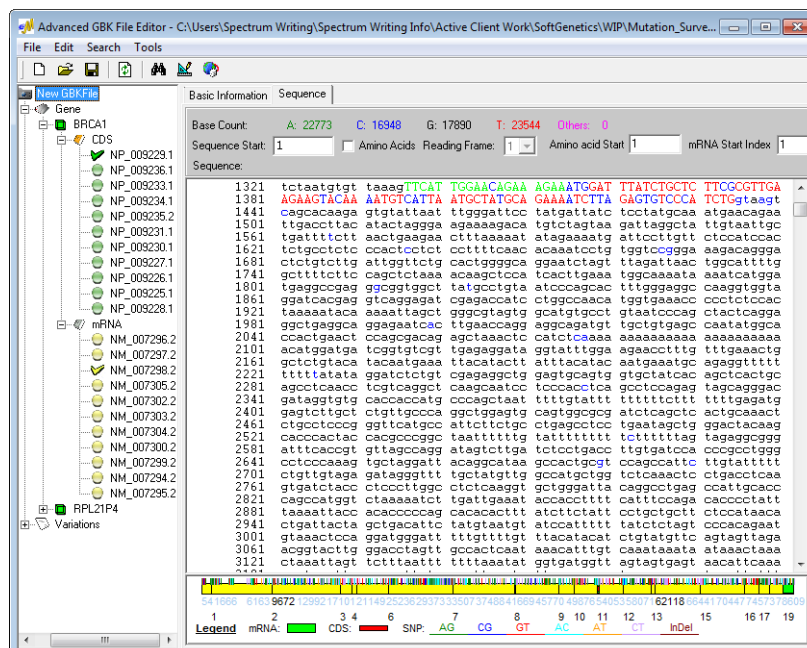




Figure 8-7: Advanced GBK File Editor window



Advanced GBK File Editor tool toolbar

Figure 8-8: Advanced GBK File Editor tool toolbar




Button	Description
	New—Clears the currently loaded GBK file and readies the editor to accept a new file.
	Open Files—Opens the Open Files dialog box which you use to load GenBank files or other files into the editor.
	Save All CDSs and all mRNAs—Saves all the CDSs and mRNAs for the loaded GenBank file in a single file.
	Refresh—Updates the display in the Advanced GBK File Editor window after you make a modification to the loaded GenBank file.
	Search in Sequence—Opens the Search in Sequence dialog box in which you can specify a nucleotide sequence for which to search. See “Advanced GBK File Editor tool functions” on page 243 .
	Adjust —Opens the Adjust File dialog box which you can specify the Sequence Start for the selected isoform. See “Advanced GBK File Editor tool functions” on page 243 .
	Search in Database—Opens the Search in Database dialog box which provides an option for accessing a SoftGenetics server and downloading a GenBank file. See “Advanced GBK File Editor tool functions” on page 243 .

Advanced GBK File Editor tool GenBank Tree File pane

The left pane in the Advanced GBK File Editor window is the GenBank Tree File pane. This pane displays all the GBK file information (protein and mRNA) in a simple tree format. Click the plus (+) and minus (-) symbols to expand and collapse the tree structure, respectively.

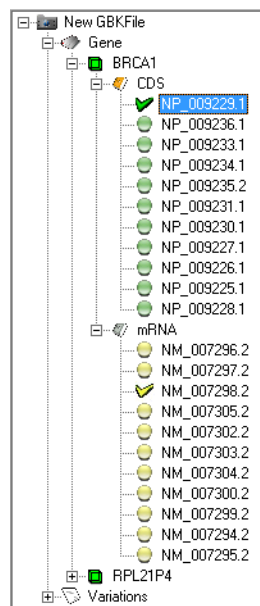Figure 8-9: Advanced GBK File Editor GenBank Tree File pane



The GenBank Tree File is interactive:

- Click an isoform in the tree structure to select it. The selected protein is marked with a green checkmark in the pane. Mutation Surveyor automatically selects the corresponding mRNA region for the CDS and marks it with a yellow checkmark.



The nucleotide sequence is the same for all isoforms. Only the mRNA and CDS locations change when you select a different isoform.


- Optionally, after you select an isoform, click File > Save As on the Advanced GBK File Editor tool main menu, and then on the dialog box that opens, indicate what information is to be saved:

Option	Description
Selected Gene Region	Saves <i>all</i> the transcripts (CDSs and mRNAs) for the gene to a new GenBank file. You can specify the name of this new GenBank file, and the location in which to save it.
Selected Gene Region; Selected mRNA Region	Saves only the transcript (CDS and mRNA) that is currently selected to a new GenBank file. You can specify the name of this new GenBank file, and the location in which to save it.

Option	Description
Seq Files for Selected CDS	An individual SEQ file is created for each exon in the currently selected CDS. The default name for each SEQ file is <GenBank filename>_exon_<n>.seq, where <n> reflects the exon number, and the files are saved in the same folder as the loaded GenBank file; for example. BRCA1_AmpID_exon_1.seq.


- Expand the Variations folder to view a list of all the recorded variations.

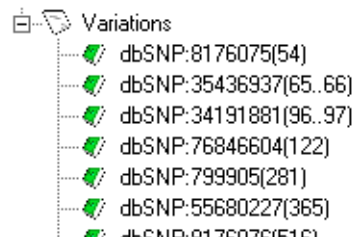Figure 8-10: Expanded Variations folder



- Double-click a variation in the Variations list to open the Variation Settings dialog box. The Variation Settings dialog box provides detailed information about the selected variation, including its alleles and its position in the gene.

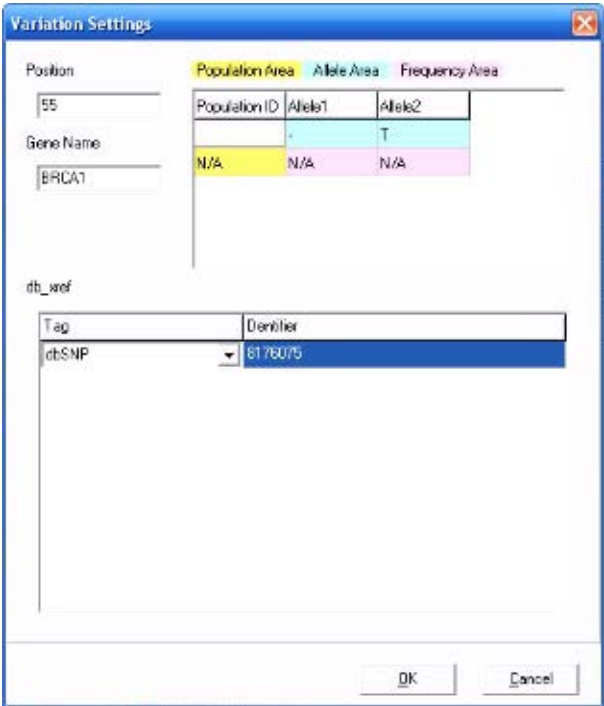Figure 8-11: Variation Settings dialog box




Functions for modifying variations, adding new variations, and adding tags and identifiers to variations are available from this dialog box. See [“Modifying and adding variations” on page 241](#).

Advanced GBK File Editor tool Tabs pane

The Tabs pane is the right pane in the Advanced GBK File Editor window. The tabs that are displayed in this pane depend on the item that is selected in the GenBank Tree File pane. If a protein or mRNA is selected in the GenBank Tree File pane, then the tabs that are displayed are the [Basic Information](#) tab and the [Sequence](#) tab. If a variation is selected in the GenBank Tree File pane, then the tabs that are displayed are the [Allele Frequency](#) tab and the [Sequence](#) tab. The [Sequence](#) tab is always the active tab when the Tabs pane first opens.

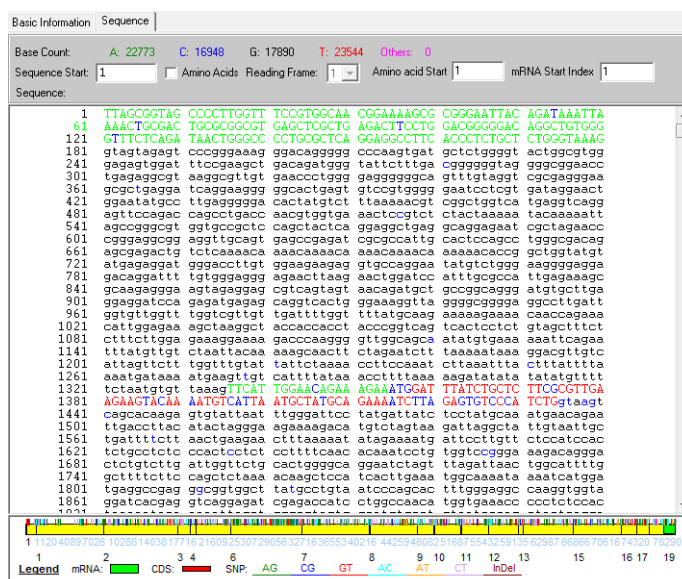Figure 8-12: Advanced GBK File Editor window Sequence tab



Sequence tab

The Sequence tab provides a visual representation of the gene. The tab is divided into three sections (upper, middle, and lower) that display different information for the gene.


- The upper section displays the *Base Count*, which is the total base count on a per nucleotide basis, with each base having a different color-code. The upper section also shows the Sequence Start, which typically has a value of one, but can be negative if promoter regions are used. You can modify the value in the Sequence Start field, or you can modify it using the Adjust option. (See [“Advanced GBK File Editor tool functions” on page 243.](#)) You can also modify the values in the Amino Acid Start or the mRNA start index fields. To view the amino acid translation for the nucleotides in the CDS, click Amino Acids.

Figure 8-13: Amino acid translation for a CDS

```
.321 tctaattgtgt taaagTTCAT TGAACAGAA AGAAATGGAT TTATCTGCTC TTCGCGTTGA
                                     M D L S A L R V E
.381 AGAAGTACAA AATGTCATTA ATGCTATGCA GAAAATCTTA GAGTGTCCTA TCTGtaagat
      E V Q N V I N A M Q K I L E C P I
```

- The middle section is a view-only textual representation of the full nucleotide sequence for the selected protein or mRNA region. Black lowercase letters indicate nucleotides in a non-coding (intron or intragenic) region. Red uppercase letters indicate nucleotides in a coding region (CDS). Green uppercase letters indicate nucleotides in an mRNA region. Blue letters indicate reported variations. (The letters are uppercase or lowercase depending on the region in which the variation occurs.) After you make a modification to the loaded GenBank file and click Refresh, Mutation Surveyor updates the display accordingly.
- The lower section is a color-coded bar chart that indicates different information about the gene. mRNA regions are shown in green and CDS regions are shown in red. Reported variations, including indels, are indicated by tick marks of different colors. The first row of numbers indicates the starting genomic base position for a CDS and the second row of numbers indicates the CDS. For example, in Figure 8-14 below, the 7th CDS has a starting genomic base position of 34,399.

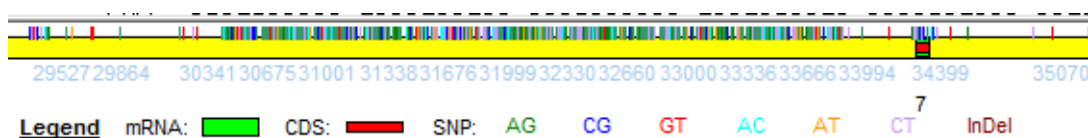Figure 8-14: CDS and genomic position numbering



The Sequence tab is interactive:

- To move the display in the middle section to the sequence for a CDS, double-click the CDS on the bar chart. Similarly, to move the display in the middle section to the sequence for an mRNA region, double-click the mRNA region on the bar chart.



You must double-click exactly the green region or the red region on the bar chart. If you click anywhere else on the bar chart, the display in the middle section does not move.

- To zoom in on a CDS region or an mRNA region in the bar chart for easier viewing, hold down the left mouse button and draw a box from the upper left hand corner of the region towards the lower right hand corner. A box is formed around the area that is being reduced for viewing.



You must make sure to include the entire CDS region or the mRNA region in the box and you cannot extend the box's upper boundary into the middle section.

- To zoom out, hold down the left mouse button and draw a box from the lower right hand corner of the region towards the upper left hand corner. (The magnification for zooming out is always 100%.)

- To open a context menu that contains options for working with the sequence text display, right-click anywhere in the sequence text display.

Figure 8-15: Sequence text display context menu

Copy	Ctrl+C
Paste	Ctrl+V
Find	Ctrl+F

Option	Description
Copy	Copies selected text in the sequence text display to your clipboard. To select text for copying, click and hold the left mouse button and drag the mouse to select the text that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.
Paste	Pastes copied text into the sequence text display.
Find	Opens the Search in Sequence dialog box. You use the options on this dialog box to search for specific text in the sequence text display. (See “Advanced GBK File Editor tool functions” on page 243.)


Basic Information tab

Mutation Surveyor pulls information directly from the loaded GenBank file and populates the fields on the Basic Information tab with this information. The information that is displayed on the tab depends on what option is selected in the GenBank Tree File pane—an isoform or an mRNA region. (See [Figure 8-18 on page 239.](#))

Figure 8-16: Advanced GBK File Editor, Basic Information tab for selected isoform

Basic Information Sequence	
Item	Value
CDS	3..54523,57756..57843,61500..61577,62078..62118,68316..68399,74334..74388,76257..76330,77748..77808,79649..79773
Region of Interest	
Amplicon ID	
External primer	
Internal primer	
CodonStart	1
product	breast cancer 1, early onset isoformBRCA1-delta9-10-11b
protein_id	NP_009236.1
note	isoform BRCA1-delta9-10-11b is encoded bytranscript variant BRCA1-delta9-10-11b: Derived byautomated computational analy
translation	MDLSALRVEEVQNVINAMQKILECPICLEIKEPVSTKCDHIFCKFCLLNQKKGPSQCPLCKNDITKRSLESTFRFSQLVEELLK
Corresponding_mRNA_Region	NM_007305.2

Field	Description
CDS	Lists the starting and ending base numbers for the first CDS, the second CDS, and so on. Within a CDS, the starting and ending base numbers are separated by two decimal points (.). CDSs are separated by commas. You can manually edit the information in the CDS field.

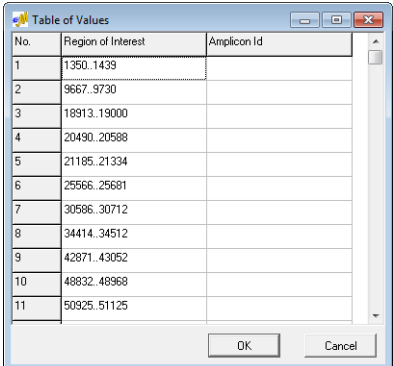Field	Description
Region of Interest/ Amplicon Id	<p>List of the ROIs defined in the CDS. Within an ROI, the starting and ending base numbers are separated by two decimal points (..). ROIs are separated by commas. Several options are available for adding and editing ROIs.</p> <ul style="list-style-type: none"> You can manually edit the information in the ROI Field. You can right-click in either the ROI or Amplicon ID field and select View of Table of Values to open the Table of Values dialog box which lists the ROIs and Amplicon IDs in table. You can modify the ROI values directly in this table, and you can right-click in this dialog box and select Insert Item to add new ROIs and Amplicon IDs. <p><i>Figure 8-17: Table of Values dialog box</i></p>  <p>Note: Amplicon ID is always an optional value; however, if your samples overlap and you are using amplicons to construct contigs, then to ensure proper alignment, the Amplicon ID should be the same as the Amplicon ID that is used in your sample filenames. (See “Use Amplicon to Construct Contig” on page 57.)</p> <ul style="list-style-type: none"> You can add ROIs that are based on the CDSs. See “Advanced GBK File Editor tool functions” on page 243.
<p>Note: Even if you modify the information for both the CDSs and the ROIs on this tab, your project’s graphical output can be limited based on the options that you have selected for Output Limits. See “Output Limits” on page 63.</p>	
External Primer Internal Primer	<p>Lists the starting and ending base numbers for the external/internal primer 5’ end and then for the external/internal primer 3’ end. For each end, the starting (From) and ending (To) base numbers are separated by two decimal points (..). The primers are separated by commas. Primer pairs are separated by semicolons.</p>

Field	Description
<ul style="list-style-type: none"> CodonStart product protein_id note translation Corresponding_mRNA_region transl_table 	Mutation Surveyor pulls the information in these fields directly from the loaded GenBank file. Mutation Surveyor deduces the corresponding mRNA region for each protein by inspecting the start and end positions of each of the exons between the protein and the mRNA. Mutation Surveyor deduces the amino acid translation relative to the nucleotides in the coding sequence.

Figure 8-18: Advanced GBK File Editor, Basic Information tab with mRNA region selected in GenBank Tree File pane

Basic Information	Sequence
Item	Value
Region	1, 180,1336, 1434,9672, 9725,18918, 18995,20495, 20583,21190, 21329,25571, 25676,34419, 34507,42876, 43047,48837
transcript_id	NM_007298.2
product	breast cancer 1, early onset, transcript variantBRCA1-delta9-11
note	Derived by automated computational analysis usinggene prediction method: BestRefseq.
Corresponding_CDS_Region	NP_009229.1

Field	Description
Region	Lists the starting and ending base numbers for the first mRNA region, the second mRNA region, and so on. Within an mRNA region, the starting and ending base number are separated by two decimal points (..). mRNA regions are separated by commas. You can manually edit the information in the Region field.
<ul style="list-style-type: none"> transcript_id product note Corresponding_CDS_region 	Mutation Surveyor pulls the information in these fields directly from the loaded GenBank file. Mutation Surveyor deduces the corresponding protein for each mRNA region by inspecting the start and end positions of each of the exons between the protein and the mRNA.

Allele Frequency tab

If a variation is selected in the GenBank Tree File pane, then the Tabs pane contains an Allele Frequency tab instead of a Basic Information tab. This tab lists all the reported variations in the GenBank file as well as any of the variations that you or other users have added to the file.




The frequency information refers to allele frequency with regards to population studies. Mutation Surveyor does not use this information and most SNPs in dbSNP do not have this information populated.

Figure 8-19: Allele Frequency tab

Basic Information		Sequence			
No	Position	Allele Number	db_xref	Frequency	Gene
1	54	2	dbSNP:8176075	N/A	BRCA1
2	56	2	dbSNP:8176077	N/A	BRCA1
3	56	0	dbSNP:8176077	N/A	BRCA1
4	60	2	dbSNP:123456	N/A	BRCA1
5	65..66	2	dbSNP:35436937	N/A	BRCA1
6	96..97	2	dbSNP:34191881	N/A	BRCA1
7	122	2	dbSNP:76846604	N/A	BRCA1
8	281	2	dbSNP:799905	N/A	BRCA1
9	365	2	dbSNP:55680227	N/A	BRCA1
10	516	2	dbSNP:8176076	N/A	BRCA1
11	1120	2	dbSNP:8176077	N/A	BRCA1
12	1221	2	dbSNP:3765640	N/A	BRCA1
13	1251..1255	2	dbSNP:56209245	N/A	BRCA1
14	1278..1280	2	dbSNP:8176078	N/A	BRCA1

To open a context menu that contains options for annotating the information in the Frequency column adding a variation, deleting a variation, and copying a variation, right-click anywhere in the Allele Frequency tab.

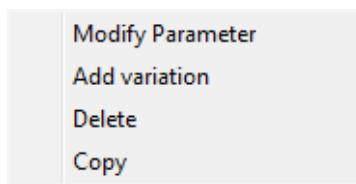Figure 8-20: Allele Frequency tab context menu



Option	Description
Modify Parameter	Opens the Variation Setting dialog box in which you can modify the selected variation. See “Modifying and adding variations” on page 241 .
Add Variation	Opens a blank Variation Setting dialog box in which you can add a new variation to the loaded GenBank file. See “Modifying and adding variations” on page 241 .

Option	Description
Delete	Deletes the selected variation line from the list of reported variations. Note: You can also delete a variation by right-clicking the variation in the GenBank Tree File pane, and then clicking Delete.
Copy	Copies the selected text in the cell to your clipboard. To copy text in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.

Modifying and adding variations

You can use functions that are available for the Allele Settings box to modify the information for a selected variation, add a variation to the loaded GenBank file, and add tags and identifiers to the variations.

Figure 8-21: Variation Settings dialog box

- To modify any of the information in the Position field or Gene Name field, double-click a displayed value to select it, and then modify the value.
- To modify the Identifier information for an existing tag, double-click the displayed value to select it, and then modify the value. You can also select a different tag for the variation on the Tag dropdown list.
- To add multiple tags for a variation, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.

- To modify the information in the Allele Frequency area for a selected variation, *triple-click* the allele that is being modified to select it, and then modify the allele as needed.

Figure 8-22: *Modifying a selected variation*

The figure shows three instances of the 'Variation Setting' dialog box, each illustrating a different type of genetic variation. Each dialog has a 'Position' field, a 'Gene Name' field, and a table with columns for 'Population ID', 'Allele1', and 'Allele2'. The 'Allele Area' is highlighted in cyan, and the 'Frequency Area' is highlighted in pink.

Position	Gene Name	Population ID	Allele1	Allele2	Type
601			A	G	Substitution
509..510			-	T	Insertion
79875..79877			ACA	-	Deletion

- To delete an allele from a variation, right-click in the Allele Frequency area, and on the context menu that opens, click Delete Allele. Allele2 is deleted. Repeat this step to delete Allele 1.
- To add a new variation:
 - Enter the Position and Gene Name for the variation.
 - Right-click in the Allele Frequency area, and on the context menu that opens, click Add Allele. Enter a value for Allele1. Repeat this step for Allele 2.
 - In the db_xref pane, select a tag from the Tag dropdown list, and then enter a value in the Identifier field. You can include both the tag and the identifier, or just one, or none. You can also include multiple tags and/or identifiers.

After you have made modifications to an existing variation, or added the values for a new variation, click OK to save your edits.

Advanced GBK File Editor tool functions

The Advanced GBK File Editor tool contains a variety of functions for working with the loaded GenBank file. You can:

- Search the loaded GenBank file for a specific sequence. See [“To search the loaded GenBank file for a specific sequence”](#) below.
- Access a GenBank database on a SoftGenetics server to download a GenBank file with the same sequence. See [“To download a GenBank file with the same sequence”](#) on page 244.
- Adjust the number of the first coding base or the sequence start. See [“To adjust the number of the first coding base or the sequence start”](#) on page 244.
- Add ROIs that are based on the CDSs. See [“To add ROIs that are based on CDSs”](#) on page 245.
- Reverse complement the currently loaded GenBank file. See [“To reverse complement the currently loaded GenBank file”](#) on page 245.

To search the loaded GenBank file for a specific sequence

1. On the Advanced GBK File Editor main menu, click Search > Search in Sequence.

The Search in Sequence dialog box opens.

Figure 8-23: Search in Sequence dialog box



2. Enter the sequence for which to search in this dialog box.
3. Optionally, to reverse complement the sequence before searching, click Complementary.

If the sequence is found, the sequence text display in the middle section is moved to the first occurrence of the sequence and the text is displayed in pink on the tab.

Figure 8-24: Located sequence in sequence text display

```

241 gagagtggat ttccgaagct gacagatggg tattctttga cggggggtag gggcggaac


```

To download a GenBank file with the same sequence

1. On the Advanced GBK File Editor main menu, click Search > Search in Database.

The Search in Database dialog box opens. This dialog box lists the entire nucleotide sequence of the currently loaded GenBank file.

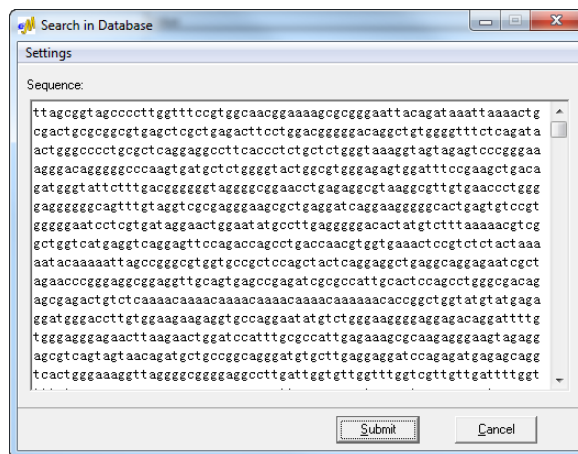Figure 8-25: Search in Database dialog box



2. Click Submit to initiate the search.

If a GenBank file is located that has the same genomic sequence as the currently loaded GenBank file, then Mutation Surveyor replaces the currently loaded GenBank file with the GenBank file from SoftGenetics's database. If a GenBank file that matches the genomic sequence of the currently loaded file *cannot* be found, then a message opens indicating this and the currently loaded GenBank file is cleared from the editor.


To adjust the number of the first coding base or the sequence start

At times, you might want to adjust the number of the first base or the first coding base in a GenBank file. For example, if a promoter region is 500 bps, then the sequence start that typically is 1 is now 501, but you still want the sequence start to be 1.

1. On the Advanced GBK File Editor main menu, click Tools > Adjust.

The Adjust File dialog box opens.

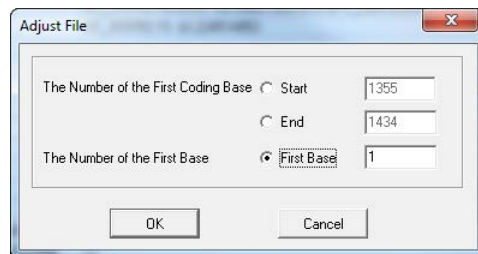Figure 8-26: Adjust File dialog box



2. Set your options as appropriate.

3. Click OK.


The Advanced GBK File Editor display is updated accordingly.


To add ROIs that are based on CDSs

1. On the Advanced GBK File Editor main menu, click Tools > Auto Create ROI.

The Create ROI for CDSs dialog box opens.

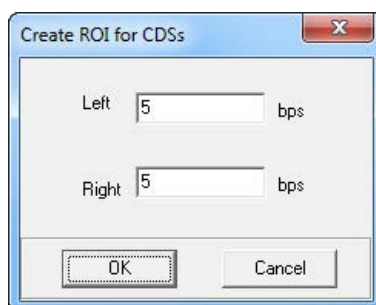Figure 8-27: Create ROI for CDSs dialog box



2. Specify the number of base pairs that are to included before (left) and after (right) each CDS.
3. Click OK.
4. Click the Refresh icon  to view these changes in the Region of Interest field on the Basic Information tab.

To reverse complement the currently loaded GenBank file

1. On the Advanced GBK File Editor main menu, click Tools > Complement.

A message opens, asking you if you want to reverse complement the currently loaded GenBank file.

2. Click Yes.

The Advanced GBK Editor display is updated accordingly, and a second message opens, asking you if you want to save the complement GenBank file.

3. Click Yes.

By default, the file name is the name of the GenBank file that was complemented and appended with the word “complement” (for example, BRCA1_NC_RC_complement.gb) and the file is saved in the same directory as the original GenBank file that was loaded. You can change either or both of these values as needed.

SEQ File Editor Tool

You [use](#) the SEQ File Editor tool to view, add, and edit CDS information in SEQ files.




You can use the SEQ File Editor tool to load and annotate FASTA (.fasta) and trace files (.scf, .ab1, and .abi) and text (.txt) files as well as .seq files.

To load a SEQ file into the SEQ File Editor tool

1. On the Mutation Surveyor main menu, click Tools > SEQ File Editor to open the SEQ File Editor window.
2. On the SEQ File Editor window main menu, click File> Open, and in the Open File dialog box, select the appropriate .seq file, and then click Open.

The .seq file is loaded into the SEQ File Editor.

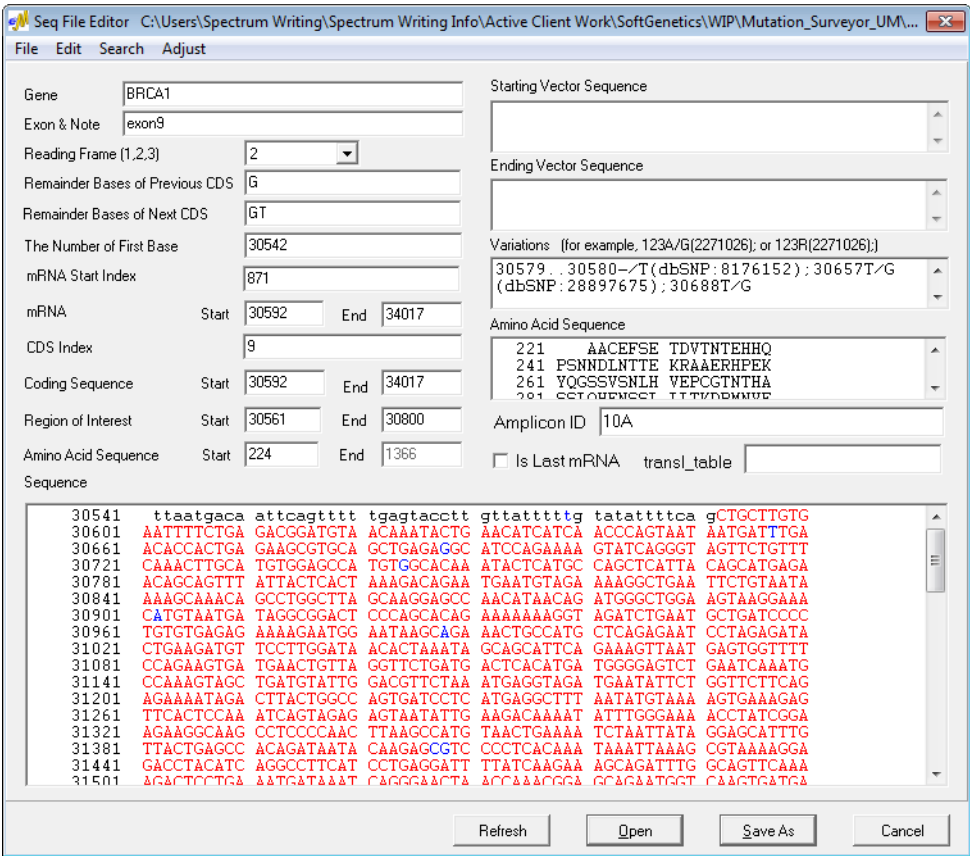Figure 8-28: SEQ File Editor window with loaded .seq file



SEQ File Editor tool layout

The SEQ File Editor window is divided into two sections—upper and lower. The upper section contains fields that are automatically populated with the information that Mutation Surveyor has pulled directly from the loaded SEQ file. You can edit much of this information to suit your working needs. The lower section is an editable textual representation of the full sequence for the single CDS. Black lowercase letters indicate nucleotides in a non-coding region. Red uppercase letters indicate nucleotides in a coding region. Blue letters indicate reported variations. (The letters are uppercase or lowercase depending on the region in which the variation occurs.) After you make a modification to the loaded SEQ file and click Refresh, Mutation Surveyor updates the display accordingly.

Field	Description
Gene	The gene name.
Exon and Note	The Exon number and any other additional information for the exon. Note: All Exon fields on all Mutation Surveyor reports are automatically populated with the value in this field, so edit this value carefully.
Reading Frame	Reading frame number for the first coding base.
Remainder Base(s) of Previous CDS	The number of bases from the previous CDS that are required to complete the first codon of the currently loaded CDS. <ul style="list-style-type: none"> • If the Reading Frame =1, then the no bases are required. The coding sequence is complete. • If the Reading Frame =2, then a single base is required from the previous CDS. • If the Reading Frame =3, then two bases are required from the previous CDS.
Remainder Base(s) of Next CDS	The number of bases from the next CDS that are required to complete the last codon of the currently loaded CDS. <ul style="list-style-type: none"> • If the Reading Frame of the next CDS=1, then the no bases are required. The coding sequence is complete. • If the Reading Frame of the next CDS=2, then two bases are required from the next CDS. • If the Reading Frame of the next CDS=3, then one base is required from the next CDS.
The Number of the First Base	The number assigned to the first nucleotide in the sequence text display. If you change this number, then numbering of all the nucleotides in the sequence text display is changed accordingly. Note: If you change the value for The Number of the First Base, the CDS numbers and the ROI numbers are not changed. To change these numbers, you must use the Adjust function. (See “SEQ File Editor tool functions” on page 250.)
mRNA Start Index	The genome position of the mRNA start is related to the mRNA numbering through the mRNA Start Index. This value is needed to determine the offset in numbering between the CDS nomenclature and the mRNA nomenclature relative to the number of bases in the untranslated region that is upstream of the CDS start.

Field	Description
mRNA	The starting and ending nucleotide numbers for the mRNA region. Note: If you modify these values, the selected sequence is read as cDNA.
CDS Index	The CDS number as pulled directly from the loaded SEQ file.
Coding Sequence	The starting and ending nucleotide numbers for the CDS. The CDS is indicated by red letters in the sequence text display.
Region of Interest	The starting and ending nucleotide numbers for the ROI defined in the CDS.
Caution: You can renumber the positions in the SEQ file (The Number of the First Base, the mRNA Start and End values, the Coding Sequence Start and End values, and the Region of Interest Start and End values) so that the numbering is relative to the CDS and not to the genomic positions (which is the numbering scheme in the GAD); however, if you do renumber the positions, then the position at the end of one sequence file might overlap the position at the beginning of the next sequence file, which can cause difficulty in your data analysis.	
Note: Even if you modify the information for both the CDSs and the ROIs on this tab, your project's graphical output can be limited based on the options that you have selected for Output Limits. See "Output Limits" on page 63 .	
Amino Acid Sequence	<p>The starting codon and ending codon numbers for the amino acid string for which the currently loaded CDS codes. If the starting and ending coding sequence positions are populated and you enter only the starting amino acid string number, and then click Refresh, Mutation Surveyor automatically calculates the ending number for the string.</p> <p>Note: To display the amino acid sequence in the standard mutation reports and the GAD, Amino Acid must be selected on the Output tab of the Project Options dialog box. See "Output tab" on page 62.</p>
<ul style="list-style-type: none"> Starting Vector Sequence Ending Vector Sequence 	<p>Trims a sample sequence according to a starting vector sequence and an ending vector sequence. To carry out a vector trim:</p> <ul style="list-style-type: none"> You must specify both the vector base sequence and the clone DNA base sequence in the Sequence field. You must specify the starting vector sequence in the Starting Vector Sequence field. You must specify the ending vector sequence in the Ending Vector Sequence field. <p>Mutation Surveyor carries out a "loose match" between the vector base sequence and the starting and ending vector sequences to determine the regions of the SEQ reference file that contain the vector sequence. When a sample trace aligns to both the vector sequence and the cDNA sequence of the SEQ reference file, Mutation Surveyor adjusts the mutation detection region inward to avoid mutation detection in the vector sequence of the sample traces.</p> <p>Note: To carry out a vector trim, Vector Trimming must be selected on the Options tab of the Project Options dialog box. See "Contig tab" on page 55. Mutation Surveyor trims the data according to the specified values and indicates the boundaries with vertical blue bars in the GAD.</p> <p>Note: You can also use the Mask Vector Sequence tool to replace base calls in the vector sequence region. See "Mask Vector Sequence Tool" on page 291.</p>

Field	Description
Variations	<p>If you modify or enter known sequence variations, the variations must adhere to a specific layout:</p> <ul style="list-style-type: none"> Substitution: <ul style="list-style-type: none"> <position><variant>/<normal>(<dbSNP_Identifier>) Example: 30565C/T(dbSNP:8176209) Insertion: <ul style="list-style-type: none"> <position_before>..<position_after>-<nucleotides>(dbSNP_identifier) Example: 30609..30610-/C(dbSNP:35466763) <position_before>..<position_after>ins<nucleotides>(<dbSNP_Identifier>) Example: 30609..30610insC(dbSNP:35466763) Deletion: <ul style="list-style-type: none"> <position_1>..<position_N><nucleotides>/-<db_SNP_Identifier>) Example: 31549..31552GCAG/-<dbSNP:84678345>) <position_1>..<position_N>del<nucleotides>/-<db_SNP_Identifier>) Example: 31549..31552delGCAG(dbSNP:84678345)
Amino Acid Sequence	The amino acid sequence that corresponds to the base sequence beginning one base after the Start base indicated for the Amino Acid Start. The string is shown using standard IUB/IUPAC amino acid codes.
Amplicon ID	Amplicon ID is always an optional value for the ROI; however, if your samples overlap and you are using amplicons to construct contigs, then to ensure proper alignment, the Amplicon Id should be the same as the Amplicon ID that is used in your sample filenames. (See "Use Amplicon to Construct Config" on page 57.)
Is Last mRNA	<p>Select this option if the SEQ file contains the last base of mRNA for the gene.</p> <p>Note: Mutation Surveyor does not automatically add a checkmark to the "Last mRNA" field for the SEQ file that contains the last base of mRNA when you create SEQ files from the Advanced GBK File Editor.</p> <p>Note: The Relative to mRNA nomenclature outputs IVS*5 for a position that is 5 bases after the mRNA, rather than IVS3123+5, indicating that the position is outside of the mRNA, and not within an intron.</p>

SEQ File Editor tool functions

The SEQ File Editor tool contains several of functions for working with the loaded SEQ file. In addition to being able to modify values directly in the editor, you can:


- Search the loaded SEQ file for a specific sequence. See [“To search the loaded SEQ file for a specific sequence”](#) below.
- Adjust the number of the first coding base or the sequence start. See [“To adjust the number of the first coding base or the sequence start”](#) on page 251.

To search the loaded SEQ file for a specific sequence

1. On the SEQ File Editor main menu, click Search > Search in Sequence.

The Search in Sequence dialog box opens.

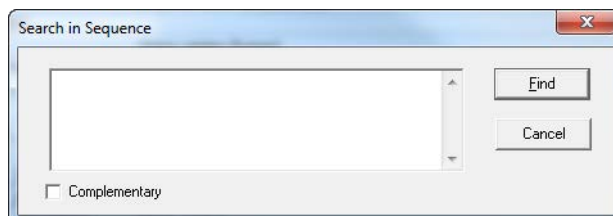Figure 8-29: Search in Sequence dialog box



2. Enter the sequence for which to search in this dialog box.
3. Optionally, to reverse complement the sequence before searching, click Complementary.

If the sequence is found, the sequence text display in the middle section is moved to the first occurrence of the sequence and the text is displayed in pink on the tab.

Figure 8-30: Located sequence in sequence text display

```
241 gagagtggat ttccgaagct gacagatggg tattctttga cgggggtag gggcggaac
```


To adjust the number of the first coding base or the sequence start

At times, you might want to adjust the number of the first base or the first coding base in a SEQ file. For example, if a promoter region is 500 bps, then the sequence start that typically is 1 is now 501, but you still want the sequence start to be 1.

1. On the Advanced GBK File Editor main menu, click Tools > Adjust.

The Adjust File dialog box opens.

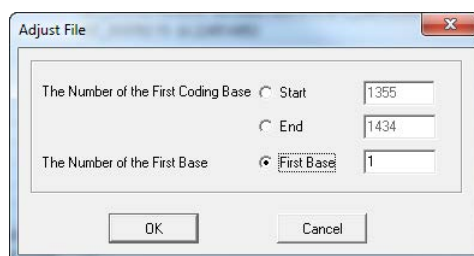Figure 8-31: Adjust File dialog box



2. Set your options as appropriate.
3. Click OK.

The SEQ File Editor display is updated accordingly.

Chapter 8
GenBank File Editors

Chapter 9

Mutation Surveyor Tools

Mutation Surveyor provides many tools for optimizing input data and analyzing results, including tools for modifying filename information to ensure the correct grouping of sample files in a project, tools for converting files from one format to another, and tools for assembling sample files and project files for comparison.

This chapter covers the following topics:

- [“Sample File Editing Tools” on page 255.](#)
- [“File Conversion Tools” on page 269.](#)
- [“Assembly Tools” on page 277.](#)
- [“Other Tools” on page 289.](#)

Chapter 9

Mutation Surveyor Tools

Section 1

Sample File Editing Tools

Section Contents

- [“Filename Editor Tool” on page 257.](#)
- [“2D Filename Match Editor tool” on page 261.](#)
- [“ND Filename Match Editor Tool” on page 264.](#)
- [“Print Header Editor Tool” on page 267.](#)

Chapter 9

Mutation Surveyor Tools

Filename Editor Tool


You use the FileName Editor tool to modify your sample filenames to ensure that your sample files will be grouped correctly in a Mutation Surveyor project. You can use the tool to [modify](#) sample filenames programmatically, or you can manually [change](#) filenames using a tab-delimited text file.



To programmatically change filenames

1. On the Mutation Surveyor main menu, click Tools > Filename Editor.

The Filename Editor window opens.

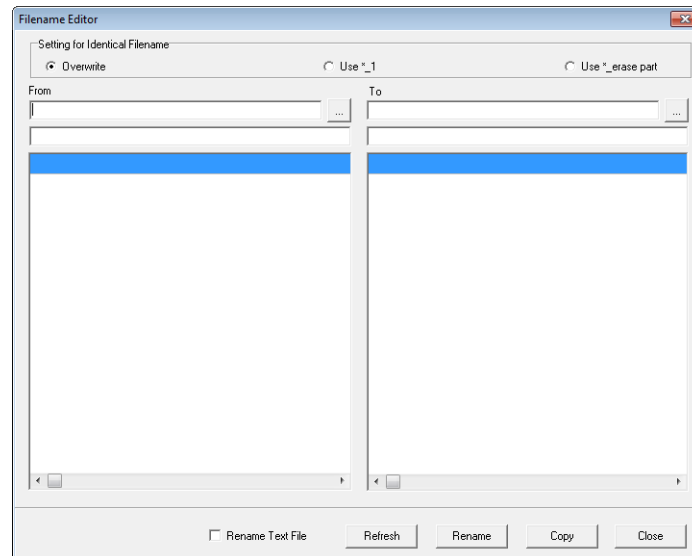Figure 9-1: *Filename Editor window*



2. Click the Browse button  next to the From field, and browse to and select the directory that contains the sample files for which you are modifying the names.
A list of all the sample files in the selected directory is displayed in the From (left) pane of the window.
3. Click the Browse button  next to the To field, and browse to and select the directory that is to contain the renamed sample files.



If you select the same directory for the From and To lists, then your original files are replaced with the edited files, so select the same directory only if you do not want to retain your original files.

4. Select the option by which to modify your existing sample filenames.

A placeholder in the form of *.* is displayed in the blank field below the From field and below the To field. The asterisk (*) in this placeholder is a wildcard character, which is a special character that represents one or more characters in a string of characters. For example, in searching, “run*” implies to search for any word that starts with “run” and that has any ending, such as running, runt, runs, rundown, and so on.

Option	Description
Overwrite	Replaces all or part of the original filename. You can replace part or all the filename with different characters, or you can replace all or part of the filename with no characters, which essentially removes the specified part of the filename.
Use*_1	Replaces all or part of the original filename and then sequentially numbers any identically named files.
Use *_erase part	Removes specified characters from the filename. The remaining characters make up the new filename.

5. In the From field, enter the part of the filename that is to be replaced or removed between the asterisks.



The list of sample files temporarily disappears from the From list as you enter the part of the filename that you want to remove. The list is displayed again after Mutation Surveyor recognizes the part of the filename that you have entered.

6. In the To field, enter the characters that are to replace the filename or part of the filename between the asterisks. If you have selected Overwrite and you simply want to remove a selected part of the filename, enter **.



Whatever value that you enter in the From field is automatically entered in the To field as well, so you must make sure to edit the value in the To field correctly.

7. Click Refresh.

The list of edited sample names is displayed in the To (right) pane of the window.

Figure 9-2: Overwrite example (characters replaced)

From	To
C:\Users\Spectrum\Writing\Spectrum\Writing Info\Active Client	C:\Users\Spectrum\Writing\Spectrum\Writing Info
Sample	*Patient*
BRCA1_012345_10A_1_F_0_Sample.scf	BRCA1_012345_10A_1_F_0_Patient.scf
BRCA1_012345_10A_1_R_0_Sample.scf	BRCA1_012345_10A_1_R_0_Patient.scf
BRCA1_012345_10A_2_F_0_Sample.scf	BRCA1_012345_10A_2_F_0_Patient.scf
BRCA1_012345_10A_2_R_0_Sample.scf	BRCA1_012345_10A_2_R_0_Patient.scf

Figure 9-3: Overwrite example (characters removed)


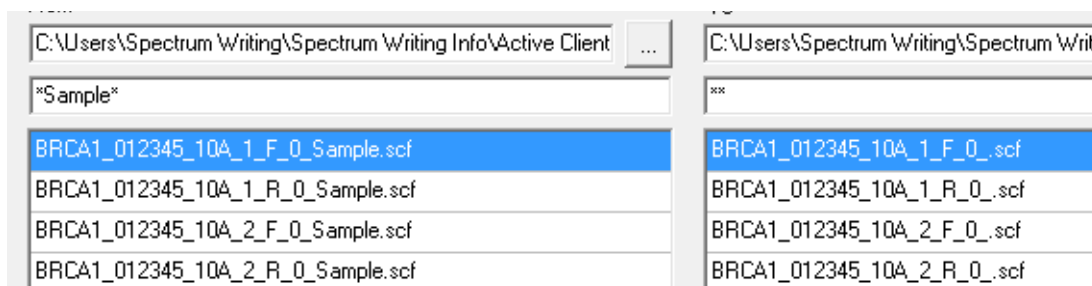


Figure 9-4: Use*_1 example

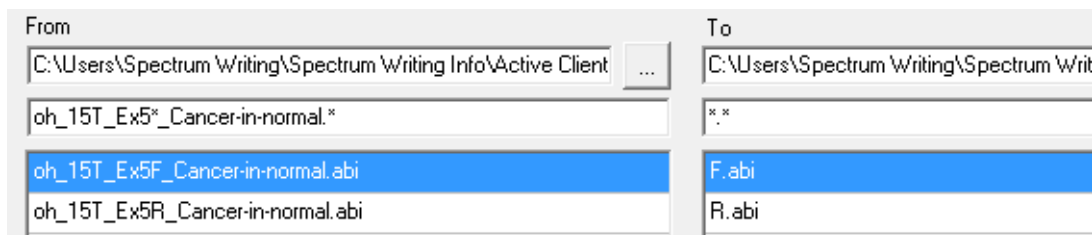


Figure 9-5: Use*_erase part example




8. Do one of the following:
 - If you selected the same directory for the From and To lists, click Rename.
 - If you selected different directories for the From and To lists, click Copy.

To manually change filenames

1. Create a tab-delimited text file that has two columns. In the first column, enter the current filenames. In the second column, enter the target filenames.
2. Save the text file in the same directory as the sample files.
3. On the Mutation Surveyor main menu, click Tools > Filename Editor.
The Filename Editor window opens.
4. Click Rename Text File.

5. Click the Browse button  next to the From field, and browse to and select the text file.

A list of all the sample files in the selected directory is displayed in the From (left) pane of the window.

6. Click the Browse button  next to the To field, and browse to and select the directory that is to contain the renamed sample files.



If you select the same directory for the From and To lists, then your original files are replaced with the edited files, so select the same directory only if you do not want to retain your original files.

7. Click Rename.

The list of edited sample names is displayed in the To (right) pane of the window.

2D Filename Match Editor tool


If your sample files are named in such a way that neither Mutation Surveyor's default project options nor Mutation Surveyor's automatic file pairing options result in your sample files being paired correctly, you should [use](#) the 2D Filename Match Editor tool to create a Grouping File. A *Grouping File* is a text file that defines how your sample files should be paired in your Mutation Surveyor project. You can point to the Grouping File in your Mutation Surveyor project, and all the sample files that are defined in the file are added to the project.

To use the 2D Filename Match Editor tool

1. On the Mutation Surveyor main menu, click Tools > 2D Filename Match Editor.

The 2D Filename Match Editor window opens.

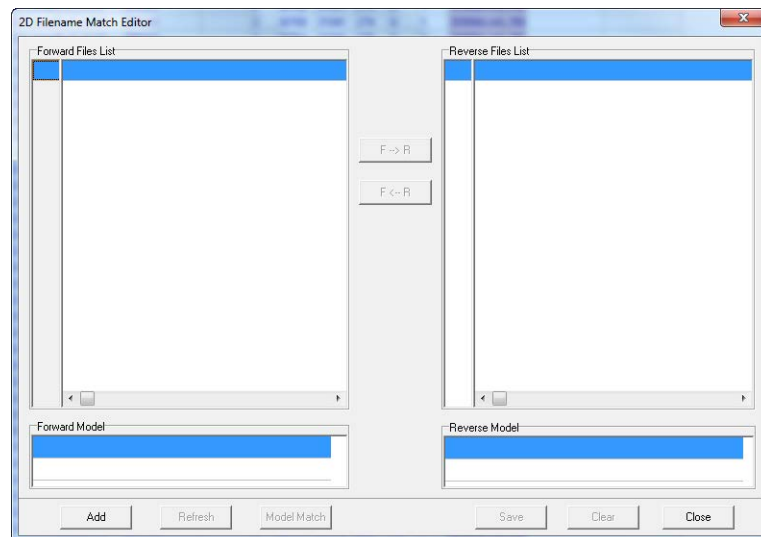Figure 9-6: 2D Filename Match Editor window




Chapter 9

Mutation Surveyor Tools

2. Click Add.

The Open Files dialog box opens.

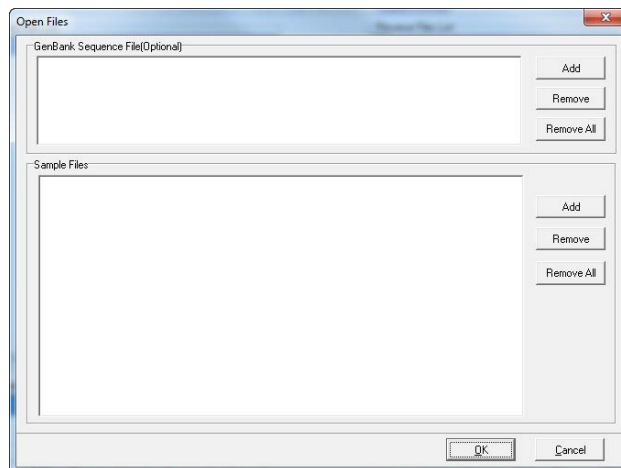Figure 9-7: Open Files dialog box



3. In the Sample Files (lower) pane, click Add, and then click Add Files or Add Directory to select the samples files that are to be grouped.




If you are adding files, CTRL-click to select multiple non-contiguous files. SHIFT-click to select multiple contiguous files. If you are adding a directory, then all the files in the directory are added in a single step.

4. Click OK.

The samples files are added to the 2D Filename Match Editor window. Mutation Surveyor initially tries to correctly group the sample files based on information in the sample file names such as *_F and *_R.

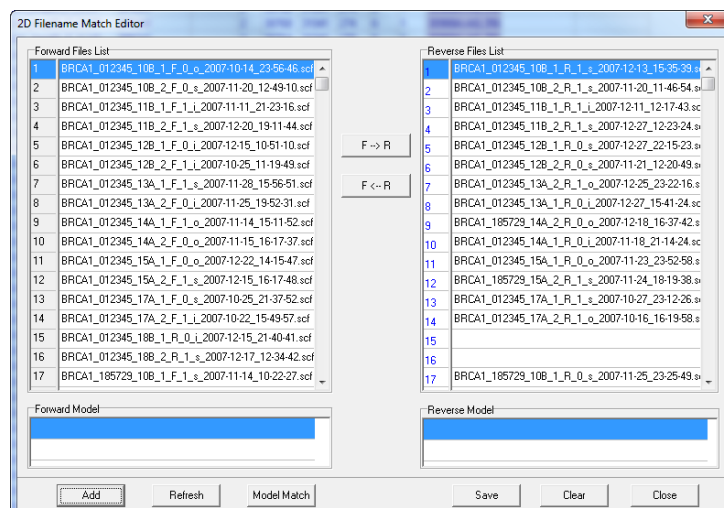Figure 9-8: 2D Filename Editor window with files paired by Mutation Surveyor




5. If the automatic pairing is correct, then continue to [Step 6](#); otherwise, continue to [Step 7](#).
6. Click Save to open the Save Result dialog box in which you can specify a name for saving the Grouping File.



By default, the Grouping File is saved as a text file in the same folder as all the sample files. You must not change this location, or the Grouping File function will not work.

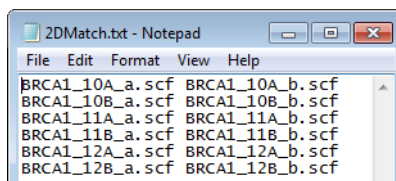Figure 9-9: 2D Grouping File example



7. If the automatic pairing is not correct, then do one or more of the following to correctly pair your sample files:
 - Add a GenBank file to assist Mutation Surveyor surveyor in aligning and pairing the sample files.
 - Select a file and use the Forward (F--> R) or Reverse (F <--R) buttons to move the sample file to the correct list. After you have moved all the necessary files, click Refresh to have Mutation Surveyor try to correctly group the sample files based on information in the sample file names such as *_F and *_R.
 - Select a file and use the Forward (F--> R) or Reverse (F <--R) buttons to move the sample file to the correct list. After you have moved all the necessary files, edit the order of the lists. (To edit the order, click a position number to select it, change the position number accordingly, and then click Refresh.)
 - Use model files to correctly pair sample files.
 - Identify two forward sample files and two reverse sample files that are appropriate to serve as “model” files.
 - Move the reverse sample files to the Reverse Files list.
 - Right-click each forward sample file and on the context menu that opens, click Copy to Model. Repeat this step for the reverse sample files.
 - Click Model Match to correctly pair the remaining sample files.
8. Click Save to open the Save As dialog box in which you can specify a name for saving the Grouping File.



By default, the Grouping File is saved as a text file in the same folder as all the sample files. You must not change this location, or the Grouping File function will not work.

ND Filename Match Editor Tool

You [use](#) the ND Filename Match Editor tool to create a Grouping file that can assist in grouping sample files correctly in the Project report and in the SGP Comparison tool. The sample files can be .scf files or .abi/ab1 files.




You do not have an option to load the ND Grouping file for the Project report after running a Mutation Surveyor project. To use the ND Grouping file for the Project report, when setting up your Mutation Surveyor project, you must select Load Grouping File on the Open files dialog box, and then click Add in the Sample Files pane to select the Grouping file. All the sample files that are listed in the Grouping file are added to the project. Although Mutation Surveyor attempts to use the information in the Grouping file to group the sample files in the project, the Grouping file's primary purpose is for the grouping of sample files in the Project report, the Project Sample Assembly subreport, and the Comparison functions.


To use the ND Filename Match Editor

1. On the Mutation Surveyor main menu, click Tools > ND Filename Match Editor.

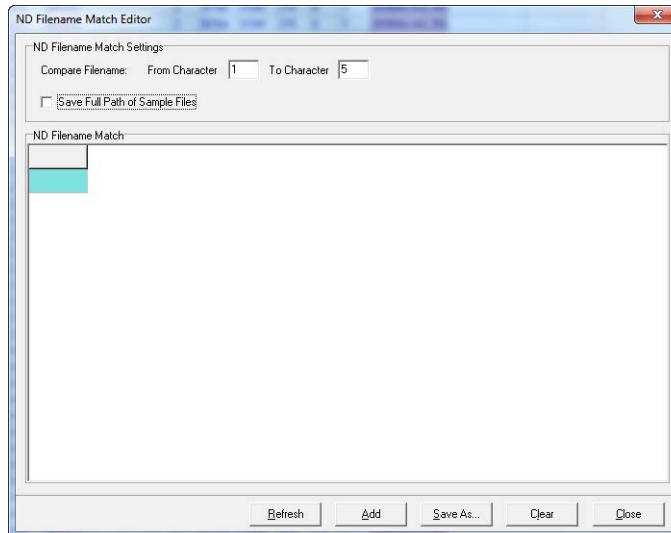Figure 9-10: ND Filename Match Editor window




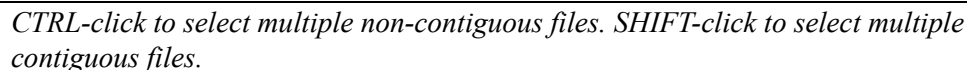2. Click Add to open the Open Sample Files dialog box and select the sample files that are to be grouped.



The sample files are displayed in the ND Filename Match Editor window. Each row in the window represents a group of sample files. By default, the sample files are first grouped by comparing the filenames from character one to character five.

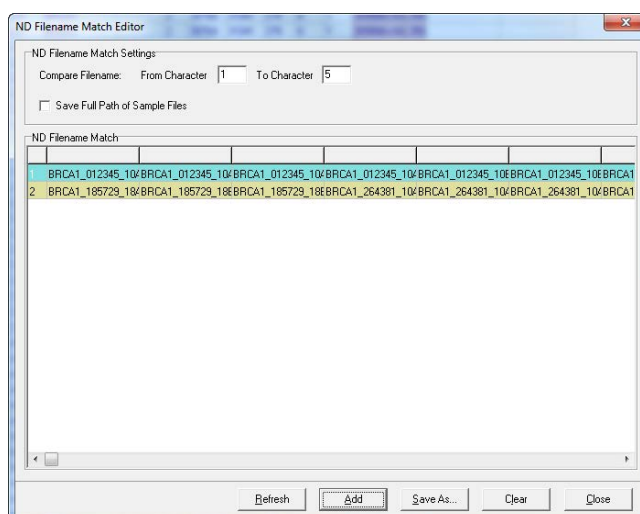
Figure 9-11: ND Filename Match Editor window with grouped sample files



3. Optionally, to modify the sample file grouping, change the character values for Compare Filename, and then click Refresh.
4. When you are satisfied with the sample grouping, do one of the following to save the Grouping file as a text file:
 - If you are saving the Grouping file in the same location as the sample files, click Save As to open the Save As dialog box in which you can specify a name and a location for saving the Grouping file as a text file.

By default, the Filename file is populated with _NDMatch.txt. You can change the name, but you cannot change the file type. The Grouping file lists each row from the display as a single line in the text file.

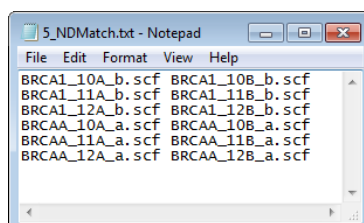Figure 9-12: ND Grouping file saved in same location as the sample files




Chapter 9

Mutation Surveyor Tools

- If you are saving the Grouping file in a *different* location than all the sample files, then click Save Full Path of Sample Files, and then click Save As to open the Save As dialog box in which you can specify a name and a location for saving the Grouping file as a text file.

By default, the Filename file is populated with _NDMatch.txt. You can change the name, but you cannot change the file type. The Grouping file lists each row from the display as a single line in the text file with the full path to the file name included.

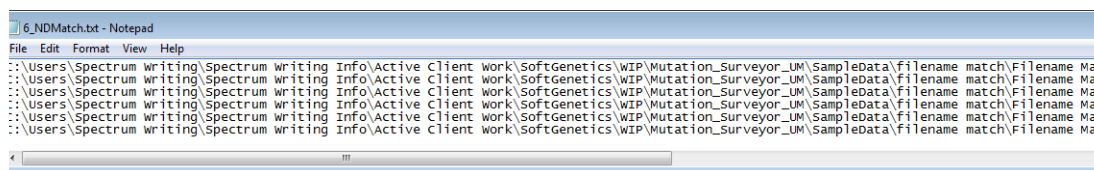Figure 9-13: ND Grouping file saved in a different location than the sample files



5. Click Close.



Even if you close, and then reopen Mutation Surveyor, the ND Filename Match Editor tool “remembers” the last settings specified for it.

Print Header Editor Tool


You **use** the Print Header Editor tool to create a custom header that can be displayed at the top of designated output files.

To use the Print Header Editor tool

1. On the Mutation Surveyor main menu, click Tools > Print Header Editor.

The Print Header Editor window opens.

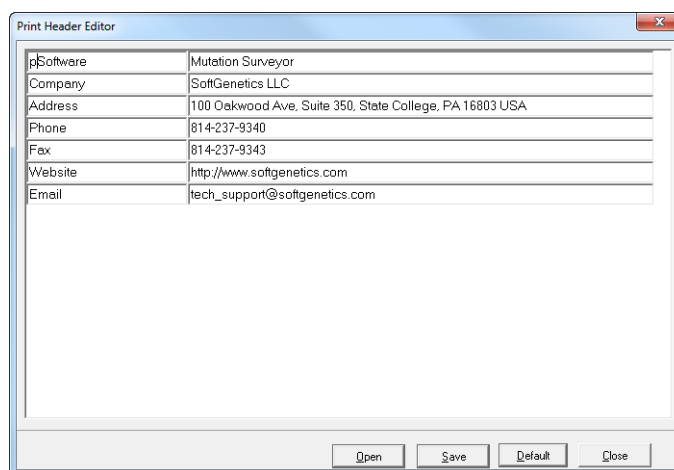Figure 9-14: *Print Header Editor window*



2. Do one of the following:
 - Modify any of the default template information in either column.
 - Click Open to select an existing custom header file to modify.



When modifying an existing custom header, click Default at any time to reset all values in both columns to their original values.

3. Optionally, add or delete rows of information as needed.
 - To delete a row from the header, right-click the row, and then click Delete a Row.
 - To insert a row into the header, right-click the row that is to be located below the inserted row, and then click Insert a Row.
 - To add a row as the last row in the header, right-click any row, and then click Add a Row.
4. Click Save to save the custom header as an .inf file.

Chapter 9

Mutation Surveyor Tools

Section 2

File Conversion Tools

Section Contents

- [Create SCF Tool on page 271](#)
- [Convert to SCF tool on page 274](#)
- [Extract SGP Tool on page 275](#)

Chapter 9

Mutation Surveyor Tools

Create SCF Tool


You **use** the Create SCF tool to create a **synthetic** or **pseudo** trace file from the sequence in a SEQ file or in a GBK file. Traces are created in both the forward and reverse direction whenever possible and are saved as .scf files.

Synthetic traces versus pseudo traces

At a minimum, you must load a GenBank file in the Create SCF tool to create SCF files. Loading sample traces is optional.


- If you do not load sample traces, then the output is one forward and one reverse trace for the first 2500 of the loaded GenBank file, or all the bases if the GenBank file contains less than 2500 bases. Each peak is 3000 RFU.
- If you load sample traces, and no match between the sample traces and the GenBank file is found, then a trace file is created for the first 2500 bases of the loaded GenBank file, or all the bases if the GenBank file contains less than 2500 bases. Each peak is 3000 RFU.
- If you load sample traces and a match between the sample traces and the GenBank file is found, then the trace option that is selected—Synthetic or Pseudo—determines the characteristics of the trace that is created.
 - If Synthetic is selected, then the resulting SCF file has a peak profile similar to that of the synthetic reference traces that Mutation Surveyor creates. The height of all four peaks at a given location are summed together and the base call and color is that of the nucleotide in the GenBank sequence at the given position.

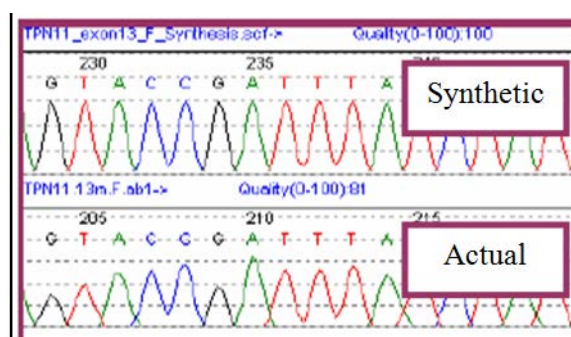Figure 9-15: Actual trace compared to a synthetic trace



- If Pseudo is selected, then the resulting SCF file has a peak profile similar to that of the original matched trace with one exception—positions where the base call differs between the loaded trace and the GenBank file are given the base call and color of the GenBank file. For example, in [Figure 9-16](#), note how the actual 179N was replaced with a pseudo 160C.

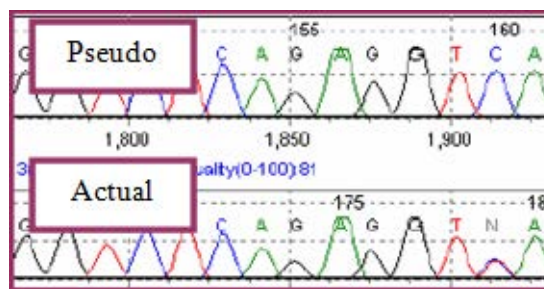Figure 9-16: Actual trace compared to pseudo trace




To use the Create SCF tool

1. On the Mutation Surveyor main menu, click Tools > Create SCF.

The Create SCF File window opens.

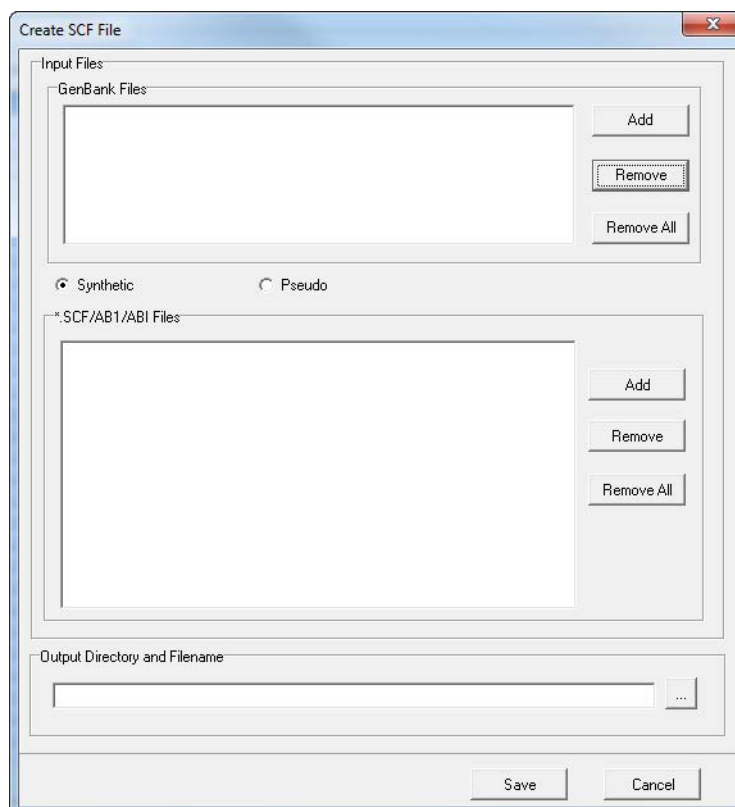
Figure 9-17: Create SCF File window



2. Leave the default selection of Synthetic as is, or click Pseudo.
3. In the GenBank files (upper) pane, click Add to open the Open Files dialog box, and select the .seq or GBK files that are to be used for creating the trace.
4. If you selected Synthetic, then continue to [Step 6](#); otherwise, continue to [Step 5](#).
5. In the *.SCF/ABI/ABI Files (lower) pane, click Add, and then click Add Files or Add Directory to select the samples files that are to be used for creating the pseudo trace.



*If you are adding files, CTRL-click to select multiple non-contiguous files.
SHIFT-click to select multiple contiguous files. If you are adding a directory, then
all the files in the directory are added in a single step.*

6. Click the Browse button  to open the Save As file dialog box in which you can specify a name and location for the saved trace file.
 - Forward and reverse synthetic trace files are saved as *_F_synthesis.scf and as *_R_synthesis.scf, respectively.
 - Forward and reverse pseudo trace files are saved as *_F_pseudo.scf and as *_R_pseudo.scf, respectively.

Convert to SCF tool


You [use](#) the Convert to SCF tool to convert an .abi/.ab1 file to an .scf file, which is a smaller file format. You can also use the Convert to SCF tool to convert a .fasta file to a .scf file; however, only the first 1200 bases in the .fasta file are converted.


To use the Convert to SCF tool

1. On the Mutation Surveyor main menu, click Tools > Convert to SCF.

The Convert to SCF File window opens.

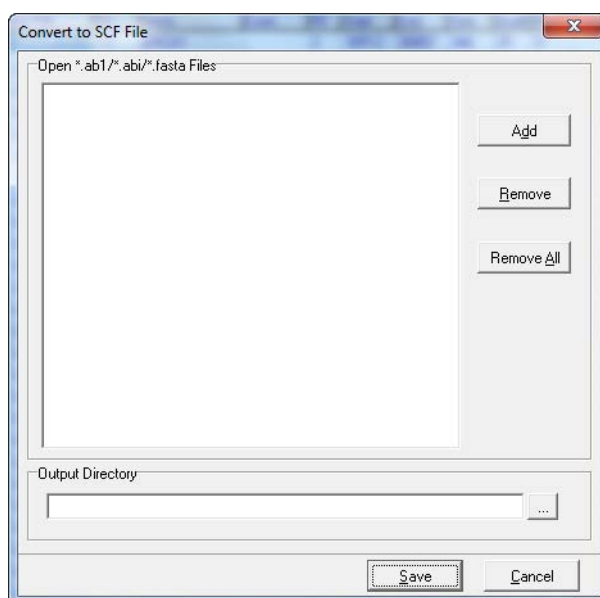Figure 9-18: Convert to SCF File window



2. Click Add to the right of the Open *.ab1/*.abi/*.fasta Files field. The Open Files dialog box opens.
Select the files that are to be converted, and then click Open to add the files to the Convert to SCF tool.
3. In the Output Directory field, leave the default location and name for the .scf file as is, or click the Browse button  to open the Browse for Folder dialog box in which you can select a different location for saving the .scf file.
4. Click Save.

Extract SGP Tool


You **use** the Extract SGP Tool to extract GenBank files, sample files, and reference files from an SGP file. The sample traces are not extracted as the originally loaded .abi/.ab1 files. Instead, they are extracted as .scf files, which is the portion of the originally loaded files that was used for analysis in the project. The reference trace is extracted whether it was a loaded trace or a generated trace. The project settings and any comments added to the project are *not* extracted.



To use the Extract SGP tool

1. On the Mutation Surveyor main menu, click Tools > Extract SGP.

The Extract SGP window opens.

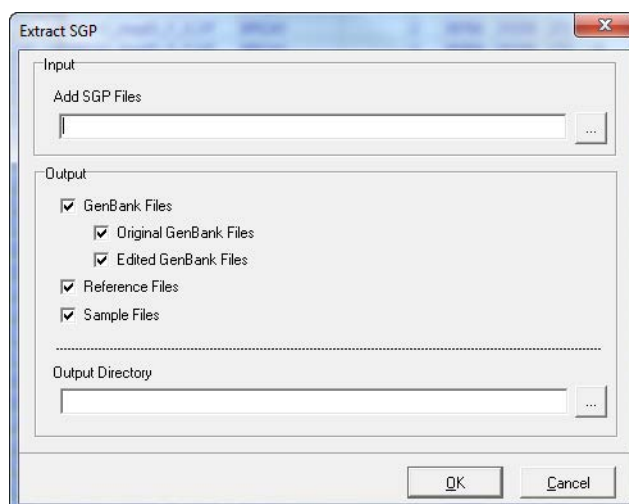Figure 9-19: *Extract SGP window*



2. In the Input pane, click the Browse button  to open the Open Mutation Project dialog box and select the mutation project files that are to be extracted.
3. In the Output pane, leave the default location and name for the extracted files as is, or click the Browse button  to open the Browse for Folder dialog box in which you can select a different location for saving the extracted files.
4. Click OK.



If you close, and then reopen Mutation Surveyor, the Extract SGP tool “remembers” the last settings specified for it.

Chapter 9

Mutation Surveyor Tools

Section 3

Assembly Tools

Section Contents

- [SGP Comparison Tool on page 279](#)
- [Mutation Assembly Tool on page 284](#)
- [Sequence Assembly Tool on page 287](#)

Chapter 9

Mutation Surveyor Tools

SGP Comparison Tool

You [use](#) the SGP Comparison tool to assemble multiple Mutation Surveyor projects and compare and contrast these projects. The SGP files to which other SGP files are being compared are the control files, or the *standard*. The SGP files that are being compared are the *unknown* files, or the sample. The report that is produced by the SGP Comparison tool is color-coded to assist in the identification of inconsistent mutations between the standard files and the unknown files. You can also use the report for somatic variant detection by comparing normal and tumor samples, comparing samples that are in the same project or multiple projects, and identifying the differences between analysts' reviews. The report is also ideal for auditing purposes.



Before you use the SGP Comparison tool, you might need to correctly group the SGP files that are being compared. See [“ND Filename Match Editor Tool” on page 264](#).

To use the SGP Comparison tool


1. On the Mutation Surveyor main menu, click Tools > SGP Comparison tool.

The SGP Comparison window opens.



If a project is already loaded in Mutation Surveyor when you open the SGP Comparison tool, then this project is automatically loaded into the tool.

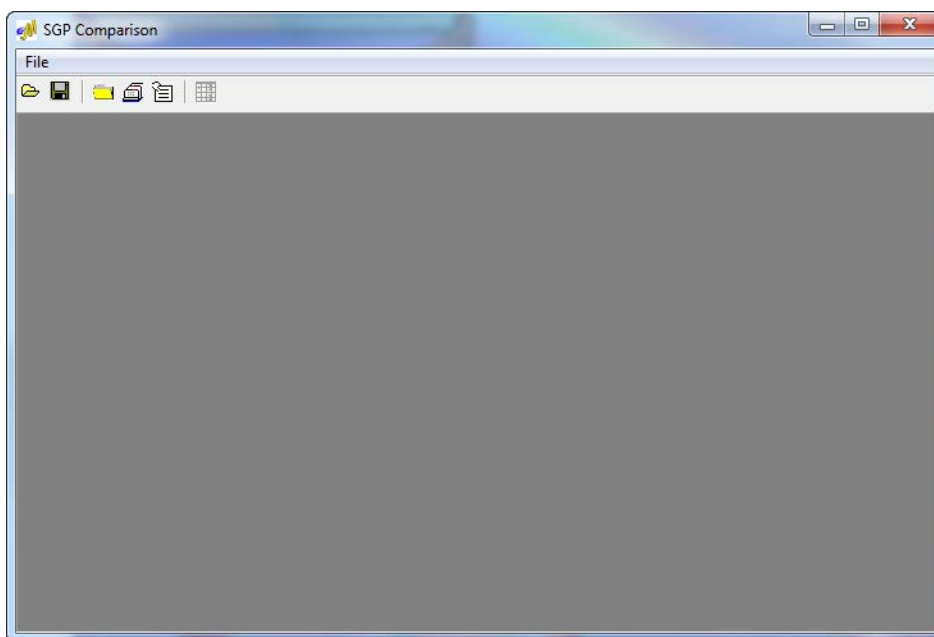
Figure 9-20: SGP Comparison window



2. On the SGP Comparison window toolbar, click the Add SGP Files button .

The Open SGP Files dialog box opens.

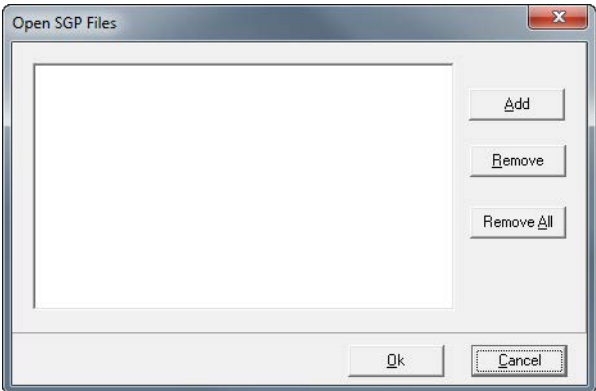Figure 9-21: Open SGP Files dialog box




3. Click Add to open the Open Files dialog box opens, select the SGP files that are to be compared, and then click OK.


The Open Files dialog box closes. The SGP files are added and displayed in the Mutation Surveyor main window.

4. On the Mutation Surveyor main menu, click Tools > SGP Comparison to bring the SGP Comparison window to the front.

All the sample files in each of the SGP files are grouped together. The separate SGP projects are highlighted in different colors.

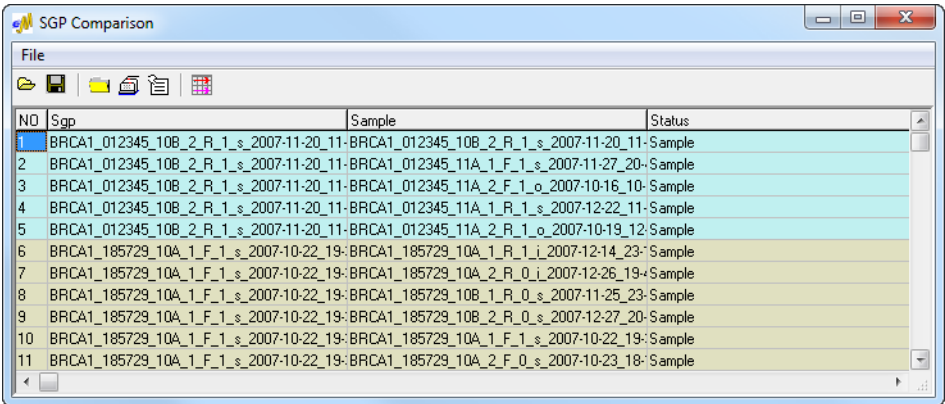
Figure 9-22: SGP Comparison window with loaded SGP files



5. To correctly group the sample files, click the Group Sample Files icon .

The Sample Grouping dialog box opens.

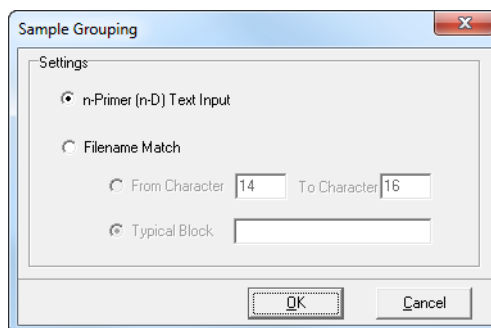Figure 9-23: Sample Grouping dialog box




6. Specify the filename substring by which Mutation Surveyor is to group all the samples files within a mutation project, and then click OK.
 - n-Primer (n-D) Text Input—Use the ND Grouping file to group the sample files. See [“ND Filename Match Editor Tool” on page 264](#).
 - Filename Match (based on one of the following):
 - From Character [] To Character []—If all the filenames in your project contain the same characters/designator in the same positions, then select this option and enter the starting and ending character positions.
 - Typical Block—If all the filenames in your project contain the same characters/designator but in different positions, then select this option and enter the value for the character/designator block.


The display in the SGP Comparison window is refreshed to reflect the groupings.

7. Click the Select Standard icon .

The Select Standard dialog box opens.

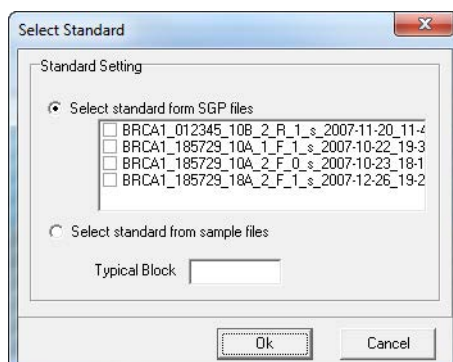Figure 9-24: Select Standard dialog box




8. Select either an entire mutation project as the standard (the default option), or click Select standard from sample files and in the Typical Block field, enter a character string that is contained in the names of all the sample files that you want to use as a standard.
9. Click OK to close the Select standard dialog box and return to the SGP Comparison window.
10. Click the SGP Comparison icon .

The SGP Comparison report opens. By default, this report shows all the mutations in all the SGP projects. The mutations are color-coded for identification.





Click Color Code link to open the Color Code dialog box which details the color coding legend used in the report.

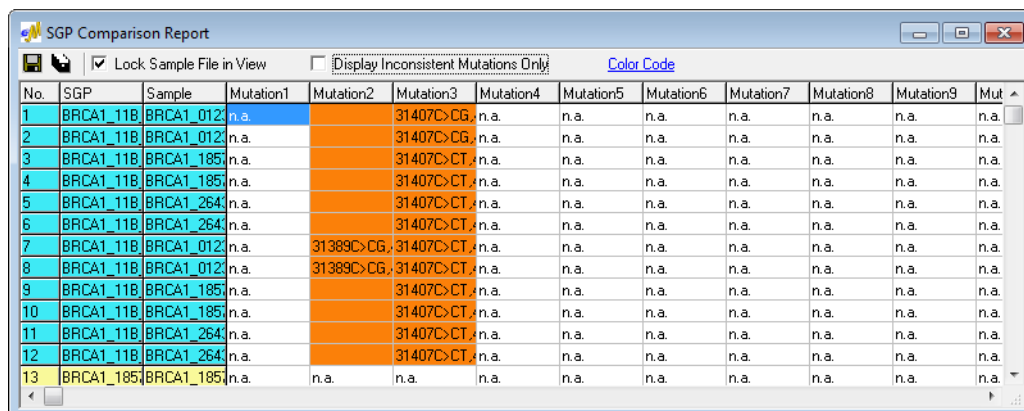Figure 9-25: SGP Comparison report



No.	SGP	Sample	Mutation1	Mutation2	Mutation3	Mutation4	Mutation5	Mutation6	Mutation7	Mutation8	Mutation9	Mut
1	BRCA1_11B	BRCA1_012	n.a.		31407C>CG	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
2	BRCA1_11B	BRCA1_012	n.a.		31407C>CG	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
3	BRCA1_11B	BRCA1_185	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
4	BRCA1_11B	BRCA1_185	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
5	BRCA1_11B	BRCA1_264	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
6	BRCA1_11B	BRCA1_264	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
7	BRCA1_11B	BRCA1_012	n.a.	31389C>CG	31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
8	BRCA1_11B	BRCA1_012	n.a.	31389C>CG	31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
9	BRCA1_11B	BRCA1_185	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
10	BRCA1_11B	BRCA1_185	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
11	BRCA1_11B	BRCA1_264	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
12	BRCA1_11B	BRCA1_264	n.a.		31407C>CT	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.
13	BRCA1_185	BRCA1_185	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.	n.a.

The report is interactive:

- To open the Color Code dialog box and view the color legend that is used in the report, click Color Code.
- To freeze the Sample File column while scrolling, leave Lock sample File in View selected.
- To view only those mutations that are inconsistent between the standard and the unknown files, click Display Inconsistent Mutations Only.
- To save the entire report, click the Save Report icon  to open the Save As dialog box in which you can specify a name and location for saving the report as well as its format (.txt, .xls/.xlsx, .xml, or .htm).
- To export only the sample files that contain inconsistent mutations as .scf files, click the Export Sample Files with Inconsistent Mutations icon  to open the Browse for Folder dialog box in which you can specify a location to which to export the files.
- To copy selected text in a cell to your clipboard, right-click and select Copy. To copy text

in a range of cells, click and hold the left mouse button and drag the mouse to select the region that you want to copy. Use standard keyboard commands or menu commands to paste the copied text into an application.

- To copy the SGP Comparison report as an image, press and hold the [Shift] key and the [Ctrl] key and then click and hold the left mouse button and draw a box around the region of the image that you want to copy. The selected region is filled with black. Right-click and select Copy to copy the image to your clipboard. Use standard keyboard commands or menu commands to paste the copied image into an application.

Mutation Assembly Tool


The Mutation Assembly Tool is especially useful for cDNA projects. You can [use](#) the tool to upload mutation project files and identify mutations in the files using cDNA nomenclature. Also, you can use the Sample Grouping options to create a consensus of all mutation calls within a single row or column for a group of sample traces. This approach is useful for merging the results of several patients that were run over multiple projects.

To use the Mutation Assembly tool

1. On the Mutation Surveyor main menu, click Tools > Mutation Assembly.

The Mutation Assembly window opens.

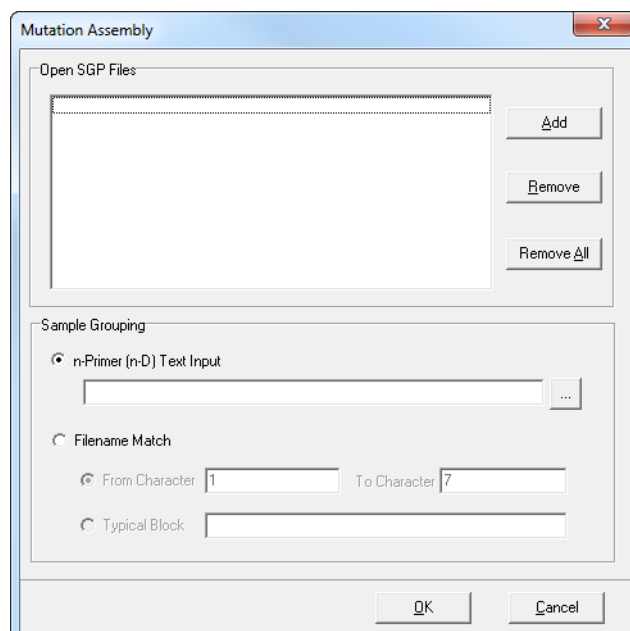Figure 9-26: Mutation Assembly window




2. Click Add to open the Open Files dialog box and select the SGP files that are to be assembled.

3. In the Sample Grouping pane, specify the filename substring by which Mutation Surveyor is to group all the samples files within the selected mutation projects, and then click OK.
 - n-Primer (n-D) Text Input—Use the ND Grouping file to group the sample files. See [“ND Filename Match Editor Tool” on page 264](#).
 - Filename Match (based on one of the following):
 - From Character [] To Character []—If all the filenames in your project contain the same characters/designator in the same positions, then select this option and enter the starting and ending character positions.
 - Typical Block—If all the filenames in your project contain the same characters/designator but in different positions, then select this option and enter the value for the character/designator block.
4. Click OK.

The Mutation Assembly report opens. This report lists all the identified mutations in the loaded mutation project files using cDNA nomenclature.

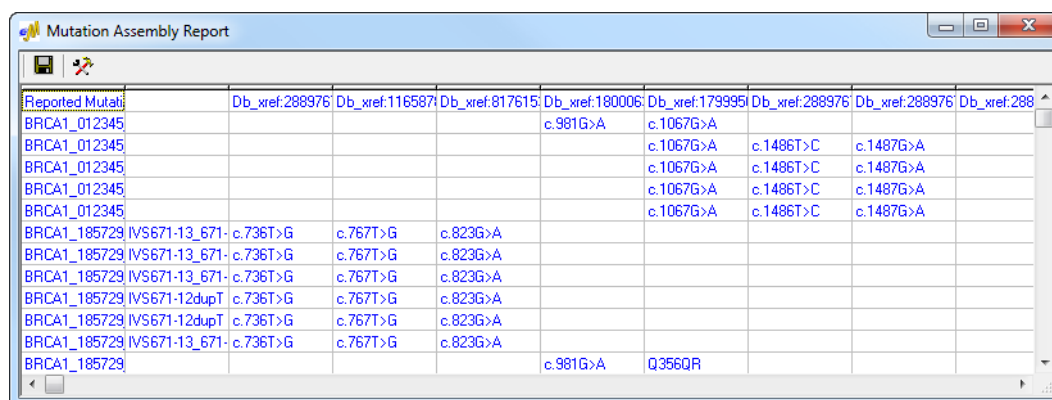Figure 9-27: Mutation Assembly report



Reported Mutati	Db_xref.288976	Db_xref.116587	Db_xref.817615	Db_xref.180006	Db_xref.179995	Db_xref.288976	Db_xref.288976	Db_xref.288
BRCA1_012345				c.981G>A	c.1067G>A			
BRCA1_012345					c.1067G>A	c.1486T>C	c.1487G>A	
BRCA1_012345					c.1067G>A	c.1486T>C	c.1487G>A	
BRCA1_012345					c.1067G>A	c.1486T>C	c.1487G>A	
BRCA1_012345					c.1067G>A	c.1486T>C	c.1487G>A	
BRCA1_185729	IVS671-13_671	c.736T>G	c.767T>G	c.823G>A				
BRCA1_185729	IVS671-13_671	c.736T>G	c.767T>G	c.823G>A				
BRCA1_185729	IVS671-13_671	c.736T>G	c.767T>G	c.823G>A				
BRCA1_185729	IVS671-12dupT	c.736T>G	c.767T>G	c.823G>A				
BRCA1_185729	IVS671-12dupT	c.736T>G	c.767T>G	c.823G>A				
BRCA1_185729	IVS671-13_671	c.736T>G	c.767T>G	c.823G>A				
BRCA1_185729				c.981G>A	Q356QR			



If you close, and then reopen Mutation Surveyor, the Mutation Assembly tool “remembers” the last settings that specified for it.

The report is interactive:



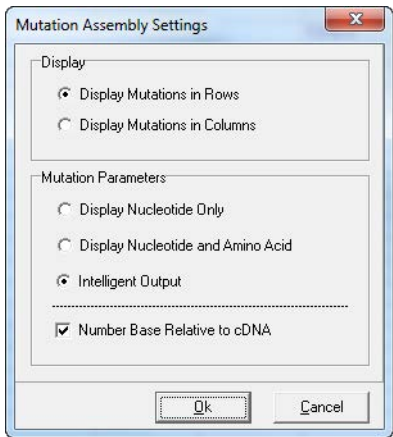
- To save the entire report as a text file, click the Save Report icon  to open the Save As dialog box in which you can specify a name and location for saving the .txt file.
- To change the information that is displayed in the report, click the Report Settings button  to open the Mutation Assembly Report Settings dialog box which contains options for filtering the information that is to be displayed in the report.

Figure 9-28: Mutation Assembly Report Settings dialog box



Option	Description
Display	<ul style="list-style-type: none">• Display Mutations in Rows—Mutations displayed as a matrix of samples versus reported mutations.• Display Mutations in Columns—Mutations displayed as a matrix of reported mutations versus samples.
Mutation Parameters	<ul style="list-style-type: none">• Display Nucleotide Only—Display only the base number and the bases involved in the mutation in the mutation code.• Display Nucleotide and Amino Acid—Display the base number, the bases involved in the mutation, and the amino acids (the normal amino acid that is affected by the mutation and the resulting amino acid) in the mutation code.• Intelligent Output—The default selection. Shows nomenclature “Relative to CDS” when the mutation call is outside of the CDS and only the resultant amino acids when the mutation call is within the CDS.
Number Base Relative to cDNA—Selected by default. The base number that is used in the mutation code is numbered relative to the cDNA region, where the beginning base for the region is Base #1, the second base in the region is Base #2, and so on. Clear this option to use the genomic base position for the base number in the mutation code.	

Sequence Assembly Tool


You **use** the Sequence Assembly to carry out de novo assembly of overlapping contigs and then save the assembled information as a .seq file. You can then use this .seq file as a GenBank file in a Mutation Surveyor project.


To use the Sequence Assembly tool

1. On the Mutation Surveyor main menu, click Tools > Sequence Assembly.

The Sequence Assembly window opens.

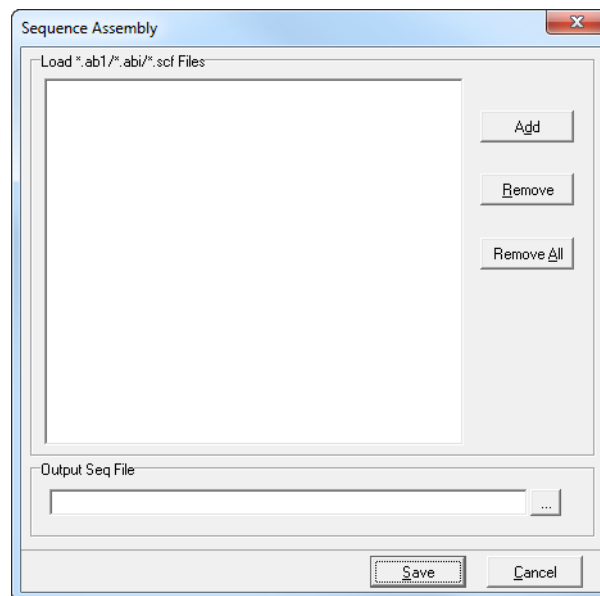Figure 9-29: Sequence Assembly window



2. Click Add to open the Open Files dialog box and select the files that are to be used for the assembly. (The files can be .scf or .abi/.ab1 files.)
3. In the Output Seq File field, leave the default location and name for the .seq file as is, or click the Browse button  next to the field to open the Save As dialog box in which you can specify a name and location for saving the .seq file.

4. Click Save.

The SEQ file is saved as one long unbroken string of nucleotides (.fasta format) in a single file.

Figure 9-30: Saved SEQ file



5. Load this file into the SEQ File Editor (see [“SEQ File Editor Tool” on page 246](#)) and save the file using the editor.

By loading and saving the SEQ file in the SEQ File Editor, the file is saved in a GenBank format, so, you can now use the file directly in Mutation Surveyor, or you can first annotate the file to identify CDSs, variations, and so on before use.

Section 4

Other Tools

Section Contents

- [Mask Vector Sequence Tool on page 291](#)
- [Output Trace Data Tool on page 293](#)
- [BasePatch Tool on page 295](#)
- [The HLA Tool on page 297](#)

Chapter 9

Mutation Surveyor Tools

Mask Vector Sequence Tool

You [use](#) the Mask Vector Sequence tool to physically modify sample files and replace base calls in the vector sequence region with N calls. These N calls are displayed as such in the GAD. Two files are created with this tool—one file with the vector region masked and one file with the insert region masked.




You can use the Mask Vector Sequence tool to replace base calls in the vector sequence region, or you can select the Vector Trim option on the Mutation Surveyor Project Settings dialog box (see [“Contig tab” on page 55](#)) and trim the sample sequence according to a starting vector sequence and an ending vector sequence that you define using the SEQ File Editor tool. (See [“SEQ File Editor Tool” on page 246](#).)

To use the Mask Vector Sequence tool

1. On the Mutation Surveyor main menu, click Tools > Mask Vector Sequence.

The Mask Vector Sequence window opens.

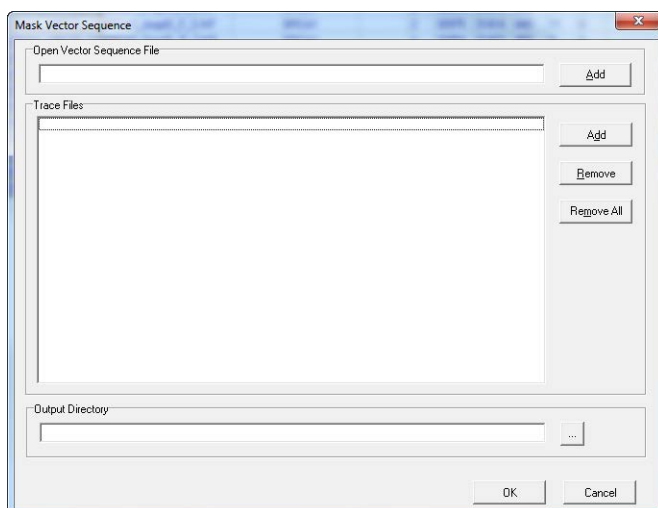Figure 9-31: Mask Vector Sequence window




2. Click Add in the Vector Sequence File pane to open the Open File dialog box and select the .seq or GBK file that contains the vector sequence.
3. Click Add in the Trace Files in the pane to open the Open File dialog box and select the trace (sample) files that are to be modified.



The tool works best when sample trace files are in the same orientation as the vector sequence.

Chapter 9

Mutation Surveyor Tools

4. In the Output Directory field, leave the default location as is, or click the Browse button  to open the Browse for Folder dialog box in which you can select a different location for saving the sample files.
5. Click OK. The output directory will now contain two copies of each sample trace file that was loaded into the tool: one copy will have the vector sequence masked (*_NN.SCF) and one copy will have the sequence of the insert masked (*_AN.SCF).
6. Open a new Mutation Surveyor project. Use the *_NN.SCF files that you saved in [Step 5](#) as the sample files and a GenBank file with the gene of interest as the GenBank file.

Output Trace Data Tool


You [use](#) the Output Trace Data tool to export information about selected traces such as frame number, base position and peak intensities to a tab-delimited text file. You can export information either by peak position or by frame.

To use the Output Trace Data tool

1. Open a Mutation Surveyor project.
2. On the Mutation Surveyor main menu, click Tools > Output Trace Data.

The Output Trace Data window opens. This window lists all the sample traces for the Mutation Surveyor project, and by default, all the traces are selected for export.

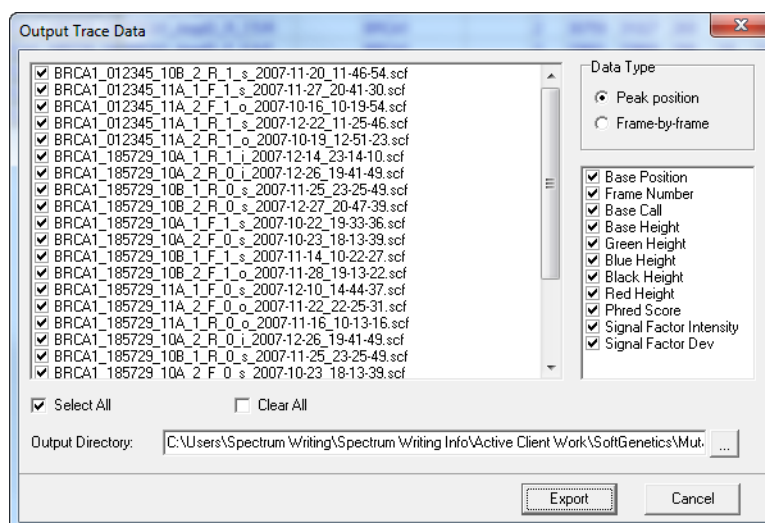Figure 9-32: Output Trace Data window



3. You can leave all sample traces selected, or you can clear the selections for the traces for which you do not want to export information.




*Click **Select All** to select all sample traces in a single step. Click **Clear All** to clear current selections in a single step.*

4. Select the way by which you want to export the data (Peak Position or Frame-by-Frame) and then *clear* the selections for the information that you do *not* want to export for each peak or each frame.




By default, all information for a data type is selected for exporting. For information about Signal Factor intensity and Signal Factor Dev for Peak Position, see [“Signal Factor Intensity and Signal Factor Dev”](#) on page 294.

5. In the Output Directory field, leave the default location and name for the text file(s) as is, or click the Browse button  to open the Browse for Folder dialog box in which you can select a different location for saving the text file(s).
6. Click Export.

The selected information is exported to a tab-delimited text file.

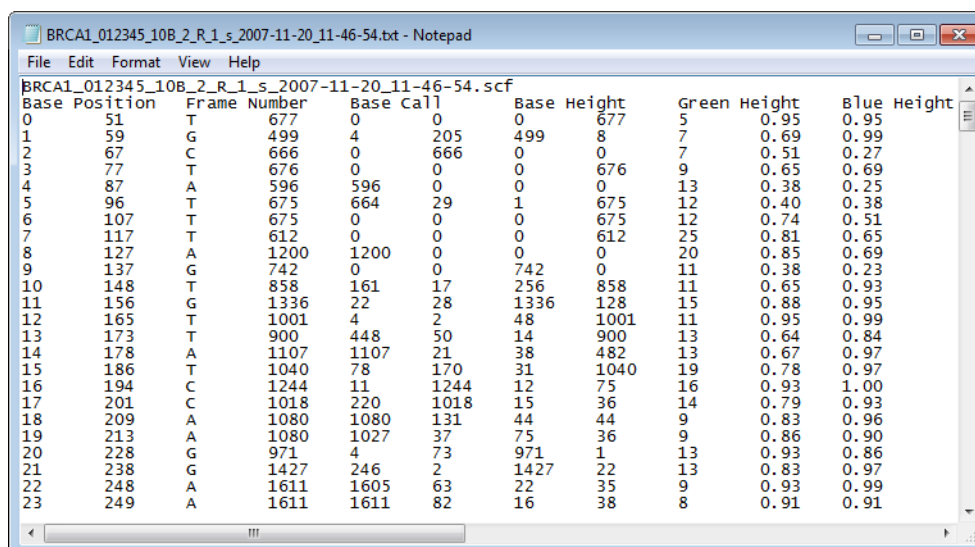Figure 9-33: Output Trace Data text file



Base Position	Frame Number	Base Call	Base Height	Green Height	Blue Height
0	51	T	677	0	677
1	59	G	499	4	205
2	67	C	666	0	666
3	77	T	676	0	676
4	87	A	596	596	0
5	96	T	675	664	29
6	107	T	675	0	675
7	117	T	612	0	612
8	127	A	1200	1200	0
9	137	G	742	0	742
10	148	T	858	161	17
11	156	G	1336	22	28
12	165	T	1001	4	2
13	173	T	900	448	50
14	178	A	1107	1107	21
15	186	T	1040	78	170
16	194	C	1244	11	1244
17	201	C	1018	220	1018
18	209	A	1080	131	44
19	213	A	1080	1027	37
20	228	G	971	4	73
21	238	G	1427	246	2
22	248	A	1611	1605	63
23	249	A	1611	1611	82

Signal Factor Intensity and Signal Factor Dev

Each base call is derived from a peak of one dye color. Secondary peaks of other dye colors might be located under this major peak. For example, if the base call for nucleotide position 100 is A:

- Signal Factor Intensity = $\text{HeightA} / (\text{HeightA} + \text{HeightT} + \text{HeightC} + \text{HeightG})$
- Signal Factor Deviation = $\text{DevA} / (\text{DevA} + \text{DevT} + \text{DevC} + \text{DevG})$, where:
 - $\text{Dev}(i) = 2.0 * \text{Height}(i) - \text{Height}(iL) - \text{Height}(iR)$ and:
 - $\text{Height}(iL)$ is the peak height 4 bases to the left, or before, position 100: $100 - 4 = \text{nucleotide position 96}$ for each peak height.
 - $\text{Height}(iR)$ is the peak height 4 bases to the right, or after, position 100: $100 + 4 = \text{nucleotide position 104}$ for each peak height.

BasePatch Tool

You **use** the BasePatch Tool to narrow and sharpen broad peaks in previously run sample traces to make the peaks more discernible in the traces, and therefore, make it easier to identify variants in the traces. You can reprocess any type of sample trace file (ABI/ABI, SCF) with the BasePatch tool.




You would consider running the BasePatch tool on sample traces for which variant identification was inconclusive or difficult because of the shapes (broadness) of their peaks.


To use the BatchPatch tool

1. On the Mutation Surveyor main menu, click Tools > BasePatch Tool.


The BasePatch Tool window opens.

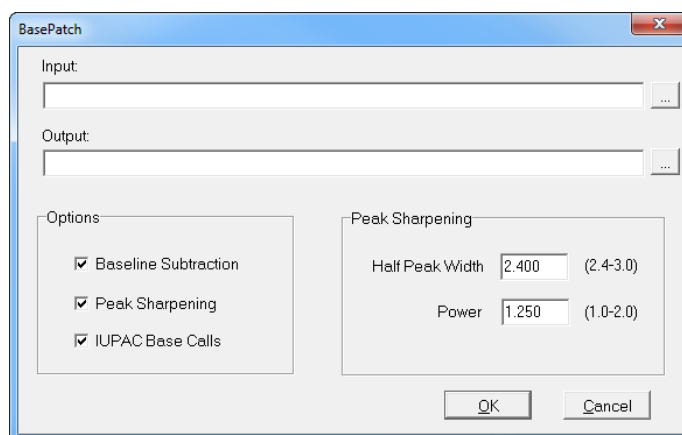Figure 9-34: BasePatch Tool window



2. Click the Browse button  next to the Input field, and browse to and select the folder that contains the sample traces that are to be reprocessed.

By default, the Output directory is set to the same as that for the Input, and the Output folder is named Processed.

3. Optionally, click the Browse button  next to the Output field, and browse to and select a different directory and/or folder in which to save the reprocessed sample traces.
4. Leave all the Options for the BasePatch tool selected.

Option	Description
Baseline Subtraction	Has a minimal affect on the reprocessing because typically the Input sample traces already have had baseline subtraction carried out.

Chapter 9
Mutation Surveyor Tools

Option	Description
Peak Sharpening	Applies the Mexican Hat Wavelet algorithm to narrow and sharpen the peaks in the sample traces. Leave the associated Peak Sharpening values (HalfPeak Width and Power) set to their default values. Note: For information about the algorithm and the Peak Sharpening values, contact tech_support@softgenetics.com for assistance.
IUPAC Base Calls	After reprocessing, the nucleotides in the sample traces are called again.

5. Click OK.

The BasePatch Tool window closes. The sample traces are reprocessed. You can now run these sample files in a Mutation Surveyor project to analyze the variants again. See [“Setting up a Single Mutation Surveyor Project” on page 70.](#)

The HLA Tool


You [use](#) the HLA Tool to analyze Human Leukocyte Antigen (HLA) data. The output includes a .dat file that can be loaded into NextGENe for additional analysis.

To use the HLA tool

1. On the Mutation Surveyor main menu, click Tools > HLA.

The HLA Tool dialog box opens.

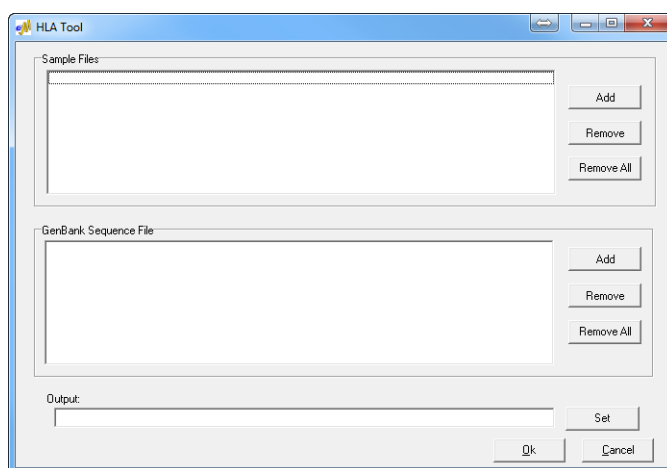Figure 9-35: HLA Tool dialog box



2. In the Samples pane, click Add to open the Open Sample File dialog box, and browse to and select the data files for analysis.
3. In the GenBank Sequence File pane, click Add to open the Open Seq/Gbk files dialog box, and browse to and select the reference GenBank file for HLA genes.



You can load multiple reference GenBank files for HLA genes.

4. In the Output field, you can leave the default value for the location of the output files as-is (the default value is the directory path for the input files), or you can click Set to select a different location.
5. Click OK.

After processing, the Mutation report is automatically displayed in the Viewing and Reporting pane of the Mutation Surveyor main window. The output includes a .dat file that can be loaded into NextGENe for additional analysis.

Chapter 10

Mutation Surveyor Specialized Applications

In addition to the specific functions for which Mutation Surveyor is designed, you can use Mutation Surveyor for some specialized applications.

This chapter covers the following topics:

- [“Methylation Detection” on page 301.](#)
- [“Mutation Quantifier Function” on page 303.](#)

Chapter 10

Mutation Surveyor Specialized Applications

Methylation Detection

A variety of options are available for analyzing methylated DNA. Mutation Surveyor is designed specifically for analyzing methylated DNA that has undergone bisulfite treatment. When you use Mutation Surveyor to analyze methylated DNA that has undergone bisulfite treatment, the following caveats apply:


- You must use a GenBank or SEQ file as the reference file.
- The GenBank or SEQ file must be loaded in the forward strand orientation. If the file is not already in this orientation, you must use the Advanced GBK File Editor tool or SEQ File Editor tool to reverse complement the sequence. See [“SEQ File Editor Tool” on page 246](#) and [“Advanced GBK File Editor Tool” on page 231](#).
- You must select Methylation on the Others tab on the Mutation Surveyor Project Settings dialog box and you must specify whether the method of detection is to be Automethylation or Set by User. Additionally, the setting to Automatically Adjust GenBank Direction, located above the Methylation section of the Others tab, must be deselected. See [“” on page 66](#).



Typically, Automethylation is the preferred option. When this option is selected, Mutation Surveyor optimizes the conditions for methylation detection which minimizes the number of mutations that are detected. By minimizing the number of detections, the analysis is simplified but its accuracy is maintained.

Bisulfite treatment of single-stranded DNA followed by PCR converts unmethylated cytosines that are outside of CpG islands to uracil. Unmethylated cytosines that are in CpG islands are also converted to uracil. Methylated cytosines are not converted. When evaluating bisulfite-treated DNA, Mutation Surveyor reads the cytosines that have been converted to uracils as thymines and methylated cytosines simply as cytosines.

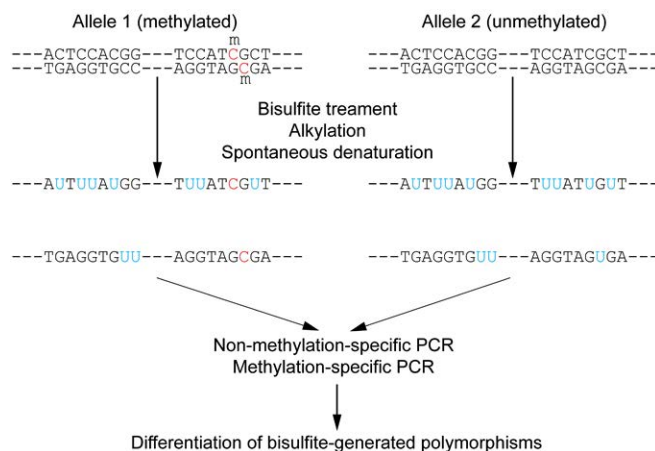Figure 10-1: Bisulfite treatment and PCR of methylated DNA



Because unmethylated sample positions in the comparison region are read as thymine, these


sample positions align to the reference file and Mutation Surveyor does not call mutations at these positions. Mutation Surveyor instead calls mutations in the comparison region only at positions where there is partial or complete methylation.



How Mutation Surveyor calls mutations for methylated DNA assumes 100% conversion of the unmethylated cytosines to uracil during bisulfite treatment. To visually evaluate and confirm this assumption, a Methylation report is available. See “Methylation report” on page 203.

Figure 10-2 is an example of a GAD for a methylation detection analysis. The top pane is the reference trace that was synthesized from the derivative GenBank file. The horizontal blue lines indicates positions where the reference file was originally a cytosine. The middle pane is the sample trace. The bottom pane is the mutation detection trace.

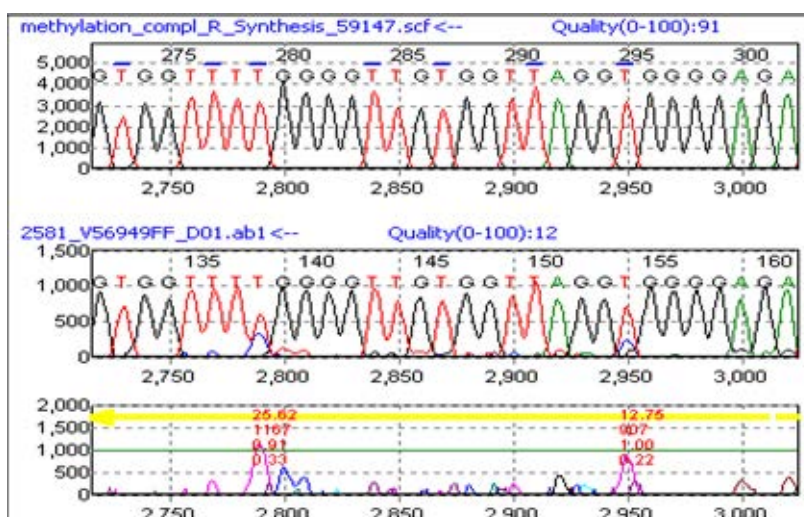Figure 10-2: GAD for methylation detection




Mutation Quantifier Function

The Mutation Quantifier function **quantifies** the allele ratios (normal allele and mutant allele) of Sanger sequencing traces at specific locations in the genome. You can use software determined standards or set your own standards for calculating the mutant peak percentage according to the following formula:


$$100 \times [\text{Mutant peak intensity} / (\text{Mutant peak intensity} + \text{Normal peak intensity})]$$

To run the Mutation Quantifier function

1. After running a Mutation Surveyor project, click on the Mutation Quantifier button  from the main toolbar.

The Mutation Quantifier dialog box opens. This dialog contains a tab for every contig in the Mutation Surveyor project.

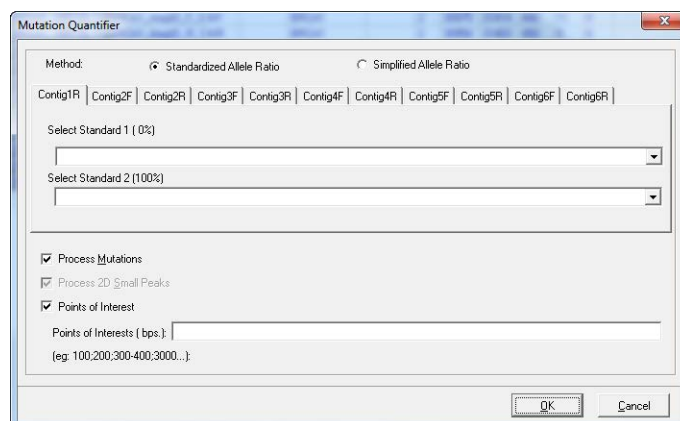Figure 10-3: Mutation Quantifier dialog box



2. Select the method by which to run the quantification—Standardized Allele Ratio or Simplified Allele Ratio.



The information that is displayed in the Mutation Quantifier report is completely different based on the method that you select. To see how two alleles vary with one sample trace, select Simplified Allele Ratio. To see how two alleles vary as compared to two Standard traces, select Standardized Allele Ratio.

3. If you selected Standardized Allele Ratio, continue to [Step 4](#); otherwise, continue to [Step 5](#).

4. For each contig in the project, do the following:
 - Standard 1—The Standard 1 dropdown list is a list of all the sample files contained in the selected contig. Standard 1 serves as one of the control files for the mutation quantification calculation. Ideally, the sample file should show a mutation peak percentage of 0%, meaning that only the normal peak is present in the sample. You can select a sample file from the Standard 1 dropdown list, or leave the Standard 1 field blank to have Mutation Surveyor select the standard file.
 - Standard 2—The Standard 2 dropdown list is a list of all the sample files contained in the selected contig. Standard 2 serves as one of the control files for the mutation quantification calculation. Ideally, the sample file should show a mutation peak percentage of 100%, meaning that the normal peak is not present in the sample. You can select a sample file from the Standard 2 dropdown list, or leave the Standard 2 field blank to have Mutation Surveyor select the standard file.



Ideally, the sample file should show a mutation peak percentage of 100% however, if necessary, you can select a heterozygous sample file that shows a mutation peak percentage of 50%.

5. Select the options for the quantification.

Option	Description
Process Mutations	Select this option to quantify substitution mutation peaks.
Process 2D Small Peaks	Select this option to quantify 2D small peaks that have been identified by Mutation Surveyor. Note: You can select this option only if Check 2D Small Peaks (Mosaic) was selected for the project. See “Check 2D Small Peaks (Mosaic)” on page 65 .
Points of Interest	Select this option to enter specific loci for quantification of mutation peaks.

6. Click OK.


The Mutation Quantifier report window opens. The information that is displayed in the report and the toolbar buttons that are available depend upon the option that you selected in [Step 2](#)—Standardized Allele Ratio or Simplified Allele Ratio. In addition, the buttons that are available on the toolbar are different. See:

- [“Mutation Quantifier report, Standardized Allele Ratio selected” on page 305.](#)
- [“Mutation Quantifier report, Simplified Allele Ratio selected” on page 307.](#)

Mutation Quantifier report, Standardized Allele Ratio selected

With Standardized Allele Ratio selected, the Mutation Quantifier report displays the percent decrease in intensity for the normal nucleotide peak and the percent increase in intensity for the mutant nucleotide peak for all identified mutations in the sample files. The report also contains a [toolbar](#) with buttons that are specific for working with the report and for changing the information that is displayed in the report.

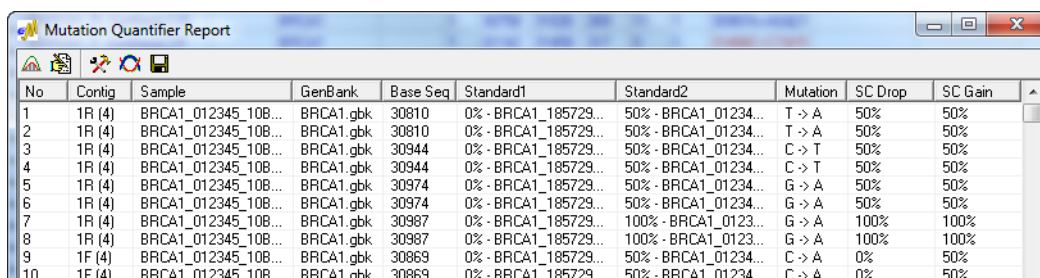Figure 10-4: Mutation Quantifier report, Standardized Allele Ratio selected



No	Contig	Sample	GenBank	Base Seq	Standard1	Standard2	Mutation	SC Drop	SC Gain
1	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30810	0% - BRCA1_185729...	50% - BRCA1_01234...	T -> A	50%	50%
2	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30810	0% - BRCA1_185729...	50% - BRCA1_01234...	T -> A	50%	50%
3	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30944	0% - BRCA1_185729...	50% - BRCA1_01234...	C -> T	50%	50%
4	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30944	0% - BRCA1_185729...	50% - BRCA1_01234...	C -> T	50%	50%
5	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30974	0% - BRCA1_185729...	50% - BRCA1_01234...	G -> A	50%	50%
6	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30974	0% - BRCA1_185729...	50% - BRCA1_01234...	G -> A	50%	50%
7	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30987	0% - BRCA1_185729...	100% - BRCA1_0123...	G -> A	100%	100%
8	1R (4)	BRCA1_012345_10B...	BRCA1.gbk	30987	0% - BRCA1_185729...	100% - BRCA1_0123...	G -> A	100%	100%
9	1F (4)	BRCA1_012345_10B...	BRCA1.gbk	30869	0% - BRCA1_185729...	50% - BRCA1_01234...	C -> A	0%	50%
10	1F (4)	BRCA1_012345_10B...	BRCA1.gbk	30869	0% - BRCA1_185729...	50% - BRCA1_01234...	C -> A	0%	50%

Column	Description
Contig	Contig number for the sample file and the direction of the sample file.
Sample	Sample filename.
GenBank	The SEQ/GBK reference file to which the sample file was compared.
Base Seq	The genomic position of the mutation in the sample file.
Standard1	The sample file used for Standard 1.
Standard2	The sample file used for Standard 2.
Mutation	The mutation shown as Normal Base > Mutant Base. Note: If the mutation is heterozygous, for example, G > GA, then only the mutant allele (in this example, A) is displayed in the Mutation Quantifier report.
SC Drop	Single Color Drop. Measures the drop in peak intensity at the normal position by normalizing the peak intensity between the reference peak from Standard 1 and the sample peak at the first four positions on both sides of the normal peak that are located greater than 1 bp away from the normal peak (referred to as the <i>local area</i>) and that have the same color as the <i>normal</i> peak.
SC Gain	Single Color Gain. Measures the increase in peak intensity at the mutant position by normalizing the peak intensity between the reference peak from Standard 2 and the sample peak at the first four positions on both sides of the variant that are located greater than 1 bp away from the mutant peak (referred to as the <i>local area</i>) and that have the same color as the <i>mutant</i> peak.
Note: Neighboring peaks at +/- 1 bps are not used for the measurement because their intensities are often affected by the mutation. Single color peaks that might contain mosaic or heterozygous mutations are also not used for the measurement. If such a peak is found in the local area, then drop or gain cannot be calculated. Note: A question mark (?) indicates inconclusive data that was the result of poor standards or poor peaks in the local area, such as mosaic or heterozygous mutations.	


The report is interactive:

- To sort the report results, click any column heading.
- To view the electropherogram for a mutation compared to the electropherograms for the standards, double-click the mutation. The mutation is marked with a red dot in the display.



In [Figure 10-5](#), the horizontal red bars were manually added to assist you with visualization. They do not normally appear in the display. For the drop in “C,” compare the Cs between Standard 1 and the sample. For the increase in “T,” compare the Ts between Standard 2 and the sample. The sample drop displays a “?” because the Cs in the local area of the sample peak are heterozygous, and reliable determination of the drop is not possible.

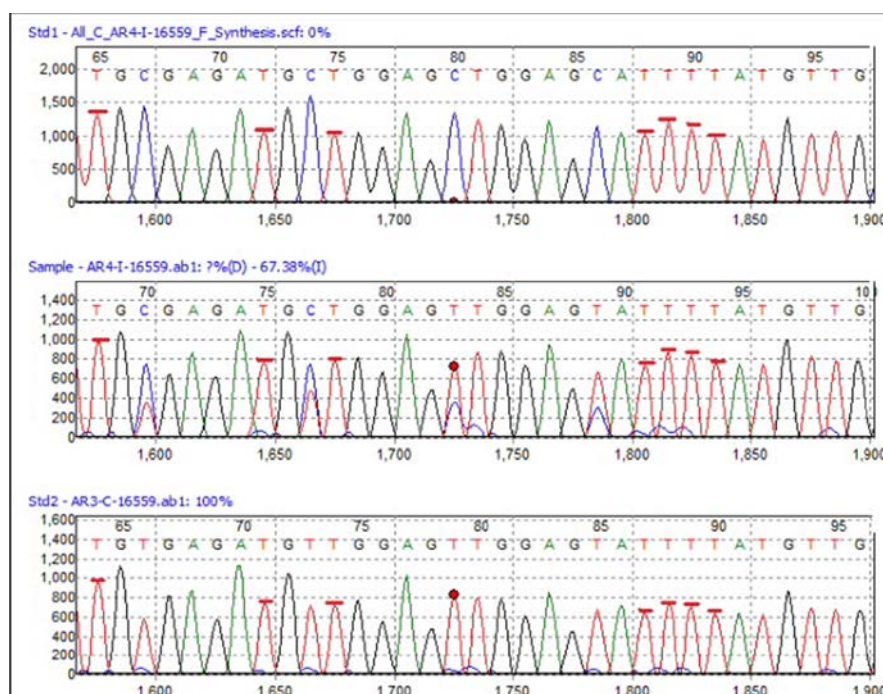
Figure 10-5: Display for mutation quantification, Standardized Allele Ratio selected



Mutation Quantifier report, Simplified Allele Ratio selected

With Simplified Allele Ratio selected, the Mutation Quantifier report displays all peak values in RFUs rather than as percentages. The report also contains a [toolbar](#) with buttons that are specific for working with the report and for changing the information that is displayed in the report.

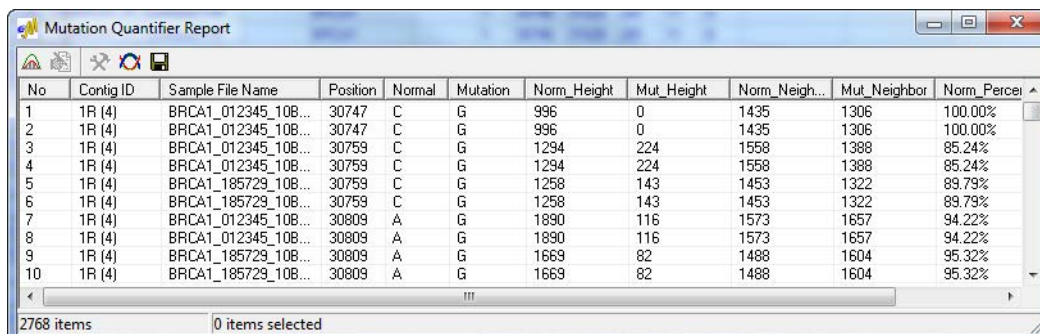Figure 10-6: Mutation Quantifier report, Simplified Allele Ratio selected



The screenshot shows a window titled "Mutation Quantifier Report" with a toolbar and a table of data. The table has 11 columns: No, Contig ID, Sample File Name, Position, Normal, Mutation, Norm_Height, Mut_Height, Norm_Neigh..., Mut_Neighbor, and Norm_Percent. The data is sorted by Norm_Percent in descending order. The first 10 rows are visible, showing mutations at positions 30747, 30759, 30759, 30759, 30759, 30809, 30809, 30809, 30809, and 30809. The status bar at the bottom indicates "2768 items" and "0 items selected".


No	Contig ID	Sample File Name	Position	Normal	Mutation	Norm_Height	Mut_Height	Norm_Neigh...	Mut_Neighbor	Norm_Percent
1	1R (4)	BRCA1_012345_108...	30747	C	G	996	0	1435	1306	100.00%
2	1R (4)	BRCA1_012345_108...	30747	C	G	996	0	1435	1306	100.00%
3	1R (4)	BRCA1_012345_108...	30759	C	G	1294	224	1558	1388	85.24%
4	1R (4)	BRCA1_012345_108...	30759	C	G	1294	224	1558	1388	85.24%
5	1R (4)	BRCA1_185729_108...	30759	C	G	1258	143	1453	1322	89.79%
6	1R (4)	BRCA1_185729_108...	30759	C	G	1258	143	1453	1322	89.79%
7	1R (4)	BRCA1_012345_108...	30809	A	G	1890	116	1573	1657	94.22%
8	1R (4)	BRCA1_012345_108...	30809	A	G	1890	116	1573	1657	94.22%
9	1R (4)	BRCA1_185729_108...	30809	A	G	1669	82	1488	1604	95.32%
10	1R (4)	BRCA1_185729_108...	30809	A	G	1669	82	1488	1604	95.32%

Column	Description
Contig	Contig number for the sample file and the direction of the sample file.
Sample	Sample filename.
Position	The genomic position of the mutation in the sample file.
Normal	The normal base as reported for the reference.
Mutation	The mutant base as reported in the GAD. Note: If the mutation is heterozygous, for example, G > GA, then only the mutant allele (in this example, A) is displayed in the Mutation Quantifier report.
Norm_Height	The peak height (in RFUs) of the normal allele.
Mut_Height	The peak height (in RFUs) of the mutant allele.
Norm_Neighbors	Average peak height (in RFUs) for the first four base positions on both sides of the normal peak that are located greater than 1 bp away from the normal peak (referred to as the <i>local area</i>) and that have same color as the <i>normal</i> peak.
Mut_Neighbor	Average peak height (in RFUs) for the first four base positions on both sides of the mutant peak that are located greater than 1 bp away from the mutant peak (referred to as the <i>local area</i>) and that have same color as the <i>mutant</i> peak.
Norm_Percent	Normal peak percentage.
Mut_Percent	Mutant peak percentage.

The report is interactive:

- To sort the report results, click any column heading *other than* Mut_Height, Norm_Neighbor, Mut_Neighbor, or Mut_Percent.
- To view the electropherogram for a mutation, double-click the mutation. The mutation is marked with a red dot in the display.

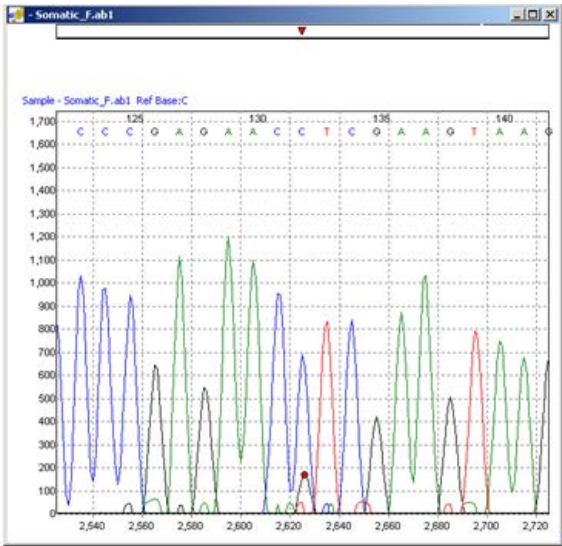Figure 10-7: Display for mutation quantification, Simplified Allele Ratio selected








Mutation Quantifier toolbar

The Mutation Quantifier report toolbar contains buttons that are specific for working with the report and for changing the information that is displayed in the report.

Figure 10-8: Mutation Quantifier report toolbar




Button	Description
	Quantify—Opens the Mutation Quantifier dialog box. See “To run the Mutation Quantifier function” on page 303.
	Edit Groups—Available only when Standardized Allele Ratio is selected. Opens the Quantification Group Editor dialog box You can use the functions on this dialog box to manually select the sample files for Standard 1 and Standard 2 and recalculate the mutation quantification based on the settings (Process Mutations, and so on) that you previously selected in the Mutation Quantifier dialog box. See “Quantification Group Editor dialog box” on page 309.

Button	Description
	Report Settings—Available only when Standardized Allele Ratio is selected. Opens the Quantifier Report Settings dialog box which contains options for filtering the information that is to be displayed in the report. See "Mutation Quantifier report settings dialog box" on page 310.
	Display Mutations with 1D/2D Settings—Toggles the GAD between the 1D and 2D views. Note: With 1D settings selected, more low confidence mutation calls might be included in reports.
	Save Report—Opens the Save Report As dialog box in which you can specify a name and location for saving the report as a .txt file.

Quantification Group Editor dialog box

You can use the functions on the Quantification Group Editor dialog box to manually select the sample files for Standard 1 and Standard 2 and recalculate the mutation quantification based on the settings (Process Mutations, and so on) that you previously selected in the Mutation Quantifier dialog box.

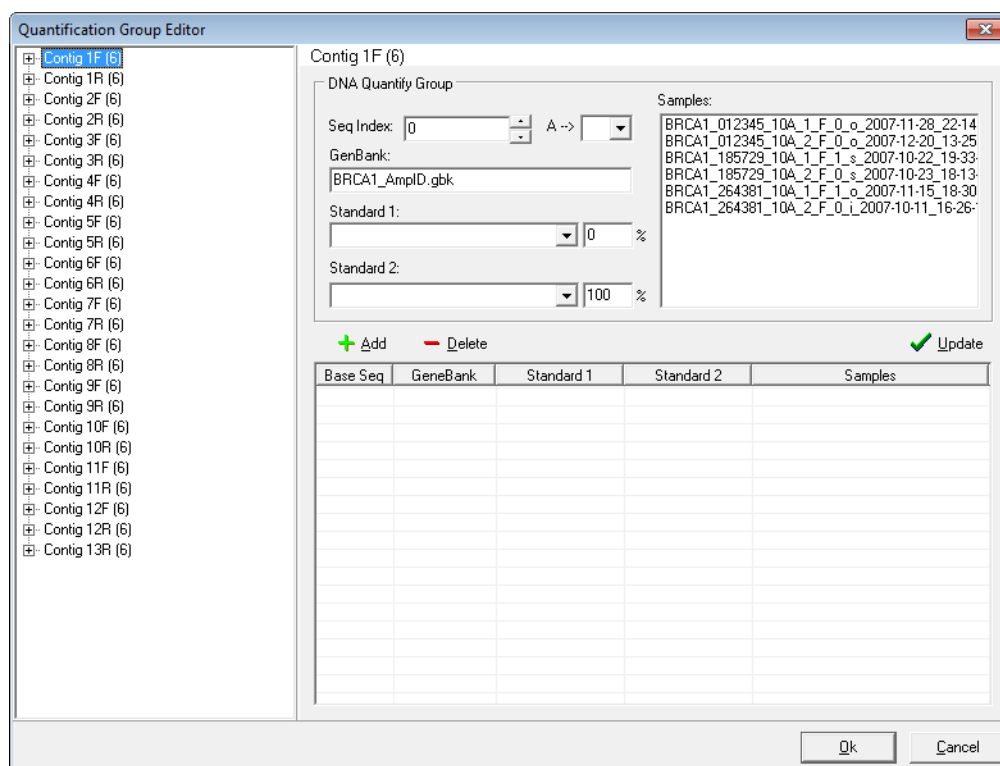Figure 10-9: Quantification Group Editor dialog box



1. In the Seq Index field, enter the contig position of a known mutation in the selected mutation project and press [Enter].

Mutation Surveyor automatically identifies the base change and in the Samples pane, displays and highlights all the sample files that contain the specific mutation. Mutation Surveyor also automatically identifies the GenBank file to which the sample files were compared and selects the sample files for Standard 1 and Standard 2.


2. Do one of the following:

If you are satisfied with Mutation Surveyor's automatic selections:	<ol style="list-style-type: none">1. Click Add to update the Mutation Quantifier table with these selections.2. Click OK to recalculate the mutation quantifications based on the settings that you previously selected in the Mutation Quantifier dialog box.
If you are not satisfied with Mutation Surveyor's automatic selections:	<ol style="list-style-type: none">1. Make any changes as needed (for example, select a different sample file for Standard 1).2. Click Update to update the Mutation Quantifier table with these selections.3. Click OK to recalculate the mutation quantifications based on the settings that you previously selected in the Mutation Quantifier dialog box.

Mutation Quantifier report settings dialog box

The Quantifier Report Settings dialog box contains options for filtering the information that is to be displayed in the report.

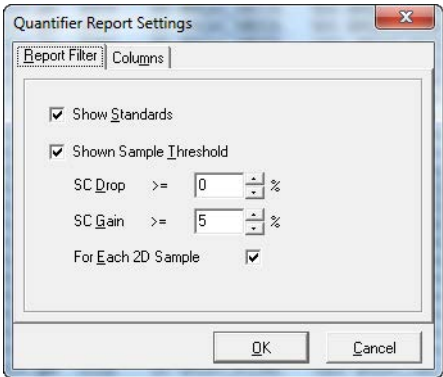Figure 10-10: Quantifier report settings dialog box



Option	Description
Report Filter tab	
Show Standards	Show the sample files used for Standard 1 and Standard 2 in the report.

Option	Description
Shown Sample Threshold	<ul style="list-style-type: none"> • SC Drop—Show only those samples with an SC Drop greater than or equal to the indicated value. • SC Gain—Show only those samples with an SC Gain greater than or equal to the indicated value. <p>Note: A sample must meet both of the specified thresholds to be displayed in the report.</p> <ul style="list-style-type: none"> • For Each 2D Sample—Selected by default. The quality of the sample pair at the same position is checked. When the SC Drop and the SC Gain of both pairs is met, then both values are displayed; otherwise, both values are hidden.
Columns tab —Select which columns are to be displayed in the Mutation Quantifier report.	

Chapter 10

Mutation Surveyor Specialized Applications

Chapter 11

Multiple Projects

Mutation Surveyor also provides the capabilities for modifying existing projects, and for working with multiple projects. For example, you might need to merge multiple projects into one project for comparison purposes, or you might need to carry out batch processing of multiple projects.

This chapter covers the following topics:

- [“Modifying Mutation Surveyor Projects” on page 315.](#)
- [“The Log File Wizard in the Mutation AutoRun Tool” on page 317.](#)
- [“The Mutation AutoRun Tool” on page 324.](#)

Chapter 11


Multiple Projects

Modifying Mutation Surveyor Projects

The Modify Mutation Projects function is available on the Mutation Surveyor main menu (File > Modify Mutation Projects) and it is also available from a button on the Project Display subreport toolbar. (See [“Project Display subreport toolbar” on page 113.](#)) The Modify Mutation Projects function provides the capabilities for modifying single or multiple projects in Mutation Surveyor. For example, you can use the Modify Mutation Projects function for the following purposes:

- To open a single project with the GenBank file that was initially used, so that you can specify sample grouping options and rerun the project.
- To select a different isoform for a single project so that you can apply the gene annotations relative to the new isoform to all mutation calls. This allows you to complete the analysis of the same project with a different isoform without needing to review your mutation calls a second time.
- To merge multiple projects into one project for comparison purposes and view the project results in a single report.

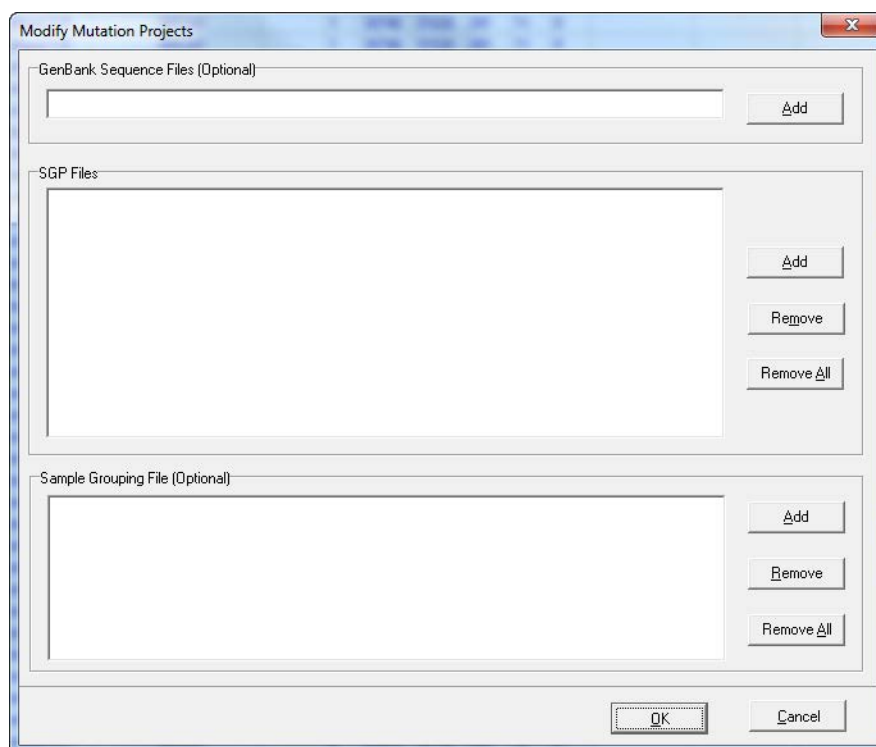Figure 11-1: Modifying Mutation Projects window



Chapter 11

Multiple Projects

Field	Description
GenBank Sequence Files (Optional)	Use the options in this pane to load a different GenBank file than what was loaded for the original project. The new GenBank file must have the same nucleotide sequence as the original GenBank file, but it can have different annotations.
SGP files	<p>Use the options in this pane to add one or more mutation projects that are being analyzed. Multiple projects are merged into a single project and the project results are contained in a single report.</p> <p>Note: If you load multiple projects and the total number of samples is ≤ 2000, then you can view the project results in the GAD and generate reports for the project. If the you load multiple projects and the total number of samples is > 2000, then you cannot view the project results in the GAD. You can only generate reports for the project.</p>
Sample Grouping File (Optional)	Use the options in this pane to load a Sample Grouping file which is a text file that defines how your sample files should be paired in your Mutation Surveyor project. See "2D Filename Match Editor tool" on page 261 .


The Log File Wizard in the Mutation AutoRun Tool

You use the Mutation AutoRun tool to carry out the batch analysis of multiple projects, where each project is referred to as a *job*. The tool scans for queued files at a user-set interval. The parameters for processing the projects are specified in a log file that you create [using](#) the Log File Wizard. You can create a job to process samples that have already been generated by your capillary electrophoresis instrument or you can create a job to process samples that have not yet been generated by your capillary electrophoresis instrument.

To open the Log File Wizard


1. On the Mutation Surveyor main menu, click Tools > Mutation AutoRun to open the Mutation AutoRun window.

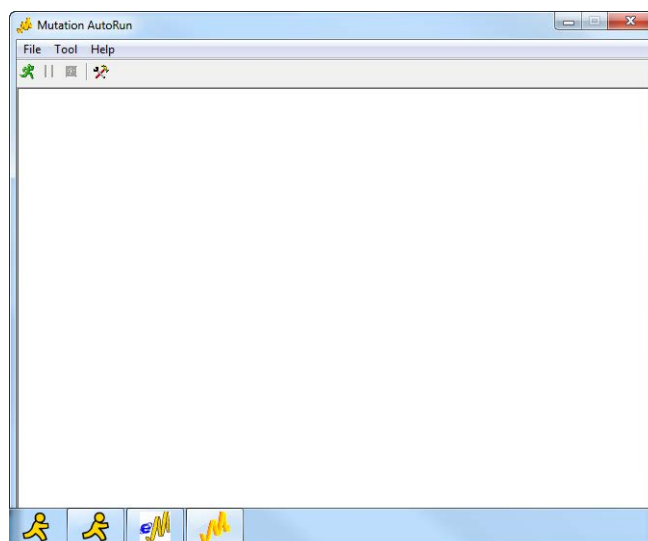Figure 11-2: Mutation AutoRun window



2. On the Mutation AutoRun menu, click Tool > Log File Wizard to open the Log File Wizard. The first page that opens is the Progress page. See [Figure 11-3 on page 318](#).

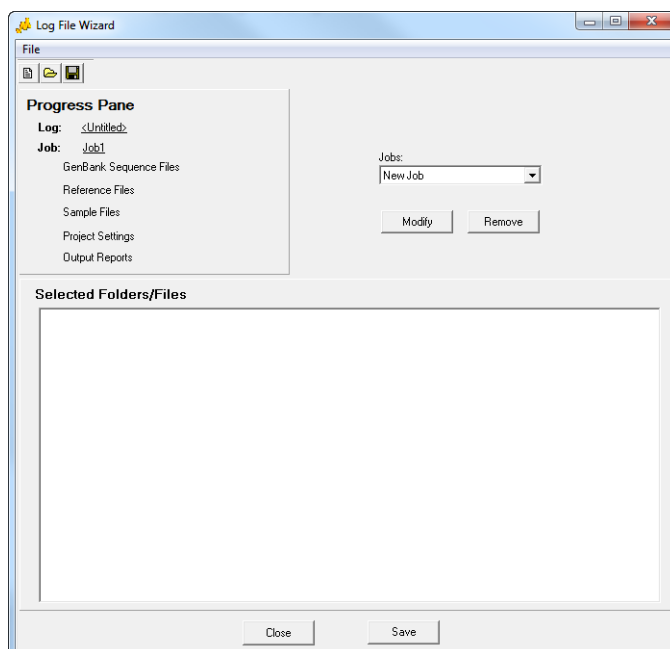Figure 11-3: Log File Wizard, Progress page



The Log File Wizard is a standard wizard consisting of multiple pages that are linked by Next and Back buttons. Each page contains instructions to guide you through the creation of a log file. After you complete the steps on a page, you click Next to move to the next page. At anytime, you can click Back as many times as needed and modify your selections for a previously completed step or steps. In addition to the standard Next and Back buttons, the Log File Wizard has page-specific links on the Progress Pane that you can double-click to open the indicated page. (These links are listed in the same order in which the pages open when you click Next.) To the right of the Progress Pane is a Jobs dropdown list of all the job files that are contained in a selected log file. Checkmarks are displayed next to the page links on the Progress Pane if the indicated information has been added for a selected job. For example, if report setting files have been added for a selected job, then a checkmark is displayed next to the Reports link.

You have two options for creating a log file to use with the Mutation AutoRun tool:


- You can set up a new log file. See [“To set up a new log file”](#) below.
- You can load a previously saved log file and modify it as needed. See [“To modify a log file”](#) on page 322.

To set up a new log file

1. On the Jobs dropdown list that is to the right of the Progress Pane, select New Job.
2. Click Modify.
3. At the prompt, enter the name for the new job.

4. Click Next.

You can now continue setting up the log file from scratch. You are prompted to enter files for a project in the following order—GenBank, Reference, and Sample. Sample files are always required for a project. GenBank files and reference files are both optional files for a job. You can add just a GenBank file, just a reference file, both files, or neither file. If you do not add either a GenBank file or a reference file, Mutation Surveyor connects to the GenBank database and searches for a GenBank file that has a sequence that matches the project's sample files. Adding GenBank files, reference files or sample files, or any other type of file (for example, custom report settings files) involves the same general series of steps:

5. Click the Browse button  on the Load Files page to open the Browse for Folder dialog box and browse to and select the folder that contains the files that are to be added to the job.

All the files that are in the selected folder are displayed in the Available Local Files pane. After you select the folder that contains the files that you want to add, the Local Path field is populated with the full directory path to the files. You can click Copy Path to copy this information to your clipboard and you can then paste it into the Local Path field in other windows if needed.



*You can create a job to process samples that have not yet been generated by your capillary electrophoresis instrument. By specifying the path where the sample files are to be located followed by *.*, all files within the folder will be processed. Because you can configure the Mutation AutoRun function to process log files at different time intervals, you can set the interval to a time that occurs after the sample files are ready for processing. For more information about delaying the start of Mutation AutoRun, see the Time option detailed in [“To use the Mutation AutoRun tool” on page 324](#).*

6. Do one of the following to add the necessary files to the job:

- Double-click each file that you want to add.
- CTRL-click to select multiple non-contiguous files or SHIFT-click to select multiple contiguous files, and then click Add > Selected.
- Click Add > All to add all the files in a single step.




When Add > All is selected, the Selected Folders/Files field contains an entry that includes the path to the folder followed by “.*.” This indicates that all contents from this folder of the appropriate file type are to be used in the analysis. If you add new files to this folder prior to running the job, the new files are also processed.*

7. After adding all the necessary files for the project, click Next.

The Input Project Settings page opens.

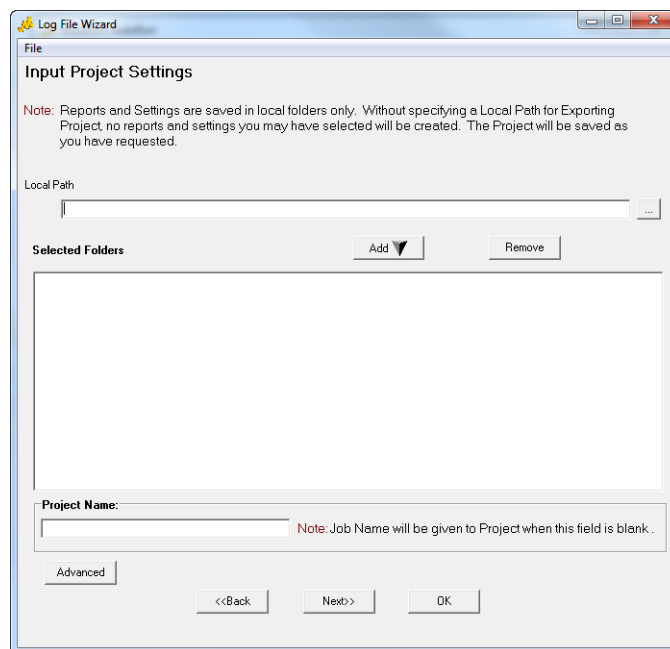Figure 11-4: Input Project Settings page





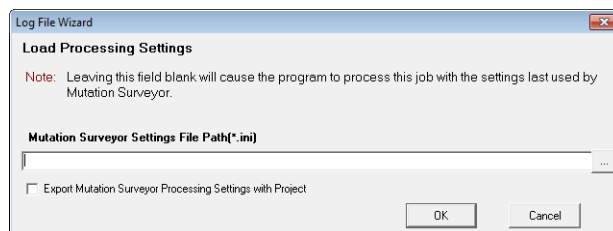
8. Click the Browse button  page to open the Browse for Folder dialog box and browse to and select the folder that is to contain the processed SGP project file, and then click Add. Any automatically generated reports are also saved to this location.
9. Optionally, you can also do one or more of the following:
 - In the Project name field, enter a name for the project.
 - Click Advanced to open the Load Process Settings dialog box, and then click the Browse button  page to open the Load Project Settings dialog box and browse to and select the settings file (.ini) file that specifies the settings that you want to use for the project.

Figure 11-5: Load Project Settings dialog box






Although specifying project settings is optional, if you do not specify any settings for the project, then Mutation Surveyor processes all the projects in the job according to the settings that were last used by Mutation Surveyor. Unless you modify the values again, all new projects are carried out according to the options that you have specified. If you load an existing project, the options revert to those set for the project. If you run a new project after loading an existing project, the new project is carried out according to the existing project settings.

- Click Export Mutation Surveyor Process Settings with Project.
10. Click OK to close the Process Settings dialog box, and then click Next.

The Input Report Information page opens.

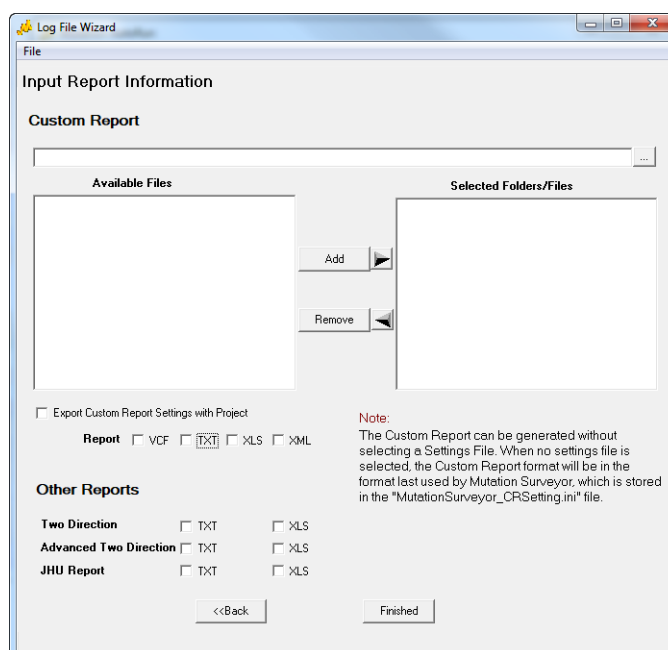Figure 11-6: Input Report Information page



11. Optionally, do one or more of the following:

- If you want to generate one or more custom reports for the project, add the appropriate custom report settings files (.ini) to the project.



If you want to create a custom report for the project, but want to use the default settings in the MutationSurveyor_Settings.ini file, then you do not need to add any settings files to the project. Instead, you can simply select the format (VCF, TXT, XLS, or XML) for the custom report.

- Click Export Custom Report Settings with Project.
- Select the format (TXT or XLS) for the other reports (Two Direction, Advanced Two Direction, or JHU) that you want to generate for the project.

12. Click Finished.

You return to the Progress Page.

13. Click Save.

The Save Log File dialog box opens.

14. Specify a name and location for the job's log file.



The log file has the extension.log and you cannot change this.

To modify a log file

When you modify a log file, you can modify the information for an existing job in the log file, you can add a new job to the log file, and you can remove a job from the log file.

1. On the Log File Wizard main menu, click File > Open.

The Open Log File dialog opens.

2. Select the log file that is to be modified, and then click Open.

You return to the Progress page. By default, the *first* job in the log file is loaded into the Log File Wizard. Checkmarks are displayed next to the page links if the information has been added for the job. For example, if report setting files have been added for the first job, then a checkmark is displayed next to the Reports link.

3. Continue to one of the following:

- [“To modify an existing job in the log file.”](#)
- [“To add a new job to a log file” on page 323.](#)
- [“To remove a job from a log file” on page 323.](#)

To modify an existing job in the log file

1. On the Jobs dropdown list that is to the right of the Progress Pane, select the job that is to be modified.

The job is loaded into the Log File Wizard. A checkmark is displayed next to a page link if the information has been added for the job. For example, if report setting files have been added for the selected job, then a checkmark is displayed next to the Reports link.

2. Modify the information as needed for the loaded job.



To modify the name for the job, click Modify on the Progress page, and at the prompt, modify the name for the job. You can use the Next and Back buttons to move through the necessary pages for modifying the job, or you can double-click a page link in the Progress Pane to open the indicated page.

3. After you have modified the information for the job, click OK on any page in the Log File Wizard.

You return to the Progress page.

4. Click Save on the Progress page to save the modifications to the job.

To add a new job to a log file

1. On the Jobs dropdown list that is to the right of the Progress Pane, select New Job.
2. Click Modify.
3. At the prompt, enter the name for the new job.
4. Click Next.
5. Add the information for the new job.
6. After you have added the information for the new job, click OK on any page in the Log File Wizard.

You return to the Progress page.

7. Click Save on the Progress page to save the new job to the log file.

To remove a job from a log file

1. On the dropdown list that is to the right of the Progress Pane, select the job that is to be deleted.
2. Click Remove.
3. Click Yes at the prompt about removing the log file.

You return to the Progress page.

4. Click Save on the Progress page to save the changes to the log file.

The Mutation AutoRun Tool

You [use](#) the Mutation AutoRun tool to carry out the batch analysis of multiple projects, where each project is referred to as a *job*. The function scans for queued files at a user-set interval. With this function, Mutation Surveyor can analyze up to 8 billion base pairs of data in eight hours.



Remember, you can also use the Mutation AutoRun tool to schedule the processing of a project where the sample files have not yet been created. See [“The Log File Wizard in the Mutation AutoRun Tool”](#) on page 317.

To use the Mutation AutoRun tool

1. On the Mutation Surveyor main menu, click Tools > Mutation AutoRun

The Mutation AutoRun window opens.


Figure 11-7: Mutation AutoRun window



2. On the Mutation AutoRun toolbar, click the Settings icon

The Mutation AutoRun Settings dialog box opens.

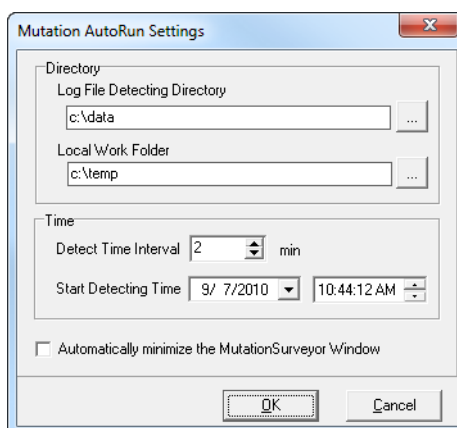Figure 11-8: Mutation AutoRun Settings dialog box



3. Specify the settings for the autorun.

Option	Description
Log File Detecting Directory	The directory in which you saved the log file. See Step 7 and Step 8 of “To set up a new log file” on page 318 in “The Log File Wizard in the Mutation AutoRun Tool” on page 317.
Local Work Folder	The temporary directory in which the analyzed data is being processed.
Time	<ul style="list-style-type: none"> • Detect Time Interval—The time interval between searches. (When Mutation Surveyor searches for log files to process.) • Start Detecting Time—The starting date and time for the search.
Minimize to Taskbar	When the Mutation AutoRun function starts, it opens Mutation Surveyor. Select this option to automatically minimize the Mutation Surveyor window after it opens to the taskbar.

4. Click OK to close the Mutation AutoRun Settings dialog box.
5. Click File > Detect.

On the specified date and time, Mutation Surveyor begins looking for unprocessed log files. It looks for these log files according to the specified time intervals, for example, every 10 minutes. When an unprocessed log file is detected, Mutation Surveyor opens the file and processes the project data according to the instructions detailed in the log file. Mutation Surveyor saves the data to the designated output folder.

Chapter 11

Multiple Projects

Chapter 12

Reviewing and Editing Mutation Projects

Mutation Surveyor automatically identifies mutation calls at positions where the variation between the reference trace and the sample trace meet or exceed detection thresholds. Mutation Surveyor also sometimes deletes mutation calls at positions that do not meet specific criteria. As a user, however, you have the option of manually adding or editing mutation calls in your mutation project, and you also have the option of overriding Mutation Surveyor's automatic actions. You can also create a custom database of known deletions that are based on the variants in a selected mutation project, and you can then use this database in other mutation projects.

This chapter covers the following topics:

- [“Adding Mutation Calls” on page 329.](#)
- [“Editing Mutation Calls” on page 335.](#)
- [“Annotating Reported Variations” on page 341.](#)
- [“Creating a User Knowledge Database for Known Deletions” on page 342.](#)
- [“Viewing the Edit History \(Audit Trail\) for a Mutation” on page 344.](#)

Chapter 12

Reviewing and Editing Mutation Projects

Adding Mutation Calls

Mutation Surveyor automatically identifies mutation calls at positions where the variation between the reference trace and the sample trace meet or exceed detection thresholds. Options are also available, however, for you to manually add mutation calls to a mutation trace. You can [add](#) substitution mutation calls, you can [add](#) homozygous indel mutation calls, and you [can](#) add heterozygous indel mutation calls.

To add a substitution

1. Open the mutation project in the GAD or in the Project Reviewer.
2. Right-click the mutation trace, and on the menu that opens, click Add Substitution.

The Add Substitution dialog box opens.

Figure 12-1: Add Substitution dialog box

The screenshot shows the 'Add Substitution' dialog box. It is divided into two main sections. The top section, 'Mutation Parameters', contains several input fields: Intensity (448), Area (1.73), Drop (0.24), Overlap (0.83), Noise (1000.00), SNRatio (1.00), Score (0.00), Frame (1265), Base# (30746), Mutation (C>AC), Comments, and Auto Comments (<<Added>>). The bottom section, 'db_xref', contains a table with two columns: 'Tag' and 'Identifier'. The table is currently empty. At the bottom of the dialog are 'OK' and 'Cancel' buttons.

3. In the Mutation field, enter the mutation, and optionally, add comments for the added mutation.



Mutation Surveyor automatically calculates the optimal values for the mutation parameters such as intensity, area, and so on for the mutation based on the position of the mutation and you should not modify these values. When you add a substitution, it can be heterozygous (for example, $A > AT$), or it can be homozygous (for example, $A > T$).

4. Optionally, to add multiple tags for the substitution, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.
5. Click OK.

The mutation trace is updated with the substitution. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.

To add a homozygous indel call

1. Open the mutation project in the GAD.
2. Press and hold the [Ctrl] key, and then click the mobility line at the position of the indel in the mutation trace.

The Add Indel dialog box opens.

Figure 12-2: Add Indel dialog box

The 'Add Indel' dialog box is shown with the following fields and values:

- Parameters:**
 - Mutation Type: Deletion (dropdown)
 - Start Index of Genbank: 134708 (text field)
 - Mutation Letters: A (text field)
 - Comments: (empty text field)
 - Auto Comments: <<Added> (text field)
- db_xref:**
 - Tag: (dropdown menu)
 - Identifier: (text field)
- Tab:** Heterozygous InDel
- Buttons:** OK, Cancel

3. Add the Indel call as appropriate.
 - If you are adding a homozygous insertion, select Insertion on the Mutation Type dropdown list, and then in the Start of GenBank Index field, enter the positions that flank the insertion as two numbers delimited by an underscore. For example, 30746_30747 means that an insertion exists between base 30746 and base 30747. The length of the insertion is determined by the Mutation Letters parameter.
 - If you are adding a homozygous deletion, then in the Start of GenBank Index field, enter the actual positions of the deletion. For example, 30746_30748 means that three bases were deleted from the sample—the base at position 30746, the base at position 30747, and the base at 30748.
4. Optionally, to add multiple tags for the indel, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.

5. Click OK.

The mutation trace is updated with the homozygous indel. The Mutation table in the GAD is updated to show this new mutation.



If you add a homozygous insertion and the bases that are inserted are identical the bases that precede them, then the mutation code is corrected in the Mutation table to reflect this duplication. See “Insertion, duplication, and deletion mutation codes” on page 131.

To add or edit a heterozygous indel (The Het-Indel tool)

You use the Heterozygous Indel Detection tool (Het-Indel tool) to add or edit heterozygous mutation calls to a mutation trace. You can open the Het-Indel tool in one of two ways:


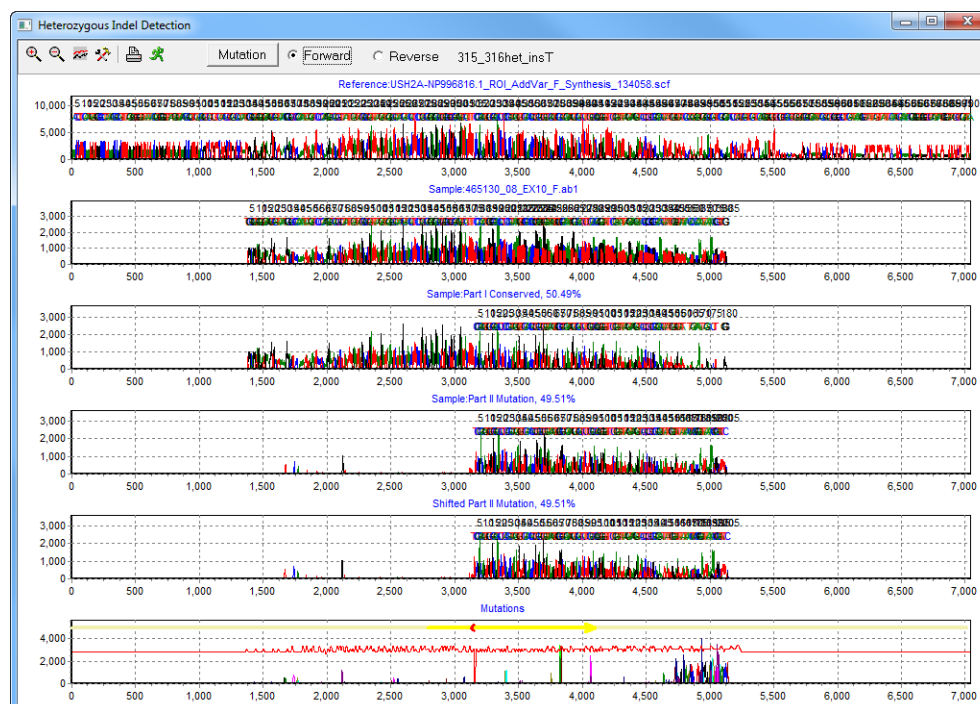
- Click the Detect Heterozygous Indels button  on the GAD toolbar.
- Press and hold the [Ctrl] key, and then click the mobility line at the position of the indel in the mutation trace in the GAD to open the Add Indel dialog box, and then on the Add Indel dialog box, click Heterozygous Indel.

Figure 12-3: Heterozygous Indel Detection window



The Heterozygous Indel Detection tool deconvolutes the mixed mutation trace into two clean traces. The *Conserved* trace shows the allele within the mixed trace that matches to the reference trace. The *Mutant* trace shows the allele within the mixed trace that does not match to the reference trace. The tool then estimates the percent contribution of conserved allele

and mutant allele. Mutation Surveyor uses several ideal peaks within the sample trace, determines the median height of the mutant allele at these locations and calculates the percent of mutant allele (M) as $[M/(M+C)]$ and the conserved allele (C) as $[C/(M+C)]$. Finally, Mutation Surveyor repositions the mutant trace so that it aligns to the reference trace and then determines whether the mutation was a heterozygous insertion or a heterozygous deletion. The indel positions are numbered relative to the positions in the reference sequence, and not to the genomic positions in the GenBank file. You can use the functions that are available from the window's toolbar for adding and editing heterozygous indel calls to the mutation trace.



When you add a heterozygous indel call to a trace or move the call upstream in the trace, the end of the comparison region is moved to the site of the heterozygous indel. As a result, all automatically added mutation calls that are downstream of the heterozygous indel are now outside of the comparison region and Mutation Surveyor deletes these calls. You can manually add mutation calls outside of the comparison region, but reliably automating these calls in these regions is not possible.

1. On the Heterozygous Indel Detection toolbar, click Mutation.

The Add/Edit Het-Indel dialog box opens.



For a complete description of all the buttons on the Heterozygous Indel Detection toolbar, see [“Heterozygous Indel Detection window toolbar”](#) on page 334.

Figure 12-4: Add/Edit Het-Indel dialog box

2. On the Mutation Type dropdown list, select the mutation type.

3. Do one of the following:
 - If you are adding a heterozygous insertion, then in the Start Index of Reference field, enter the reference positions that flank the insertion as two numbers delimited by an underscore. For example, 135_136 means that an insertion exists between base 135 and 136 of the reference. The length of the insertion is determined by the Mutation Letters parameter.
 - If you are adding a heterozygous deletion, then in the Start Index of Reference field, enter the actual positions of the deletion. For example, 235_237 means that three bases were deleted from the sample—the base at reference position 235, the base at reference position 236, and the base at reference position 237.
4. Optionally, to add multiple tags for the het-indel, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.
5. Click OK.

Three results are possible:

- The mutation trace is updated with the insertion or deletion. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.



If you add a heterozygous insertion and the bases that are inserted are identical to the bases that precede them, then the mutation code is corrected in the Mutation table to reflect this duplication. See [“Insertion, duplication, and deletion mutation codes” on page 131](#).







- If Mutation Surveyor cannot deconvolute the mixed trace after you have added an insertion or deletion call because of poor trace quality, or because of the complexity of the call, then a warning message opens, indicating that Mutation Surveyor cannot display the mutation call as you have specified and asking you if you want to keep it. Click Yes to save the mutation call to the trace. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.
- A message opens indicating that Mutation Surveyor was able to detect a different mutation call than the one specified and asking if you want to use Mutation Surveyor’s call. Click No to save the mutation call as you have specified it.

Heterozygous Indel Detection window toolbar

The Heterozygous Indel Detection window toolbar contains buttons that are specific for working with the display and for changing the information that is displayed.

Figure 12-5: Heterozygous Indel Detection window toolbar



Option/Button	Description
	Zoom In—Zooms in on a region of the report for viewing. Note: You can also manually zoom in on a region for viewing. Click and hold the left mouse button and draw a box from the upper left hand corner of the pane towards the lower right hand corner. A box is formed around the area that is being reduced for viewing.
	Zoom Out—Zooms out from a region of the report for viewing. Note: You can also manually zoom out from a region for viewing. Click and hold the left mouse button and draw a box from the lower right hand corner of the pane towards the upper left hand corner. The magnification for zooming out is always 100%.
	Show Lines—Provides options for showing all nucleotides in the traces and mutation traces, hiding all nucleotides in the traces and mutation traces, or showing only selected nucleotides in the traces and mutation traces.
	Settings—Opens a menu that contains selections for the layout of the display and for navigating the display. <ul style="list-style-type: none"> • Y auto—Clear this option to turn off the Y-axis auto-scaling for each of the traces. • Mouse Zoom—Clear this option to turn off the mouse zooming functions. (You can use only the Zoom In and Zoom Out buttons to change the zoom level of the display.) • Show Grid—Clear this option to remove the grid from the traces.
	Print Preview—Opens the Preview window, which displays a preview of what is to be printed for the Heterozygous Indel display. You can use the options on the Preview window's toolbar to adjust variables such as the color scale and the page layout before you print the display. Note: Before you click the Print Preview button, make sure to adjust the zoom level so that the area of the display that you want to print is displayed in the Preview window.
	Auto Identify the Start—Runs a more rigorous detection algorithm for deconvoluting the mixed trace and aligning the mutant trace to the reference trace. Note: Typically, you run this algorithm if you think that Mutation Surveyor made an incorrect het-indel call, or that het-indel call is missing.
Mutation	Opens the Add/Edit Het-Indel dialog box which contains options for adding or editing a het-indel call in a mutation trace. See “To add or edit a heterozygous indel (The Het-Indel tool)” on page 331 or “To edit heterozygous indels” on page 337 .
Forward/ Reverse	Indicates the sample trace that is currently displayed in the window.

Editing Mutation Calls

When you edit a mutation call, you can keep the mutation call type the same at the position, or you can change the call type at the position. If a mutation call of one type was made at a given position and you want to have a call of another type at the position, then you should first delete the original mutation call at the position, and then add the different type of mutation call at the position. You can [edit](#) substitution calls, you can [edit](#) homozygous indels, you can [edit](#) heterozygous indels, and you can [edit](#) Del/Ins.

To edit substitutions

You can edit a substitution call from the GAD or from the Project Reviewer.

1. Right-click the call in the Mutation table in the GAD, or in the Project Reviewer, and on the context menu that opens, click Edit.

The Edit Substitution dialog box opens.

Figure 12-6: Edit Substitution dialog box

2. In the Mutation field, enter the mutation, and optionally, add comments for the modified mutation.



Mutation Surveyor automatically calculates the optimal values for the mutation parameters such as intensity, area, and so on for the mutation based on the position of the mutation and you should not modify these values. When you add a substitution, it can be heterozygous (for example, A > AT), or it can be homozygous (for example, A > T).

3. Optionally, to add multiple tags for the substitution, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.
4. Click OK.

The mutation trace is updated with the substitution. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.

To edit homozygous indels

You can edit a homozygous indel call from the GAD or from the Project Reviewer.

1. Right-click the call in the Mutation table in the GAD or in the Project Reviewer, and on the context menu that opens, click Edit.

The Edit Indel dialog box opens.

Figure 12-7: Edit Indel dialog box

db_xref	
Tag	Identifier

2. Edit the indel call as appropriate.
 - If you are editing a homozygous insertion, then in the Start of GenBank Index field, enter the positions that flank the insertion as two numbers delimited by an underscore. For example, 30746_30747 means that an insertion exists between base 30746 and base 30747. The length of the insertion is determined by the Mutation Letters parameter.
 - If you are editing a homozygous deletion, then in the Start of GenBank Index field, enter the actual positions of the deletion. For example, 30746_30748 means that three bases were deleted from the sample—the base at position 30746, the base at position 30747, and the base at 30748.
3. Optionally, to add multiple tags for the homozygous indel, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.

4. Click OK.

The mutation trace is updated with the homozygous indel. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.



If you add a homozygous insertion and the bases that are inserted are identical to the bases that precede them, then the mutation code is corrected in the Mutation table to reflect this duplication. See “Insertion, duplication, and deletion mutation codes” on page 131.

To edit heterozygous indels

You use the Heterozygous Indel Detection tool (Het-Indel tool) to add or edit heterozygous mutation calls to a mutation trace.


1. Do one of the following to open the Het-Indel tool:
 - Double-click the het-indel call in the Mutation report to open the Heterozygous Indel Detection tool.
 - Click the Detect Heterozygous Indels button  on the GAD toolbar.
 - Press and hold the [Ctrl] key, and then click the mobility line at the position of the indel in the mutation trace in the GAD to open the Add Indel dialog box, and then on the Add Indel dialog box, click Heterozygous Indel.


Figure 12-8: Heterozygous Detection window



2. On the Heterozygous Indel Detection toolbar, click Mutation.

The Add/Edit Het-Indel dialog box opens.

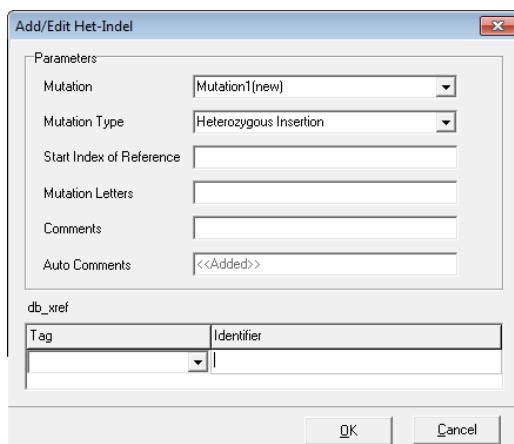Figure 12-9: Add/Edit Het-Indel dialog box



3. On the Mutation Type dropdown list, select the mutation type.
4. Do one of the following:
 - If you are adding a heterozygous insertion, then in the Start Index of Reference field, enter the positions of the reference that flank the insertion as two numbers delimited by an underscore. For example, 135_136 means that an insertion exists between base 135 and 136 of the reference. The length of the insertion is determined by the Mutation Letters parameter.
 - If you are adding a heterozygous deletion, then in the Start Index of Reference field, enter the reference positions of the deletion. For example, 235_237 means that three bases were deleted from the sample—the base at reference position 235, the base at reference position 236, and the base at reference position 237.
5. Optionally, to add multiple tags for the heterozygous indel, right-click anywhere in the db_xref pane, and on the context menu that opens, click Add Line. A new blank line for a tag and identifier is added to the pane.
6. Click OK.

Three results are possible:

- The mutation trace is updated with the insertion or deletion. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.



If you add a heterozygous insertion and the bases that are inserted are identical to the bases that precede them, then the mutation code is corrected in the Mutation table to reflect this duplication. See [“Insertion, duplication, and deletion mutation codes”](#) on page 131.

- If Mutation Surveyor cannot deconvolute the mixed trace after you have added an insertion or deletion call because of poor trace quality, or because of the complexity of the call, then a warning message opens, indicating that Mutation Surveyor cannot display the mutation call as you have specified and asking you if you want to keep it. Click Yes to save the mutation call to the trace. The Mutation table in the GAD and in the Project Reviewer are also updated to show this new mutation.
- A message opens indicating that Mutation Surveyor was able to detect a different mutation call than the one specified and asking if you want to use Mutation Surveyor's call. Click No to save the mutation call as you have specified it.


To edit Del/Ins

A *Del/Ins* is a combination of a deletion and an insertion at the same location. In many cases, you must edit such a mutation call to get the correct mutation call output. The following is a summary of the approaches that SoftGenetics recommends for adding and editing homozygous and heterozygous Del/Ins.

To edit a homozygous Del/Ins

1. You might initially need to delete substitutions that were falsely identified at these positions.
2. If no homozygous Indel was identified, then add the homozygous deletion as described in [“To add a homozygous indel call” on page 330](#).
3. At the same starting position, add the homozygous insertion as described in [“To add a homozygous indel call” on page 330](#).

To edit a heterozygous Del/Ins

1. Do one of the following to pen the Heterozygous Indel Detection tool.
 - Double-click the het-indel call in the Mutation report to open the Heterozygous Detection tool.
 - Click the Detect Heterozygous Indels button  on the GAD toolbar.
 - Press and hold the [Ctrl] key, and then click the mobility line at the position of the indel in the mutation trace in the GAD to open the Add Indel dialog box, and then on the Add Indel dialog box, click Heterozygous Indel.
2. Edit Mutation1 to the heterozygous deletion relative to reference trace positions.

3. Add Mutation2 as the heterozygous insertion relative to reference trace positions.

Two results are possible:

- The mutation trace is updated with the insertion or deletion. The Mutation table in the GAD and in the Project Reviewer are also updated to show these modified mutations.



If you add a heterozygous insertion and the bases that are inserted are identical the bases that precede them, then the mutation code is corrected in the Mutation table to reflect this duplication. See [“Insertion, duplication, and deletion mutation codes” on page 131](#).

- If Mutation Surveyor cannot deconvolute the mixed trace after you have edited an insertion or deletion call because of poor trace quality, or because of the complexity of the call, then one of two warning messages opens:
 - A warning message opens indicating that Mutation Surveyor cannot display the mutation call as you have specified and asking you if you want to keep it. Click Yes to save the mutation call to the trace. The Mutation table in the GAD and in the Project Reviewer are also updated to show these modified mutations.
 - A warning message opens asking you if you want to use the mutation call that is automatically detected at the position. Click No to save the mutation calls as you have specified them. The Mutation table in the GAD and in the Project Reviewer are also updated to show these modified mutations.

Annotating Reported Variations

You can [annotate](#) *reported* variations in a sample trace from the GAD or from the Project Reviewer report. Mutation Surveyor automatically modifies the GenBank file that is stored for the selected project with these annotations.

To annotate a variation

1. Right-click the call in the Mutation table in the GAD, or in the Project Reviewer report, and on the context menu that opens, click Edit.



Remember, a reported variant has a purple background in the Mutation table in the GAD and Project Reviewer report.

An Edit (mutation type) dialog box opens. The type of dialog box that opens depends on the type of mutation call that was selected for annotation (Substitution or Indel) but both dialog boxes have the lower db_xref pane. The pane is automatically populated with all the reported annotations for the position.

Figure 12-10: db_xref pane

Tag	Identifier
dbSNP	28897676

2. Optionally, to modify an existing annotation, select a new tag for it on the tag dropdown list.
3. To add a new annotation, right-click in the pane, and on the context menu that opens, click Add a Line.

A new blank line is added for the annotation.

4. On the tag dropdown list, select a tag for the new annotation.
5. In the Identifier field, enter the identifier for the new annotation.



If you add an annotation in error, or want to delete an existing annotation, right-click the annotation in the db_xref pane, and on the context menu that opens, click Delete a Line.

6. Click OK.

Mutation Surveyor automatically adds the new annotations to the GenBank file that is stored for the selected project.



You can use the Extract SGP tool to export the updated GenBank file. See [“Extract SGP Tool”](#) on page 275.

Creating a User Knowledge Database for Known Deletions

Mutation Surveyor automatically identifies mutation calls at positions where the variation between the reference trace and the sample trace meet or exceed detection thresholds. If, however, you know that a mutation is being called as the result of a false positive or a sequencing artifact, and the GenBank file that was used for the project contains chromosome information, then you can delete the called mutation and [add](#) this deletion to a User Knowledge database. For this User Knowledge database, a false positive and an artifact are defined as follows:

- A false positive is dependent on the chromosome position. As a result, this same variant is considered to be a false positive for all subsequent analyses.
- An artifact is dependent on not only the chromosome position, but also its relative location in the trace. For example, in one trace, a variant is located at the 5' end of the trace, but in another trace, the same variant is located in the middle of the trace. As a result, if this variant is marked as an artifact in the trace where it was found at the 5' end, the variant is not considered to be an artifact in the trace where it was located in the middle.



For an artifact, the User Knowledge Database stores a range of positions, relative to location within a trace. This is more ideal for mutation calls that routinely occur at the beginning or end of an amplicon as a mutation call that is centered in an amplicon is more than likely to be a real mutation.


You can use this custom database of known deletions in other projects. If a trace in a project has a mutation called at a position that has been identified as a false positive or artifact in the User Knowledge database, then Mutation Surveyor automatically flags the position, which streamlines the process of identifying variations of clinical significance for the project.

To create a custom variant database for known deletions

1. Right-click the call in the Mutation table in the GAD or in the Project Reviewer, and on the context menu that opens, click Delete Mutation and Add to Database.

The Delete Mutation dialog box opens.

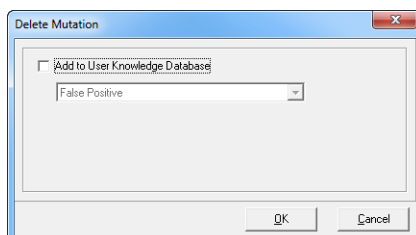Figure 12-11: Delete Mutation dialog box





The Delete Mutation and Add to Database option is available only if the requirements of an imported whole human genome reference and the inclusion of a GenBank file with chromosome coordinate in the project are met.


2. Do one of the following:
 - To delete the mutation without adding it to the User Knowledge database, click OK.
 - To add the mutation to the User Knowledge database, select Add to User Knowledge Database, and then complete the remaining steps in this procedure.
3. Leave the reason for the deletion set to the default value of False Positive, or select Artifact.
4. Click OK.

The Delete Mutation dialog box closes. The deleted mutation (position and reason) are saved to the User Knowledge database. If a trace in another project has a mutation called at this exact position, then Mutation Surveyor automatically flags this position.

Viewing the Edit History (Audit Trail) for a Mutation

Analysis options are available for a mutation in a variety of locations in Mutation Surveyor, including various reports and displays and the context menus for the Mutation table and the GAD. Any edit action (addition, substitution, confirmation, or deletion) that you carry out for a mutation is displayed in the Actions column for the mutation in the Mutation table and this action is also automatically added to the audit trail for the mutation. An option is available from all these locations to view the edit history for a mutation. To view the edit history for a mutation, right-click the mutation (for example, in the Mutation table), and on the context menu that opens, click View Edit History to open the Edit History dialog box.

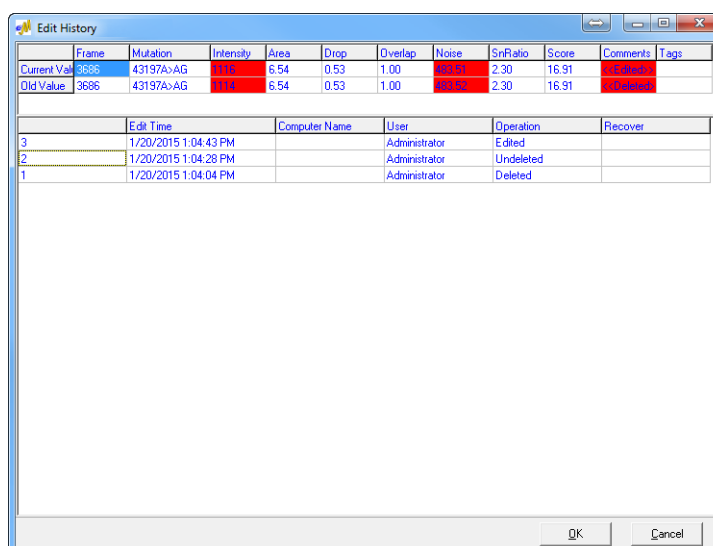Figure 12-12: Edit History dialog box



The lower half of the Edit History dialog box displays all the edit operations that have been carried for the selected mutation. The date and time and the username for the user who carried out the edit is displayed for each edit. When you select an edit entry in the lower pane, a selected series of old and new values is displayed in the upper half of the dialog box. If the edit resulted in a change for a mutation value, then the old and new values are highlighted in red.

Index

Numerics

2 Directions tab on the Mutation Project Settings dialog box 65

A

activity log file

viewing for your Mutation Surveyor users 43

Advanced GBK File Editor tool .. 231

functions 243

GenBank Tree File pane..... 233

Tabs pane..... 235

toolbar 232

Advanced Two Direction report . 140

context menu 144

toolbar 142

viewing and analysis options .. 144

allele ratios

quantifying for Sanger sequencing traces 303

Amino Acid frame in the Graphical Analysis Display 79

amino acid text comparison function in the Project Sample Assembly subreport 154

assumptions for the manual 16

audit trail

see edit history

B

BasePatch tool 295

Bentley report 196

context menu 198

report settings 198

toolbar 197

viewing and analysis options .. 197

Browser pane

Mutation Surveyor main window... 27

C

CDS information

adding and modifying for a SEQ file in the SEQ File Editor tool . 246

Clinical report..... 116

color codes

for mutation codes in the Mutation report..... 130

Color tab for a Custom report 176

Confidence mark

for Custom report 173

for JHU report 187

consensus sequence text function in the Project Sample Assembly subreport..... 151

context menu

Advanced Two Direction report 144

Bentley report..... 198

Emory report 194

Full Gene report..... 220

HGVS report 161

JHU report..... 188

Methylation report 206

MSM_Song report..... 198

Mutation report..... 132

PrettyBase report..... 202

Project report 155

Project Reviewer report 101

Two Direction report..... 137

Contig tab on the Mutation Project Settings dialog box..... 55

conventions used

in Mutation Surveyor 31

in the manual..... 15

Copyright..... 3

Custom report

Color tab..... 176

Confidence mark for..... 173

creating a template for 164

defined 164

Display tab 172

Filters tab 170

Format tab..... 166

Mutation tab 173

Nomenclature tab..... 177

Other tab 179

Customer reports

Bentley report..... 196

Emory report 193

Full Gene report 215

FullBase report..... 222

JHU report..... 185

Methylation report 203

Graphic Display of Methylations report..... 205

MSM_Song report..... 196

PrettyBase report 200

Tassel report 190

Vertex report 207

Vertex2 report 209

Vertex3 report 212

D

Database tab

on the Mutation Project Settings
dialog box..... 69

Del/Ins

defined 339
editing for a mutation trace 339

Display tab for a Custom report . 172

Display tab on the Mutation Project
Settings dialog box 64

E

edit history

viewing for a mutation 344

Emory report 193

context menu 194

toolbar 193

viewing and analysis options .. 194

Export GAD function

defined 94

F

Filters tab for a Custom report ... 170

Format tab for a Custom report.. 166

Full Gene report..... 215

context menu 220

Format report settings 216

Header report settings 217

Output report settings 217

toolbar 219

viewing and analysis options .. 220

FullBase report 222

layout 223

report settings 222

toolbar 223

viewing and analysis options... 224

functions

in the Advanced GBK File Editor
tool 243

in the SEQ File Editor tool..... 250

G

GAD

see Graphical Analysis Display

GAD report..... 114

toolbar 115

GenBank file

adding and modifying variations
for in the Advanced GBK File
Editor tool 241

as a project file type in Mutation
Surveyor..... 50

automatically downloading for a
project 51

downloading from NCBI 227

editing and annotating with the
Advanced GBK File Editor tool 231

overview of 227

GenBank Tree File pane

Advanced GBK File Editor tool 233

genomic terms

conventions used for in Mutation
Surveyor..... 24

Graphic Display of Methylations
report..... 205

Graphic Display report 108

layout..... 109

toolbar 110

Graphical Analysis Display

annotating reported variations in
its mutation traces 341

defined..... 77

exporting information for..... 94

mutation peak identification..... 82

Mutation table..... 83

mutation trace legend in 80

navigation and viewing options for
90

opening..... 77

toolbar for 84

Graphical Analysis Display
components

Amino Acid frame 79

Nucleotide Sequence frame 78

Sample Traces frame 79

Tracks frame 79

graphical output

conventions used for in Mutation
Surveyor 25

Graphical reports

Clinical report 116

GAD report 114

Graphic Display report..... 108

Project Display subreport..... 111

Nucleotide Text Display report 120

Project Display subreport 111

Project Reviewer report..... 99

group

defined..... 38

adding..... 38

deleting..... 38

editing..... 38

H

- heterozygous indel
 - adding to a mutation trace in the comparison region 331
 - adding to a mutation trace outside the comparison region 331
 - editing for a mutation trace 337
- Heterozygous Indel Detection tool... 331
- HGVS report 157
 - context menu 161
 - report settings 157
 - toolbar 160
 - viewing and analysis options .. 161
- homozygous indel
 - adding to a mutation trace 330
 - editing for a mutation trace 336

I

- Input tab on the Mutation Project Settings dialog box 54
- installing
 - Mutation Surveyor..... 22

J

- JHU report 185
 - Confidence mark for 187
 - context menu 188
 - report settings 187
 - toolbar 186
 - viewing and analysis options .. 187

L

- layout
 - FullBase report..... 223
 - Graphic Display report 109
 - Nucleotide Text Display report 121
 - Project Reviewer report 99
 - SEQ File Editor tool 247
 - Two Direction report..... 134
 - Vertex3 report 213

M

- main menu
 - Mutation Surveyor main window... 26
- main window for Mutation Surveyor. 26
- methyated DNA
 - analyzing..... 301
- Methylation report 203
 - context menu 206
 - report settings 204
 - toolbar 204
- mobility line
 - defined 82
- Modify Mutation Projects function
 - defined 315
 - using 315
- MSM_Song report..... 196
 - context menu 198
 - report settings 198
 - toolbar 197
 - viewing and analysis options .. 197
- mutation
 - adding to the User Knowledge Database when deleted 342
 - viewing the edit history for..... 344
- Mutation AutoRun tool
 - Log File Wizard in 317
 - using to analyze multiple projects . 324
- mutation calls
 - adding to a mutation trace..... 329
 - editing for a mutation trace..... 335
- mutation code
 - defined 129
- mutation codes
 - color codes for in the Mutation report..... 130
 - Two Direction report..... 136
- mutation peak identification in the Graphical Analysis Display 82
- Mutation Project Settings dialog box
 - 2 Directions tab 65
 - Contig tab..... 55
 - Database tab..... 69
 - Display tab 64
 - Input tab 54
 - Mutation tab 59
 - Others tab 67
 - Output tab 62
- Mutation Quantifier function
 - defined 303
 - running 303
- Mutation report..... 127
 - context menu..... 132
 - report codes 129
 - report fields 128
 - settings

Index

- variation tracks..... 87
- viewing and analysis options .. 132
- Mutation Surveyor
 - conventions
 - genomic terms 24
 - mutation terms 24
 - reports and graphical output ... 25
 - installing 22
 - main window 26
 - Browser pane..... 27
 - main menu 26
 - title bar 26
 - toolbar 27
 - project file types 50
 - setting up a project in..... 70
 - starting 23
 - system requirements for 21
- Mutation tab for a Custom report 173
- Mutation tab on the Mutation Project Settings dialog box 59
- Mutation table in the Graphical Analysis Display..... 83
- mutation terms
 - conventions used for in Mutation Surveyor..... 24
- mutation trace
 - adding a heterozygous indel to in the comparison region 331
 - adding a heterozygous indel to outside the comparison region 331
 - adding a homozygous indel to 330
 - adding a substitution to 329
 - adding mutation calls to 329
 - editing a Del/Ins for 339
 - editing a heterozygous indel for ... 337
 - editing a homozygous indel for 336
 - editing a substitution for 335
 - editing the annotations in 341
 - editing the mutation calls for ... 335
 - reprocessing with the BasePatch tool 295
- mutation trace legend in the Graphical Analysis Display 80
- mutation_server_project_settings 52
- N**
 - Nomenclature tab for a Custom report..... 177
 - Nucleotide Sequence frame in the Graphical Analysis Display 78
 - Nucleotide Text Display report... 120
 - layout..... 121
 - settings..... 120
 - toolbar 122
- O**
 - organization of the manual..... 16
 - Other tab for a Custom report 179
 - Others tab
 - on the Mutation Project Settings dialog box..... 67
 - Output tab on the Mutation Project Settings dialog box..... 62
 - overview
 - GenBank file..... 227
- P**
 - PrettyBase report..... 200
 - context menu..... 202
 - report settings..... 201
 - toolbar 200
 - viewing and analysis options.. 201
 - project
 - analyzing multiple projects using the Mutation AutoRun tool..... 324
 - automatically downloading a GenBank file for..... 51
 - modifying single or multiple 315
 - modifying the settings for 52
 - setting up in Mutation Surveyor 70
 - Project Display subreport 111
 - report settings..... 111
 - toolbar 113
 - project file types
 - GenBank file..... 50
 - reference file..... 50
 - sample file 50
 - Project report..... 146
 - context menu..... 155
 - report settings..... 146
 - toolbar 148
 - viewing and analysis options.. 155
 - Project Reviewer report
 - annotating reported variations in its mutation traces 341
 - context menu..... 101
 - layout..... 99
 - report settings..... 104
 - toolbar 102
 - Project Sample Assembly subreport 150
 - amino acid text comparison function in 154

consensus sequence text function
in 151

sequence text comparison
function in 152

Project Viewer report 99

R

reference file

- as a project file type in Mutation
Surveyor 50

report codes

- Mutation report 129

report fields

- Mutation report 128

report settings

- Bentley report 198
- Full Gene report 216, 217
- FullBase report 222
- HGVS report 157
- JHU report 187
- Methylation report 204
- MSM_Song report 198
- Nucleotide Text Display report 120
- PrettyBase report 201
- Project Display subreport 111
- Project report 146
- Project Reviewer report 104
- Two Direction report 136
- Verrtex3 report 212
- Vertex2 report 209

reports

- see Graphical reports, Standard
reports, Custom report, or
Customer reports
- conventions used for in Mutation
Surveyor 25

S

sample file

- as a project file type in Mutation
Surveyor 50

Sample Traces frame in the
Graphical Analysis Display 79

- legend for 80
- mobility line in 82
- mutation peak detection in 82

Sanger sequencing traces

- quantifying the allele ratios for 303

SEQ file

- adding and modifying CDS
information for in the SEQ File
Editor tool 246

SEQ File Editor tool 246

- functions 250
- layout 247

sequence text comparison function
in the Project Sample Assembly
subreport 152

settings

- Full Gene report 217

special information about the
manual 15

Standard reports

- Advanced Two Direction report
140
- HGVS report 157
- Mutation report 127
- Project report 146
- Project Sample Assembly
subreport 150
- Two Direction report 134

substitution

- adding to a mutation trace 329

editing for a mutation trace 335

system requirements for Mutation
Surveyor 21

T

Tabs pane

- Advanced GBK File Editor tool 235

Tassel report 190

- toolbar 191
- viewing and analysis options... 192

template

- creating for a Custom report ... 164

title bar

- Mutation Surveyor main window...
26

toolbar

- Advanced GBK File Editor tool 232
- Advanced Two Direction report
142
- Bentley report 197
- Emory report 193
- Full Gene report 219
- FullBase report 223
- GAD report 115
- Graphic Display report 110
- Graphical Analysis Display 84
- HGVS report 160
- JHU report 186
- Methylation report 204
- MSM_Song report 197
- Mutation Surveyor main window...
27
- Nucleotide Text Display report 122
- PrettyBase report 200
- Project Display subreport 113

Index

Project report	148	user management		Project report.....	155
Project Reviewer report	102	configuring.....	31	Tassel report	192
Tassel report.....	191	defined	31	Two Direction report.....	136
Two Direction report.....	135	turning off	36	Vertex2 report.....	211
Vertex2 report	211	turning on	34	Vertex3 report.....	214
Vertex3 report	214	using the manual.....	15		
tools		V			
Advanced GBK File Editor	231	variations			
BasePatch.....	295	adding and modifying in the			
Heterozygous Indel Detection		Advanced GBK File Editor tool	241		
(Het_Indel)	331	annotating for a mutation trace in			
Mutation AutoRun	324	the GAD or from the Project			
Log File Wizard.....	317	Reviewer report.....	341		
SEQ File Editor	246	Vertex report	207		
Tracks frame in the Graphical		Vertex2 report	209		
Analysis Display.....	79	report settings	209		
Two Direction report.....	134	toolbar	211		
context menu	137	viewing and analysis options...	211		
layout	134	Vertex3 report	212		
mutation codes.....	136	layout.....	213		
report settings	136	report settings	212		
toolbar	135	toolbar	214		
viewing and analysis options ..	136	viewing and analysis options...	214		
U		viewing and analysis options			
user		Advanced Two Direction report			
defined	38	144			
adding	43	Emory report	194		
deleting	43	Full Gene report	220		
editing	43	FullBase report.....	224		
viewing the activity for in a log file.		Graphical Analysis Display.....	90		
43		HGVS report.....	161		
User Knowledge Database		JHU report.....	187		
creating for known mutation		Mutation report	132		
deletions.....	342	PrettyBase report	201		

Glossary

comparison region

A region where the sample traces align to the reference traces, and all the mutation project *analysis* settings have been met, and therefore, mutations, with the exception of heterozygous indels, can be called in this region. The comparison region is marked by two vertical blue bars in the mutation trace.

